


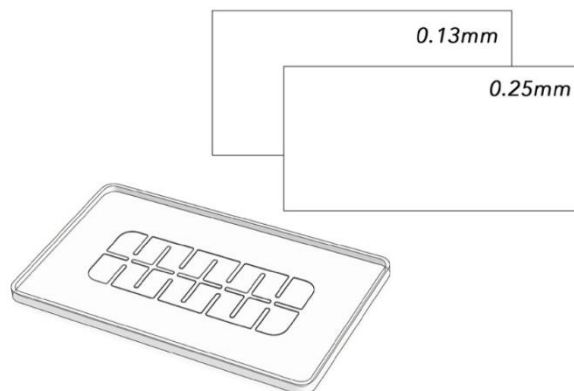

Using the Ward's LiveSlide®

A Ward's LiveSlide® set contains 3 pieces, the base slide and a set of coverslips, all of which are reusable. The 2 coverslips are different thicknesses. One is very thin (0.13mm), perfect for high magnification. The second is a bit thicker (0.25mm) for all applications.

Note: Each coverslip has a thin plastic film on each side for protection, one green and one clear. Remove both films by either peeling, or using a small piece of tape on the corner and lifting. Discard the films.



Using a Transfer Pipette, place the desired number of specimen droplets in the wells on the base slide.

Then gently place the coverslip over the culture well



Place the Ward's LiveSlide® on a microscope stage or under a digital microscope, and enjoy exploring the magnified world!

Order Yours Today!

Order #	Description
470313-232	5-pack of LiveSlides®
470309-714	Bulk box of 50 LiveSlides®

