

Quick Guide

Auto *lite* Software Installation

ECN: 710-2072

- 1 Start PC**

➡ Log onto PC as an Administrator and shut all running software down

2 Insert software source

➡ Insert the software source, unzip if the installation set is zipped.

3 Support folder



Support



Framework



SetupFramework.exe

➡ Open the Support folder

➡ Open the Framework folder

➡ Run SetupFramework.exe and follow the on-screen instructions, performing a system restart if update requested


4 Install SQL Server



Support



Database

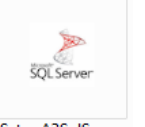

SetupA3SqlServer.exe

➡ Open the Support folder

➡ Open the Database folder

➡ Run SetupA3SqlServer.exe (see SQL Installation within the manual) and on completion perform a PC restart

For single user PC: Choose SQL LocalDB
For multiuser PC: Choose SQL Server

5

Run Set-up



- ➔ Run Setup.exe once again and follow the on-screen instructions (see *Auto lite* Full Installation Guide for details within Installation -> Manuals-> Quick Guides)

6

Connect the Auto *lite*


- ➔ Connect the power cord to the Auto *lite* and plug the unit in, and connect the USB cable from the Auto *lite* to a spare USB port in the PC

7


Adding users to SQL server

The following steps need to be carried out when logged in as an Administrator.

When Auto *lite* is installed users need to be a member of one or more of these groups.

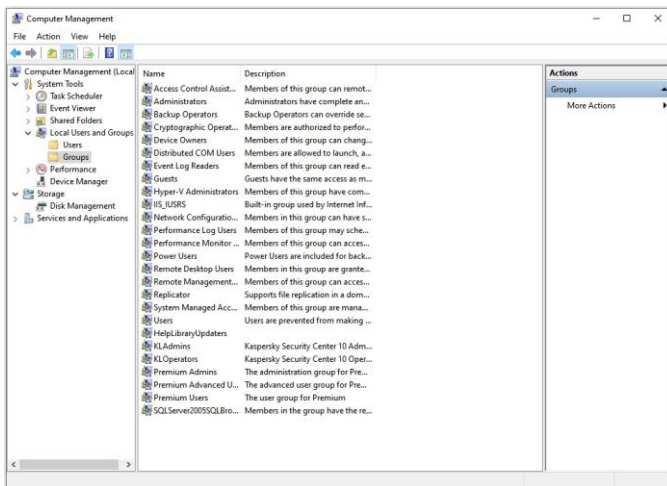


- ➔ From the Windows Control Panel select Administrative Tools then Computer Management
- ➔ Expand (+) local users and groups then double click on Groups
- ➔ Right click on Auto *lite* User and select add to group

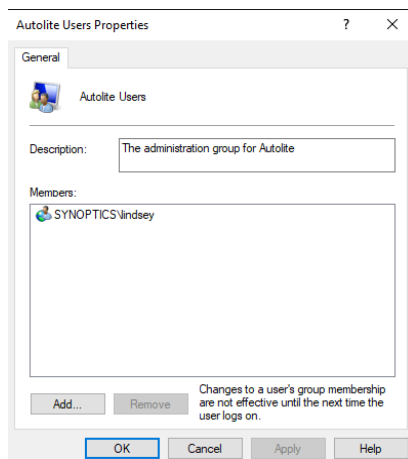


- ➔ Click Add

8

Adding users to SQL server (cont.)

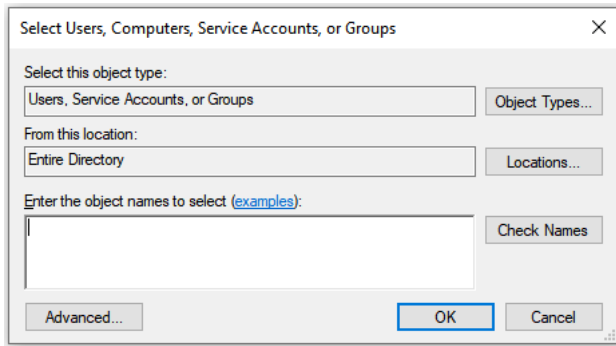

➔ To add a single user add their Windows log in name

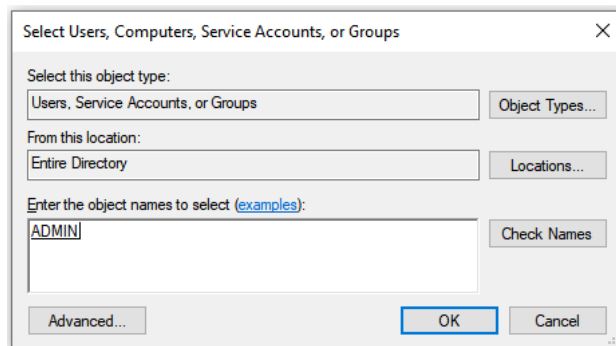

➔ Click Check Names

➔ Click OK

➔ Close computer management window

Auto lite is now ready to be used