Eutech Testr 35 multi-parameter pocket testers

Instruction Manual

PCTEST35, PCSTEST35
Revision Date: 13-11-2015

Language Index	
English	Page 2 - 5
Deutsch	Page 6 - 9
Français	Page 10 - 13
Nederlands	Page 14 - 16
Español	Page 17 - 18

MULTI-PARAMETER TESTER 35

pH/ Conductivity/ TDS/ Salinity/ Temperature

CONDITIONING by immersing the sensor in tap water for 30 minutes before use. DO NOT use de-ionized water.

SAVING THE SET UP

After any manual changes have been done during set up, you are required to save the setting.

1. Press both ▲ and ▼ keys at the same time and release.
2. Press again until you see **SA YES**
3. Use either ▲ OR ▼ key to toggle between YES and NO. Selecting YES will save your changes/ calibrations.
NOTE: If you did NOT make changes or recalibrate, this screen will NOT appear.

SYSTEM SETUP

1. Turn the tester off.
2. Press and hold ENT key and press ON/OFF to enter the Setup mode.
3. Release both keys, the display will show **SET UP**, followed by **PARA**
4. Press ▲ or ▼ to select **SYSt**
5. Press ENT to select
6. Use the ▲ or ▼ keys to go through the menu
7. Press ENT to select one of the following options:
 - a. **Unit rSt** Resets meter calibration to factory defaults
 - i. Press ENT to select
 - ii. Press ▲ or ▼ to select **PH rSt** or **EC rSt**
 - iii. Press ENT to select
 - iv. Press ▲ or ▼ to select **CAL rSt** (calibration reset) or **FCt rSt** (reset all to factory defaults)
 - v. Repeat if necessary for pH or EC.
 - b. **SEt A.OFF** for turning on or off the Auto-Off function (meter shuts down after 8,5 minutes)
 - i. Press ENT to select
 - ii. Press ▲ or ▼ to select **YES** or **NO**
 - iii. Press ENT to confirm
 - c. **SEt t.C** for setting the temperature coefficient
 - i. Press ENT to confirm
 - ii. Press ▲ or ▼ to adjust the value
 - iii. Press ENT to confirm
 - d. **SEt ATC** for turning on or off the Automatic Temperature Compensation
 - i. Press ENT to select
 - ii. Press ▲ or ▼ to select **YES** or **NO**
 - iii. Press ENT to confirm
 - e. **SEt °C°F** for setting the unit of measurement for temperature
 - i. Press ENT to select
 - ii. Press ▲ or ▼ to select **°C** or **°F**
 - iii. Press ENT to confirm

8. Save changes as described at "SAVING THE SET UP".

PARAMETER SETUP

1. Turn the tester off.
2. Press and hold ENT key and press ON/OFF to enter the Setup mode.
3. Release both keys, the display will show **SET UP**, followed by **PARA**
4. Press ▲ or ▼ to select **PARA**
5. Press ENT to select
6. Use the ▲ or ▼ to select PH, COND, TDS or SALT
7. Press ENT to select:
 - a. **PH** settings, buffers and calibration points
 - i. Press ▲ or ▼ to select **USA** or **NIST** buffers
 - ii. Press ENT to confirm
 - iii. Press ▲ or ▼ to select **3-Pt** or **5-Pt** (3 or 5 point calibration)
 - iv. Press ENT to confirm and go back to menu
 - b. **COND** settings, auto / manual calibration and calibration points
 - i. Press ▲ or ▼ to select **YES A.CAL** or **nO A.CAL** (auto calibration on or off)
 - ii. Press ENT to confirm
 - iii. Press ▲ or ▼ to select **YES SPC** (Single Point Calibration) or **nO SPC** (Multi Point Calibration)
 - iv. Press ENT to confirm and go back to menu
 - c. **TDS** settings for adjusting the TDS factor
 - i. Press ▲ or ▼ to change the value **0.67 FACt**
 - ii. Press ENT to confirm and go back to menu
 - d. **SALT** settings, measure in ppt or % (PEr Unit)
 - i. Press ▲ or ▼ to change the value **PPt Unit** or **PEr Unit**
 - ii. Press ENT to confirm and go back to menu
8. Save changes as described at "SAVING THE SET UP".


CALIBRATION

1. Turn the tester on.
2. Press MODE to select the desired parameter (pH, Cond, TDS or Salt)
3. Press CAL.
4. Place the tester in the calibration solution.
5. Wait for the value to stabilize.
 - a. **pH Calibration & Automatic Conductivity calibration:**
 - i. Press ENT to confirm
 - ii. Place the tester in the next calibration solution
 - iii. Repeat steps 5 to 5b for additional points or CAL to save when calibration is done
 - b. **Manual Conductivity Calibration / TDS Calibration and Salt Calibration:**
 - i. Press ▲ or ▼ to enter the correct value
 - ii. Press ENT to confirm, tester goes back in measurement mode

TEMPERATURE Calibration

This is only done once in a long while, calibration for temperature will remain accurate unless used under severe conditions.

1. Press ON/OFF Place a standard thermometer in the sample for reference.
2. Place tester into the same sample.
3. Allow some time for sensor to stabilize.
4. Press MODE to select the pH mode.
5. Press CAL.
6. Press HOLD for 5 seconds and wait till screen changes.
7. Use ▲ / ▼ keys to adjust until you reach the correct temperature of your sample according to the reference thermometer.
8. Press ENT to confirm.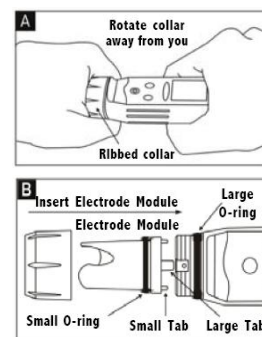

HOLD Function: Freezes the display.


1. Press HOLD to freeze the measurement. 'HOLD' will be displayed.
2. Press HOLD again to release.

MEASUREMENT

1. Press ON/OFF to switch the meter on.
2. Dip the electrode 2 to 3cm into the sample. Stir and let the reading stabilize.
3. Press MODE to view the desired parameter

CHANGING BATTERIES: (performance can be improved when 1 bar battery life remaining)

1. Open battery compartment lid. Remove old batteries noting polarity (flat side of battery is +).
2. Replace with (4) fresh ones. LR44 or A76 batteries

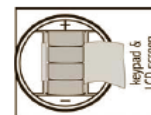CHANGING THE SENSOR

Electrode Replacement: necessary when calibration fails or when readings are erroneous in new calibration standards with hydrated electrode sensor and fresh batteries.

1. Grip the ribbed collar with electrode facing you. Twist the collar counter clockwise (Fig A). Save the ribbed tester collar and O-ring.
2. Pull the electrode off the tester.
3. Align the 4 tabs on the new module with the 4 slots on the tester (Fig B).
4. Gently push the new module onto the slots.
5. Push the smaller O-ring fully onto the electrode.
6. Push the collar over the module and twist it clockwise.
7. Note: Recalibrate your tester after electrode replacement.

Warranty

The waterproof testers are warranted to be free from manufacturing defects for 1 year and the electrode module is warranted for 6 months, unless otherwise stated. If repair, adjustment or replacement is necessary and has not been

the result of abuse or misuse within the time period specified, please return the tester – freight prepaid – and correction will be made without charge. Out of warranty products will be repaired on a charge basis.

Return of Items

Authorization must be obtained from your distributor before returning items for any reason. When applying for authorization, please include information regarding the reason the item(s) are to be returned.

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