



DWK LIFE SCIENCES
DURAN WHEATON KIMBLE

DURAN® AUTOCLAVE BOTTLE AND PUSH-ON CAP SYSTEM

NEW

The safer,
faster solution
for laboratory
sterilization



DWK Life Sciences SALES PARTNER

 **vwr**™
part of avantor

order on
[vwr.com](https://www.vwr.com)

DURAN®

AUTOCLAVE BOTTLE AND PUSH-ON CAP SYSTEM

Steam sterilization – a proven method for sterilizing liquids

Steam sterilization is one of the most reliable and widely adopted methods for ensuring **liquid sterility**. Typically performed in an **autoclave**, the process works by exposing the material directly to steam at the required temperature and pressure for a defined period.

When autoclaving liquids in glass bottles, it is essential that screw closures are only loosely applied. If bottles are sealed tightly, pressure can build up, potentially leading to breakage

or even explosive failure. **Safe autoclaving** of sealed bottles is only possible if the autoclave is equipped with **support pressure compensation**.

Alternatives to screw closures include vented caps and ported caps that can accommodate small sterile-venting filters. While effective, such closures may pose challenges for validation in GLP processes and should be carefully evaluated before use.

DURAN® Autoclave Bottle and Cap – safe, simple, and reliable

The **DURAN® Autoclave Bottle** and **Push-on Cap System** provides a convenient and secure solution for steam sterilization of liquid media, eliminating the need for pre- or post-autoclaving adjustments.

Available in three sizes, the DURAN® Autoclave Bottle is paired with a self-venting, push-on cap, ensuring safe and efficient


sterilization. Manufactured from **DURAN® borosilicate 3.3 glass**, these bottles offer exceptional thermal shock resistance, making them a trusted choice for applications where repeatable steam sterilization is essential. The accompanying Autoclave Push-on Cap is molded from a high-quality natural rubber formulation, delivering both reliability and peace of mind.

Suitability of bottle closures for steam sterilization




How the DURAN® Autoclave Push-on Cap works

The Autoclave Push-on Cap fits securely over the specially designed rim of the DURAN® Autoclave Bottle. During the autoclave cycle, it allows safe venting of steam and air. As the bottle cools, a vacuum forms and the cap automatically seals, with the concave center providing a clear visual check of integrity. For easy access, the push-on cap can be removed via the side tab, minimizing the risk of contact with sterile contents.

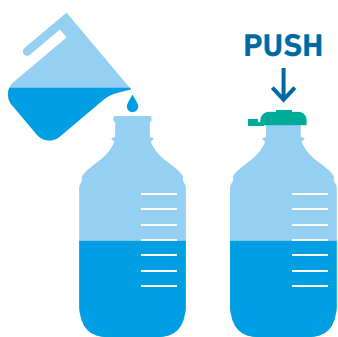


Preparation

Fill the bottle with liquid media up to 75% of its capacity, then push on the cap.

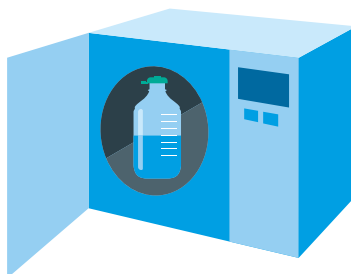

Sterilization

Place the bottle in the autoclave. During the cycle, steam is vented automatically through the valve of the cap.

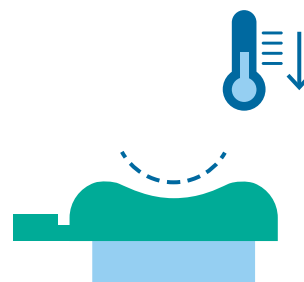

Post-Cycle

As the bottle cools, the valve closes, creating a vacuum inside. The center of the cap becomes concave, providing a visual indication of the hermetic seal.

Applications


- Steam sterilization of liquid media and reagents
- Media preparation for anaerobic microbiology
- Sterilizing animal drinking water in gnotobiology facilities



Specially designed bottle neck fits **self-venting Push-on Cap** for safe sterilization

Push-on natural rubber cap for simple, secure, and easy sealing


Autoclavable at 121 °C, for **reliable steam sterilization**

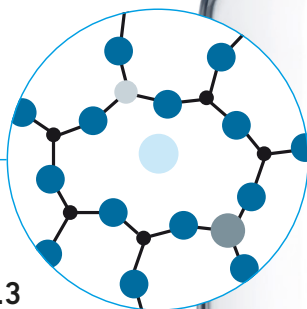

Made from **DURAN® Borosilicate Glass 3.3** with outstanding thermal and chemical resistance

Lot-specific **Retrace Code** for full manufacturing traceability and certification



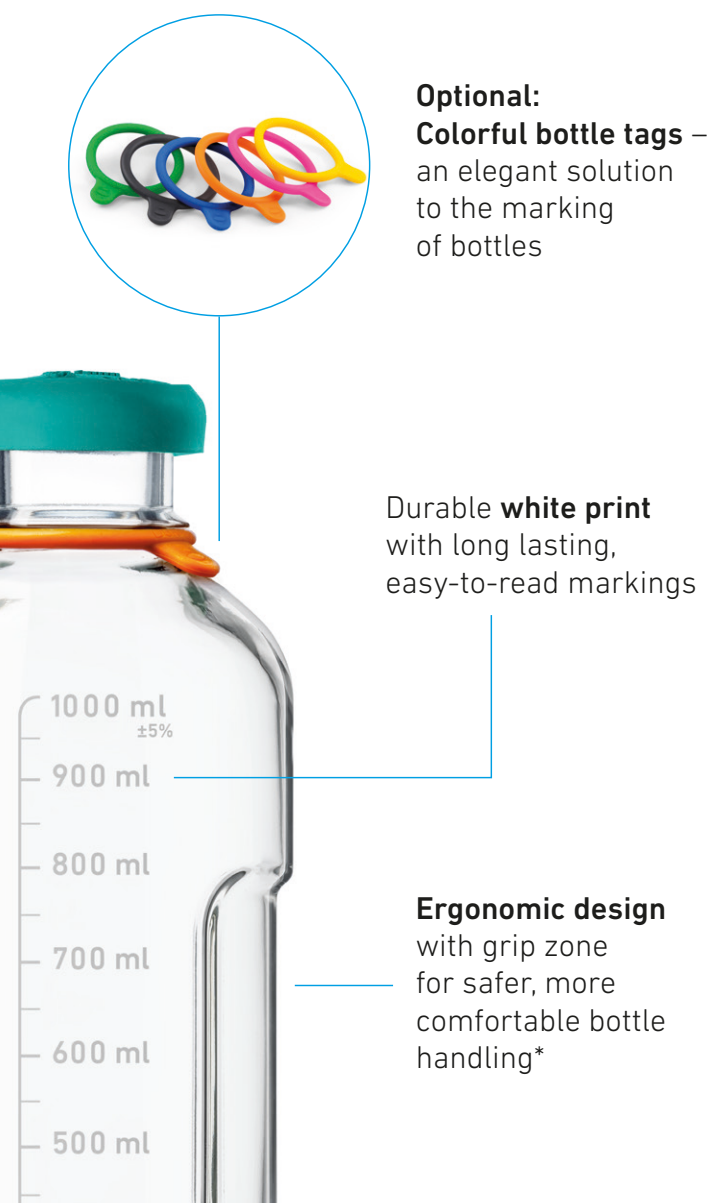

Optional:
Colorful bottle tags –
an elegant solution
to the marking
of bottles

Durable **white print**
with long lasting,
easy-to-read markings

Ergonomic design
with grip zone
for safer, more
comfortable bottle
handling*

The DURAN® Autoclave Bottle and Push-on Cap system benefits:

- **Minimized Sterility Risk**
Prevents contamination by minimizing manual interventions.
- **Enhanced User Safety**
Reduces operator risk by removing manual closure steps during post-autoclave handling.
- **Easy Identification**
Labels and color-coded bottle tags ensure clear, organized identification in the lab.
- **Optimized Space and Ergonomics**
Slim design and ergonomic grip zones improve space utilization and handling comfort.*
- **Time Efficiency**
Eliminates manual cap tightening, saving time and improving autoclaving efficiency.
- **Streamlined Ordering and Inventory**
Available in matched quantities for easier ordering and inventory management.

Ordering Information

Description	Pack Format	Cat. No.
DURAN® Autoclave Bottle with Push-on NR caps 500 ml	4 bottles & 4 caps	79060-108
DURAN® Autoclave Bottle with Push-on NR caps 1000 ml	4 bottles & 4 caps	79060-112
DURAN® Autoclave Bottle with Push-on NR caps 2000 ml	1 bottle & 4 caps	79060-110
DURAN® Autoclave Push-on NR Cap (Replacement caps)	4 caps	76459-164
DURAN® Autoclave Push-on NR Cap (Replacement caps) Bulk pack	100 caps	79060-114
DURAN® YOUTILITY® GL 45 Bottle Tag, Orange	20	76539-390
DURAN® YOUTILITY® GL 45 Bottle Tag, Yellow	20	76264-238
DURAN® YOUTILITY® GL 45 Bottle Tag, Green	20	76264-240
DURAN® YOUTILITY® GL 45 Bottle Tag, Blue	20	76264-242
DURAN® YOUTILITY® GL 45 Bottle Tag, Purple	20	76264-244
DURAN® YOUTILITY® GL 45 Bottle Tag, Black	20	76264-246
DURAN® YOUTILITY® Blank Laboratory Labels	100	76264-248



Excellence in your hands

DWK Life Sciences SALES PARTNER



[vwr.com](https://www.vwr.com)

Any prices, product, and/or services details are current when published and subject to change. Offers subject to product availability. Void where prohibited by law or company policy. Any claims concerning products manufactured by third parties are the sole claims of the manufacturer and not those of Avantor or its affiliates. All products are subject to the terms and conditions of sale of Avantor or its affiliates, including the applicable limitations of liability contained therein. The information herein should not be considered a warranty of any kind. Please carefully review any product information sheets and safety information for the product. Trademarks are owned by Avantor Inc. or its affiliates, unless otherwise noted. | Visit [vwr.com](https://www.vwr.com) to view our privacy policy and additional disclaimers. © 2026 Avantor Inc. All rights reserved.

0826 Lit. No. 040100W