

Corning® HYPERStack® Cell Culture Vessels

Closed Systems for High Yield Cell Growth


Corning's High Yield PERFORMANCE (HYPER) Platform

The HYPERStack cell culture vessel combines the best of two Corning products: the Corning CellSTACK® and the Corning HYPERFlask® cell culture vessels.

The utilization of the gas-permeable film technology provided in the spatial footprint of the CellSTACK vessel allows the HYPERStack platform to be among the most efficient, scalable cell culture vessels for adherent cell culture available today.


Features and Benefits

- More cells – provides up to 5X the growth surface area of a traditional cell culture vessel of comparable footprint
- Closed system – no open fluid manipulations
- Scalable product – multiple size offerings support scale-up and scale-out
- Ergonomic design – easier manipulation with handling equipment and accessories
- Fixed media volume – 0.2 mL/cm² fills vessel for less volumetric waste

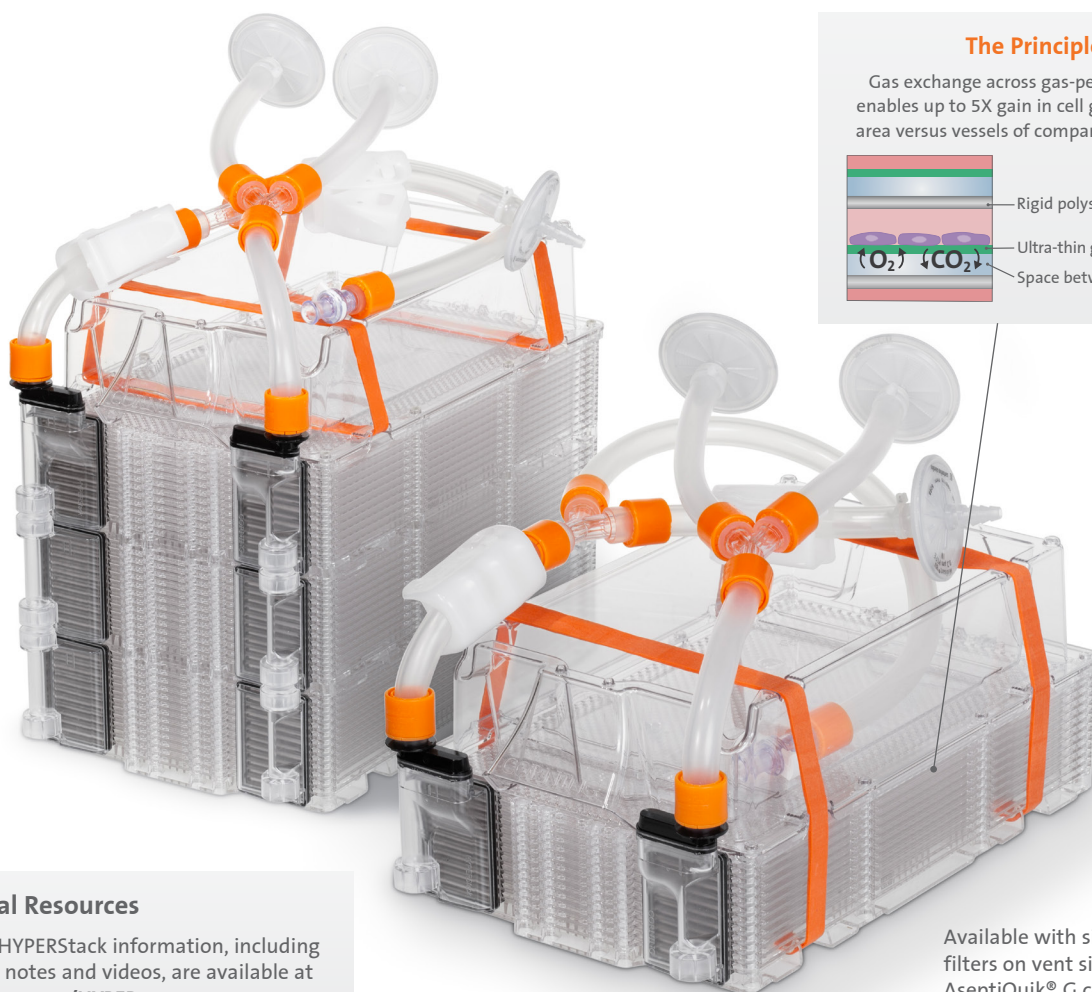

The Principle

Gas exchange across gas-permeable film enables up to 5X gain in cell growth surface area versus vessels of comparable footprint.

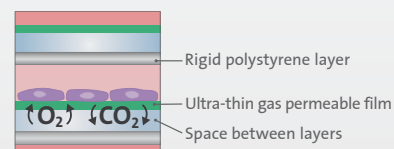


Additional Resources

Additional HYPERStack information, including application notes and videos, are available at www.corning.com/HYPER.

Available with single or dual filters on vent side; MPC or AsepticQuik® G connectors on fill side

Corning® HYPERStack® Cell Culture Vessels Improve Scale-up Efficiency

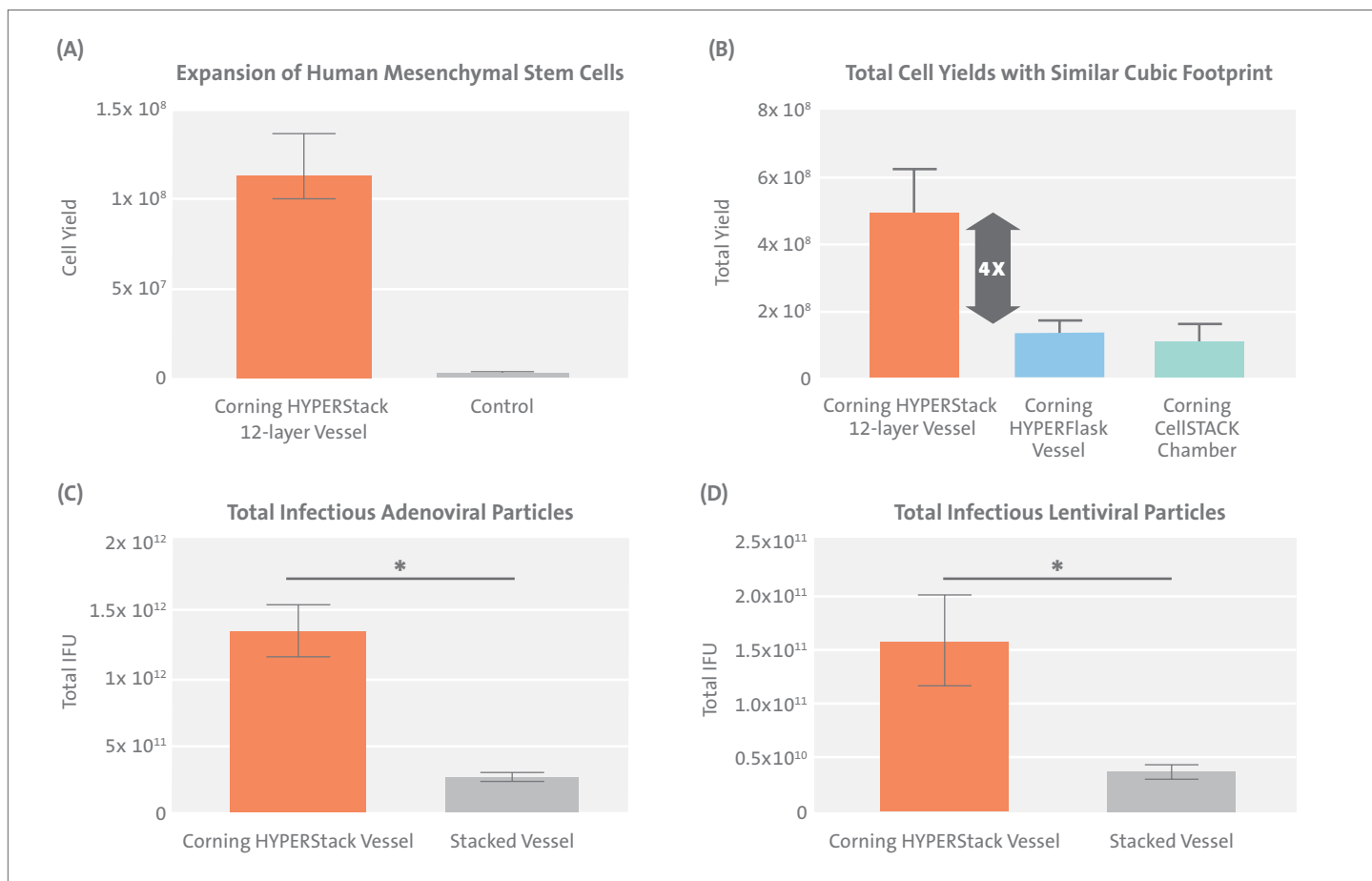

(A) Final expansion of Human Mesenchymal Stem Cells in Corning HYPERStack 12-layer or control T-flask cell culture vessels with Corning CellBIND® surface treatment using hMSC medium. (B) Total Corning HYPERStack vessel yields of Vero cell cultures were 4X greater than yields from Corning CellSTACK® vessels with similar cubic footprint. (C) Corning HYPERStack vessels generated a significantly higher amount of total infectious adenoviral particles versus stacked vessels of similar cubic footprint. Total infectious adenoviral particles were calculated based on titers and the volume of each fraction (cells and medium). (D) Corning HYPERStack vessels generated a significantly higher amount of total infectious lentiviral particles compared to stacked vessels of similar cubic footprint. *Paired t-test, p < 0.05, N=3.

The Properties



Corning HYPERStack 12-layer

Surface	6,000 cm ²
Volume	1.31L
Filled weight (approx.)	2.2 kg
Dimensions (L x W)	13.42 x 8.15 in. (342 x 207 mm)
Vessel Height	2.8 in. (71 mm)
Height with Accessory Tray	5.5 in. (140 mm)



Corning HYPERStack 36-layer

Surface	18,000 cm ²
Volume	3.92L
Filled weight (approx.)	6.6 kg
Dimensions (L x W)	13.42 x 8.15 in. (342 x 207 mm)
Height with Accessory Tray	10.97 in. (278 mm)



Ordering Information

Corning® HYPERStack® Cell Culture Vessels

VWR Cat. No.	Corning Cat. No.	Description	Growth Area (cm ²)	Qty/Pk	Qty/Cs
76463-392	20012	HYPERStack 12-layer vessel, CellBIND® surface, single vent filter, sterile	6,000	1	4
76463-396	20036	HYPERStack 36-layer vessel, CellBIND surface, single vent filter, sterile	18,000	1	2
76463-394	20013	HYPERStack 12-layer vessel, non-treated, single vent filter, sterile	6,000	1	4
76463-398	20037	HYPERStack 36-layer vessel, non-treated, single vent filter, sterile	18,000	1	2
77888-540	24012	HYPERStack 12-layer vessel, CellBIND surface, dual vent filter, sterile	6,000	1	4
77888-544	24036	HYPERStack 36-layer vessel, CellBIND surface, dual vent filter, sterile	18,000	1	2
77888-542	24013	HYPERStack 12-layer vessel, non-treated, dual vent filter, sterile	6,000	1	4
77888-546	24037	HYPERStack 36-layer vessel, non-treated, dual vent filter, sterile	18,000	1	2
77889-608	23012	HYPERStack 12-layer vessel, CellBIND surface, dual vent filter, AseptiQuik® G connector, sterile	6,000	1	4
77889-612	23036	HYPERStack 36-layer vessel, CellBIND surface, dual vent filter, AseptiQuik® G connector, sterile	18,000	1	2
77889-610	23013	HYPERStack 12-layer vessel, non-treated, dual vent filter, AseptiQuik® G connector, sterile	6,000	1	4
77889-614	23037	HYPERStack 36-layer vessel, non-treated, dual vent filter, AseptiQuik® G connector, sterile	18,000	1	2

Accessories

VWR Cat. No.	Corning Cat. No.	Description	Qty/Pk	Qty/Cs
89204-710	431518	2L Erlenmeyer flask with dip tube with 0.2 µm vent, male MPC, chemically resistant, heat sealable flexible tubing, 1/4" ID, 3/8" OD	1	3
10790-810	11501	5L Erlenmeyer flask with dip tube with 0.2 µm filter, male MPC chemically resistant, heat sealable flexible tubing, 1/4" ID, 3/8" OD	1	2
89184-632	10043	Disposable tubing set for use with 850 cm ² polystyrene roller bottle, 3/8" ID x 1/2" OD, chemically resistant, heat sealable, thermoplastic elastomer tubing, 0.2 µm filter, MPC	1	2
89184-640	431644	850 cm ² polystyrene bottle, easy grip cap, non-treated, sterile	1	40
76478-956	10047	HYPERStack Nest accessory	1	1
78068-514	11370	HYPERStack manifold, MPC connectors, 5-link, polypropylene, standard tubing lengths, sterile	1	2
78068-516*	11380*	HYPERStack manifold, MPC connectors, 5-link, polypropylene, extended tubing lengths, sterile	1	2
78068-518	11630	HYPERStack manifold, AseptiQuik® G connectors, 5-link, polypropylene, standard tubing lengths, sterile	1	2
78068-512	11640	HYPERStack manifold, AseptiQuik® G connectors, 5-link, polypropylene, extended tubing lengths, sterile	1	2

*Made to order. Please contact your local Account Representative for more information.

Warranty/Disclaimer: Unless otherwise specified, all products are for research use or general laboratory use only.* Not intended for use in diagnostic or therapeutic procedures. Not for use in humans. These products are not intended to mitigate the presence of microorganisms on surfaces or in the environment, where such organisms can be deleterious to humans or the environment. Corning Life Sciences makes no claims regarding the performance of these products for clinical or diagnostic applications. *For a listing of US medical devices, regulatory classifications or specific information on claims, visit www.corning.com/resources.

Corning's products are not specifically designed and tested for diagnostic testing. Many Corning products, though not specific for diagnostic testing, can be used in the workflow and preparation of the test at the customers discretion. Customers may use these products to support their claims. We cannot make any claims or statements that our products are approved for diagnostic testing either directly or indirectly. The customer is responsible for any testing, validation, and/or regulatory submissions that may be required to support the safety and efficacy of their intended application.

CORNING

For additional product or technical information, visit vwr.com/corning, call 1.800.932.5000, or contact your Avantor representative.



avantorsciences.com

Any prices, product, and/or services details are current when published and subject to change. Offers subject to product availability. Void where prohibited by law or company policy. Any claims concerning products manufactured by third parties are the sole claims of the manufacturer and not those of Avantor or its affiliates. All products are subject to the terms and conditions of sale of Avantor or its affiliates, including the applicable limitations of liability contained therein. The information herein should not be considered a warranty of any kind. Please carefully review any product information sheets and safety information for the product. Trademarks are owned by Avantor, Inc. or its affiliates, unless otherwise noted. | Visit vwr.com/avantorsciences.com to view our privacy policy and additional disclaimers. © 2025 Avantor, Inc. All rights reserved.

AseptiQuik® is a registered trademark of Colder Products Company.

The information contained within is accurate as of the date of publication and subject to change without notice.

For a listing of trademarks, visit www.corning.com/clstrademarks. All other trademarks are the property of their respective owners.

© 2021-2025 Corning Incorporated. All rights reserved. 10/25 CLS-CC-057 REV4 VWR

1025 Lit. No. 030478WREV