

SPECIFICATIONS

Range: –50 to 150 °C
Battery: 357A button cell
alkaline
or
357 silver-oxide (Cat No. 1039)

Features a convenient stand and hook that is specifically designed to be used in a dishwasher. Waterproof unit is designed for rough handling.

OPERATION

Press the ON/OFF button once to turn on the unit. The unit will turn on and begin reading temperature in Celsius. To switch between Fahrenheit/Celsius, press the °C/°F button. Insert unit into the desired location to begin measuring temperature.

MAXIMUM/MINIMUM

To view the Maximum and Minimum temperature the thermometer has achieved since the unit has been turned on, press the MAX/MIN button to view the Minimum temperature. Press MAX/MIN button again within 3 seconds to view the maximum temperature.

Note: Turning off the unit will reset the Maximum and Minimum memory.

HOLD FUNCTION

To “freeze” a temperature displayed on the LCD, press the HOLD button once to continuously show the current temperature. “HOLD” will flash on the display to indicate that you are currently using the HOLD function. To exit HOLD, press the HOLD button, and the current temperature will be displayed.

ALL OPERATIONAL DIFFICULTIES

If this thermometer does not function properly for any reason, replace the battery with a new high-quality battery (see “Battery Replacement” section). Low battery power can occasionally cause any number of “apparent” operational difficulties. Replacing the battery with a new fresh battery will solve most difficulties.

BATTERY REPLACEMENT

Erratic readings, a faint display, no display, or low battery icon  are all indications that the battery must be replaced. To replace the battery, with the unit facing forward, unscrew the metal housing cover counterclockwise until the front half is removed. Carefully remove the faceplate, and then gently pry the electronics component out of its metal casing. With the unit removed from its casing, turn it over to expose six screws. Using a Phillips® #000 screwdriver or equivalent, remove the six screws, gently remove the plastic casing, and then remove the exposed battery. Replace the battery with a new 1.5 volt 357 silver oxide battery (Replacement Cat. No. 1039). Make certain the positive (+) side is facing out. Replace the plastic cap onto the unit and screw the six screws back into the electronics component. Return component to its metal casing, replace the faceplate, and screw on the metal housing cover by turning it clockwise.

Technical service

Web Resources

Visit the VWR's website at www.vwr.com for:

- Français, Español, Deutsch, Italiano, Português version
- Complete technical service contact information
- Access to VWR's Online Catalogue, and information about accessories and related products
- Additional product information and special offers

Contact us for information or technical assistance contact your local VWR representative or visit. www.vwr.com.

Warranty

VWR International warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of purchase. If a defect is present, VWR will, at its option, repair, replace, or refund the purchase price of this product at no charge to you, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. For your

protection, items being returned must be insured against possible damage or loss. This warranty shall be limited to the replacement of defective products. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

**GO TO VWR.COM FOR THE LATEST NEWS,
SPECIAL OFFERS AND DETAILS OF YOUR LOCAL
VWR DISTRIBUTOR**

Cat. No. 89204-738 / 620-9100

©2015 92-4203-60 Rev. 1 011115



**TRACEABLE[®]
DISHWASHER
THERMOMETER
INSTRUCTIONS**