


Operating manual

VWR[®] MU 6100 L - Multi Meter

	EU cat. no	NA cat. no
MU 6100 L S1	665-0309	76460-494
MU 6100 L S2	665-0310	76460-496

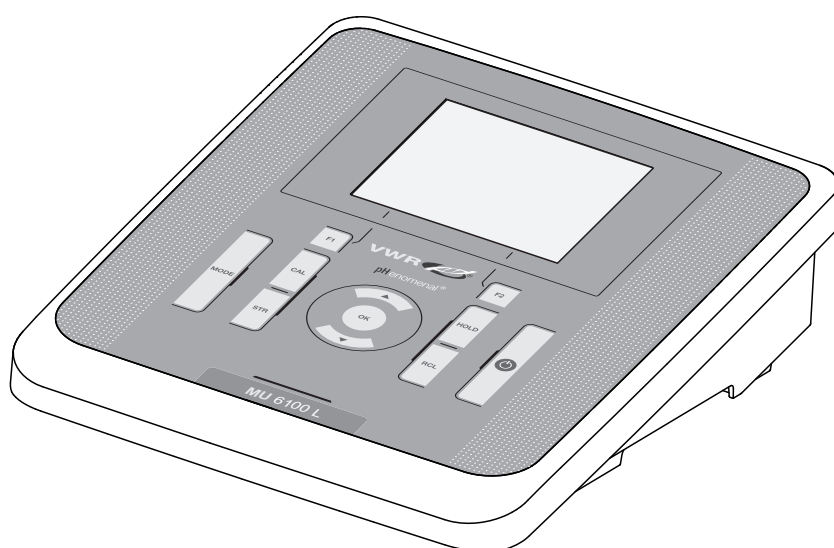

CE

UK
CA

Legal Address of Manufacturer:

Europe

VWR International BV
Researchpark Haasrode 2020
Geldenaaksebaan 464
B-3001 Leuven
+32 16 385011
<http://be.vwr.com>

United States

VWR International LLC
100 Matsonford Rd
Radnor, PA 19087
+1 800-932-5000
www.vwr.com

Importer to UK:

VWR International Ltd
Hunter Boulevard, Magna Park
Lutterworth, Leicestershire, LE17 4XN
<http://uk.vwr.com>

Country of origin: Germany.

MU 6100 L - Contents

1	Safety	7
1.1	Safety information	7
1.1.1	Safety information in the operating manual	7
1.1.2	Safety signs on the meter	7
1.1.3	Further documents providing safety information	7
1.2	Safe operation	8
1.2.1	Authorized use	8
1.2.2	Requirements for safe operation	8
1.2.3	Unauthorized use	8
2	Overview	9
2.1	MU 6100 L meter	9
2.2	Sensors	9
3	Technical data	10
3.1	Measuring ranges, resolution, accuracy	10
3.1.1	pH/ORP	10
3.1.2	Oxi	11
3.1.3	Cond.	12
3.2	General data	13
4	Commissioning	15
4.1	Scope of delivery	15
4.2	Power supply	15
4.3	Initial commissioning	15
4.3.1	Inserting the batteries	16
4.3.2	Connect the power pack	16
4.3.3	Stand	17
5	Operation	18
5.1	General operating principles	18
5.1.1	Keypad	18
5.1.2	Display	19
5.1.3	Status information (meter)	19
5.1.4	Connectors	20
5.2	Switching on the meter	20
5.3	Switching off	21
5.4	Navigation	21
5.4.1	Operating modes	21

5.4.2	Measured value display	22
5.4.3	Menus and dialogs	22
5.4.4	Elements in menus and dialogs	22
5.4.5	Navigation example 1: Setting the language	23
5.4.6	Example 2 on navigation: Setting the date and time	25
5.5	Channel display	27
5.5.1	Display of several sensors in the measuring mode	27
6	pH value	28
6.1	Measuring	28
6.1.1	Measuring the pH value	28
6.1.2	Measuring the temperature	29
6.2	Calibration	30
6.2.1	Why calibrate?	30
6.2.2	When do you have to calibrate?	30
6.2.3	Automatic calibration (AutoCal)	30
6.2.4	Manual calibration (AnyCal)	33
6.2.5	Calibration points	36
6.2.6	Calibration data	37
7	ORP voltage	39
7.1	Measuring	39
7.1.1	Measuring the ORP	39
7.1.2	Measuring the temperature	40
7.2	ORP calibration	40
8	Ion concentration	41
8.1	Measuring	41
8.1.1	Measuring the ion concentration	41
8.1.2	Measuring the temperature	42
8.2	Calibration	43
8.2.1	Why calibrate?	43
8.2.2	When to calibrate?	43
8.2.3	Calibration (ISE Cal)	43
8.2.4	Calibration standards	46
8.2.5	Calibration data	46
9	Dissolved oxygen	49
9.1	Measuring	49
9.1.1	Measuring D.O.	49
9.1.2	Measuring the temperature	50
9.2	Calibration	51
9.2.1	Why calibrate?	51
9.2.2	When to calibrate?	51

9.2.3	Calibration procedure	51
9.2.4	Calibration in water vapor-saturated air (air calibration vessel)	51
9.2.5	Calibrating with a comparison measurement (OxiComp)	52
9.2.6	Calibration data	54
10	Conductivity	56
10.1	Measuring	56
10.1.1	Measuring the conductivity.	56
10.1.2	Measuring the temperature	57
10.2	Calibration	58
10.2.1	Why calibrate?	58
10.2.2	When to calibrate?	58
10.2.3	Determining the cell constant (calibration in control standard)	58
10.2.4	Calibration data	59
11	Settings	61
11.1	Measurement settings	61
11.1.1	Settings for pH measurements.	61
11.1.2	Buffer sets for calibration	62
11.1.3	Calibration interval	64
11.1.4	Settings for ORP measurements	64
11.1.5	Settings for D.O. sensors	65
11.1.6	Calibration interval	66
11.1.7	Settings for conductivity measuring cells	67
11.1.8	Calibration interval	69
11.1.9	Temperature compensation	69
11.1.10	Setting the TDS factor	70
11.2	Sensor-independent settings	71
11.2.1	<i>System</i>	71
11.2.2	<i>Data storage</i>	72
11.2.3	<i>Automatic Stability control</i>	72
11.3	Reset.	72
11.3.1	Resetting the measurement settings.	72
11.3.2	Resetting the system settings	74
12	Data memory	75
12.1	Manual storage	75
12.2	Automatic data storage at intervals	76
12.3	Measurement data memory	78
12.3.1	Editing the measurement data memory	78
12.3.2	Erasing the measurement data memory	80
12.3.3	Measurement dataset	80
12.3.4	Memory locations	80

13	Transmitting data (USB interface)	81
13.1	Options for data transmission	81
13.2	Connecting a PC	82
13.3	MultiLab Importer	82
14	Maintenance, cleaning, disposal, accessories	83
14.1	Maintenance	83
14.1.1	General maintenance activities	83
14.1.2	Replacing the batteries	83
14.2	Cleaning	84
14.3	Packing	84
14.4	Disposal	84
14.5	Accessories	85
14.5.1	General information	85
14.5.2	pH / ORP	86
14.5.3	ISE	88
14.5.4	Conductivity	88
14.5.5	D.O.	89
15	What to do if...	90
15.1	pH/ORP	90
15.2	Dissolved oxygen	91
15.3	Conductivity	93
15.4	General information	94
16	Firmware update	95
17	Glossary	96
18	Index	99
19	Technical service	108
20	Warranty	108
21	Compliance with local laws and regulations	108

1 Safety

1.1 Safety information

1.1.1 Safety information in the operating manual

This operating manual provides important information on the safe operation of the meter. Read this operating manual thoroughly and make yourself familiar with the meter before putting it into operation or working with it. The operating manual must be kept in the vicinity of the meter so you can always find the information you need.

Important safety instructions are highlighted in this operating manual. They are indicated by the warning symbol (triangle) in the left column. The signal word (e.g. "CAUTION") indicates the level of danger:

**WARNING**

indicates a possibly dangerous situation that can lead to serious (irreversible) injury or death if the safety instruction is not followed.

**CAUTION**

indicates a possibly dangerous situation that can lead to slight (reversible) injury if the safety instruction is not followed.

NOTE

indicates a possibly dangerous situation where goods might be damaged if the actions mentioned are not taken.

1.1.2 Safety signs on the meter

Note all labels, information signs and safety symbols on the meter and in the battery compartment. A warning symbol (triangle) without text refers to safety information in this operating manual.

1.1.3 Further documents providing safety information

The following documents provide additional information, which you should observe for your safety when working with the measuring system:

- Operating manuals of sensors and other accessories
- Safety datasheets of calibration or maintenance accessories (such as buffer solutions, electrolyte solutions, etc.)

1.2 Safe operation

1.2.1 Authorized use

This meter is authorized exclusively for pH, ISE, ORP, dissolved oxygen and conductivity measurements in a laboratory environment.

Only the operation and running of the meter according to the instructions and technical specifications given in this operating manual is authorized (see section 3 TECHNICAL DATA, page 10).

Any other use is considered unauthorized.

1.2.2 Requirements for safe operation

Note the following points for safe operation:

- The meter may only be operated according to the authorized use specified above.
- The meter may only be supplied with power by the energy sources mentioned in this operating manual.
- The meter may only be operated under the environmental conditions mentioned in this operating manual.
- The meter may only be opened if this is explicitly described in this operating manual (example: Inserting the batteries).

1.2.3 Unauthorized use

The meter must not be put into operation if:

- it is visibly damaged (e.g. after being transported)
- it was stored under adverse conditions for a lengthy period of time (storing conditions, see section 3 TECHNICAL DATA, page 10).

2 Overview

2.1 MU 6100 L meter

The MU 6100 L meter enables you to perform measurements (pH, U, conductivity, dissolved oxygen)) quickly and reliably.

The MU 6100 L provides the maximum degree of operating comfort, reliability and measuring certainty for all applications.

2.2 Sensors

A measuring system ready to measure consists of the MU 6100 L meter and a suitable sensor.

Suitable sensors are pH electrodes, ORP electrodes, D.O. sensors and conductivity measuring cells.

3 Technical data

3.1 Measuring ranges, resolution, accuracy

3.1.1 pH/ORP

Measuring ranges, resolution

Variable	Measuring range	Resolution
pH	-2.0 ... +20.0	0.1
	-2.00 ... +20.00	0.01
	- 2.000 ... + 19.999	0.001
U [mV]	-2500 ... +2500	1
	-1200.0 ... +1200.	0.1
T [°C]	-5.0 ... +105.0	0.1
T [°F]	23.0 ... +221.0	0.1

Manual temperature input

Variable	Range	Increment
T _{manual} [°C]	-25 ... +130	1
T _{manual} [°F]	-13 ... +266	1

Accuracy (± 1 digit)

Variable	Accuracy	Temperature of the test sample
pH / range *		
-2.0 ... +20.0	± 0.1	+15 °C ... +35 °C
-2.00 ... +20.00	± 0.01	+15 °C ... +35 °C
- 2.000 ... + 19.999	± 0.005	+15 °C ... +35 °C
U [mV] / range		
-2500 ... +2500	± 1	+15 °C ... +35 °C
-1200.0 ... +1200.	± 0.3	+15 °C ... +35 °C
T [°C] / temperature sensor		
NTC 30	± 0.1	
PT 1000	± 0.1	

* when measuring in a range of ± 2 pH around a calibration point



The accuracy values specified here apply exclusively to the meter. The accuracy of the electrodes and buffer solutions has to be taken into account additionally.

3.1.2 Oxi

Measuring ranges, resolution (depending on the sensor)

Variable	Measuring range	Resolution
Concentration [mg/l]	0 ... 20.00 (0 ... 20.0) 20.0 ... 90.0 (20 ... 90)	0.01 (0.1) 0.1 (1)
Saturation [%]	0 ... 200.0 (0 ... 600) 0 ... 600	0.1 (1) 1
D.O. partial pressure [mbar]	0 ... 200.0 (0 ... 1250) 0 ... 1250	0.1 (1) 1
T [°C]	0 ... 50,0	0.1

Accuracy (± 1 digit)

Variable	Accuracy
Concentration [mg/l]	± 0.5 % of measured value at ambient temperature of +5 °C ... +30 °C
Saturation [%]	± 0.5% of measured value when measuring in the range of ± 10 K around the calibration temperature
D.O. partial pressure [mbar]	± 0.5 % of measured value at ambient temperature of +5 °C ... +30 °C
T [°C] / temperature sensor	
NTC 30	± 0.1
PT 1000	± 0.1

Correction functions

Temperature compensation	Accuracy better than 2 % at 0 ... +40 °C
Salinity correction	0 ... 70.0 SAL
Air pressure correction	Automatic through integrated pressure sensor in the range of 500 ... 1100 mbar



The accuracy values specified here apply exclusively to the meter. The accuracy of the D.O. sensors has also to be taken into account.

3.1.3 Cond

Measuring ranges,
resolution
(depending on the sen-
sor)

Variable	Measuring range	Resolution
χ [μS/cm]	0.000 ... 1.999* 0.00 ... 19.99** 0.0 ... 199.9 200 ... 1999	0.001 0.01 0.1 1
χ [mS/cm]	2.00 ... 19.99 20.0 ... 199.9 200 ... 1000	0.01 0.1 1
ρ (Resistivity) [Ohm*cm]	1.000 ... 1.999 2.00 ... 19.99 20.0 ... 199.9 200 ... 1999	0.001 0.01 0.1 1
ρ (Resistivity) [kOhm*cm]	2.00 ... 19.99 20.0 ... 199.9 200 ... 1999	0.01 0.1 1
ρ (Resistivity) [MOhm*cm]	2.00 ... 19.99** 20.0 ... 199.9*	0.01 0.1
SAL	0.0 ... 70.0 according to the IOT table	0.1
TDS	0 ... 1999 mg/l 2.00 ... 19.99 g/l 20.0 ... 199.9 g/l	1 0.01 0.1
T [°C]	-5.0 ... +105.0	0.1
T [°F]	+23.0 ... +221.0	0.1

* only possible with cells of the cell constant, 0.010 cm⁻¹

** only possible with cells of the cell constant, 0.010 cm⁻¹ or 0.090 ... 0.110 cm⁻¹

Cell constants

Cell constant C	Values
Can be calibrated in the ranges	0.450 ... 0.500 cm ⁻¹ 0.800 ... 0.880 cm ⁻¹
Adjustable	0.090 ... 0.110 cm ⁻¹ 0.010 cm ⁻¹ (fixed) 0,250 ... 1,000 cm ⁻¹

Reference temperature

Reference temperature	Values
Adjustable	20 °C (Tref20) 25 °C (Tref25)

Accuracy (± 1 digit)

Variable	Accuracy	Temperature of the test sample
χ and ρ / temperature compensation		
None (Off)	± 0.5 %	
Nonlinear (nLF)	± 0.5 %	0 °C ... +35 °C according to EN 27 888
	± 0.5 %	+35 °C ... +50 °C enhanced nLF function
Linear (lin)	± 0.5 %	+10 °C ... +75 °C
SAL / range		
0.0 ... 42.0	± 0.1	+5 °C ... +25 °C
	± 0.2	+25 °C ... +30 °C
TDS [mg/l]		
	± 0.5 %	
T [°C] / temperature sensor		
NTC 30	± 0.1	
PT 1000	± 0.1	



The accuracy values specified here apply exclusively to the meter. The accuracy of the measuring cell has also to be taken into account.



Further data are given in the documentation of your sensor.

3.2 General data

Dimensions

ca. 240 x 190 x 80 mm

Weight

Approx. 1.0 kg

Mechanical structure

Type of protection IP 43

Electrical safety

Protective class III

Test certificates

CE

Ambient conditions

Storage -25 °C ... +65 °C

Operation -10 °C ... +55 °C

Admissible relative humidity
 Yearly mean: < 75 %
 30 days/year: 95 %
 Other days: 85 %

Power supply

Batteries	4 x 1.5 V alkali-manganese batteries, type AA
Rechargeable batteries	4 x 1.2 V NiMH rechargeable batteries, type AA (no charging function)
Operational life	Up to 1000 h without / 150 h with illumination
Power pack (charging device)	Input: 100 ... 240 V ~ / 50 ... 60 Hz / 0.5 A Output: 9 V = / 1,1 A Connection max. overvoltage category II Primary plugs contained in the scope of delivery: Euro, US, UK and Australian.

pH sensor input

Input resistance	$> 5 \cdot 10^{12} \text{ ohm}$
Input current	$< 1 \cdot 10^{-12} \text{ A}$

USB interface

Type	USB 1.1 USB B (device), data output
Baud rate	Adjustable: 1200, 2400, 4800, 9600, 19200 Baud
Data bits	8
Stop bits	2
Parity	None
Handshake	RTS/CTS
Cable length	Max. 3 m

Guidelines and norms used

EMC	EU directive 2014/30/EU EN 61326-1 FCC Class A
Meter safety	EU directive 2014/35/EU EN 61010-1
IP protection class	EN 60529
RoHS	EU directive 2011/65/EU

4 Commissioning

4.1 Scope of delivery

- MeterMU 6100 L
- 4 batteries 1.5 V Mignon type AA
- Power pack
- Stand
- Stand holder
- Short instructions
- CD-ROM with
 - USB drivers
 - detailed operating manual
 - Software MultiLab Importer

4.2 Power supply

The MU 6100 L is supplied with power in the following ways:


- Mains operation with the supplied power pack
- Battery operation (4 x alkaline manganese batteries, type AA)
- USB operation via a connected USB-B cable

4.3 Initial commissioning

Perform the following activities:

- Insert the supplied batteries
- For mains operation: Connect the power pack
- If necessary, mount the stand
- Switch on the meter
(see section 5.2 SWITCHING ON THE METER, page 20)
- Set the date and time
(see section 5.4.6 EXAMPLE 2 ON NAVIGATION: SETTING THE DATE AND TIME, page 25)

4.3.1 Inserting the batteries

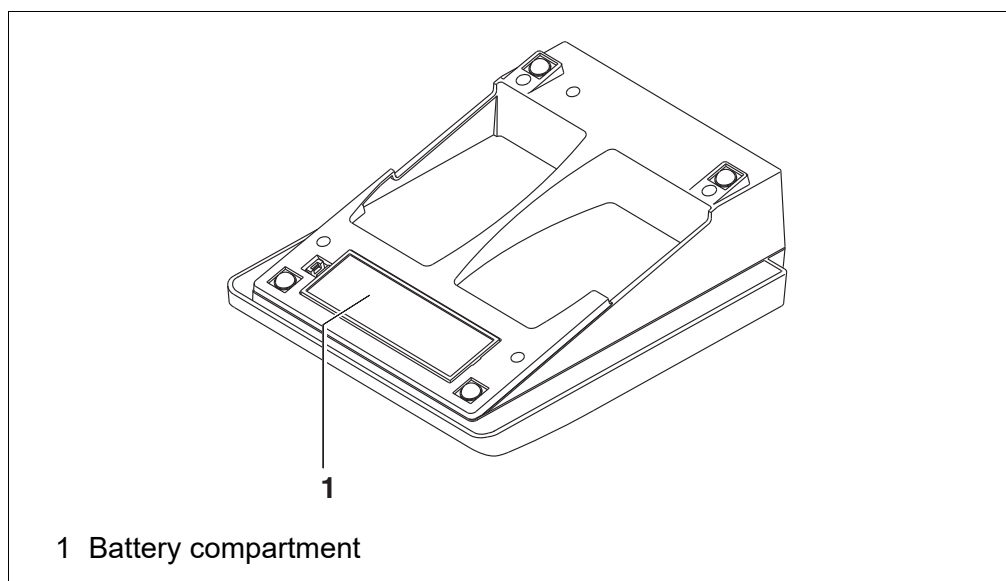

1. Open the battery compartment (1) on the underside of the meter.

**CAUTION**

Make sure that the poles of the batteries are positioned correctly.

The $\pm$ signs on the batteries must correspond to the $\pm$ signs in the battery compartment.



You can operate the meter either with normal batteries or with rechargeable batteries (Ni-MH). In order to charge the batteries, an external charging device is required.

2. Place four batteries (type Mignon AA) in the battery compartment.
3. Close the battery compartment.
4. Set the date and time
(see section 5.4.6 EXAMPLE 2 ON NAVIGATION: SETTING THE DATE AND TIME, page 25).

4.3.2 Connect the power pack

**CAUTION**

The line voltage at the operating site must lie within the input voltage range of the original power pack (see section 3 TECHNICAL DATA, page 10).



CAUTION
Use original power packs only.


1. Connect the plug of the power pack to the socket for the power pack on the MU 6100 L.
2. Connect the original power pack to an easily accessible power outlet.

4.3.3 Stand

The stand base can be mounted at the right side of the meter.

1. Screw the stand base to the underside of the meter.
2. Insert the stand rod in the stand base.

An arrangement of the meter with the stand may look as follows:

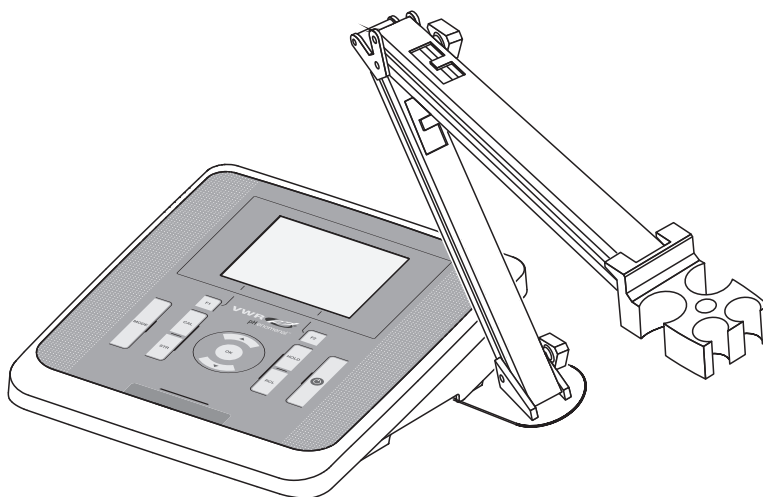

5 Operation

5.1 General operating principles


5.1.1 Keypad

In this operating manual, keys are indicated by brackets <.> .

The key symbol (e.g. **<OK>**) generally indicates a short keystroke (under 2 sec) in this operating manual. A long keystroke (approx. 2 sec) is indicated by the underscore behind the key symbol (e.g. **<OK__>**).

<F1> : <F1__> : <F2> : <F2__> :	Softkeys providing situation dependent functions, e.g.: <F2> /[USB output]: Outputs data to the USB interface <F2__> /[USB output]: Configures the automatic data output to the USB interface
<On/Off> :	Switches the meter on or off
<MODE> :	Selects the measured parameter / Quits the settings
<CAL> : <CAL__> :	Calls up the calibration procedure Displays the calibration data
<STR> : <STR__> :	Saves a measured value manually Opens the menu for the automatic save function
<RCL> : <RCL__> :	Displays the manually stored measured values Displays the automatically stored measured values
<▲><▼> : <▲__><▼__> :	Menu control, navigation Increments, decrements values Increments, decrements values continuously
<OK> : <OK__> :	Opens the menu for measurement settings / confirms entries Opens the menu for system settings
<HOLD>	Freezes the measured value (HOLD function)


5.1.2 Display

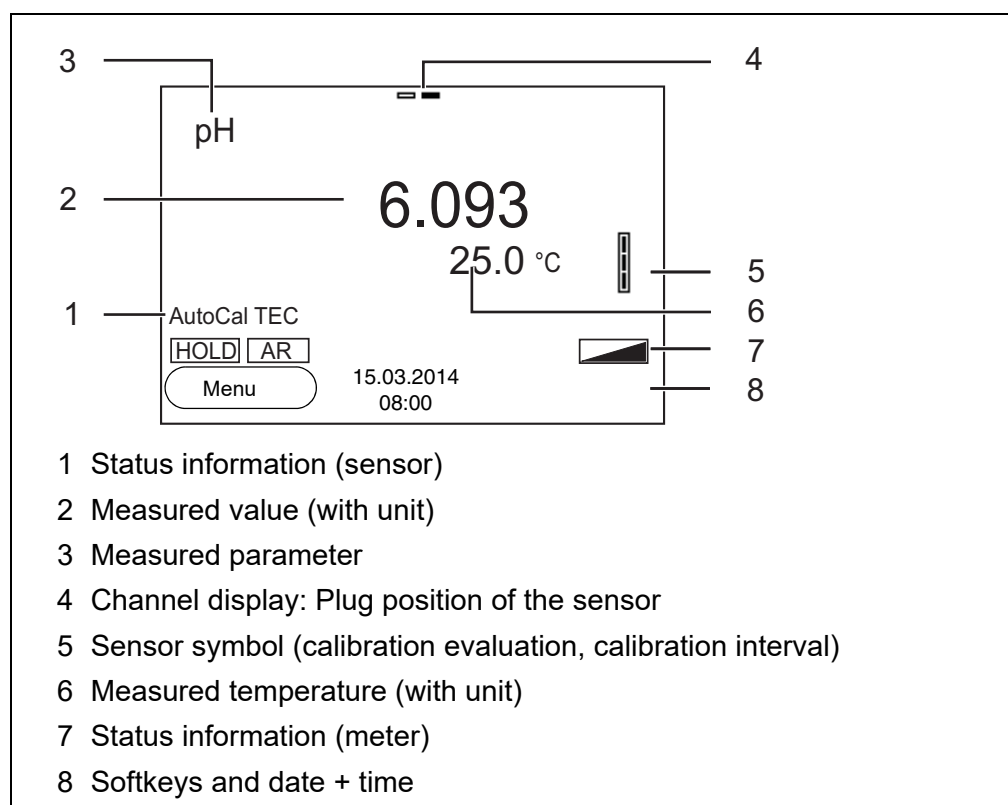

5.1.3 Status information (meter)

AR	Stability control (AutoRead) is active (calibration)
HOLD	Measured value is frozen (<HOLD> key)
	Batteries are almost empty
	Data are automatically output to the USB-B interface at intervals

5.1.4 Connectors

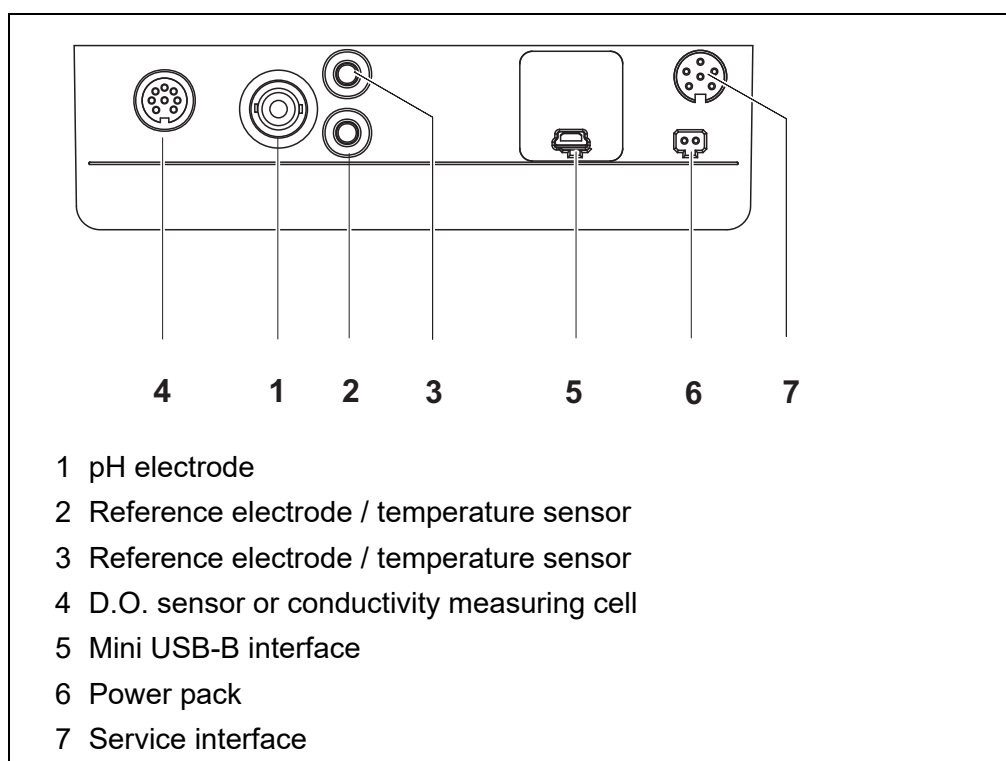


CAUTION

Only connect sensors to the meter that cannot return any voltages or currents that are not allowed (> SELV and > current circuit with current limiting).
 Almost all customary sensors fulfill these conditions.

5.2 Switching on the meter

1. Switch the meter on with **<On/Off>**.
 The meter performs a self-test.
 The display shows the manufacturer's logo while the self-test is being performed.
 The measured value display appears.

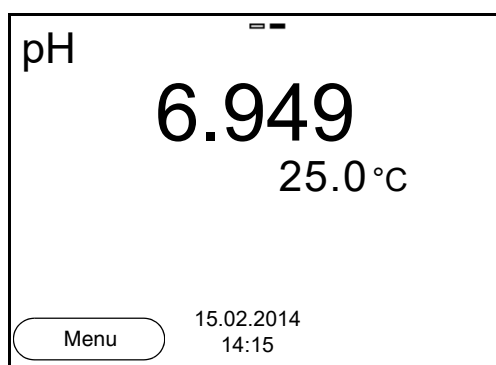

5.3 Switching off

1. Switch the printer off with **<On/Off>**.

Automatic shut-off function

The instrument has an automatic shut-off function in order to save the batteries (see section 11.2.1 SYSTEM, page 71). The automatic shut-off function switches off the meter if no key is pressed for an adjustable period.

The automatic shut-off function is not active

- if the communication cable is connected
- if the power pack is connected
- if the *Automatic data storage* function is active, or with automatic data transmission

Display illumination

The meter automatically switches off the display illumination if no key is pressed for 30 seconds. The illumination is switched on with the next keystroke again.

You can also generally switch the display illumination on or off (see section 11.2.1 SYSTEM, page 71).

5.4 Navigation

The principles of navigation in menus and dialogs are explained in the following sections.

5.4.1 Operating modes

The instrument has the following operating modes:

Operating mode	Explanation
Measuring	The measurement data of the connected sensor are shown in the measured value display
Calibration	The course of a calibration with calibration information, functions and settings is displayed
Storing in memory	The meter stores measuring data automatically or manually
Transmitting data	The meter transmits measuring data and calibration records to a USB-B interface automatically or manually.
Setting	The system menu or a sensor menu with submenus, settings and functions is displayed

Only those displays and functions are available in the active operating mode that are currently being required.

5.4.2 Measured value display

In the measured value display, open the setting menus with **<OK>**. The current functions of the softkeys are shown on the display.

- Use **<OK>** (short pressure) to open the menu for calibration and measurement settings for the displayed measured parameter.
- Use **<OK__ >** (long keystroke (approx. 2 s) to open the *Storage & config* menu with the sensor-independent settings.

Use the keys of the keypad to carry out further functions such as storage or calibration (see section 5.1.1 KEYPAD, page 18). These functions are not available in other operating situations.

5.4.3 Menus and dialogs


The menus for settings and dialogs in procedures contain further subelements.

- To select a subelement, use the **<▲><▼>** keys. The current selection is displayed with a frame.
- To make further settings, switch to the next higher menu level with **<F1>**[Back].
- Use **<MODE>** to return to the measured value display.

5.4.4 Elements in menus and dialogs

- Submenus

The name of the submenu is displayed at the upper edge of the frame. Submenus are opened by confirming with **<OK>**. Example:

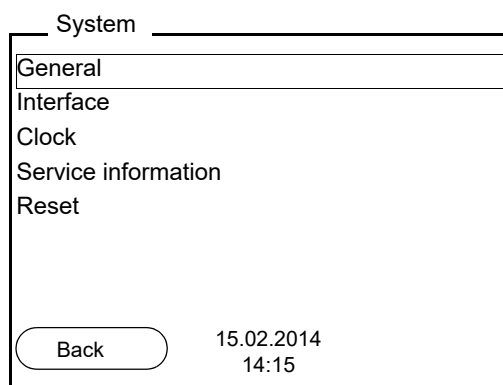

- Settings

Settings are indicated by a colon. The current setting is displayed on the right-hand side. The setting mode is opened with **<OK>**. Subsequently, the setting can be changed with **<▲><▼>** and **<OK>**. Example:

General	
Language:	Deutsch
Beep:	Off
Illumination:	On
Contrast:	50 %
Switchoff time:	1 h
Temperature unit:	°C
Stability control:	On
Back 15.02.2014 14:15	

- **Functions**

Functions are designated by the name of the function. They are immediately carried out by confirming with **<OK>**.

Example: Display the *Calibration record* function.

pH	
Calibration record	
Calibration data storage	
Buffer:	TEC
One point calibration:	Yes
Calibration interval:	7 d
Unit for slope:	mV/pH
<i>i</i> 4.00 7.00 10.00 (20 °C)	
Back 15.02.2014 14:15	


- **Messages**

Information is marked by the **i** symbol. It cannot be selected. Example:


pH	
Calibration record	
Calibration data storage	
Buffer:	TEC
One point calibration:	Yes
Calibration interval:	7 d
Unit for slope:	mV/pH
<i>i</i> 4.00 7.00 10.00 (20 °C)	
Back 15.02.2014 14:15	

5.4.5 Navigation example 1: Setting the language


1. Press the **<On/Off>** key.
The measured value display appears.
The instrument is in the measuring mode.

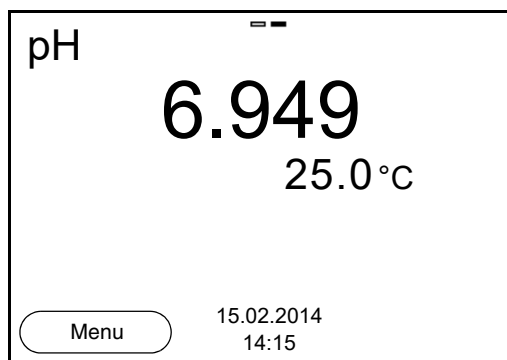


2. Open the *Storage & config* menu with **<OK>**. The instrument is in the setting mode.

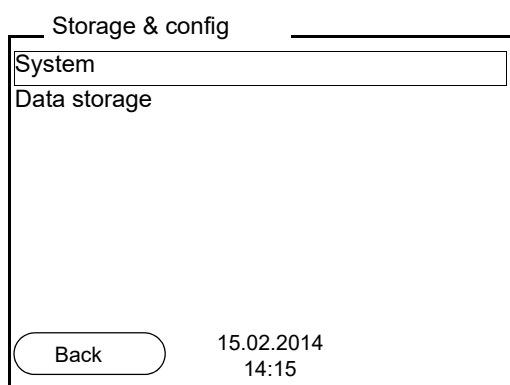


3. Select the *System* submenu with **<▲><▼>**. The current selection is displayed with a frame.
4. Open the *System* submenu with **<OK>**.

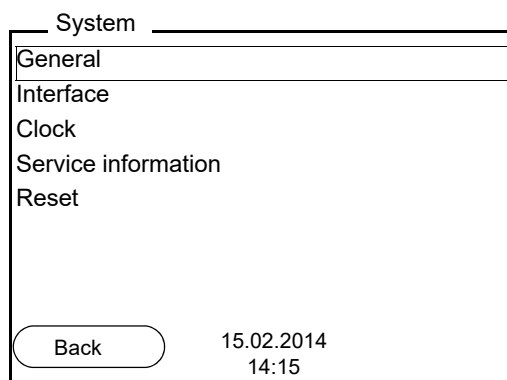

5. Select the *General* submenu with **<▲><▼>**. The current selection is displayed with a frame.
6. Open the *General* submenu with **<OK>**.

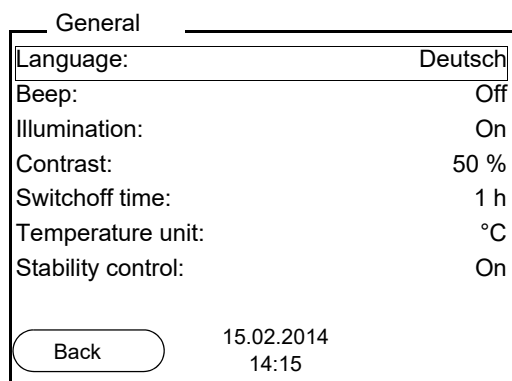

7. Open the setting mode for the *Language* with **<OK>**.

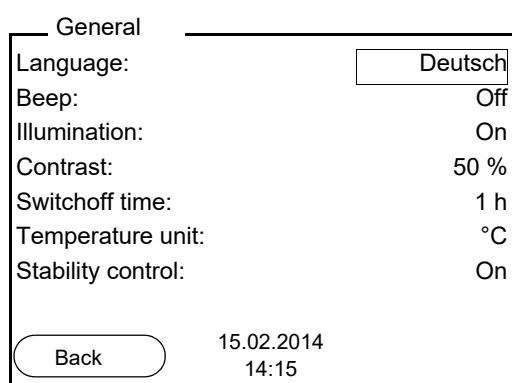

8. Select the required language with **<▲><▼>**.
9. Confirm the setting with **<OK>**.
The meter switches to the measuring mode.
The selected language is active.

5.4.6 Example 2 on navigation: Setting the date and time

The meter has a clock with a date function. The date and time are shown in the measured value display.

When storing measured values and calibrating, the current date and time are automatically stored as well.

The correct setting of the date and time and date format is important for the following functions and displays:

- Current date and time
- Calibration date
- Identification of stored measured values.

Therefore, check the time at regular intervals.



After a fall of the supply voltage (empty batteries), the date and time are reset.

The date format can be switched from the display of day, month, year (*dd.mm.yyyy*) to the display of month, day, year (*mm/dd/yyyy* or *mm.dd.yyyy*).


1. In the measured value display:
Open the *Storage & config* menu with **<OK>**.
The instrument is in the setting mode.
2. Select and confirm the *System / Clock* menu with **<▲><▼>** and **<OK>**.
The setting menu for the date and time opens up.
3. Select and confirm the *Time* menu with **<▲><▼>** and **<OK>**.
The hours are highlighted.

Clock	
Date format:	dd.mm.yyyy
Date:	15.02.2014
Time:	14:15:25
Back 15.02.2014 14:15	

4. Change and confirm the setting with **<▲><▼>** and **<OK>**.
The minutes are highlighted.
5. Change and confirm the setting with **<▲><▼>** and **<OK>**.
The seconds are highlighted.
6. Change and confirm the setting with **<▲><▼>** and **<OK>**.
The time is set.
7. If necessary, set the *Date* and *Date format*. The setting is made similarly to that of the time.
8. To make further settings, switch to the next higher menu level with **[Back]<F1>**.
or
Switch to the measured value display with **<MODE>**.
The instrument is in the measuring mode.

5.5 Channel display

The MU 6100 L manages the connected sensors and displays which sensor is plugged to which connection.

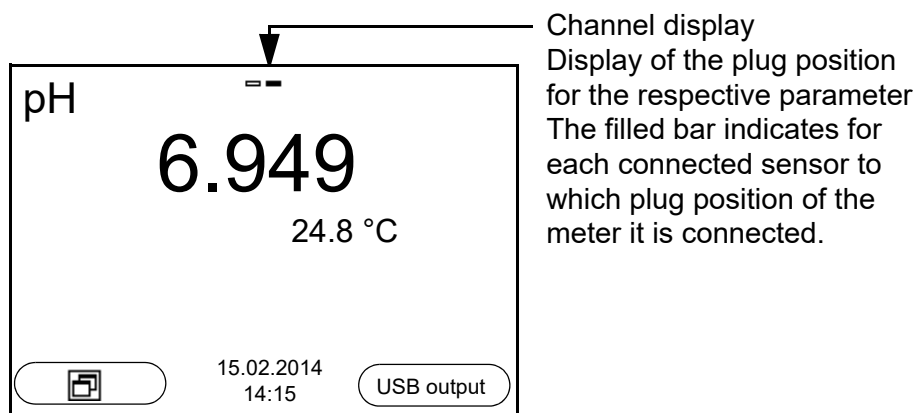


5.5.1 Display of several sensors in the measuring mode

The measured values of the connected sensors can be displayed in the following ways:

- Clear display of all connected sensors
- Detailed display of one sensor

With the softkey you can very easily switch between the two display types. The suitable softkey is displayed depending on the operating situation.

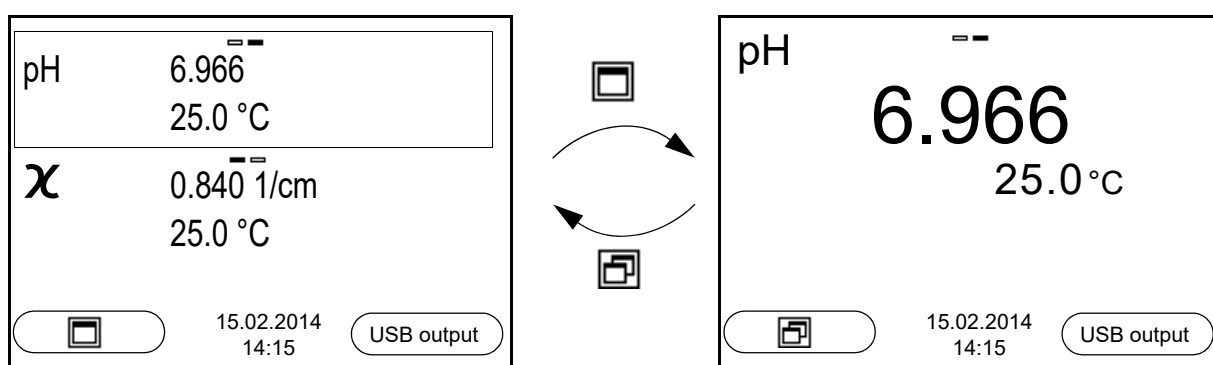

6 pH value

6.1 Measuring

6.1.1 Measuring the pH value


NOTE

When connecting an grounded PC/printer, measurements cannot be performed in grounded media as the values would be incorrect. The USB interface is not galvanically isolated.



To ensure the high measurement accuracy of the measuring system, always measure with a calibrated electrode (see section 6.2 CALIBRATION, page 30).

1. Connect the pH electrode to the meter.
2. If necessary, select the measured parameter with **<MODE>**.
3. When measuring without temperature sensor:
 - Temper the test sample, or measure the current temperature.
 - Enter the temperature value in the menu.
4. Immerse the pH electrode in the test sample.
The measured value is checked for stability (automatic stability control).
The display of the measured parameter flashes.
5. Wait for a stable measured value.
The display of the measured parameter no longer flashes.

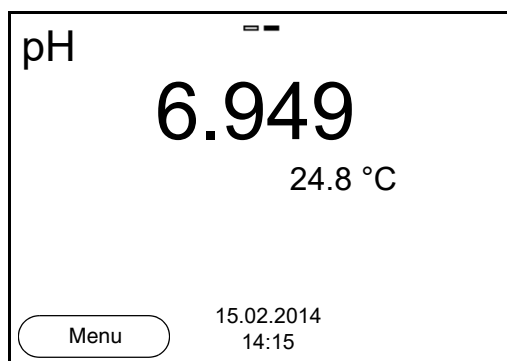

Stability control (AutoRead)

The stability control function (*AutoRead*) continually checks the stability of the measurement signal. The stability has a considerable impact on the reproducibility of measured values (see section 11.2.3 AUTOMATIC STABILITY CONTROL, page 72).

Criteria for a stable measured value

The *Stability control* function checks whether the measured values are stable

within the monitored time interval.

Measured parameter	Time interval	Stability in the time interval
pH value	15 seconds	Δ : better than 0.01 pH
Temperature	15 seconds	Δ : better than 0.5 °C

The minimum duration until a measured value is assessed as stable is the monitored time interval. The actual duration is mostly longer.

Freezes the measured value (HOLD function)

With the HOLD function, you can freeze the current measured value. The displayed measured value stops changing until you switch the HOLD function off.

1. Freeze the measured value with **<HOLD>**.
The [HOLD] status indicator is displayed.
2. Release the frozen measured value again with **<HOLD>**.
The HOLD function is switched off.
The [HOLD] status display disappears.

6.1.2 Measuring the temperature

For reproducible pH measurements, it is essential to measure the temperature of the test sample.

You have the following options to measure the temperature:

- Automatic measurement of the temperature with the temperature sensor (NTC30 or Pt1000) integrated in the sensor.
- Manual determination and input of the temperature.

The measuring instrument recognizes whether a suitable sensor is connected and automatically switches on the temperature measurement.

The display of the temperature indicates the active temperature measuring mode:

Temperature sensor	Resolution of the temp. display	Temp. measurement
yes	0.1 °C	Automatic with temperature sensor
-	1 °C	Manual

If you wish to measure (or calibrate) without temperature sensor, proceed as follows:

1. Measure the current temperature of the test sample.
2. In the **<OK>/pH/Man. temperature** menu, set the temperature value with **<▲><▼>**.

6.2 Calibration

6.2.1 Why calibrate?

pH electrodes age. This changes the zero point (asymmetry) and slope of the pH electrode. As a result, an inexact measured value is displayed. Calibration determines and stores the current values of the zero point and slope of the electrode.

Thus, you should calibrate at regular intervals.

6.2.2 When do you have to calibrate?

- After connecting a sensor
- Routinely within the framework of the company quality assurance
- When the calibration interval has expired

6.2.3 Automatic calibration (AutoCal)

Make sure that in the sensor menu, *Buffer* menu, the buffer set is correctly selected (see 11.1.1 SETTINGS FOR PH MEASUREMENTS, PAGE 61).


Use any one to five buffer solutions of the selected buffer set in ascending or descending order.

Below, calibration with Technical buffers (*TEC*) is described. When other buffer sets are used, other nominal buffer values are displayed. Apart from that, the procedure is identical.




If single-point calibration was set in the menu, the calibration procedure is automatically finished with the measurement of buffer solution 1 and the calibration record is displayed.

1. Connect the pH electrode to the meter.
2. When measuring without temperature sensor:
 - Temper the buffer solution, or measure the current temperature.
 - Enter the temperature value in the menu.
3. In the measured value display, select the measured parameter pH or mV with **<MODE>**.
4. Start the calibration with **<CAL>**.
The calibration display for the first buffer appears (voltage display).

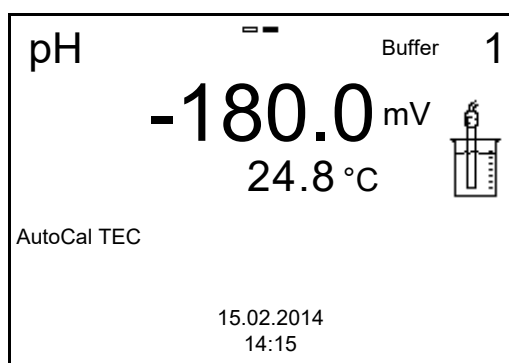

5. Thoroughly rinse the electrode with deionized water.
6. Immerse the electrode in the first buffer solution.
7. When measuring without temperature sensor:
 - Temper the buffer solution, or measure the current temperature.
 - Enter the temperature value in the menu.
8. Start the measurement with **<OK>**.
The measured value is checked for stability (stability control). The [AR] status indicator is displayed. A progress bar is displayed and the display of the measured parameter flashes.

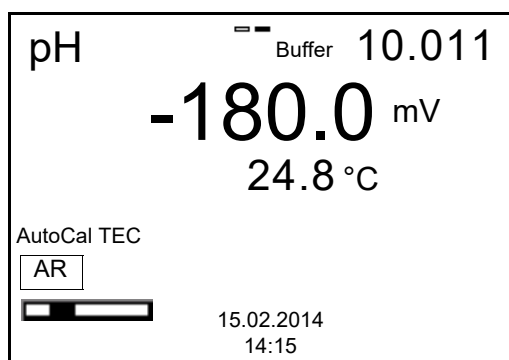

9. Wait for the end of the measurement with stability control or accept the calibration value with **<OK>**.
The calibration display for the next buffer appears (voltage display).
10. If necessary, finish the calibration procedure as a single-point calibration with **<MODE>**.
The calibration record is displayed.




For **single-point calibration**, the instrument uses the Nernst slope (-59.2 mV/pH at 25 °C) and determines the zero point of the electrode.

Continuing with two-point calibration

11. Thoroughly rinse the electrode with deionized water.
12. Immerse the electrode in the second buffer solution.


13. When measuring without temperature sensor:
 - Temper the buffer solution, or measure the current temperature.
 - Enter the temperature value in the menu.
14. Start the measurement with **<OK>**.
 The measured value is checked for stability (stability control). The [AR] status indicator is displayed. A progress bar is displayed and the display of the measured parameter flashes.

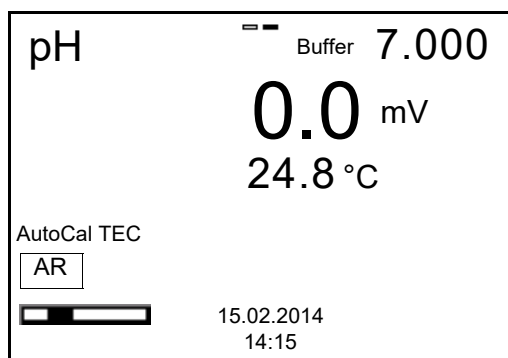

15. Wait for the measurement with stability control to be completed or terminate the stability control and take over the calibration value with **<OK>**.
 The calibration display for the next buffer appears (voltage display).
16. If necessary, finish the calibration procedure as a two-point calibration with **<MODE>**.
 The calibration record is displayed.

Continuing with three- to five-point calibration

17. Thoroughly rinse the electrode with deionized water.
18. Immerse the electrode in the next buffer solution.
19. When measuring without temperature sensor:
 - Temper the buffer solution, or measure the current temperature.
 - Enter the temperature value in the menu.
20. Start the measurement with **<OK>**.
 The measured value is checked for stability (stability control). The [AR] status indicator is displayed. A progress bar is displayed and the display of the measured parameter flashes.

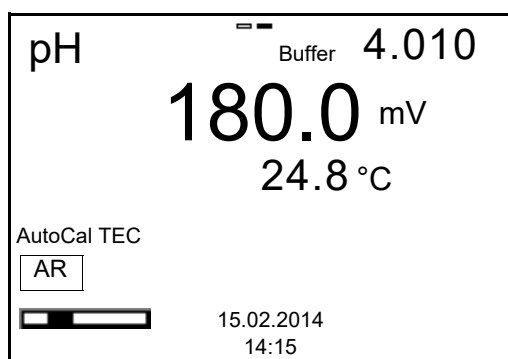

21. Wait for the measurement with stability control to be completed or terminate the stability control and take over the calibration value with **<OK>**. The calibration display for the next buffer appears (voltage display).
22. If necessary, use **<MODE>** to finish calibration or Continue calibrating using the next buffer with **<OK>**.



Calibration is automatically completed after the last buffer of a buffer set has been measured. Then the calibration record is displayed.

The calibration line is determined by linear regression.

6.2.4 Manual calibration (AnyCal)


Make sure that in the sensor menu, *Buffer* menu, the AnyCal buffer set is correctly selected (see section 11.1.1 SETTINGS FOR PH MEASUREMENTS, page 61).

Use any one to five buffer solutions in ascending or descending order.




If single-point calibration was set in the menu, the calibration procedure is automatically finished with the measurement of buffer solution 1 and the calibration record is displayed.

1. Connect the pH electrode to the meter.
The pH measuring window is displayed.
2. When measuring without temperature sensor:
 - Temper the buffer solution, or measure the current temperature.
 - Enter the temperature value in the menu.
3. In the measured value display, select the measured parameter pH or mV with **<MODE>**.
4. Start the calibration with **<CAL>**.
The calibration display appears (voltage display).

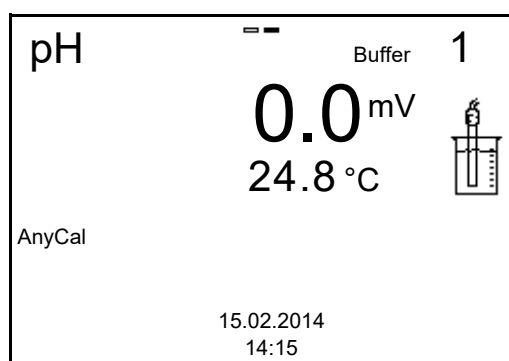


5. Thoroughly rinse the electrode with deionized water.
6. Immerse the electrode in the first buffer solution.

7. When measuring without temperature sensor:
 - Temper the buffer solution, or measure the current temperature.
 - Enter the temperature value in the menu.
8. Start the measurement with **<OK>**.
 The measured value is checked for stability (stability control). The [AR] status indicator is displayed. A progress bar is displayed and the display of the measured parameter flashes.

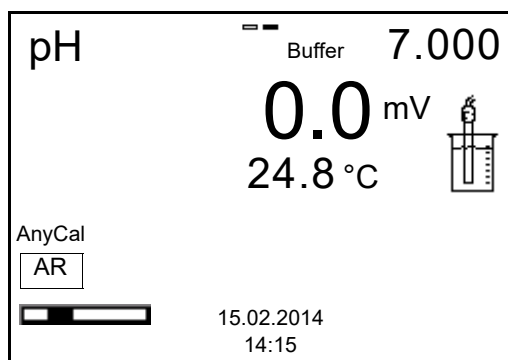

9. Wait for the end of the measurement with stability control or accept the calibration value with **<OK>**.
 The calibration display for the setting of the nominal buffer value appears.

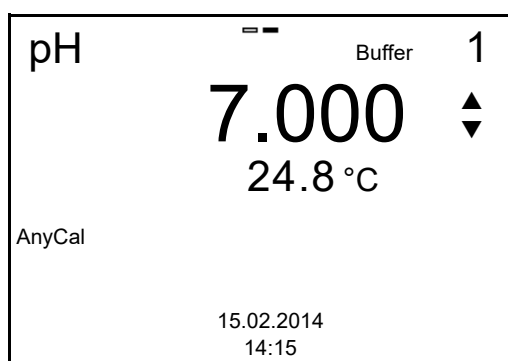


10. Set the nominal buffer value for the measured temperature with **<▲ ><▼ >**.
11. Accept the set calibration value with **<OK>**.
 The calibration display for the next buffer appears (voltage display).
12. If necessary, finish the calibration procedure as a single-point calibration with **<MODE>**.
 The calibration record is displayed.



For **single-point calibration**, the instrument uses the Nernst slope (-59.2 mV/pH at 25 °C) and determines the zero point of the electrode.

Continuing with two-point calibration


13. Thoroughly rinse the electrode with deionized water.
14. Immerse the electrode in the second buffer solution.
15. When measuring without temperature sensor:
 - Temper the buffer solution, or measure the current temperature.
 - Enter the temperature value in the menu.
16. Start the measurement with **<OK>**.
The measured value is checked for stability (stability control). The [AR] status indicator is displayed. A progress bar is displayed and the display of the measured parameter flashes.
17. Wait for the measurement with stability control to be completed or terminate the stability control and take over the calibration value with **<OK>**.
The calibration display for the setting of the nominal buffer value appears.

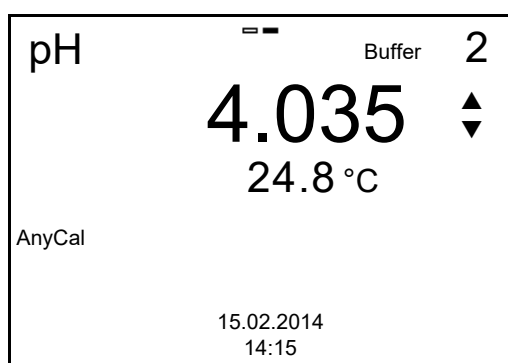

18. Set the nominal buffer value for the measured temperature with **<▲ ><▼ >**.
19. Accept the set calibration value with **<OK>**.
The calibration display for the next buffer appears (voltage display).
20. Finish the calibration procedure as a two-point calibration with **<MODE>**.
The calibration record is displayed.

Continuing with three- to five-point calibration

21. Thoroughly rinse the electrode with deionized water.
22. Immerse the electrode in the next buffer solution.
23. When measuring without temperature sensor:
 - Temper the buffer solution, or measure the current temperature.
 - Enter the temperature value in the menu.
24. Start the measurement with **<OK>**.
The measured value is checked for stability (stability control). The [AR] status indicator is displayed. A progress bar is displayed and the display of the measured parameter flashes.
25. Wait for the measurement with stability control to be completed or terminate the stability control and take over the calibration value with **<OK>**.
The calibration display for the setting of the nominal buffer value appears.

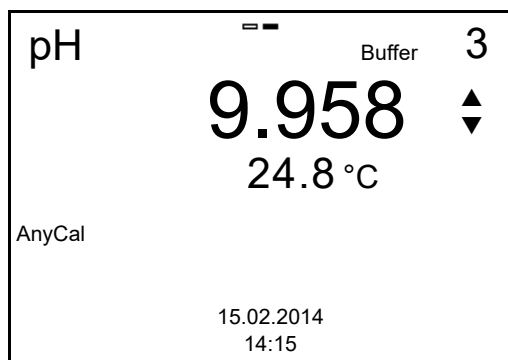

26. Set the nominal buffer value for the measured temperature with **<▲ ><▼ >**.
27. Accept the set calibration value with **<OK>**.
The calibration display for the next buffer appears (voltage display).
28. Use **<MODE>** to finish calibration or
Continue calibrating using the next buffer with **<OK>**.



After the fifth buffer has been measured the calibration is automatically finished. Then the calibration record is displayed.

The calibration line is determined by linear regression.

6.2.5 Calibration points

Calibration can be performed using one to five buffer solutions in any order (single-point to five-point calibration). The meter determines the following values and calculates the calibration line as follows:

	Determined values	Displayed calibration data
1-point	<i>Asy</i>	● Zero point = <i>Asy</i> ● Slope = Nernst slope (-59.2 mV/pH at 25 °C)
2-point	<i>Asy</i> <i>Slp.</i>	● Zero point = <i>Asy</i> ● Slope = <i>Slp.</i>
3-point to 5-point	<i>Asy</i> <i>Slp.</i>	● Zero point = <i>Asy</i> ● Slope = <i>Slp.</i> The calibration line is calculated by linear regression.



You can display the slope in the units, mV/pH or % (see section 11.1.1 SETTINGS FOR PH MEASUREMENTS, page 61).

6.2.6 Calibration data



The calibration record is automatically transmitted to the interface after calibrating.

Displays the calibration data

The calibration record of the last calibration is to be found under the menu item, **<OK>** / *Calibration* / *Calibration record*. To open it, press the **<CAL__>** key in the measured value display.

Subsequently, you can transmit the displayed calibration data to the interface, e.g. to a PC, with the **<F2>**/[USB output] key.

Displaying the calibration data memory

The calibration records of the last calibrations are to be found in the menu, **<OK>** / *Calibration* / *Calibration data storage*.

Menu item	Setting/function	Explanation
<i>Calibration</i> / <i>Calibration data storage</i> / <i>Display</i>	-	Displays the calibration record. Further options: ● Scroll through the calibration records with <▲>/<▼>. ● Output the displayed calibration record to the interface with <F2>/[USB output]. ● Quit the display with <F1>/[Back] or <OK>. ● Switch directly to the measured value display with <MODE>.
<i>Calibration</i> / <i>Calibration data storage</i> / <i>Output to USB</i>	-	Outputs the calibration records to the interface.

Calibration evaluation

After calibrating, the meter automatically evaluates the calibration. The zero point and slope are evaluated separately. The worse evaluation of both is taken into account. The evaluation appears on the display and in the calibration

record.

Display	Calibration record	Zero point [mV]	Slope [mV/pH]
	+++	-15 ... +15	-60.5 ... -58.0
	++	-20 ... <-15 or >+15 ... +20	>-58.0 ... -57.0
	+	-25 ... <-20 or >+20 ... +25	-61.0 ... <-60.5 or >-57.0 ... -56.0
	-	-30 ... <-25 or >+25 ... +30	-62.0 ... <-61.0 or >-56.0 ... -50.0
Clean the electrode according to the electrode operating manual			
Error	Error	<-30 oder >+30	<-62.0 oder >-50.0
Error elimination (see section 15 WHAT TO DO IF..., page 90)			

Calibration record (example)

```

MU 6100 L
Ser. no. 11292113

CALIBRATIONpH
15.02.2014 15:55

AutoCal TEC
Buffer 1          4.01
Buffer 2          7.00
Buffer 3          10.01
Voltage 1         184.0 mV
Voltage 2         3.0 mV
Voltage 3        -177.0 mV
Temperature 1     24.0 °C
Temperature 2     24.0 °C
Temperature 3     24.0 °C
Slope            -60.2 mV/pH
Asymmetry        4.0 mV
Sensor           +++

etc...
```

7 ORP voltage


7.1 Measuring

7.1.1 Measuring the ORP

NOTE

When connecting an grounded PC/printer, measurements cannot be performed in grounded media as the values would be incorrect. The USB interface is not galvanically isolated.

1. Connect the ORP electrode to the meter.
2. If necessary, select the U (mV) display with **<MODE>**.
3. When measuring without temperature sensor:
 - Temper the test sample, or measure the current temperature.
 - Enter the temperature value in the menu.
4. Rinse the ORP electrode and immerse it in the test sample. The measured value is checked for stability (automatic stability control). The display of the measured parameter flashes.
5. Wait for a stable measured value. The display of the measured parameter no longer flashes.

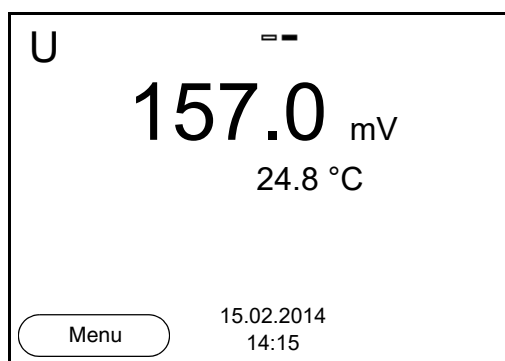

Stability control (AutoRead)

The stability control function (*AutoRead*) continually checks the stability of the measurement signal. The stability has a considerable impact on the reproducibility of measured values (see section 11.2.3 AUTOMATIC STABILITY CONTROL, page 72).

Criteria for a stable measured value

The *Stability control* function checks whether the measured values are stable within the monitored time interval.

Measured parameter	Time interval	Stability in the time interval
ORP	15 seconds	Δ : better than 0.3 mV
Temperature	15 seconds	Δ : better than 0.5 °C

The minimum duration until a measured value is assessed as stable is the monitored time interval. The actual duration is mostly longer.

Freezes the measured value (HOLD function)

With the HOLD function, you can freeze the current measured value. The displayed measured value stops changing until you switch the HOLD function off.

1. Freeze the measured value with **<HOLD>**.
The [HOLD] status indicator is displayed.
2. Release the frozen measured value again with **<HOLD>** or **<MODE>**.
The HOLD function is switched off.
The [HOLD] status display disappears.

7.1.2 Measuring the temperature

For reproducible ORP measurements, it is essential to measure the temperature of the test sample.

You have the following options to measure the temperature:

- Automatic measurement of the temperature by the temperature sensor (NTC30 or Pt1000) integrated in electrode.
- Manual determination and input of the temperature.

The measuring instrument recognizes whether a suitable sensor is connected and automatically switches on the temperature measurement.

The display of the temperature indicates the active temperature measuring mode:

Temperature sensor	Resolution of the temp. display	Temp. measurement
yes	0.1 °C	Automatic with temperature sensor
-	1 °C	Manual

If you wish to measure without temperature sensor, proceed as follows:

1. Measure the current temperature of the test sample.
2. Set the temperature value with **<▲><▼>**.
or
In the **<OK>/U/Man. temperature** menu, set the temperature value with **<▲><▼>**.

7.2 ORP calibration

ORP electrodes are not calibrated. You can, however, check ORP electrodes by measuring the ORP of a test solution and comparing the value with the nominal value.

8 Ion concentration

8.1 Measuring

8.1.1 Measuring the ion concentration

NOTE

When connecting an grounded PC/printer, measurements cannot be performed in grounded media as the values would be incorrect. The USB interface is not galvanically isolated.



Incorrect calibration of ion sensitive electrodes will result in incorrect measured values. Calibrate regularly before measuring.




For precise ISE measurements the temperature difference between measurement and calibration should not be greater than 2 K. Therefore, adjust the temperature of the standard and measuring solutions accordingly. If the temperature difference is greater the *[TempErr]* warning appears in the measured value display.

1. Connect the ISE electrode to the meter.
The pH/U measuring window is displayed.
2. If necessary, select the ISE display (unit, mg/l) with **<MODE>**.
3. When measuring without temperature sensor:
 - Temper the test sample, or measure the current temperature.
4. Calibrate or check the meter with the electrode.



While no valid calibration is available, e.g. in the delivery condition, "Error" appears in the measured value display.

5. Immerse the electrode in the test sample.

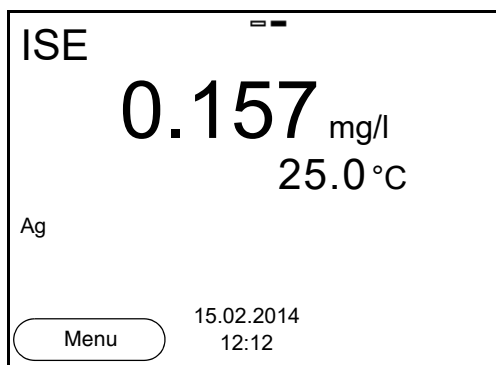

Stability control (AutoRead)

The stability control function (*AutoRead*) continually checks the stability of the measurement signal. The stability has a considerable impact on the reproducibility of measured values (see section 11.2.3 AUTOMATIC STABILITY CONTROL,

page 72).

Criteria The AutoRead criteria affect the reproducibility of the measured values. The following criteria can be adjusted:

- *high*: highest reproducibility
- *medium*: medium reproducibility
- *low*: lowest reproducibility



Increasing reproducibility also causes the response time to increase until a measured value is evaluated as stable.

Freezes the measured value (HOLD function)

With the HOLD function, you can freeze the current measured value. The displayed measured value stops changing until you switch the HOLD function off.

1. Freeze the measured value with **<HOLD>**.
The [HOLD] status indicator is displayed.
2. Release the frozen measured value again with **<HOLD>**.
The HOLD function is switched off.
The [HOLD] status display disappears.

8.1.2 Measuring the temperature

For reproducible ion-selective measurements, it is essential to measure the temperature of the test sample.

You have the following options to measure the temperature:

- Measurement by an external temperature sensor.
- Manual determination and input of the temperature.

The measuring instrument recognizes whether a suitable sensor is connected and automatically switches on the temperature measurement.

The display of the temperature indicates the active temperature measuring mode:

Temperature sensor	Resolution of the temp. display	Temp. measurement
yes	0.1 °C	Automatic with temperature sensor
-	1 °C	Manual

If you wish to measure (or calibrate) without temperature sensor, proceed as follows:

1. Measure the current temperature of the test sample.

2. When measuring without temperature sensor:
 - Temper the test sample, or measure the current temperature.

8.2 Calibration

8.2.1 Why calibrate?

Ion-selective electrodes age and are temperature-dependent. This changes the slope. As a result, an inexact measured value is displayed. Calibration determines the calibration line of the electrode and stores this value in the meter.

Thus, you should calibrate before each measurement (if possible), and at regular intervals.


8.2.2 When to calibrate?

- Before any measurement if possible
- After connecting another ISE electrode
- When the sensor symbol flashes, e.g. after a voltage interruption (empty batteries)


8.2.3 Calibration (ISE Cal)

ISE Cal is the conventional **two-point to seven-point calibration procedure** that uses 2 to 7 freely selectable standard solutions. The concentration expected in the measurement determines the concentration of the calibration standards.


1. Connect the ISE electrode to the meter.
The pH/U/ISE measuring window is displayed.
2. Keep the standard solutions ready.
3. When measuring without temperature sensor:
 - Temper the test sample, or measure the current temperature.
4. In the measured value display, select the ISE measuring window with **<▲ >** **<▼ >** and **<MODE>**.
5. If necessary, change the unit of the measurement result and calibration standards in the *ISE setup/Unit* menu.
6. Start the calibration with **<CAL>**.
The calibration display appears.

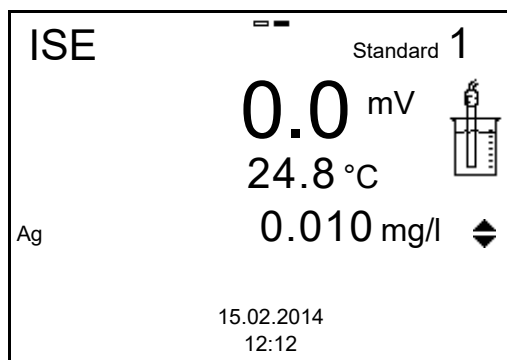

7. Thoroughly rinse the electrode with distilled water.
8. Immerse the electrode in standard solution 1.
9. When calibrating without temperature sensor:
 - Measure the temperature of the standard solution using a thermometer.
 - Use **<F2>/[↑]** to select the setting of the temperature.
 - Use **<▲ > <▼ >** to set the temperature.
 - Use **<F2>/[↑]** to select the setting of the concentration.
10. Set the concentration of the standard solution with **<▲ > <▼ >** and press **<OK>**.
 The standard solution is measured.
 The measured value is checked for stability (AutoRead).

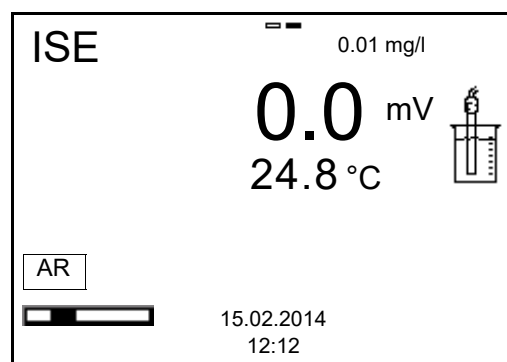

11. Wait for the end of the AutoRead measurement or accept the calibration value with **<OK>**.
 The calibration display for the next standard solution appears.

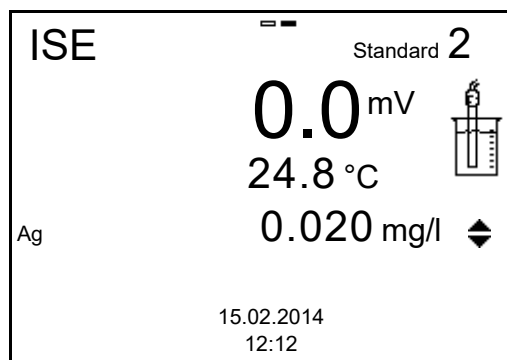


Continuing with two-point calibration

12. Thoroughly rinse the electrode with distilled water.
13. Immerse the electrode in standard solution 2.
14. When calibrating without temperature sensor:
 - Measure the temperature of the standard solution using a thermometer.
 - Use <F2>/[$\updownarrow$] to select the setting of the temperature.
 - Use <▲ > <▼ > to set the temperature.
 - Use <F2>/[$\updownarrow$] to select the setting of the concentration.
15. Set the concentration of the standard solution with <▲ > <▼ > and press <OK>.


The standard solution is measured.

The measured value is checked for stability (AutoRead).

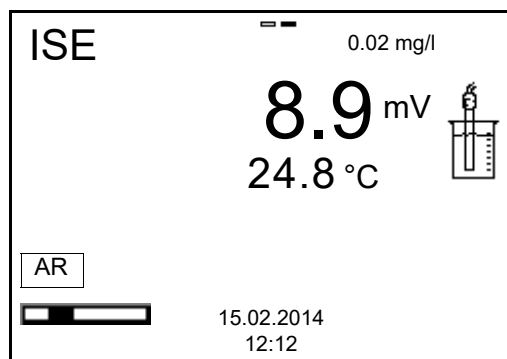

16. Wait for the end of the AutoRead measurement or accept the calibration value with <OK>.

The calibration display for the next standard solution appears.

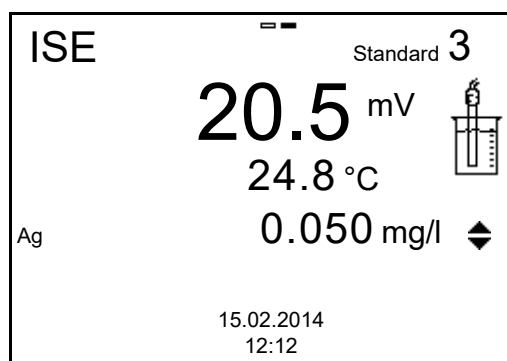

17. Press **<OK>** to continue with three-point calibration.
or
Finish the calibration procedure as a two-point calibration with **<MODE>**.
The new calibration values are displayed.

Continuing with three- to seven-point calibration

Repeat the steps 12 to 17 in the same way with the third and further standard solutions as necessary. The new calibration values are displayed after the last calibration step was completed.



Based on the calibration data, the calibration curve is determined in sections, according to the Nernst equation modified by Nikolski.

8.2.4 Calibration standards

Use two to seven different standard solutions. The standard solutions have to be selected in either increasing or decreasing order.



Select the unit of the standard solution and measurement result in the *ISE setup/Unit* menu.

Standard solution (Std 1 - 7)	Values
Unit [mg/l]	0.010 ... 500.000
Unit [mol/l]	0.100 ... 5.000 µmol/l 10.00 ... 5.000 mmol/l
Unit [mg/kg]	0.010 ... 500.000
Unit [ppm]	0.010 ... 500.000
Unit [%]	0.001 ... 50.000



The measurement precision is also dependent on the selected standard solutions. Therefore, the selected standard solutions should cover the value range expected of the following concentration measurement.

If the measured electrode potential is outside the calibrated range, the *[ISEErr]* warning is displayed.

8.2.5 Calibration data

Displays the calibration data

The calibration record of the last calibration is to be found under the menu item, **<OK>** / *Calibration* / *Calibration record*. To open it, press the **<CAL__>** key in the measured value display.

Subsequently, you can transmit the displayed calibration data to the interface, e.g. to a PC, with the **<F2>/[USB output]** key.

Displaying the calibration data memory

The calibration records of the last calibrations are available in the menu, **<OK> / Calibration / Calibration data storage**.

Menu item	Setting/function	Explanation
<i>Calibration / Calibration data storage / Display</i>	-	Displays the calibration record. Further options: ● Scroll through the calibration records with <▲><▼>. ● Output the displayed calibration record to the interface with <F2>/[USB output]. ● Quit the display with <F1>/[Back] or <OK>. ● Switch directly to the measured value display with <MODE>.
<i>Calibration / Calibration data storage / Output to USB</i>	-	Outputs the calibration records to the interface.

Calibration evaluation

After calibrating, the meter automatically evaluates the calibration.

Display	Calibration record	Magnitude of the slope [mV]
	+++	50.0 ... 70.0 or 25.0 ... 35.0
<i>Error</i> Error elimination (see section 15 WHAT TO DO IF..., page 90)	<i>Error</i>	< 50 or > 70 or < 25 or > 35

Calibration record (example)

```

MU 6100 L
Ser. no. 12345678

CALIBRATIONISE
18.01.2013 08:09:10

Standard 1      0.010 mg/l
Standard 2      0.020 mg/l
Voltage 1       38.5 mV
Voltage 2       58.0 mV
Temperature 1   24.0 °C
Temperature 2   24.0 °C
Ion type        Ag
Slope           64.7 mV
Sensor          +++
  
```


9 Dissolved oxygen


9.1 Measuring

9.1.1 Measuring D.O.

NOTE

When connecting an grounded PC/printer, measurements cannot be performed in grounded media as the values would be incorrect. The USB interface is not galvanically isolated.

1. Connect the D.O. sensor to the meter.
The D.O. measuring screen is displayed.
2. If necessary, select the measured parameter with **<MODE>**.
3. Immerse the D.O. sensor in the test sample.

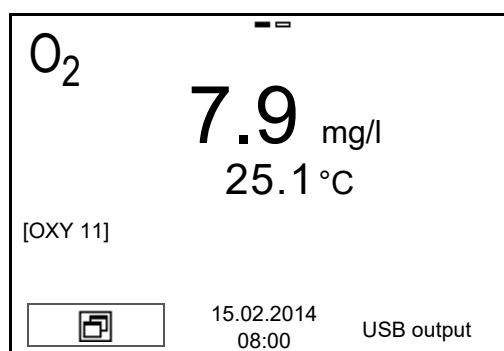

4. If necessary, select the measured parameter with **<MODE>**.
 - D.O. concentration [mg/l]
 - D.O. saturation [%]
 - D.O. partial pressure [mbar].

Salinity correction

When measuring the concentration of solutions with a salt content of more than 1 g/l, a salinity correction is required. For this, you have to measure and input the salinity of the measured medium first. When the salinity correction is switched on, the **[SAL]** indicator is displayed in the measuring window.



You can switch the salinity correction on or off and enter the salinity in the menu for calibration and measurement settings (see section 11.1.5 SETTINGS FOR D.O. SENSORS, page 65).

Stability control (AutoRead)

The stability control function (*AutoRead*) continually checks the stability of the measurement signal. The stability has a considerable impact on the reproducibility of measured values (see section 11.2.3 AUTOMATIC STABILITY CONTROL,

page 72).

Criteria for a stable measured value

The *Stability control* function checks whether the measured values are stable within the monitored time interval.

Measured parameter	Time interval	Stability in the time interval
D.O. concentration	20 seconds	Δ : better than 0.05 mg/l
D.O. saturation	20 seconds	Δ : better than 0.6 %
D.O. partial pressure	20 seconds	Δ : Better than 1.2 mbar
Temperature	15 seconds	Δ : better than 0.5 °C

The minimum duration until a measured value is assessed as stable is the monitored time interval. The actual duration is mostly longer.

Freezes the measured value (HOLD function)

With the HOLD function, you can freeze the current measured value. The displayed measured value stops changing until you switch the HOLD function off.

1. Freeze the measured value with **<HOLD>**.
The [HOLD] status indicator is displayed.
2. Release the frozen measured value again with **<HOLD>**.
The HOLD function is switched off.
The [HOLD] status display disappears.

9.1.2 Measuring the temperature

For reproducible D.O. measurements, it is essential to measure the temperature of the test sample.

The temperature is automatically measured by the temperature sensor (NTC30 or Pt1000) integrated in the sensor.



9.2 Calibration

9.2.1 Why calibrate?

D.O. sensors age. This changes the slope of the D.O. sensor. Calibration determines the current slope of the sensor and stores this value in the instrument.

9.2.2 When to calibrate?

- After connecting a sensor
- Routinely within the framework of the company quality assurance
- When the calibration interval has expired

9.2.3 Calibration procedure

The MU 6100 L provides 2 calibration procedures:

- Calibration in water vapor-saturated air.
Use an OxiCal[®] air calibration vessel for the calibration.
- Calibration via a comparison measurement (e.g. Winkler titration according to DIN EN 25813 or ISO 5813). At the same time, the relative slope is adapted to the comparison measurement by a correction factor. When the correction multiplier is active, the *[Factor]* indicator appears in the measuring window.

9.2.4 Calibration in water vapor-saturated air (air calibration vessel)

For this calibration procedure, the *Comparison meas.* setting must be set to *Off* in the *Calibration* menu.


Proceed as follows to calibrate the instrument:

1. Connect the D.O. sensor to the meter.
2. Put the D.O. sensor into the air calibration vessel.



The sponge in the air calibration vessel must be moist (not wet). Leave the sensor in the air calibration vessel for a time long enough to adjust.

3. Start the calibration with **<CAL>**.
The last calibration data (relative slope) is displayed.

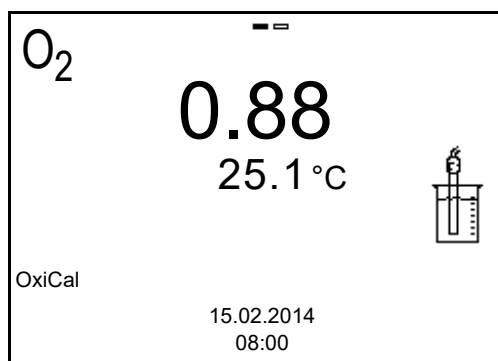

4. Start the measurement with **<OK>**.
The measured value is checked for stability (stability control).
The [AR] status indicator is displayed. A progress bar is displayed and the display of the measured parameter flashes.
5. Wait for the end of the measurement with stability control or accept the calibration value with **<OK>**.
The calibration record is displayed and output to the interface.
6. Using **<F1>**/[Continue] or **<OK>**, switch to the measured value display.
The meter displays a D.O. saturation of approx. 101.7 % in the air calibration vessel.


9.2.5 Calibrating with a comparison measurement (OxiComp)

For this calibration procedure, the *Comparison meas.* setting must be set to *On* in the *Calibration* menu.




Before calibrating via a comparison measurement, the sensor should be calibrated in the air calibration vessel.

1. Connect the D.O. sensor to the meter.
2. Immerse the D.O. sensor in the reference solution.
3. Start the calibration with **<CAL>**.

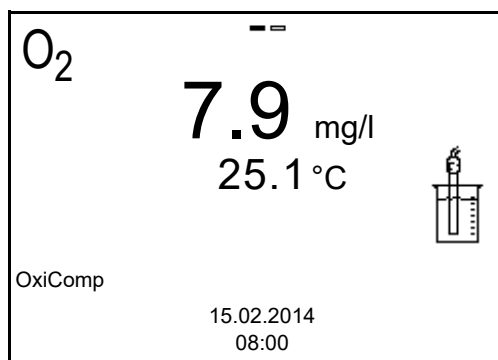

4. Start the measurement with **<OK>**.
The measured value is checked for stability (stability control).
The [AR] status indicator is displayed. The measured parameter flashes.
5. Wait for the end of the measurement with stability control or accept the calibration value with **<OK>**.
The factor that was set last is displayed.

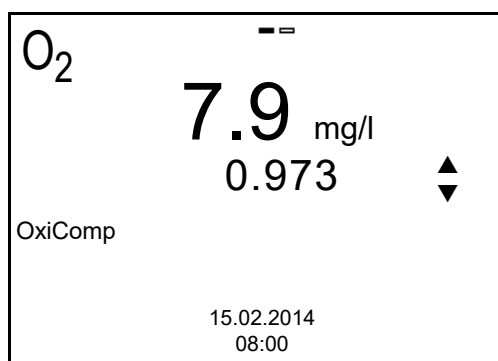

6. Using **<▲ > <▼ >**, set the correction factor to adjust the displayed concentration value to the nominal value (value of the comparison measurement). Subsequently, accept the correction factor with **<OK>**.
The meter switches to the measured value display.
The status display *[Factor]* is active.

9.2.6 Calibration data



The calibration record is automatically transmitted to the interface after calibrating.

Displays the calibration data

The calibration record of the last calibration is to be found under the menu item, **<OK>** / *Calibration* / / *Calibration record*. To open it, press the **<CAL__>** key in the measured value display.

Subsequently, you can transmit the displayed calibration data to the interface, e.g. to a PC, with the **<F2>**[*USB output*] key.

Displaying the calibration data memory

The calibration records of the last calibration procedures are available in the menus, **<OK>**/Calibration / / Calibration data storage and **<OK>**/Storage & config / Data storage/ / Calibration data storage.

Menu item	Setting/function	Explanation
<i>Calibration / Calibration data storage / Display</i> or <i>Data storage / Calibration data storage / Display</i>	-	Displays the calibration record. Further options: ● Scroll through the calibration records with <▲><▼>. ● Output the displayed calibration record to the interface with <F2>[<i>USB output</i>]. ● Quit the display with <F1>/ [<i>Back</i>] or <OK>. ● Switch directly to the measured value display with <MODE>.
<i>Calibration / Calibration data storage / Output to USB</i> or <i>Data storage / Calibration data storage / Output to USB</i>	-	Outputs the calibration records to the interface.

Calibration evaluation

After calibration, the meter automatically evaluates the current status of the calibration. The evaluation appears on the display and in the calibration record.

Display	Calibration record	Relative slope
	+++	$S = 0.8 \dots 1.25$
	++	$S = 0.7 \dots 0.8$
	+	$S = 0.6 \dots 0.7$
Error Error elimination (see section 15 WHAT TO DO IF..., page 90)	Error	$S < 0.6$ or $S > 1.25$

Calibration record (example)

```

MU 6100 L
Ser. no. 11292113

CALIBRATION OXY 11
15.02.2014 16:13:33

S = 0.88 25.0 °C
Sensor +++

```

10 Conductivity

10.1 Measuring

10.1.1 Measuring the conductivity

NOTE


When connecting an grounded PC/printer, measurements cannot be performed in grounded media as the values would be incorrect. The USB interface is not galvanically isolated.

1. Connect the conductivity measuring cell to the measuring instrument. The conductivity measuring screen is displayed.
2. If necessary, select the measured parameter with **<MODE>**.
3. Check the *Measuring cell* settings and cell constant for the connected conductivity measuring cell. If necessary, correct the settings.



The selection of the measuring cell and the setting of the cell constant is done in the measurement settings menu for conductivity (see section 11.1.7 SETTINGS FOR CONDUCTIVITY MEASURING CELLS, page 67). The cell constant to be set must either be taken from the operating manual of the measuring cell or is printed on the measuring cell.

4. Immerse the conductivity measuring cell in the test sample.

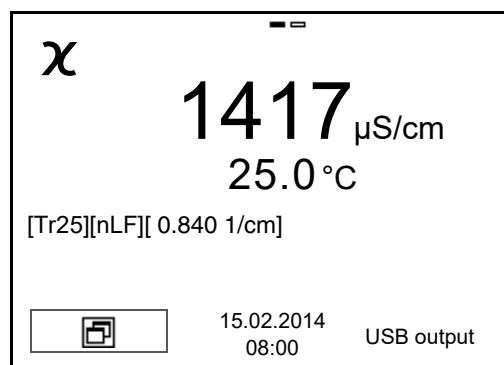

5. If necessary, select the measured parameter with **<MODE>**.
 - Conductivity [$\mu\text{S}/\text{cm}$] / [mS/cm]
 - Resistivity [$\Omega \cdot \text{cm}$] / [$\text{k}\Omega \cdot \text{cm}$] / [$\text{M}\Omega \cdot \text{cm}$]
 - Salinity SaL []
 - Total dissolved solids TDS [mg/l] / [g/l]



The factor to calculate the total dissolved solids is set to 1.00 in the factory. You can adjust this factor to meet your requirements in the range 0.40 ... 1.00. The factor is set in the *Measurement* menu for the parameter, TDS.

Stability control (AutoRead)

The stability control function (*AutoRead*) continually checks the stability of the measurement signal. The stability has a considerable impact on the reproducibility of measured values (see section 11.2.3 AUTOMATIC STABILITY CONTROL, page 72).

Criteria for a stable measured value

The *Stability control* function checks whether the measured values are stable within the monitored time interval.

Measured parameter	Time interval	Stability in the time interval
Conductivity χ	10 seconds	Δ : better than 1.0% of measured value
Temperature	15 seconds	Δ : better than 0.5 °C

The minimum duration until a measured value is assessed as stable is the monitored time interval. The actual duration is mostly longer.

Freezes the measured value (HOLD function)

With the HOLD function, you can freeze the current measured value. The displayed measured value stops changing until you switch the HOLD function off.

1. Freeze the measured value with **<HOLD>**.
The [HOLD] status indicator is displayed.
2. Release the frozen measured value again with **<HOLD>**.
The HOLD function is switched off.
The [HOLD] status display disappears.

10.1.2 Measuring the temperature

For reproducible conductivity measurements, it is essential to measure the temperature of the test sample.

The temperature is automatically measured by the temperature sensor (NTC30 or Pt1000) integrated in the sensor.



The conductivity sensor CO 11 has an integrated temperature sensor.

10.2 Calibration

10.2.1 Why calibrate?

Aging slightly changes the cell constant, e.g. due to coatings. As a result, an inexact measured value is displayed. The original characteristics of the cell can often be restored by cleaning the cell. Calibration determines the current value of the cell constant and stores this value in the meter.

Thus, you should calibrate at regular intervals (we recommend: every 6 months).

10.2.2 When to calibrate?

- After connecting a sensor
- Routinely within the framework of the company quality assurance
- When the cleaning interval has expired

10.2.3 Determining the cell constant (calibration in control standard)

You can determine the actual cell constant of the conductivity measuring cell by calibrating with the control standard in the following ranges:


- $0.800 \dots 0.880 \text{ cm}^{-1}$
(e.g. CO 11 with a nominal cell constant of 0.840 cm^{-1})
- $0.450 \dots 0.500 \text{ cm}^{-1}$
(e.g. conductivity measuring cells with a nominal cell constant of around 0.475 cm^{-1})

The cell constant is determined in the control standard, 0.01 mol/l KCl . Cell constants outside the ranges quoted above cannot be calibrated.

In the default condition, the calibrated cell constant of the meter is set to 0.840 cm^{-1} (conductivity measuring cell CO 11).

For this calibration procedure, the *Measuring cell* setting must be set to cal in the *Measurement* menu. Proceed as follows to determine the cell constant:

1. Connect the conductivity measuring cell to the measuring instrument.
2. In the measured value display, select the conductivity parameter with **<MODE>**.
3. Start the calibration with **<CAL>**.
The cell constant that was calibrated last is displayed.

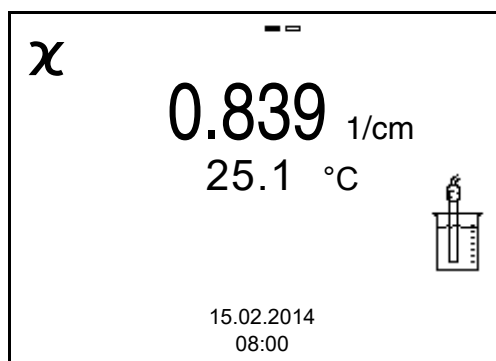

4. Immerse the conductivity measuring cell in the control standard solution, 0.01 mol/l KCl.
5. Start the measurement with **<OK>**.
The [AR] status indicator is displayed. A progress bar is displayed and the display of the measured parameter flashes.
6. Wait for the end of the measurement with stability control or accept the calibration value with **<OK>**.
The calibration record is displayed and output to the interface.
7. Using **<F1>**/[Continue] or **<OK>**, switch to the measured value display.

10.2.4 Calibration data



The calibration record is automatically transmitted to the interface after calibrating.

Displays the calibration data

The calibration record of the last calibration is to be found under the menu item, **<OK>** / Calibration / Calibration record. To open it, press the **<CAL__>** key in the measured value display.

Subsequently, you can transmit the displayed calibration data to the interface, e.g. to a PC, with the **<F2>**[USB output] key.

Displaying the calibration data memory

The calibration records of the last calibration procedures are available in the **<OK>**/Calibration / Calibration data storage and **<OK>**/Storage & config / Data storage / Calibration data storage menu.

Menu item	Setting/function	Explanation
<i>Calibration / Calibration data storage / Display</i> or <i>Data storage / Calibration data storage / Display</i>	-	Displays the calibration record. Further options: ● Scroll through the calibration records with <▲>/<▼>. ● Output the displayed calibration record to the interface with <F2>/[USB output]. ● Quit the display with <F1>/[Back] or <OK>. ● Switch directly to the measured value display with <MODE>.
<i>Calibration / Calibration data storage / Output to USB</i> or <i>Data storage / Calibration data storage / Output to USB</i>	-	Outputs the calibration records to the interface.

Calibration evaluation

After calibration, the meter automatically evaluates the current status of the calibration. The evaluation appears on the display and in the calibration record.

Display	Calibration record	Cell constant [cm^{-1}]
	+++	within the ranges $0.450 \dots 0.500 \text{ cm}^{-1}$ or $0.800 \dots 0.880 \text{ cm}^{-1}$
Error Error elimination (see section 15 WHAT TO DO IF..., page 90)	Error	outside the ranges $0.450 \dots 0.500 \text{ cm}^{-1}$ or $0.800 \dots 0.880 \text{ cm}^{-1}$

Calibration record (example)

```

MU 6100 L
Ser. no. 11292113

CALIBRATION Cond
15.02.2014 16:13:33

Cell constant      0.841 1/cm      25.0 °C
Sensor            +++

```

11 Settings

11.1 Measurement settings

11.1.1 Settings for pH measurements

The settings for pH measurements are made in the menu for calibration and measurement settings of the pH/ORP measurement. To open the settings, display the required measured parameter in the measured value display and press the **<OK>** key. After completing the settings, switch to the measured value display with **<MODE>**.

In the following table, only those settings are listed that concern the pH measurement.

Default settings are printed in **bold**.

Menu item	Possible setting	Explanation
<i>Calibration / Calibration record</i>	-	Displays the calibration record of the last calibration.
<i>Calibration / Calibration data storage / Display</i>	-	Displays the calibration record. Further options: ● Scroll through the calibration records with <▲><▼>. ● Output the displayed calibration record to the interface with <F2>/[USB output]. ● Output all calibration records to the interface with <F2__>/[USB output]. ● Quit the display with <F1>/[Back] or <OK>. ● Switch directly to the measured value display with <MODE>.
<i>Calibration / Calibration data storage / Output to USB</i>	-	Outputs the calibration records to the interface.
<i>Calibration / Buffer</i>	TEC AnyCal NIST/DIN TEC 2 ...	Buffer sets to be used for pH calibration (see section 6.2 CALIBRATION, page 30).
<i>Calibration / One point calibration</i>	Yes No	Quick calibration with 1 buffer

Menu item	Possible setting	Explanation
<i>Calibration / Serial number (sensor)</i>	-	Entry of the serial number of the connected sensor. The serial number is output in the calibration record. ● Change the contents of the highlighted position with <▲ ><▼ >. ● Go to the next position with <F2>/[▶]. ● When the serial number has been completely entered, confirm with <OK>.
<i>Calibration / Calibration interval</i>	1 ... 7 ... 999 d	Calibration interval for the pH electrode (in days). The meter reminds you to calibrate regularly by the flashing sensor symbol in the measuring screen.
<i>Calibration / Unit for slope</i>	mV/pH %	Unit of the slope. The % display refers to the Nernst slope of -59.2 mV/pH (100 x determined slope/Nernst slope).
<i>Alternative temperature</i>	On Off	Takes the temperature value from the 2nd sensor.
<i>Man. temperature</i>	-25 ... +25 ... +130 °C	Entry of the manually determined temperature. For measurements without temperature sensor only.
<i>Resolution pH</i>	0.001 0.01 0.1	Resolution of the pH display
<i>Reset</i>	-	Resets all sensor settings to the delivery condition (see section 11.3.1 RESETTING THE MEASUREMENT SETTINGS, page 72).

11.1.2 Buffer sets for calibration

You can use the buffer sets quoted in the table for an automatic calibration. The pH values are valid for the specified temperature values. The temperature dependence of the pH values is taken into consideration during calibration.

No.	Buffer set *	pH values	at
1	TEC EU Technical buffers EU	2.00 4.00 7.00 10.00	20 °C
2	TEC US Technical buffers US		
3	AnyCal	Any	Any
4	NIST/DIN DIN buffers according to DIN 19266 and NIST Traceable Buffers	1.679 4.006 6.865 9.180 12.454	25 °C
5	TEC 2 Technical buffers	4.010 7.000 10.011	25 °C
6	Merck 1 *	4.000 7.000 9.000	20°C
7	Merck 2 *	1.000 6.000 8.000 13.000	20°C
8	Merck 3 *	4.660 6.880 9.220	20°C
9	Merck 4 *	2.000 4.000 7.000 10.000	20°C
10	Merck 5 *	4.010 7.000 10.000	25 °C

* Brand names or trade names are trademarks of their respective owners protected by law.



The buffers are selected in the menu, pH / **<OK>** / *Calibration / Buffer* (see 11.1.1 SETTINGS FOR PH MEASUREMENTS, PAGE 61).

11.1.3 Calibration interval

The calibration evaluation is displayed as a sensor symbol.

The sensor symbol flashes after the adjusted calibration interval has expired. It is still possible to measure.



To ensure the high measuring accuracy of the measuring system, calibrate after the calibration interval has expired.

Setting the calibration interval

The calibration interval is set to 7 days in the factory. You can change the interval (1 ... 999 days):

1. Open the menu for measurement settings with **<OK>**.
2. In the *Calibration / Calibration interval* menu, set the calibration interval with **<▲><▼>**.
3. Confirm the setting with **<OK>**.
4. Quit the menu with **<MODE>**.

11.1.4 Settings for ORP measurements

The settings for ORP measurements are made in the menu for calibration and measurement settings of the pH/ORP measurement. To open the settings, display the required measured parameter in the measured value display and press the **<OK>** key. After completing the settings, switch to the measured value display with **<MODE>**.

In the following table, only those settings are listed that influence the ORP measurement.

Default settings are printed in **bold**.

Menu item	Possible setting	Explanation
<i>Man. temperature</i>	-25 ... +25 ... +130 °C	Entry of the manually determined temperature. For measurements without temperature sensor only.
<i>Resolution mV</i>	0.1 <i>1</i>	Resolution of the mV display
<i>Reset</i>	-	Resets all sensor settings to the delivery condition (see section 11.3.1 RESETTING THE MEASUREMENT SETTINGS, page 72).

11.1.5 Settings for D.O. sensors

The settings are available in the menu for measurement and calibration settings. To open the settings, display the required measured parameter in the measured value display and press the **<OK>** key. After completing the settings, switch to the measured value display with **<MODE>**.

Default settings are printed in **bold**.

Menu item	Possible setting	Explanation
<i>Calibration / Calibration record</i>	-	Displays the calibration record of the last calibration.
<i>Calibration / Calibration data storage / Display</i>	-	Displays the calibration record. Further options: ● Scroll through the calibration records with <▲><▼>. ● Output the displayed calibration record to the interface with <F2>/[USB output]. ● Output all calibration records to the interface with <F2__>/[USB output]. ● Quit the display with <F1>/[Back] or <OK>. ● Switch directly to the measured value display with <MODE>.
<i>Calibration / Calibration data storage / Output to USB</i>	-	Outputs the calibration records to the interface.
<i>Calibration / Serial number (sensor)</i>	-	Entry of the serial number of the connected sensor. The serial number is output in the calibration record. ● Change the contents of the highlighted position with <▲ ><▼ >. ● Go to the next position with <F2>/[►]. ● When the serial number has been completely entered, confirm with <OK>.
<i>Calibration / Calibration interval</i>	1 ... 14 ... 999 d	<i>Calibration interval</i> for the D.O. sensor (in days). The meter reminds you to calibrate regularly by the flashing sensor symbol in the measuring screen.

Menu item	Possible setting	Explanation
<i>Calibration / Comparison meas.</i>	<i>On</i> <i>Off</i>	Enables to adjust the measured value with the aid of a comparison measurement, e.g. Winkler titration. For details, see section 9.2 CALIBRATION, page 51.
<i>Sal correction</i>	<i>On</i> <i>Off</i>	Manual salt content correction for concentration measurements.
<i>Salinity</i>	0.0 ... 70.0	Salinity or salinity equivalent for the salt content correction. This function is only available for concentration measurements if the manual salt content correction is switched on.
<i>Reset</i>	-	Resets all sensor settings to the delivery condition (see section 11.3.1 RESETTING THE MEASUREMENT SETTINGS, page 72).

11.1.6 Calibration interval

The calibration evaluation is displayed as a sensor symbol.

The sensor symbol flashes after the adjusted calibration interval has expired. It is still possible to measure.



To ensure the high measuring accuracy of the measuring system, calibrate after the calibration interval has expired.

Setting the calibration interval

The calibration interval is set to 14 days in the factory. You can change the interval (1 ... 999 days):

1. Open the menu for measurement settings with **<OK>**.
2. In the *Calibration / Calibration interval* menu, set the calibration interval with **<▲><▼>**.
3. Confirm the setting with **<OK>**.
4. Quit the menu with **<MODE>**.

11.1.7 Settings for conductivity measuring cells

The settings are made in the *Measurement* menu for the measured parameter, conductivity. To open the settings, display the required measured parameter in the measured value display and press the **<OK>** key. After completing the settings, switch to the measured value display with **<MODE>**.

Default settings are printed in **bold**.

Menu item	Possible setting	Explanation
<i>Calibration / Calibration record</i>	-	Displays the calibration record of the last calibration.
<i>Calibration / Calibration data storage / Display</i>	-	Displays the calibration record. Further options: ● Scroll through the calibration records with <▲><▼>. ● Output the displayed calibration record to the interface with <F2>/ [USB output]. ● Output all calibration records to the interface with <F2__>/ [USB output]. ● Quit the display with <F1>/ [Back] or <OK>. ● Switch directly to the measured value display with <MODE>.
<i>Calibration / Calibration data storage / Output to USB</i>	-	Outputs the calibration records to the interface.
<i>Calibration / Serial number (sensor)</i>	-	Entry of the serial number of the connected sensor. The serial number is output in the calibration record. ● Change the contents of the highlighted position with <▲ ><▼ >. ● Go to the next position with <F2>/ [►]. ● When the serial number has been completely entered, confirm with <OK>.
<i>Calibration / Calibration interval</i>	1 ... 150 ... 999 d	<i>Calibration interval</i> for the measuring cell (in days). The meter reminds you to calibrate regularly by the flashing sensor symbol in the measuring screen.

Menu item	Possible setting	Explanation
Measurement / Measuring cell / Type	Cal <i>C 0.100</i> <i>C 0.010</i> <i>man</i>	<i>Measuring cell used</i> Measuring cells whose cell constant is determined by calibration in the KCL control standard solution. Calibration ranges: 0.450 ... 0.500 cm⁻¹ and 0.800 ... 0.880 cm⁻¹. The currently valid cell constant is displayed in the status line. Measuring cell C 0.100, nominal cell constant 0.100 cm⁻¹. The cell constant can be adjusted in the range 0.090 ... 0.110 cm⁻¹. Measuring cell C 0.010, nominal cell constant 0.010 cm⁻¹. The cell constant is permanently set. Any measuring cells with freely adjustable cell constants.
Measurement / Measuring cell / Cell const. man	0,250 ...	Display and setting option of the cell constant of any measuring cells (man).
Measurement / Measuring cell / Cell const. LR325/ 01	0.090 ... 0.100 ... 0.110 cm ⁻¹	Display and setting options for the cell constant of the measuring cell C 0.100.
Measurement / Temp. comp. (TC) / Method	nLF <i>Lin</i> <i>Off</i>	Procedure for temperature compensation (see section 11.1.9 TEMPERATURE COMPENSATION, page 69). This setting is only available for the measured parameters, χ and ρ .
Measurement / Temp. comp. (TC) / Linear coeff.	0.000 ... 2.000 ... 10.000 %/K	Coefficient of the linear temperature compensation. This setting is only available when the linear temperature compensation is set.

Menu item	Possible setting	Explanation
<i>Measurement / Temp. comp. (TC) /Reference temp.</i>	20 °C 25 °C	Reference temperature This setting is only available for the measured parameters, χ and ρ .
<i>Measurement / TDS factor</i>	0.40 ... 1.00	Multiplier for TDS value
<i>Reset</i>	-	Resets all sensor settings to the delivery condition (see section 11.3.1 RESETTING THE MEASUREMENT SETTINGS, page 72).

11.1.8 Calibration interval

The calibration evaluation is displayed as a sensor symbol.

The sensor symbol flashes after the adjusted calibration interval has expired. It is still possible to measure.



To ensure the high measuring accuracy of the measuring system, calibrate after the calibration interval has expired.

Setting the calibration interval

The calibration interval is set to 150 days in the factory. You can change the interval (1 ... 999 days):

1. Open the menu for measurement settings with **<OK>**.
2. In the *Calibration / Calibration interval* menu, set the calibration interval with **<▲><▼>**.
3. Confirm the setting with **<OK>**.
4. Quit the menu with **<MODE>**.

11.1.9 Temperature compensation

The calculation of the temperature compensation is based on the preset reference temperature, 20 °C or 25 °C. It appears on the display as Tr20 or Tr25.

You can select one of the following temperature compensation methods:

- **Nonlinear temperature compensation (nLF)** according to EN 27 888
- **Linear temperature compensation (lin)** with adjustable coefficient in the range 0.000 ... 10.000 %/K
- **No temperature compensation (off)**



The reference temperature and temperature compensation are set in the *Measurement* menu for the parameter, conductivity (see section 11.1.7 SETTINGS FOR CONDUCTIVITY MEASURING CELLS, page 67).

Application tips Set the temperature compensation suitable for your test sample:

Test sample	Temperature compensation	Display
Natural water (ground water, surface water, drinking water)	<i>nLF</i> according to EN 27 888	<i>nLF</i>
Ultrapure water	<i>nLF</i> according to EN 27 888	<i>nLF</i>
Other aqueous solutions	<i>Lin</i> Set linear temperature coefficient 0.000 ... 10.000 %/K	<i>Lin</i>
Salinity (seawater)	Automatic <i>nLF</i> according to IOT (International Oceanographic Tables)	<i>Sal, nLF</i>

11.1.10 Setting the TDS factor

The factor to calculate the total dissolved solids is set to 1.00 in the delivery condition.

You can adjust this factor to meet your requirements in the range 0.40 ... 1.00.

The factor is set in the menu for the parameter TDS (*Measurement / TDS factor*).

11.2 Sensor-independent settings

11.2.1 System

To open the *Storage & config* menu, press the **<OK>** key in the measured value display. After completing the settings, switch to the measured value display with **<MODE>**.

Default settings are printed in **bold**.

Menu item	Possible setting	Explanation
<i>System / General / Language</i>	<i>Deutsch</i> English <i>(more)</i>	Selects the menu language
<i>System / General / Beep</i>	On Off	Switches on/off the beep on keystroke
<i>System / General / Illumination</i>	Auto On Off	Switches the display illumination on/off
<i>System / General / Contrast</i>	0 ... 50 ... 100 %	Changes the display contrast
<i>System / General / Switchoff time</i>	10 min ... 1h ... 24 h	Adjusts the shut-off time
<i>System / General / Temperature unit</i>	°C °F	Temperature unit, degrees Celsius or degrees Fahrenheit. All temperature values are displayed with the selected unit.
<i>System / General / Stability control</i>	On Off	Switches on or off the automatic stability control during measurement
<i>System / Interface / Baud rate</i>	1200, 2400, 4800 , 9600, 19200	Baud rate of the data interface
<i>System / Interface / Output format</i>	ASCII CSV	Output format for data transmission (see section 13 TRANSMITTING DATA (USB INTERFACE), page 81)
<i>System / Interface / Decimal separator</i>	Dot (xx.x) Comma (xx,x)	Decimal separator
<i>System / Interface / Output header</i>		Output of a header for <i>Output format: CSV</i>

Menu item	Possible setting	Explanation
<i>System / Clock</i>	<i>Date format</i> <i>Datum</i> <i>Time</i>	Time and date settings (see section 5.4.6 EXAMPLE 2 ON NAVIGATION: SETTING THE DATE AND TIME, page 25)
<i>System / Service information</i>		Hardware version and software version of the meter are displayed.
<i>System / Reset</i>	-	Resets the system settings to the delivery condition (see section 11.3.2 RESETTING THE SYSTEM SETTINGS, page 74).

11.2.2 Data storage

This menu contains all functions to display, edit and erase stored measured values and calibration records (see section 12 DATA MEMORY, page 75).

11.2.3 Automatic Stability control

The automatic *Stability control* (AutoRead) function continuously checks the stability of the measurement signal. The stability has a considerable impact on the reproducibility of measured values.

You can activate or switch off the automatic *Stability control* function (see section 11.2.1 SYSTEM, page 71).

The measured parameter flashes on the display

- as soon as the measured value is outside the stability range
- when the automatic *Stability control* is switched off.

11.3 Reset

You can reset (initialize) all sensor settings and sensor-independent settings separately from each other.

11.3.1 Resetting the measurement settings



The calibration data are reset to the default settings together with the measuring parameters. Recalibrate after performing a reset.

The following settings for pH measurements are reset to the default settings

with the *Reset* function:

pH

Setting	Default settings
<i>Buffer</i>	<i>TEC</i>
<i>Calibration interval</i>	7 d
<i>Unit for slope</i>	mV/pH
<i>Measured parameter</i>	pH
Resolution pH	0.001
Resolution mV	0.1
Asymmetry	0 mV
Slope	-59.2 mV
<i>Man. temperature</i>	25 °C
<i>One point calibration</i>	No

The sensor settings are reset under the *Reset* menu item in the menu for calibration and measurement settings. To open it in the measured value display, press the **<OK>** (or **<F1>/[Menu]**) key.

Oxi

The calibration data are reset to the default settings together with the measuring parameters. Recalibrate after performing a reset.

The following settings for D.O. measurements are reset to the default settings with the *Reset* function:

Setting	Default settings
Kal.-Intervall	14 d
Comparison meas.	Off
<i>Measured parameter</i>	D.O. concentration
<i>Relative slope (S_{Rel})</i>	1.00
<i>Salinity (value)</i>	0,0
<i>Salinity (function)</i>	Off

The sensor settings are reset under the *Reset* menu item in the menu for calibration and measurement settings. To open it in the measured value display, press the **<OK>** (or **<F1>/[Menu]**) key.

Cond

The following settings for conductivity measurements are reset to the default

settings with the *Reset* function:

Setting	Default settings
<i>Kal.-Intervall</i>	150 d
<i>Measured parameter</i>	χ
<i>Cell constant (C)</i> <i>(calibrated)</i>	0.475 cm ⁻¹ or 0.840 cm ⁻¹ (nominal cell constant of the conductivity measuring cell last calibrated)
<i>Cell constant (C)</i> <i>(set)</i>	0.840 1/cm
<i>Temperature compensation</i>	nLF
<i>Reference temperature</i>	25 °C
<i>Temperature coefficient (TC) of the linear temperature compensation</i>	2.000 %/K
<i>TDS multiplier</i>	1.00

The sensor settings are reset under the *Reset* menu item in the menu for calibration and measurement settings. To open it in the measured value display, press the **<OK>** (or **<F1>/[Menu]**) key.

11.3.2 Resetting the system settings

The following system settings can be reset to the default status:

Setting	Default settings
<i>Language</i>	English
<i>Beep</i>	On
<i>Baud rate</i>	4800 Baud
<i>Output format</i>	ASCII
<i>Decimal separator</i>	Dot (xx.x)
<i>Contrast</i>	50 %
<i>Illumination</i>	Auto
<i>Switchoff time</i>	1 h
<i>Stability control</i>	On
<i>Temperature unit</i>	°C

The system settings are reset in the menu, *Storage & config / System / Reset*. To open the *Storage & config* menu in the measured value display, press the **<OK__ >** (or **<F1__>/[Menu]**) key.

12 Data memory

You can transmit measured values (datasets) to the data memory:

- Manual data memory (see section 12.1 MANUAL STORAGE, page 75)
- Automatic storage at intervals (see section 12.2 AUTOMATIC DATA STORAGE AT INTERVALS, page 76)



With each data storage process, the current datasets of the sensors indicated on the display are transmitted to the interface at the same time.

12.1 Manual storage

You can transmit a measurement dataset to the data memory as follows. With each data storage process, the current datasets of the sensors indicated on the display are transmitted to the interface at the same time.

1. Press the **<STR>** key shortly.
The menu for manual data storage appears.

Manual data storage 4 von 200

15.02.2014 11:24:16
 pH 7.0 24.8 °C AR +++
 O2 7.80 mg/l 24.8 °C AR +++

ID number: 1

Continue

Back
15.02.2014
14:15

2. If necessary, change and confirm the ID number (1 ... 10000) with **<▲><▼>** and **<OK>**.
The dataset is stored. The meter switches to the measured value display.

If the memory is full The following window appears if all 200 storage locations are occupied:

A warning dialog box titled "Warning" with the text "Data storage full. Erase?". It contains two radio buttons labeled "Yes" and "No", with "No" selected. At the bottom left is a "Back" button, and at the bottom right is the date and time "15.02.2014 14:15".

You have the following options:

- To erase the entire memory, confirm *Yes*.
- To cancel the storing process and switch to the measured value display, confirm *No*. Then you can e.g. store the recorded data to a PC (see section 12.3.1 EDITING THE MEASUREMENT DATA MEMORY, page 78) and subsequently erase the data memory (see section 12.3.2 ERASING THE MEASUREMENT DATA MEMORY, page 80).

12.2 Automatic data storage at intervals

The storing interval (*Interval*) determines the time interval between automatic data storing processes. With each data storage process, the current datasets of the sensors indicated on the display are transmitted to the interface at the same time.

Configuring the automatic memory function

1. Press the **<STR__>** key.
The menu for automatic data storing appears.

A configuration menu titled "Automatic data storage". It displays the following settings: ID number 1, Interval 30 s, and Duration 180 min. A "Continue" button is shown. Below it is a graphical display of memory usage with a progress bar labeled "0d03h00min" and a "Back" button. The date and time "15.02.2014 14:15" are displayed. Three numbered arrows point to specific elements: 1 points to the "Duration" field, 2 points to the "Continue" button, and 3 points to the graphical display area.


1 Specified entire storing duration
2 Max. available storing duration
3 Graphical display of the memory usage

Settings You can configure the automatic data storing function with the following settings:

Menu item	Possible setting	Explanation
<i>ID number</i>	<i>1 ... 10000</i>	ID number for the dataset series.
<i>Interval</i>	<i>1 s, 5 s, 10 s, 30 s, 1 min, 5 min, 10 min, 15 min, 30 min, 60 min</i>	Storing interval. The lower limit of the storing interval can be restricted by the number of free memory locations. The upper limit is restricted by the storing duration.
<i>Duration</i>	<i>1 min ... x min</i>	Storing duration. Specifies after which time the automatic data storing should be terminated. The lower limit of the storing duration is restricted by the storing interval. The upper limit is restricted by the number of free memory locations.

Starting the automatic storing function

To start the automatic storing function, select *Continue* with **<▲>****<▼>** and confirm with **<OK>**. The meter switches to the measured value display.

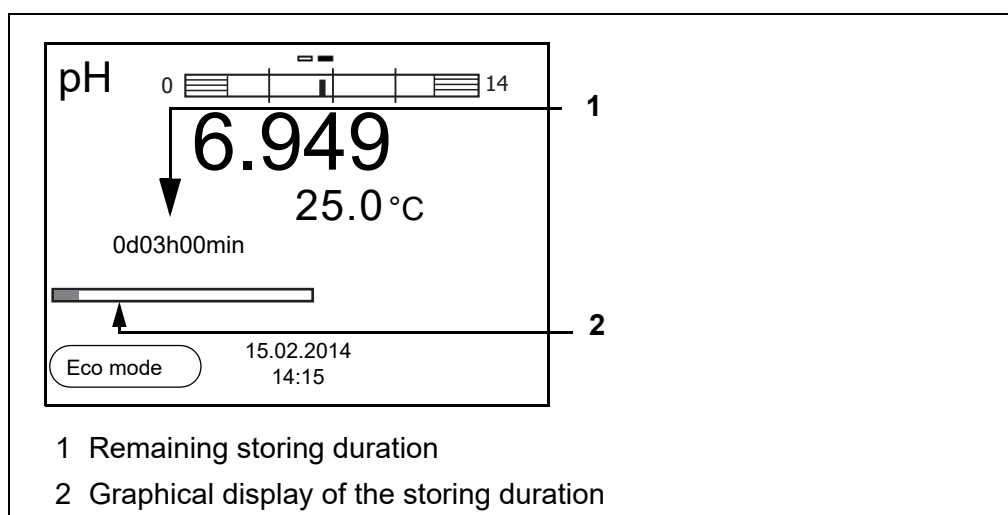

The active automatic storage function can be recognized by the progress bar in the function display. The progress bar indicates the remaining storage duration.



If the automatic storage function is activated, only the following keys are active: Softkeys, **<MODE>**, **<STR__>** and **<On/Off>**. The other keys and the automatic shut-off function are deactivated.

Energy saving mode ([Eco mode])

If the automatic storing function is active, the meter provides an energy saving mode ([Eco mode]) to avoid unnecessary energy consumption. The energy

saving mode switches off functions of the meter that are not required for the automatic storage of measurement data (such as the display). By pressing any key the energy saving mode is switched off again.

Terminating the automatic memory function prematurely

Proceed as follows to switch off the automatic data storing function before the adjusted storing duration has expired:

1. Press the **<STR__>** key.
The following window appears.

A rectangular dialog box with a black border. At the top left, the word "Warning" is displayed. Below it, the text "Stop automatic storage?" is centered. Underneath, there are two options: "Yes" and "No", each with a corresponding radio button. The "No" option is currently selected. At the bottom left, there is a rounded rectangular button labeled "Back". To the right of the button, the date and time "15.02.2014 14:15" are displayed.

2. Using **<▲><▼>**, select Yes and confirm with **<OK>**.
The meter switches to the measured value display.
The automatic data storing function is terminated.

12.3 Measurement data memory

12.3.1 Editing the measurement data memory

The contents of the manual or automatic measurement data memory can be shown on the display and output to the interface.

Each of the measurement data memories has a function to erase the entire contents.

Editing the data memory

The memory is edited in the menu, *Storage & config/ Data storage*. To open the *Storage & config* menu, press the **<OK__>** key in the measured value display.

Open the manual or automatic storage directly with the **<RCL>** or **<RCL__>** key.



The settings are explained here using the manual data memory as an example. The same settings and functions are available for the automatic data memory.

Settings

Menu item	Setting/ function	Explanation
<i>Data storage / Manual data storage / Display</i>	-	Displays all measurement datasets page by page. Further options: ● Scroll through the datasets with <▲><▼>. ● Output the displayed dataset to the interface with <F2>/[USB output]. ● Quit the display with <F1>/[Back].
<i>Data storage / Manual data storage / Erase</i>	-	Erases the entire manual measurement data memory. All calibration data remain stored when this action is performed.
<i>Data storage / Manual data storage / Output to USB</i>	-	Outputs all stored measurement data to the interface.

Display presentation of a dataset

Manual data storage	3 of 64	◆
15.02.2014 11:24:16	ID number: 2	
pH 7.0 24.8 °C AR +++		
O2 7.80 mg/l 24.8 °C AR +++		
Back	15.02.2014 14:15	

Representation of a dataset (USB output)

15.02.2014 09:56:20
MU 6100 L
Ser. no. 08502113
ID number 2
pH1 6.012 24.8 °C, AR, S: +++
O2 7.80 24.8 °C, AR, S: +++
15.02.2014 10:56:20
MU 6100 L
Ser. no. 08502113
ID number 2
pH1 6.012 24.8 °C, AR, S: +++
O2 7.80 24.8 °C, AR, S: +++

Quitting the display

To quit the display of stored measurement datasets, you have the following

options:

- Switch directly to the measured value display with **<MODE>**.
- Quit the display and move to the next higher menu level with **<F1>/[Back]**.

12.3.2 Erasing the measurement data memory

Erasing the measurement data memory (see section 12.3.1 EDITING THE MEASUREMENT DATA MEMORY, page 78).

12.3.3 Measurement dataset

A complete dataset consists of:

- ID number
- Date/time
- Measured values of the connected sensors
- Measured temperature value of the connected sensors or manually set temperature
- AutoRead info: The *AR* indicator appears with the measured value if the AutoRead criterion was met while storing (stable measured value). Otherwise, there is no *AR* indicator.
- Calibration evaluation: +++, ++, +, -, or no evaluation

12.3.4 Memory locations

The MU 6100 L meter has two measurement data memories. The measured values recorded either manually or automatic are stored separately in individual measurement data memories.

Data memory	Maximum number of datasets
<i>Manual data storage</i>	200
<i>Automatic data storage</i>	5000

13 Transmitting data (USB interface)

13.1 Options for data transmission

Via the USB interface you can transmit data to a PC. The following table shows which data are transmitted to the interface in which way:

Data	Control	Operation / description
Current data-sets of the sensors indicated on the display	Manual	● With <F2>[/USB output]. ● Simultaneously with every manual data storage process (see section 12.1 MANUAL STORAGE, page 75).
	Automatic, at intervals	● With <F2__>[/USB output]. Then you can set the transmission interval. ● Simultaneously with every automatic data storing process (see section 12.2 AUTOMATIC DATA STORAGE AT INTERVALS, page 76).
Stored measured values	Manual	● Displayed dataset with <F2>[/USB output] after calling up from the memory. ● All datasets with the <i>Output to USB</i> function. (see section 12.3.1 EDITING THE MEASUREMENT DATA MEMORY, page 78).
Calibration records	Manual	● Calibration record with <F2>[/USB output] (see section 6.2.6 CALIBRATION DATA, page 37).
	Automatic	● At the end of a calibration procedure.



The following rule applies: With the exception of the menus, shortly pressing the **<F2>[/USB output]** key generally outputs the display contents to the interface (displayed measured values, measurement datasets, calibration records).

13.2 Connecting a PC

Connect the MU 6100 L to the PC via the USB interface.

NOTE

When connecting an grounded PC/printer, measurements cannot be performed in grounded media as the values would be incorrect. The USB interface is not galvanically isolated.

Installation of the USB driver on the PC

System requirements of the PC for installation of the USB driver:

1. Insert the supplied installation CD in the CD drive of your PC.
2. Install the driver from the CD.
Follow the Windows installation instructions as necessary.
3. Connect the MU 6100 L to the PC via the USB interface.
The meter is listed as a virtual COM interface among the connections in the Windows instrument manager.

13.3 MultiLab Importer

With the aid of the MultiLab Importer software, you can record and evaluate measurement data with a PC.



More detailed information can be found in the MultiLab Importer operating manual.

14 Maintenance, cleaning, disposal, accessories

14.1 Maintenance


14.1.1 General maintenance activities

The only maintenance activity required is replacing the batteries.



See the relevant operating manuals of the electrodes for instructions on maintenance.

14.1.2 Replacing the batteries

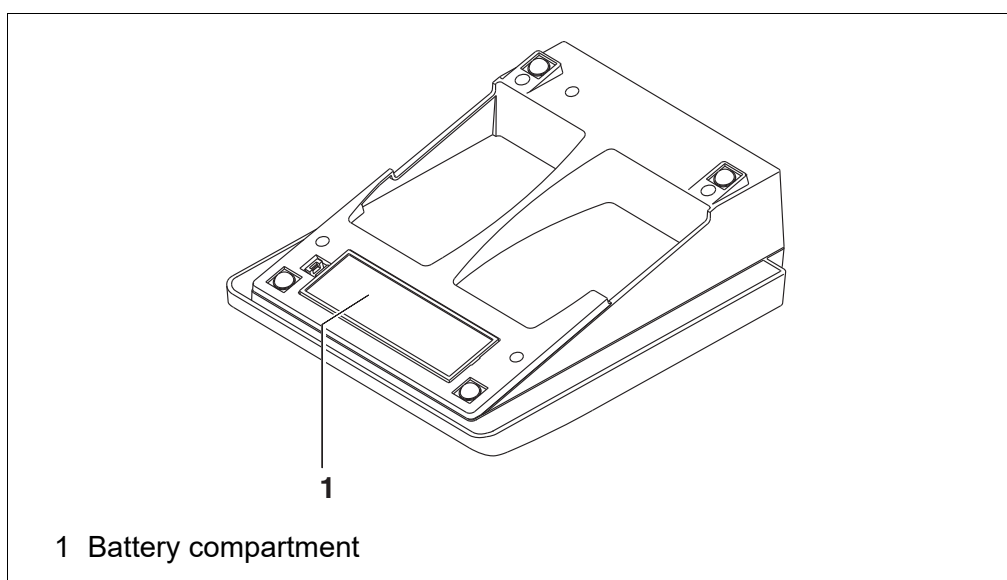

1. Open the battery compartment (1) on the underside of the meter.
2. Remove the batteries from the battery compartment.



CAUTION

Make sure that the poles of the batteries are positioned correctly.

The $\pm$ signs on the batteries must correspond to the $\pm$ signs in the battery compartment.



You can operate the meter either with normal batteries or with rechargeable batteries (Ni-MH). In order to charge the batteries, an external charging device is required.

3. Place four batteries (type Mignon AA) in the battery compartment.
4. Close the battery compartment.
5. Set the date and time
(see section 5.4.6 EXAMPLE 2 ON NAVIGATION: SETTING THE DATE AND TIME, page 25).



When the batteries are nearly discharged, the  status indicator is displayed.



Dispose of used batteries according to the local regulations of your country.

End users within the European Union are obligated to return used batteries (even ecologically compatible ones) to a collection point set up for recycling purposes.

Batteries are marked with the crossed-out waste container symbol. Therefore, they may not be disposed with the domestic waste.

14.2 Cleaning

Occasionally wipe the outside of the measuring instrument with a damp, lint-free cloth. Disinfect the housing with isopropanol as required.



CAUTION

The housing is made of synthetic material (ABS). Thus, avoid contact with acetone or similar detergents that contain solvents. Remove any splashes immediately.

14.3 Packing

This meter is sent out in a protective transport packing. We recommend: Keep the packing material. The original packing protects the meter against damage during transport.



14.4 Disposal

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste.

Instead it's your responsibility to correctly dispose of your equipment at lifecycle-end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

Thank you

14.5 Accessories

14.5.1 General information

Accessories	Order no. (Catalogue No.)	
	EU	NA
Storage tube	662-1167	89236-592
COMMUNICATION KIT (CD-ROM, USB cable, Manual)	662-1225	76470-840
Articulated stand for 4 electrodes and temperature sensor	662-1169	-
Storage vessel	662-1248	-

14.5.2 pH / ORP

Electrodes (pH)	Order no. (Catalogue No.)	
	EU	NA
pHenomenal 110 PH ELECTRODE PHENOMENAL GEL EPOXY BNC	662-1156	76460-452
pHenomenal 111 PH ELECTRODE PHENOMENAL 3IN1 GEL 1M BNC	662-1157	76460-454
pHenomenal 111-3 PH ELECTRODE PHENOMENAL 3IN1 GEL 3M BNC	662-1158	-
pHenomenal 220 PH ELECTRODE PHENOMENAL NACHF. GLAS BNC	662-1159	76460-456
pHenomenal 221 PH ELECTRODE PHENOMENAL 3IN1 GLAS 1M	662-1161	76460-458
pHenomenal 221-3 PH ELECTRODE PHENOMENAL 3IN1 GLAS 3M	662-1162	-
pHenomenal MIC 220 PH ELECTRODE PHENOMENAL MICRO NACHF.BNC	662-1163	-
pHenomenal SPEAR 220 PH ELECTRODE PHENOMENAL KOMB. SPEER TYP	662-1164	-
pHenomenal LS 221 PH ELEKTRODE PHENOMENAL 3IN1 GLAS 1M	662-1247	76460-462
GENERAL PH EKTRODE GEL EPOXY BNC SJ113	662-1382	76460-468
GENERAL PH ELEKTRODE GEL EPOXY BNC DJ113	662-1385	76460-470
SURFACE PH ELEKTRODE GEL EPOXY BNC SF113	662-1388	76460-472
SEMI MICRO PH ELEKTRODE GEL GLASS SM123	662-1391	76460-474
GENERAL PH ELEKTRODE KCL GLASS BNC SJ223	662-1395	76460-476
SURFACE PH ELEKTRODE KCL GLASS BNC SF223	662-1398	76460-478

Electrodes (pH)	Order no. (Catalogue No.)	
	EU	NA
SEMI MICRO PH ELEKTRODE KCL GLASS SM223	662-1402	76460-480
SPEAR PH ELEKTRODE KCL GLASS BNC SP223	662-1405	76460-482

Electrodes (ORP)	Order no. (Catalogue No.)	
	EU	NA
pHenomenal ORP 220 REDOX ELECTRODE PHENOMENAL KOMB. 1M BNC	662-1165	76460-460
REDOX ELEKTRODE GEL EPOXY BNC RD113	662-1408	76460-484
REDOX ELEKTRODE GEL GLASS BNC RD223	662-1409	76460-486

Solutions	EU	NA
Buffer pH 4 AVS TITRINORM, 100 ml	32095.184	-
Buffer pH 7 AVS TITRINORM, 100 ml	32096.187	-
Buffer pH 10 AVS TITRINORM, 100 ml	32040.185	-
Buffer NIST pH 4.01, 30 x 30 ml	1.99001.0001	-
Buffer NIST pH 7, 30 x 30 ml	1.99002.0001	-
Buffer NIST pH 10, 30 x 30 ml	1.99004.0001	-
Buffer pH 1.68, 500 mL	-	BDH5006-500ML
Buffer pH 4, 500 mL	-	BDH5024-500ML
Buffer pH 7, 500 mL	-	BDH5052-500ML
Buffer pH 10, 500 mL	-	BDH5078-500ML
Buffer pH 12.45, 500 mL	-	BDH5096-500ML
Buffer NIST pH 4.006, 500 mL	-	BDH5018-500ML
Buffer NIST pH 6.865, 500 mL	-	BDH5040-500ML
Buffer NIST pH 9.18, 500 mL	-	BDH5066-500ML
Buffer NIST pH 10.012, 500 mL	-	BDH5072-500ML
Storage Solution (3 moles/l KCl), 100 ml	83605.180	BDH7296-0
Cleaning Solution Pepsine/Hydrochloric acid, 100 ml	83603.180	89207-612

14.5.4 Conductivity

Measuring cells	Order no. (Catalogue No.)	
	EU	NA
pHenomenal CO 11 CONDUCTIVITY SENSOR PHENOMENAL 1,5M 8PIN	663-0147	76470-822

Solutions	EU	NA
KCl 0.01 mol/L: 1.413 mS/cm, 100 mL	83607.180	-
KCl 0.01 mol/L: 1.413 mS/cm, 500 mL	83607.290	89236-544
KCl 0.1 mol/L: 12.8 mS/cm, 500 mL	83608.260	89236-546

14.5.5 D.O.

Measuring cells	Order no. (Catalogue No.)	
	EU	NA
pHenomenal OXY 11 OXYGEN SENSOR PHENOMENAL 3M 8 PIN	664-0042	76470-820
Accessories	Order no. (Catalogue No.)	
	EU	NA
MAINTENANCE KIT OXYGEN	664-0049	76460-466

15 What to do if...

15.1 pH/ORP



More information and instructions on cleaning and exchange of sensors are given in the documentation of your sensor.

Error message *OFL, UFL*

The measured value is outside the measuring range.

Cause	Remedy
Electrode:	
– Air bubble in front of the junction	– Remove air bubble
– Air in the junction	– Extract air or moisten junction
– Cable broken	– Replace the electrode
– Gel electrolyte dried out	– Replace the electrode
– The measured value is outside the measuring range of the meter	– Use a suitable electrode

Error message, *Error*

Cause	Remedy
Electrode:	
– The values determined for zero point and slope of the electrode are outside the allowed limits.	– Recalibrate
– Junction contaminated	– Clean the junction
– Electrode broken	– Replace the electrode
Buffer solutions:	
– Incorrect buffer solutions	– Change calibration procedure
– Buffer solutions too old	– Use only once. Note the shelf life
– Buffer solutions depleted	– Change solutions

No stable measured value	Cause	Remedy
	Electrode:	
	– Junction contaminated	– Clean the junction
	– Membrane contaminated	– Clean membrane
	Test sample:	
	– pH value not stable	– Measure with air excluded if necessary
	– Temperature not stable	– Temper if necessary
	Electrode + test sample:	
	– Conductivity too low	– Use a suitable electrode
	– Temperature too high	– Use a suitable electrode
	– Organic liquids	– Use a suitable electrode

Obviously incorrect measured values

Cause	Remedy
Electrode:	
– Electrode unsuitable	– Use a suitable electrode
– Temperature difference between buffer and test sample too great	– Adjust temperature of buffer or sample solutions
– Measurement procedure not suitable	– Follow special procedure

15.2 Dissolved oxygen



More information and instructions on cleaning and exchange of sensors are given in the documentation of your sensor.

Error message OFL, UFL

The measured value is outside the measuring range.

Cause	Remedy
– The measured value is outside the measuring range of the meter	– Use a suitable D.O. sensor

Error message, <i>Error</i>	Cause	Remedy
	– D.O. sensor contaminated – Electrolyte depleted	– Clean the D.O. sensor – Change the electrolyte and membrane cap – If necessary, replace the D.O. sensor

Error message <i>Leak</i>	Cause	Remedy
	– Membrane cap not screwed on tight enough – Hole in the membrane	– Screw membrane head tighter – Replace and refill the membrane cap

15.3 Conductivity



More information and instructions on cleaning and exchange of sensors are given in the documentation of your sensor.

Error message
OFL, UFL

The measured value is outside the measuring range.

Cause

- The measured value is outside the measuring range of the meter

Remedy

- Use a suitable measuring cell

Error message,
Error

Cause

- Measuring cell contaminated
- Calibration solution not suitable

Remedy

- Clean cell and replace it if necessary
- Check the calibration solutions

15.4 General information

Sensor symbol flashes	Cause – Calibration interval expired	Remedy – Recalibrate the measuring system
Display 	Cause – Batteries almost empty	Remedy – Replace the batteries (see section 14.1 MAINTENANCE, page 83)
Meter does not react to keystroke	Cause – Operating condition undefined or EMC load unallowed	Remedy – Processor reset: Press the <OK> and <On/Off> key simultaneously
You want to know which software version is in the meter	Cause – E.g., a question by the service department	Remedy – Switch on the meter. Open the menu, <OK__ > / <i>Storage & config / System / Service information</i>. The instrument data are displayed.

16 Firmware update

Available firmware updates are provided on the Internet.

With the firmware update program and a PC you can update the firmware of the MU 6100 L to the newest version.

For the update you have to connect the meter to a PC.

For the update via the USB interface, the following is required:

- a free USB interface (virtual COM port) on the PC
- the driver for the USB interface (on the enclosed CD-ROM)
- the USB cable (included in the scope of delivery of the MU 6100 L).

1. Install the downloaded firmware update on a PC.
An update folder is created in the Windows start menu.
If an update folder already exists for the meter (or meter type), the new data is displayed there.
2. In the windows start menu, open the update folder and start the firmware update program.
3. Using the USB interface cable, connect the MU 6100 L to a USB interface (virtual COM port) of the PC.
4. Switch on the MU 6100 L.
5. In the firmware update program, start the update process with OK.
6. Follow the instructions of the firmware update program.
During the programming process, a corresponding message and a progress bar (in %) are displayed.
The programming process takes approx. three minutes. A terminatory message is displayed after a successful programming process. The firmware update is completed.
7. Disconnect the MU 6100 L from the PC.
The MU 6100 L is ready for operation again.

After switching the meter off and on you can check whether the meter has taken over the new software version (see YOU WANT TO KNOW WHICH SOFTWARE VERSION IS IN THE METER, PAGE 94).

17 Glossary

pH/ORP/ISE

Asymmetry	see zero point
Electromotive force of an electrode	The electromotive force U of the electrode is the measurable electromotive force of an electrode in a solution. It equals the sum of all the galvanic voltages of the electrode. Its dependency on the pH results in the electrode function, which is characterized by the parameters, slope and zero point.
Junction	The junction is a porous body in the housing wall of reference electrodes or electrolyte bridges. It arranges the electrical contact between two solutions and makes the electrolyte exchange more difficult. The expression, junction, is also used for ground or junction-less transitions.
ORP	The ORP is caused by oxidizing or reducing substances dissolved in water if these substances become effective on an electrode surface (e.g. a gold or platinum surface).
pH value	The pH value is a measure of the acidic or basic effect of an aqueous solution. It corresponds to the negative decadic logarithm of the molal hydrogen ions activity divided by the unit of the molality. The practical pH value is the value of a pH measurement.
Potentiometry	Name of a measuring technique. The signal (depending on the measured parameter) of the electrode is the electrical potential. The electrical current remains constant.
Slope	The slope of a linear calibration function.
Zero point	The zero point of a pH electrode is the pH value at which the electromotive force of the pH electrode at a specified temperature is zero. Normally, this is at 25 °C.

Dissolved oxygen

D.O. partial pressure	Pressure caused by the oxygen in a gas mixture or liquid.
D.O. saturation	Short name for the relative D.O. saturation.
OxiCal®	Name for a procedure to calibrate D.O. measuring systems in water vapor-saturated air.
Salinity	The absolute salinity S_A of seawater corresponds to the relationship of the mass of dissolved salts to the mass of the solution (in g/Kg). In practice, this dimension cannot be measured directly. Therefore, the practical salinity according to IOT is used for oceanographic monitoring. It is determined by measuring the electrical conductivity.
Salt content	General designation for the quantity of salt dissolved in water.
Slope (relative)	The relative slope expresses the relation of the slope value to the value of a theoretical reference sensor of the same construction type.

Conductivity

Cell constant C	Characteristic quantity of a conductivity measuring cell, depending on the geometry.
Conductivity	Short form of the expression, specific electrical conductivity. It corresponds to the reciprocal value of the resistivity. It is a measured value of the ability of a substance to conduct an electric current. In water analysis, the electrical conductivity is a dimension for the ionized substances in a solution.
Reference temperature	Fixed temperature value to compare temperature-dependent measured values. For conductivity measurements, the measured value is converted to a conductivity value at a reference temperature of 20 °C or 25 °C.
Resistance	Short name for the specific electrolytic resistance. It corresponds to the reciprocal value of the electrical conductivity.
Salinity	The absolute salinity S_A of seawater corresponds to the relationship of the mass of dissolved salts to the mass of the solution (in g/Kg). In practice, this dimension cannot be measured directly. Therefore, the practical salinity according to IOT is used for oceanographic monitoring. It is determined by measuring the electrical conductivity.
Salt content	General designation for the quantity of salt dissolved in water.
Temperature coefficient	Value of the slope α of a linear temperature function. $\mathcal{R}_{T_{\text{Ref}}} = \mathcal{R}_{\text{Meas}} * \frac{1}{1 + \alpha * (T - T_{\text{Ref}})}$
Temperature compensation	Name of a function that considers the temperature influence on the measurement and converts it accordingly. Depending on the measured parameter to be determined, the temperature compensation functions in different ways. For conductimetric measurements, the measured value is converted to a defined reference temperature. For potentiometric measurements, the slope value is adjusted to the temperature of the test sample but the measured value is not converted.

General topics

Adjusting	To manipulate a measuring system so that the relevant value (e.g. the displayed value) differs as little as possible from the correct value or a value that is regarded as correct, or that the difference remains within the tolerance.
AutoRange	Name of the automatic selection of the measuring range.
Calibration	Comparing the value from a measuring system (e.g. the displayed value) to the correct value or a value that is regarded as correct. Often, this expression is also used when the measuring system is adjusted at the same time (see adjusting).

Channel	A channel is a display indication that corresponds to a physical connection on the meter.
Measured parameter	The measured parameter is the physical dimension determined by measuring, e.g. pH, conductivity or D.O. concentration.
Measured value	The measured value is the special value of a measured parameter to be determined. It is given as a combination of the numerical value and unit (e.g. 3 m; 0.5 s; 5.2 A; 373.15 K).
Molality	Molality is the quantity (in Mol) of a dissolved substance in 1000 g solvent.
Reset	Restoring the original condition of all settings of a measuring system.
Resolution	Smallest difference between two measured values that can be displayed by a meter.
Stability control (AutoRead)	Function to control the measured value stability.
Standard solution	The standard solution is a solution where the measured value is known by definition. It is used to calibrate a measuring system.
Temperature function	Name of a mathematical function expressing the temperature behavior of a test sample, a sensor or part of a sensor.
Test sample	Designation of the test sample ready to be measured. Normally, a test sample is made by processing the original sample. The test sample and original sample are identical if the test sample was not processed.

18 Index

A

Air calibration beaker	51
Automatic switch-off function	21
AutoRead	
ORP	39, 41, 49, 57
pH	28

B

Battery compartment	16, 83
---------------------------	--------

C

Calibration	
ISE	43
pH	30, 40
Calibration evaluation	54, 60
ISE	47
pH	37
Calibration interval	64, 66, 69
Calibration points	
pH	36
Comparison measurement (D.O.)	51
Connecting a PC	82
Connectors	20
Control standard	58

D

Dataset	80
Date and time	25
Default settings	
Measured parameter	72, 73
System settings	74
Display	19

E

Energy saving mode	77
--------------------------	----

F

Firmware update	95
-----------------------	----

I

Initial commissioning	15
Initialize	72
Interval for calibration	64, 66, 69

K

Keys	18
------------	----

M

Measured value display	22
------------------------------	----

Measurement accuracy	64, 66, 69
Measurement data memory	
Edit	78
Erase	78
Measurement dataset	80
Measuring	49, 56
ISE	41
ORP	39
pH	28
Menus (navigation)	22
Messages	23

P

pH buffer sets	62
Print	81

R

Reset	72
-------------	----

S

Scope of delivery	15
Single-point calibration	
pH	31, 34
Slope	
ISE	43
pH	30
Slope, relative	51
Stability control	
Automatic	72
Storing in memory	75
Automatic	76
Manual	75
Storing interval	76

T

TDS factor	70
Temperature compensation	69
Temperature measurement	50
ISE	42
pH	29, 40
Three-point calibration	
ISE	46
pH	32, 35
Transmitting data	81
Automatic	81
Manual	81
Transmitting measured values	81
Two-point calibration	
ISE	45
pH	31, 35

Z

Zero point of pH electrode 30

19 Technical service

Web resources

Visit the VWR website at www.vwr.com for:

- Complete technical service contact information
- Access to VWR's Online Catalogue, and information about accessories and related products
- Additional product information and special offers

Contact us:

For information or technical assistance contact your local VWR representative or visit www.vwr.com.

20 Warranty

VWR warrants that this product will be free from defects in material and workmanship for a period of three (3) years from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

21 Compliance with local laws and regulations

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorizations necessary to run or use the product in its local environment. VWR will not be held liable for any related omission or for not obtaining the required approval or authorization, unless any refusal is due to a defect of the product.

Local VWR offices in Europe and Asia Pacific

Austria

VWR International GmbH
Graumannsgasse 7
1150 Vienna
Tel.: +43 01 97 002 0
info.at@vwr.com

Belgium

VWR International BV
Researchpark Haasrode 2020
Geldenaaksebaan 464
3001 Leuven
Tel.: +32 016 385 011
vwr.be@vwr.com

Canada

VWR International
2360 Argentia Road
Mississauga, Ontario L5N 5Z7
Tel.: +1 800 932 5000
Canada_Orders@vwr.com

China

VWR (Shanghai) Co., Ltd
Bld.No.1, No.3728 Jinke Rd,
Pudong New District
Shanghai 200123- China
Tel.: +400 821 8006
info_china@vwr.com

Czech Republic

VWR International s. r. o.
Veetee Business Park
Pražská 442
CZ - 281 67 Strábrná Skalice
Tel.: +420 321 570 321
info.cz@vwr.com

Denmark

VWR International A/S
Tobaksvejen 21
2860 Søborg
Tel.: +45 43 86 87 88
info.dk@vwr.com

Finland

VWR International Oy
Valimotie 9
00380 Helsinki
Tel.: +358 09 80 45 51
info.fi@vwr.com

France

VWR International S.A.S.
Le Périgares – Bâtiment B
201, rue Carnot
94126 Fontenay-sous-Bois cedex
Tel.: 0 825 02 30 30* (national)
Tel.: +33 (0) 1 45 14 85 00 (international)
info.fr@vwr.com
* 0,18 € TTC/ min + prix appel

Germany

VWR International GmbH
Hilpertstraße 20a
D - 64295 Darmstadt
Tel.: 0800 702 00 07* (national)
Tel.: +49 (0) 6151 3972 0 (international)
info.de@vwr.com
*Freecall

Hungary

VWR International Kft.
Simon László u. 4.
4034 Debrecen
Tel.: +36 (52) 521-130
info.hu@vwr.com

India

Avantor Performance Materials India Limited
17th Floor, Building No. 5, Tower C
DLF Cyber City Phase – III
Gurgaon - 122002, Haryana
Tel.: +91-1244-65-6700
help@avantorinc.com

Ireland

VWR International Ltd
International (Northern Ireland) Ltd
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: +353 01 88 22 222
sales.ie@vwr.com

Italy

VWR International S.r.l.
Via San Giusto 85
20153 Milano (MI)
Tel.: +39 02-3320311
info.it@vwr.com

Korea

VWR International ~
17, Daehak 4-ro, Yeongtong-gu,
Suwon-si, Gyeonggi-do
Tel.: +82 31 645 7256
saleskorea@avantorsciences.com

The Netherlands

VWR International B.V.
Postbus 8198
1005 AD Amsterdam
Tel.: +31 020 4808 400
info.nl@vwr.com

Mexico

VWR International, S.de R.L. de C.V.
Km. 14.5 Carretera
Tlalnepantla-Cuautitlán
Col. Lechería
Tultitlán Edo. de México
CP 54940
Tel.: +52 (55) 5005 0100
vwrmx@vwr.com

Middle East & Africa

VWR International FZ-LLC
DSP Laboratory Complex
125, Floor 01
Dubai, United Arab Emirates
Tel.: +971 4 5573271
info.mea@vwr.com

Norway

VWR International AS
Brynsalleen 4
0667 Oslo
Tel.: +47 22 90 00 00
info.no@vwr.com

Poland

VWR International Sp. z o.o.
Limbowa 5
80-175 Gdansk
Tel.: +48 58 32 38 200
info.pl@vwr.com

Portugal

VWR International –
Material de Laboratório, Lda
Centro Empresarial de Alfragide
Rua da Indústria, nº 6
2610-088 Amadora
Tel.: +351 21 3600 770
info.pt@vwr.com

Singapore

VWR Singapore Pte Ltd
18 Gul Drive
Singapore 629468
Tel: +65 6505 0760
sales.sg@vwr.com

Spain

VWR International Eurolab S.L.U.
C/ Tecnología 5-17
A-7 Llinars Park
08450 - Llinars del Vallès
Barcelona
Tel.: +34 902 222 897
info.es@vwr.com

Sweden

VWR International AB
Fagerstagatan 18a
163 94 Stockholm
Tel.: +46 (0) 8 621 34 00
kundservice.se@vwr.com

Switzerland

VWR International GmbH
Lerzenstrasse 16/18
8953 Dietikon
Tel.: +41 (0) 44 745 13 13
info.ch@vwr.com

UK

VWR International Ltd
Customer Service Centre
Hunter Boulevard - Magna Park
Lutterworth
Leicestershire
LE17 4XN
Tel.: +44 (0) 800 22 33 44
uksales@vwr.com

United States

VWR International, LLC
100 Matsonford Road
Building One Suite 200
Radnor, PA 19087
Tel.: +1 800 932 5000
VWRCustomerService@vwr.com