Instruction manual

VWR® Rocking Platform Shakers & 3-D Rotator Wavers Incubating Rocking Platform Shakers & Incubating 3-D Rotator Wavers

EU cat. no

Rocker Platform Shaker: 444-0756

3-D Rotator Waver: 444-0759

Incubating Rocker Platform Shaker: 444-0762

Incubating 3-D Rotator Waver: 444-0765

EN	- English		1
FR	- Français		16
ES	- Español		30
DE	- Deutsch		46
IT	- Italiano		61
PT	- Português		76
BG	- Български		91
CZ	- Čeština		95
DK	- Dansk		99
EE	- Eesti		103
FI	- Suomi		107
GR	- Ελληνικά		111
HR	- Hrvatska		115
HU	- Magyar		119
IS	- Íslenska		123
LT	- Lietuvių		127
LV	- Latviešu		133
NL	- Nederlands		137
NO	- Norsk		141
PL	- Polski		145
RO	- Română		149
SI	- Slovenščina		153
SK	- Slovenčina		157
SV	- Svenska		161

Legal address of Manufacturer

Europe

VWR International bv,
Researchpark Haasrode 2020
Geldenaaksebaan 464 B-3001
Leuven
+3216385011
be.vwr.com

UK Importer:

VWR International Ltd
Hunter Boulevard, Magna Park
Lutterworth, Leicestershire, LE17 4XN
uk.vwr.com

Country of Origin

United States

Table of Contents

Safety information	4
Package Contents	4
Installation	4
Intended Use	4
Maintenance & Servicing	5
Environmental Conditions	5
Specifications	
<i>Rocking Platform</i>	6
<i>3-D Rotator Waver</i>	6
<i>Incubating Rocking Platform Shaker</i>	6
<i>Incubating 3-D Rotator Waver</i>	6
Control Panel	
<i>Rocker/Waver</i>	7
<i>Incubating Rocker/Waver</i>	8
Operating Instructions	
Rocker/Waver	
<i>Getting Ready</i>	9
<i>Setting Tilt</i>	9
<i>Setting Speed</i>	9
<i>Setting Time</i>	9
<i>Setting Timed Mode</i>	9
<i>Turning Off Unit</i>	10
Incubating Rocker/Waver	
<i>Getting Ready</i>	10
<i>Set Temperature</i>	10
<i>Caution Hot Indicator</i>	10
<i>Setting Tilt</i>	10
<i>Setting Speed</i>	10
<i>Setting Time</i>	11
<i>Setting Timed Mode</i>	11
<i>Turning Off Unit</i>	11
Troubleshooting	12
Accessories	12
Technical Service	13
Warranty	13
Disposal	14
Distributor Contact Information	15

Safety Instructions

Please read the entire instruction manual before operating the Rocker/Waver.

WARNING! DO NOT use the Rocker/Waver in a hazardous atmosphere or with hazardous materials for which the unit was not designed. Also, the user should be aware that the protection provided by the equipment may be impaired if used with accessories not provided or recommended by the manufacturer or used in a manner not specified by the manufacturer.

Always operate unit on a level surface for best performance and maximum safety.

DO NOT lift unit by the platform tray or lid.

CAUTION! To avoid electrical shock, completely cut off power to the unit by disconnecting the power cord from the unit or unplug from the wall outlet. Disconnect unit from the power supply prior to maintenance and servicing.

Spills should be removed promptly. Biohazard spills should be cleaned up using approved laboratory procedures. Solvent spills are a fire hazard. Stop the unit immediately, and **DO NOT** operate until cleanup is complete and vapors have dissipated (motor brushes arc and may ignite flammable vapors).

DO NOT immerse the unit for cleaning.

DO NOT operate the unit if it shows signs of electrical or mechanical damage.

Earth Ground - Protective Conductor Terminal

Alternating Current

Package Contents

*Rocking Platform Shaker or 3-D Rotator Waver or
Incubating Rocker Platform Shaker or Incubating 3-D Rotator Waver*

Description	Quantity
234cm detachable power cords (EU and UK) CH adapter	3
Non-skid rubber mat (not included with incubating units)	1

Installation

Upon receiving the VWR Rocker/Waver, check to ensure that no damage has occurred in shipment. It is important that any damage that occurred in transport is detected at the time of unpacking. If you do find such damage the carrier must be notified immediately.

After unpacking, place the Rocker/Waver on a level bench or table, away from explosive vapors. Ensure that the surface on which the unit is placed will withstand typical heat produced by the unit and place the unit a minimum of six (6) inches from vertical surfaces. Always place the unit on a sturdy work surface.

The Rocker/Waver is supplied with a power cord that is inserted into the IEC connector on the back of the unit first, then it can be plugged into a properly grounded outlet. The 230V unit plugs into a 230 volt, 50/60 Hz source.

Intended use

These Rockers and Wavers are intended for general laboratory use. For research use only. Not intended for any animal or human therapeutic or diagnostic use.

Maintenance & Servicing

The Rocker/Waver is built for long, trouble-free, dependable service. No lubrication or other technical user maintenance is required. It needs no user maintenance beyond keeping the surfaces clean. The unit should be given the care normally required for any electrical appliance. Avoid wetting or unnecessary exposure to fumes. Spills should be removed promptly. Do not use a cleaning agent or solvent on the front panel which is abrasive or harmful to plastics, nor one which is flammable. Always ensure the power is disconnected from the unit prior to any cleaning. If the unit ever requires service, contact your VWR representative.

Environmental Conditions

Open Air Rockers and Wavers

Operating Conditions: Indoor use only.

Temperature*: For use in CO₂ environments, incubators or cold rooms ranging from -10 to 60°C

Humidity: 80% relative humidity, non-condensing

Altitude: 0 to 2000 M above sea level

Non-Operating Storage:

Temperature*: -20 to 65°C

Humidity: maximum 80% relative humidity, non-condensing

Installation Category II and Pollution Degree 2 in accordance with IEC 664.

* **Avoid cold starts:** Unit is not designed to start after being in a cold room environment. Bring unit into cold room from a room temperature environment, operate and remove unit from cold room as soon as operation is complete.

Incubating Rockers and Wavers

Operating Conditions: Indoor use only.

Temperature: 5 to 40°C

Humidity: 80% relative humidity, non-condensing

Altitude: 0 to 2000 M above sea level

Non-Operating Storage:

Temperature: -20 to 65°C

Humidity: maximum 80% relative humidity, non-condensing

Installation Category II and Pollution Degree 2 in accordance with IEC 664

Product Specifications

Platform Rocker

Overall dimensions (L x W x H): 43,2 x 27,9 x 12,7cm

Platform dimensions (L x W): 32,4 x 25,4cm

Electrical (50/60 Hz): 230 volts AC, 5 amps, 25 watts

Fuses: 5mm x 20mm, 5 amp quick acting

Speed range: 1 to 50rpm*

Speed accuracy: ±1rpm

Tilt angle: 0 to 15°*

Weight capacity: 4,5kg**

Timer: digital, 1 second to 9999 minutes

(increased in 1 second increments)

Controls: see page 7

Ship weight: 7kg

* Maximum speed/tilt angle may vary with heavy or unbalanced loads.

** Centered on tray.

3-D Rotator Waver

Overall dimensions (L x W x H): 43,2 x 27,9 x 15,2cm

Platform dimensions (L x W): 29,9 x 22,2cm

Electrical (50/60 Hz): 230 volts AC, 5 amps, 25 watts

Fuses: 5mm x 20mm, 5 amp quick acting

Speed range: 1 to 30rpm*

Speed accuracy: ±1rpm

Tilt angle: 0 to 20°*

Weight capacity: 2,3kg**

Timer: digital, 1 second to 9999 minutes

(increased in 1 second increments)

Controls: see page 7

Ship weight: 7.3kg

* Maximum speed/tilt angle may vary with heavy or unbalanced loads.

** Centered on tray.

Incubating Rocker

Overall dimensions (L x W x H): 43,2 x 27,9 x 26,7cm

Platform dimensions (L x W): 25,4 x 19,1cm

Electrical (50/60 Hz): 230 volts AC, 5 amps, 450 watts

Fuses: 5mm x 20mm, 5 amp quick acting

Temperature range: ambient +5° to 65°C

Temperature uniformity: +/-0,5°C at 37°C

Speed range: 1 to 50rpm*

Speed accuracy: ±1rpm

Tilt angle: 0 to 15°*

Weight capacity: 4,5kg**

Timer: digital, 1 second to 9999 minutes

(increased in 1 second increments)

Controls: see page 8

Ship weight: 10kg

* Maximum speed/tilt angle may vary with heavy or unbalanced loads.

** Centered on tray.

Incubating Waver

Overall dimensions (L x W x H): 43,2 x 27,9 x 26,7cm

Platform dimensions (L x W): 23,5 x 18,4cm

Electrical (50/60 Hz): 230 volts AC, 5 amps, 450 watts

Fuses: 5mm x 20mm, 5 amp quick acting

Temperature range: ambient +5° to 65°C

Temperature uniformity: +/-0,5°C at 37°C

Speed range: 1 to 30rpm*

Speed accuracy: ±1rpm

Tilt angle: 0 to 20°*

Weight capacity: 2,3kg**

Timer: digital, 1 second to 9999 minutes

(increased in 1 second increments)


Controls: see page 8

Ship weight: 10kg

* Maximum speed/tilt angle may vary with heavy or unbalanced loads.

** Centered on tray.


Control Panel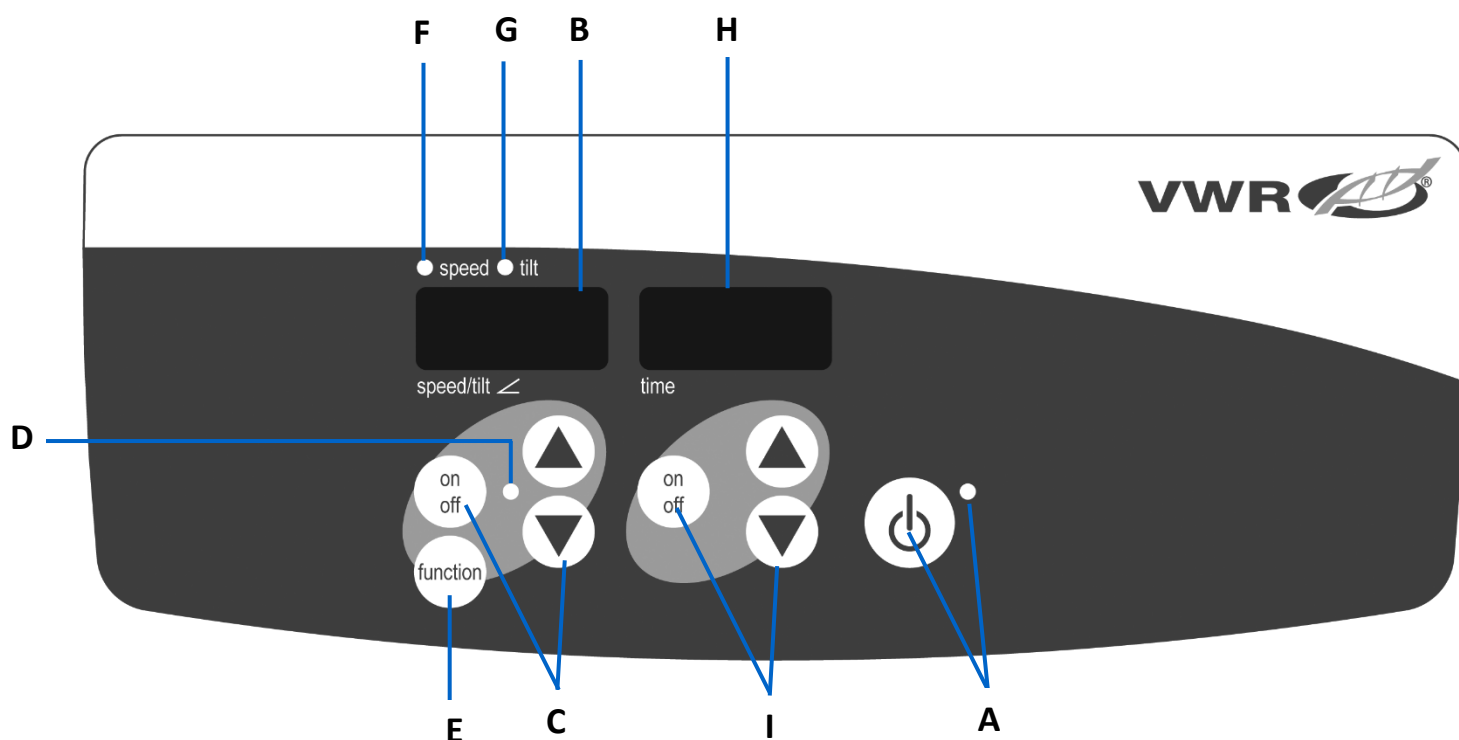

Rocker/Waver:

The front panel of the Rocker/Waver contains all the controls and displays needed to operate the unit.

- A. Standby button/standby indicator light:** The standby indicator light will illuminate when the unit is plugged in. The unit will be in standby mode. Press the standby button to activate the speed/tilt and time functions. The standby indicator light will shut off and the speed/tilt display and time display will illuminate. Press the standby button again and the unit will once again be in standby mode.
- B. Speed/tilt display:** Displays the speed and tilt of the unit.
- C.** Up/down arrows for set-point control. On/off button starts/stops the rocking/waving function.
- D.** The speed/tilt indicator light will be illuminated when the unit is rocking/waving.
- E. Function button:** Press to choose the function you are setting speed or tilt.
- F. Speed indicator light:** Illuminates when the speed is displayed.
- G. Tilt indicator light:** Illuminates when the tilt angle is displayed.
- H. Time display:** Displays accumulated time (continuous mode) or how much time is remaining (timed mode).
- I.** Up/down arrows for set-point control. On/off button starts/stops the time function. The display range is from 0 to 9999 minutes in one (1) second increments. The display will indicate minutes and seconds until the timer reaches 99 minutes and 59 seconds (99:59), then the display will automatically display minutes up to 9999.

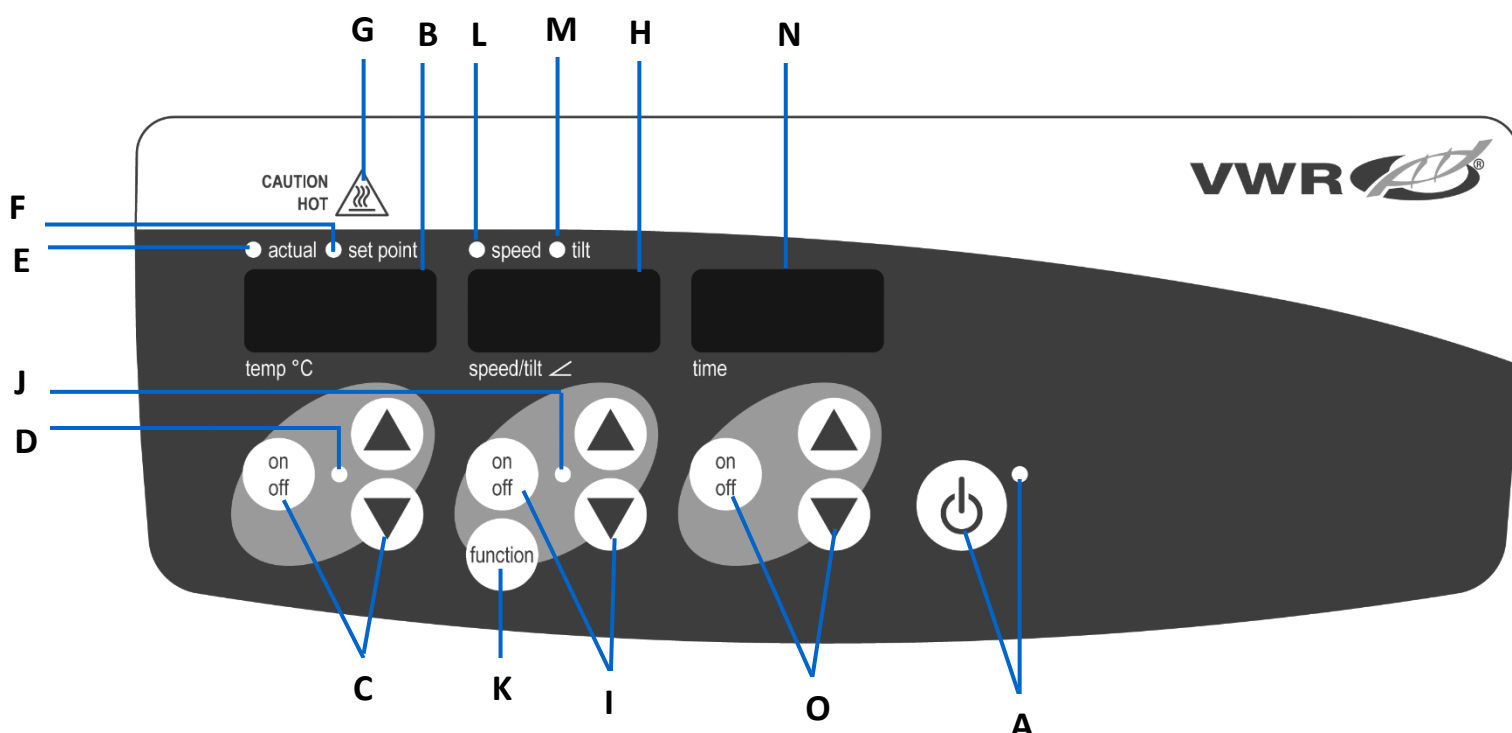Incubating Rocker/Waver:

The front panel of the Incubating Rocker/Waver contains all the controls and displays needed to operate the unit.

- A. Standby button/standby indicator light:** The standby indicator light will illuminate when the unit is plugged in. The unit will be in standby mode. Press the standby button to activate the temperature, speed/tilt, and time functions. The standby indicator light will shut off and the temperature, speed/tilt and time displays will illuminate. Press the standby button again and the unit will once again be in standby mode.
- B. Temperature display:** Displays the actual/set-point temperatures in conjunction with the actual/set-point indicator lights.
- C.** Up/down arrows for set-point control. On/off button starts/stops the heat function.
- D.** The heat indicator light will be illuminated when the unit is heating.
- E. Actual indicator light:** Illuminates when the temperature displayed is the actual temperature of the air in the chamber.
- F. Set-point indicator light:** Illuminates when the set-point temperature is displayed.
- G. Caution hot indicator light:** Illuminates when the air temperature of the chamber is above 40°C (104°F).
- H. Speed/tilt display:** Displays the speed and tilt of the unit.
- I.** Up/down arrows for set-point control. On/off button starts/stops the rocking/waving function.
- J.** The speed/tilt indicator light will be illuminated when the unit is rocking/waving.
- K. Function button:** Press to choose the function you are setting, speed, or tilt.
- L. Speed indicator light:** Illuminates when the speed is displayed.
- M. Tilt indicator light:** Illuminates when the tilt angle is displayed.
- N. Time display:** Displays accumulated time (continuous mode) or how much time is remaining (timed mode).
- O.** Up/down arrows for set-point control. On/off button starts/stops the time function. The display range is from 0 to 9999 minutes in one (1) second increments. The display will indicate minutes and seconds until the timer reaches 99 minutes and 59 seconds (99:59), then the display will automatically display minutes up to 9999.

Operating Instructions

Rocker/Waver

The VWR Rocker/Wavers are used for gentle mixing of laboratory samples. These units have been designed for the speed/tilt and time functions to work independently of one another. The speed/tilt can be re-set without re-setting the timer and the timer can be stopped and started without interrupting the rocking/waving functions.

1. Getting ready:

- a. Plug the power cord into a properly grounded outlet. The standby indicator light will illuminate, verifying power to the unit. Manually move the platform tray to a near horizontal "home" position before powering unit.
- b. Press the standby button to move the unit from standby mode. The standby indicator light will shut off and the speed/tilt and time displays will illuminate and display the previously used settings.

2. Setting tilt: Electronic tilt enables the user to adjust the angle of tray movement while the unit is rocking/waving or while the unit is stopped.

- a. Press the function button below the speed/tilt display until the tilt indicator light illuminates. You are now ready to set the tilt angle.
- b. Press the up/down arrows below the speed/tilt display until you reach the desired angle. When you release the arrow button, the display will blink off and then on indicating the new set angle has been accepted. The unit will complete one rotation at the previously set angle, then smoothly change to the newly programmed angle. The tilt indicator light will flash until the transition to the new tilt angle is complete.

3. Setting speed:

- a. Press the function button below the speed/tilt display until the speed indicator light illuminates. You are now ready to set the speed.
- b. Press the up/down arrows below the speed/tilt display until you reach the desired speed. When you release the arrow button, the display will blink off and then on indicating the new set speed has been accepted.
- c. Press the on/off button to start the rocking/waving function. The indicator light below the speed/tilt display will flash until the transition to the set speed is complete, then the light will stay illuminated indicating the rocking/waving function is in use.
- d. To stop rocking/waving function, press the on/off button below the speed/tilt display. The unit will complete one full rotation then stop in the horizontal "home" position. The speed indicator light will flash until the rotation is complete. When the cycle is complete, the unit will automatically move to standby mode and the standby indicator light will illuminate.

Operating Tips

When the unit is running at slow speeds or high angles, making large changes to the tilt angle or speed may take several minutes to complete. The quickest way to make large changes to the speed or tilt angle is to stop the unit, change the speed or tilt angle, then restart by pressing the on/off button below the speed/tilt display.

4. Setting time to zero (0:00) and continuous mode: Accumulated time.

- a. Press and hold the on/off button below the time display. After three (3) seconds, the display will indicate the previous set time.
- b. Simultaneously press both the up and down arrows, the display will indicate zero (0:00). The unit time is now set to zero (0:00) minutes. Alternately, you can use the up/down arrows to get to zero (0:00).
- c. Press the on/off button below the time display. The display will indicate the accumulated time. The up/down arrows will become inactive. To stop timer, press the on/off button again. **IMPORTANT:** This will **NOT** interrupt the rocking/waving function. Press the on/off button below the speed/tilt display to interrupt the rocking/waving function.
- d. To re-set, press and hold the on/off button below the time display. After three (3) seconds, the display will indicate the previous set time, which was zero (0:00).

5. Setting timed mode: Programmed time.

- a. Press the up/down arrows below the time display until you reach the desired time.
- b. Start this function by pressing the on/off button below the time display. The unit will run for the selected time, the up/down arrows will become inactive while the timer is running. The unit will stop rocking/waving when time display reaches zero (0:00). Four (4) audible beeps will indicate the time down function is complete. The time display will default back to the set time. To repeat for the same time, simply depress the on/off button again.
- c. To interrupt an automatic timing cycle before it is completed, press the on/off button below the time display. The time display will flash off and on to indicate the time function is on "hold". **IMPORTANT:** This will **NOT** interrupt the rocking/waving function. Press the on/off button below the speed/tilt display to interrupt the rocking/waving function. Restart the timer by pressing the on/off button below the time display. Unit will continue counting down to zero (0:00). When the display reaches zero (0:00), you will hear the four (4) audible beeps that indicate the time down function is complete and the rocking/waving function will cease.

6. Turning unit off:

- a. To turn the unit off, press the standby button. The speed/tilt and time displays will be blank, the standby indicator light will illuminate. The Rocker/Waver should be kept in standby mode when not in use. To completely cut off power to the unit, disconnect the power cord from the unit or un-plug from the wall outlet.

Operating Tips

Built-in memory maintains the last used speed/tilt and time settings during a power interruption.

A built-in program will shut power off to the motor if the platform tray is prevented from rocking/waving, or the unit is overloaded beyond its recommended weight capacity.

Incubating Rocker/ Wavers

The VWR Incubating Rocker/Wavers are used to prepare samples for testing. These units have been designed for the temperature, speed/tilt, and time functions to work independently of one another. The temperature and speed/tilt can be re-set without re-setting the timer, and the timer can be stopped and started without interrupting the heating or rocking/waving functions.

1. Getting ready:

- a. Plug the power cord into a properly grounded outlet. The standby indicator light will illuminate, verifying power to the unit. Manually move the platform tray to a near horizontal "home" position before powering unit.
- b. Press the standby button to move the unit from standby mode. The standby indicator light will shut off and the temperature, speed/tilt and time displays will illuminate and display the previously used settings.

2. Set temperature:

- a. Press the up/down arrows below the temperature display until you reach the desired temperature. When you release the arrow button, the display will blink off and then on indicating the new set temperature has been accepted.
- b. Press the on/off button to start the heating function. The indicator light below the temperature display will illuminate to indicate the heating function is in use. You will hear five (5) audible beeps that indicate set temperature has been reached.
- c. Set-point temperature adjustments can be made without interrupting heating using the up/down arrows below the temperature display. After the change has been made and you release the button, the display will blink off and then on indicating the new set temperature has been accepted.
- d. To stop heating, press the on/off button below the temperature display, the heat indicator light will turn off.

Caution Hot Indicator:

The caution hot indicator light warns that the temperature of the air in the chamber is above 40°C. The light will illuminate and remain on when the temperature of the air in the chamber reaches approximately 40°C. When the heat is turned off, the caution hot indicator light will stay illuminated until the temperature of the air in the chamber is less than 40°C.

3. Setting tilt: Electronic tilt enables the user to adjust the angle of tray movement while the unit is rocking/waving or while the unit is stopped.

- a. Press the function button below the speed/tilt display until the tilt indicator light illuminates. You are now ready to set the tilt angle.
- b. Press the up/down arrows below the speed/tilt display until you reach the desired angle. When you release the arrow button, the display will blink off and then on indicating the new set angle has been accepted. The unit will complete one rotation at the previously set angle, then smoothly change to the newly programmed angle. The tilt indicator light will flash until the transition to the new tilt angle is complete.

4. Setting speed:

- a. Press the function button below the speed/tilt display until the speed indicator light illuminates. You are now ready to set the speed.
- b. Press the up/down arrows below the speed/tilt display until you reach the desired speed. When you release the arrow button, the display will blink off and then on indicating the new set speed has been accepted.
- c. Press the on/off button to start the rocking/waving function. The indicator light below the speed/tilt display will flash until the transition to the set speed is complete, then the light will stay illuminated indicating the rocking/waving function is in use.

Operating Tips

When the unit is running at slow speeds or high angles, making large changes to the tilt angle or speed may take several minutes to complete. The quickest way to make large changes to the speed or tilt angle is to stop the unit, change the speed or tilt angle, then restart by pressing the on/off button below the speed/tilt display.

5. Setting time to zero (0:00) and continuous mode: Accumulated time.

- a. Press and hold the on/off button below the time display. After three (3) seconds, the display will indicate the previous set time.
- b. Simultaneously press both the up and down arrows; the display will indicate zero (0:00). The unit time is now set to zero (0:00) minutes. Alternately, you can use the up/down arrows to get to zero (0:00).
- c. Press the on/off button below the time display. The display will indicate the accumulated time. The up/down arrows will become inactive. To stop timer, press the on/off button again. **IMPORTANT:** This will **NOT** interrupt the rocking/waving function. Press the on/off button below the speed/tilt display to interrupt the rocking/waving function.
- d. To re-set, press and hold the on/off button below the time display. After three (3) seconds, the display will indicate the previous set time, which was zero (0:00).

6. Setting timed mode: Programmed time.

- a. Press the up/down arrows below the time display until you reach the desired time.
- b. Start this function by pressing the on/off button below the time display. The unit will run for the selected time, the up/down arrows will become inactive while the timer is running. The unit will stop rocking/waving when time display reaches zero (0:00). Four (4) audible beeps will indicate the time down function is complete. The time display will default back to the set time. To repeat for the same time, simply depress the on/off button again.
- c. To interrupt an automatic timing cycle before it is completed, press the on/off button below the time display. The time display will flash off and on to indicate the time function is on "hold". **IMPORTANT:** This will **NOT** interrupt the rocking/waving function. Press the on/off button below the speed/tilt display to interrupt the rocking/waving function. Restart the timer by pressing the on/off button below the time display, unit will continue counting down to zero (0:00). When the display reaches zero (0:00), you will hear the four (4) audible beeps that indicate the time down function is complete and the rocking/waving function will cease.

7. Turning unit off:

- a. To turn unit off, press the standby button. The temperature, speed/tilt and time displays will be blank, the standby indicator light will illuminate. The Incubating Rocker/Waver should be kept in standby mode when not in use. To completely cut off power to the unit, disconnect the power cord from the unit or un-plug from the wall outlet.

Operating Tips

Built-in memory maintains the last used temperature, speed/tilt, and time settings during a power interruption. A built-in program will shut power off to the motor if the platform tray is prevented from rocking/waving, or the unit is overloaded beyond its recommended weight capacity.

Troubleshooting

Review the information in the table below to troubleshoot operating problems. During operation, any rattling or ticking sounds may indicate a loose screw on the platform tray, tray attachment or accessory. All accessories should be sufficiently tightened in place before starting the unit.

Problem	Cause	Solution
E1	RTD open or temperature over 100°C (<i>Incubating Units</i>)	This error cannot be fixed by the end user. Unplug the unit and please contact your VWR representative for repair.
E2	RTD shorted or temperature below 0°C (<i>Incubating Units</i>)	This error cannot be fixed by the end user. Unplug the unit and please contact your VWR representative for repair.
E3	Mechanical obstruction Ceased motor Drive belt broken	Press the standby button to clear this error. If the E3 error persists, along with grinding, knocking, or rubbing noises, un-plug the unit and contact your VWR representative for repairs. In the event of an errant E3 error, such as someone inadvertently touching the platform tray while the unit is running, the unit will automatically go to the horizontal "home" position and restart. In the event of an obstruction left under the tray, which would cause the unit to continually restart, the unit will attempt to auto home and restart four (4) times and then stop running, displaying an E3 error. This error can be cleared by the user by pressing the standby button
E4	Maximum load exceeded	Press the standby button to clear this error. Be sure the load is within the maximum load specification before restarting the unit. If problem persists, unplug the unit and please contact your VWR representative for repair.

Accessories

Dimpled Mat

Designed to hold centrifuge tubes, vials, culture tubes, and micro-tubes securely in place. Mat can be easily removed for cleaning and transporting tubes from bench to tray.


Description	Dimensions	Cat. No.
Dimpled Mat for Rocking Platform Shaker	32,4 x 25,4 cm	444-0768
Dimpled Mat for 3-D Rotator Waver	29,9 x 22,2 cm	444-0770

Stacking Tray (*not for use with Incubating Rockers or Wavers*)

Easily attaches to the lower platform tray to add a second tier for higher capacity applications. Second tier tray mounts 8,9 cm above the lower tray.

Description	Cat. No.
Stacking Tray for Rocking Platform Shaker	444-0772
Stacking Tray for 3-D Rotator Waver	444-0773

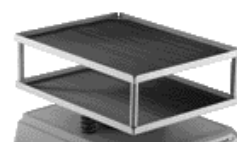Technical service

Web Resources

Visit the VWR website at www.vwr.com for:

- Complete technical service contact information
- Access to the VWR Online Catalogue, and information about accessories and related products
- Additional product information and special offers

Contact us For information or technical assistance contact your local VWR representative or visit. www.vwr.com.

Warranty


VWR warrants that this product will be free from defects in material and workmanship for a period of five (5) years from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non-performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

Compliance with local laws and regulations

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorisations necessary to run or use the Product in its local environment. VWR will not be held liable for any related omission or for not obtaining the required approval or authorisation, unless any refusal is due to a defect of the product.

Equipment disposal

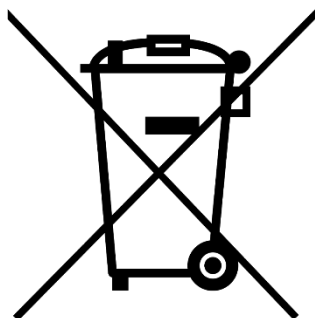This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste.

Instead it's your responsibility to correctly dispose of your equipment at lifecycle -end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

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India

VWR Lab Products Private Limited
No.139. BDA Industrial Suburb,
6th Main, Tumkur Road, Peenya Post,
Bangalore, India – 560058
Tel.: +91 80 28078400
Email: vwr_india@vwr.com

Ireland

VWR International Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: +353 01 88 22 222
Email: sales.ie@vwr.com

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