
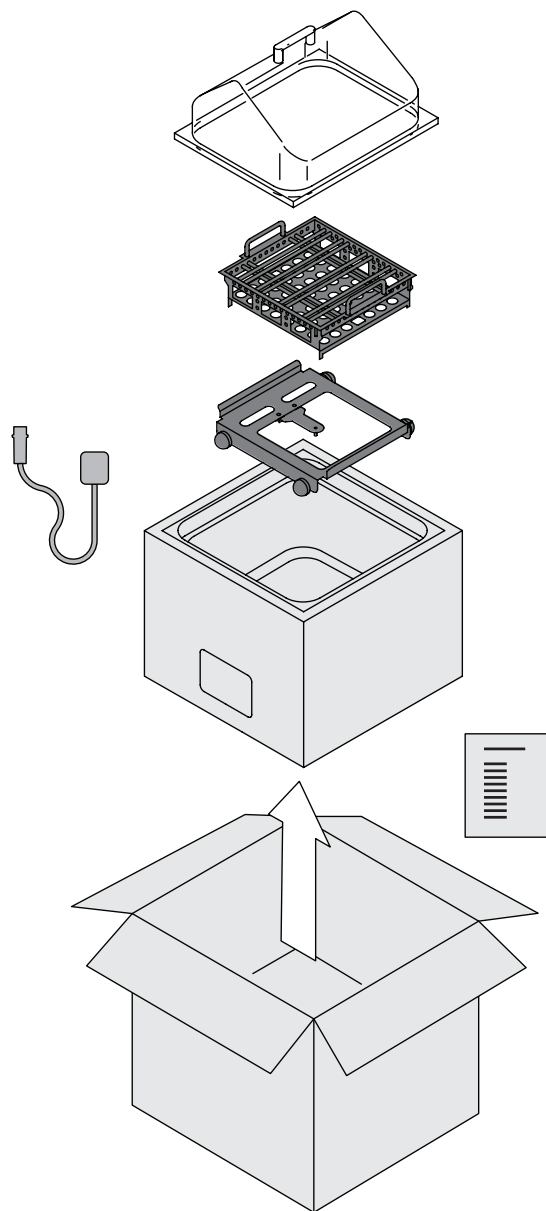
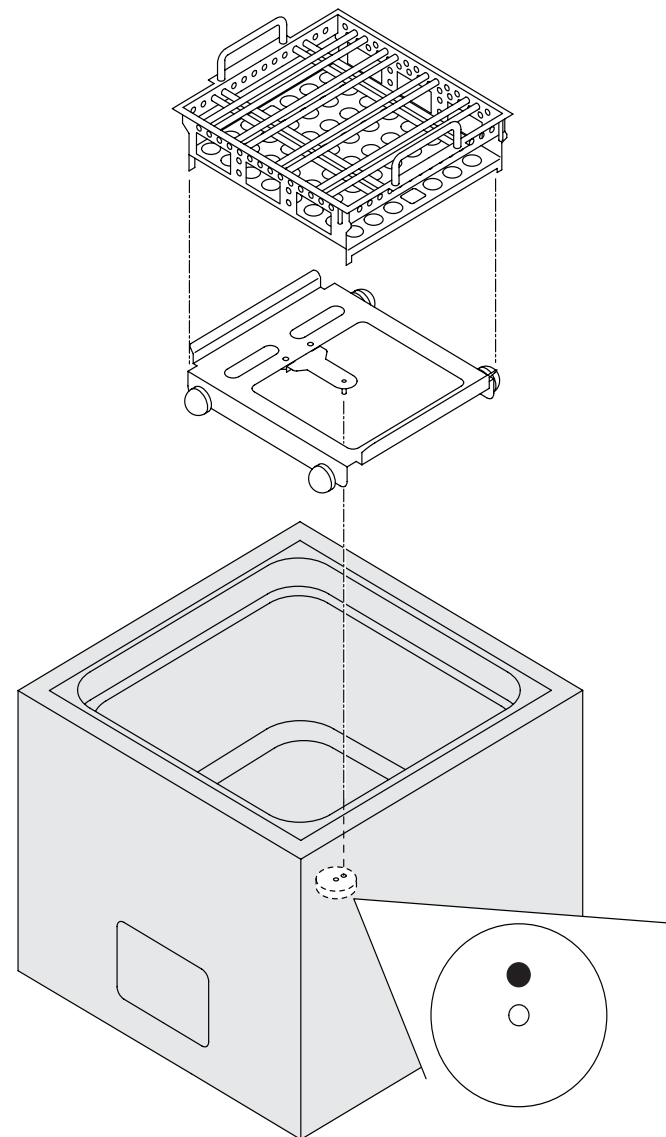
1


2

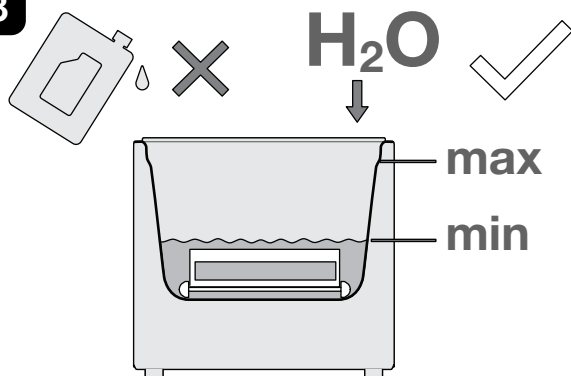
3

4

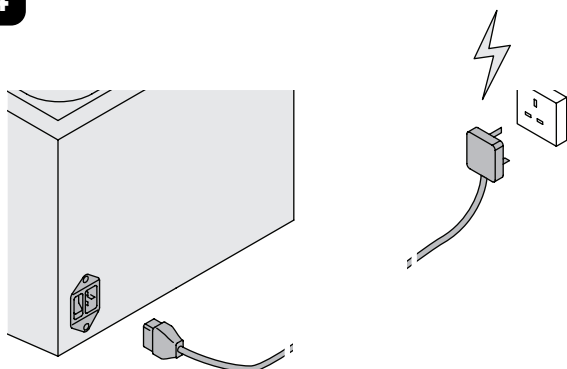
5

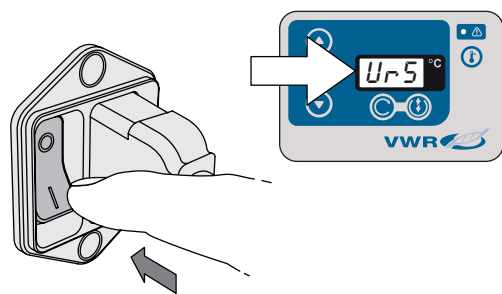
6

°C

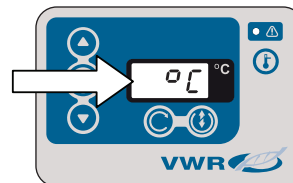
SET


SET

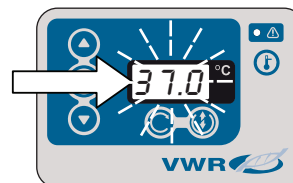SET



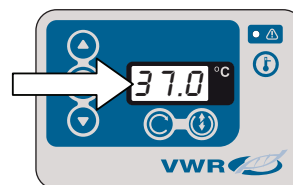
°C

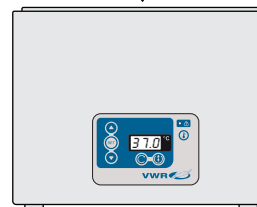
7

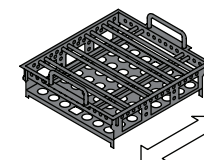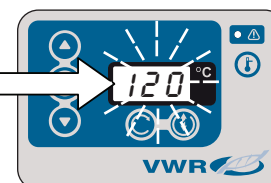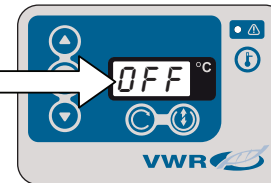
rpm

stop



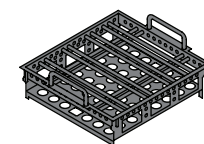
start
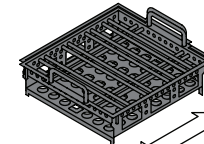
>60°C

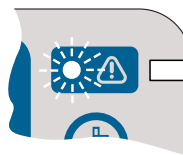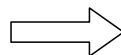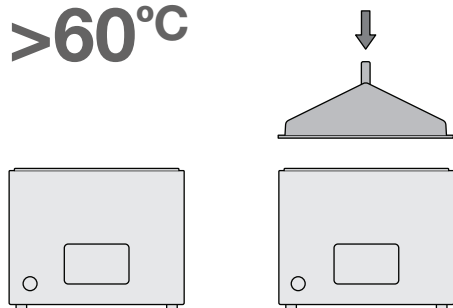Specification

Specification

Mains supply: 220 – 240V @ 50/60Hz

Earth connection required

Ambient operating temp: + 5°C to 40°C

Ambient storage temp: – 20°C to + 60°C

Current Consumption





Model	Size	Current
462 – 0493	12 L	3.6A
462 – 0494	18 L	6.2A

Legal Address of Manufacturer Europe

VWR International bvba
Researchpark Haasrode 2020
Geldenaaksebaan 464
B-3001 Leuven
+ 32 16 385011
<http://be.vwr.com>

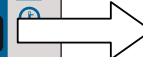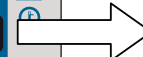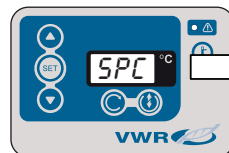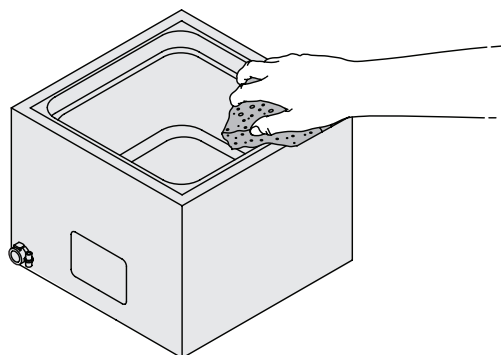H₂O ✓

Service

If service is required, switch off the unit and contact your local VWR representative or visit vwr.com.

Online resources
www.vwr.com

Safety information

EN Original instructions

FR Notice originale

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This product is a general purpose thermostatically controlled water bath designed for indoor laboratory use by or supervised by a professional user.

Read the whole of the instructions. Safety may be impaired if they are not followed.

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Surfaces and water can be hot during and after use.

This bath is only intended for use with water. Use of other fluids may invalidate warranty and present a risk of fire or explosion.

Place on a stable flat surface to reduce the risk of accidental spillage.

No user serviceable parts. Risk of electric shock after disassembly or operation with covers removed.

Not for use in environments with a risk of flammable or explosive gases.

To be operated within the limits listed below.

Only use the mains cord provided. Ensure that the mains plug and the switch are easily accessible.

Use only with a mains supply which includes a ground connection.

A clearance of >10cm around the bath is required to ensure adequate air flow.

Take care when operating after a power interruption. Shaking will restart at the previously set speed.

Take care when operating with vessels in the tray. Accidental water spillage can occur when shaking is started. Reduce the water level as required.

If a potentially hazardous liquid is spilt onto the equipment, disconnect it from the power supply and have it checked by a competent person. It is the user's responsibility to carry out appropriate decontamination if hazardous material is spilt on the equipment.

There are accessible moving parts in the bath. Observe safe working practises when operating a shaking bath. Users should not interfere with moving parts when in use and prevent accidental contact from items such as fingers, hair or clothing.

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