

VWR® Gel Documentation system Basic Image Capture - Quick Guide

European Catalogue number: 1730-1470

1

Turn on

Select



Turn on the camera (ensure the battery is charged)



The Canon control software launches automatically when you turn on the camera. If the software does not start automatically click on Start/All Programmes/Canonutilities/EOS Utility/EOS Utility

2

Capturing an image

Select

Live View shoot. ...



Select Live View shoot to open live window




Ensure the system has been installed correctly, as per the Installation Quick Guide

3

Settings

Select

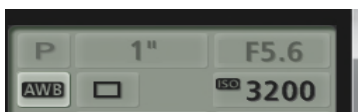

Select




The AWB button and from the White Balance menu select Cloudy from the menu

4

Focus setting

Select

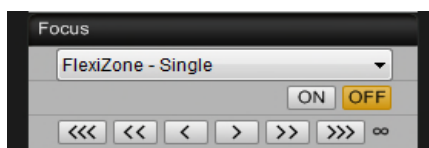

From the Remote Live View window

To adjust zoom turn the lens of the camera



To adjust focus use the arrows

5

Grab image

Press



From the EOS 4000D window

The camera will set an exposure time relevant to the light levels coming from the sample



Then turn off the camera

6

Save and export

- ➔ Images are saved as JPEG files
- ➔ Export a TIFF or BMP

CONTACT VWR:
For further details , please consult the camera Manual
Visit www.vwr.com to find your local VWR contact



BIC1016