Multilanguage short instruction manual & Troubleshooting

VWR[®] DISSOLEVD OXYGEN METER DO-230

EU Catalogue Number: 664-0236

Version: 1.0
Issued: 01, December, 2020

LANGUAGES


English.....	3
Français.....	5
Deutsch.....	7
Italiano.....	9
Español.....	11
Português	13
Hrvatski.....	15
Český.....	17
Dansk.....	19
Nederlandse.....	21
Suomen.....	23
Magyar.....	25
Latviešu.....	27
Lietuvių.....	29
Norsk.....	31
Polski.....	33
Română.....	35
Slovák	37
Slovenski.....	39
Svenska.....	41

FOR MORE DETAILS, PLEASE CONSULT THE FULL MANUAL

English


The DO-230 oxygen meter allows you to measure oxygen in water in % or mg / l, oxygen in air in % and temperature. Turn the device on with the  key. Other operating activities are performed using the touch screen.

Sockets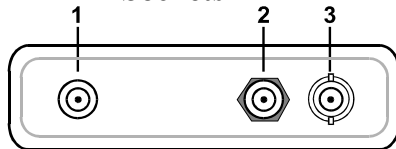


1	Power adapter socket
2	Temperature sensor socket
3	Oxygen sensor socket

Measurement screen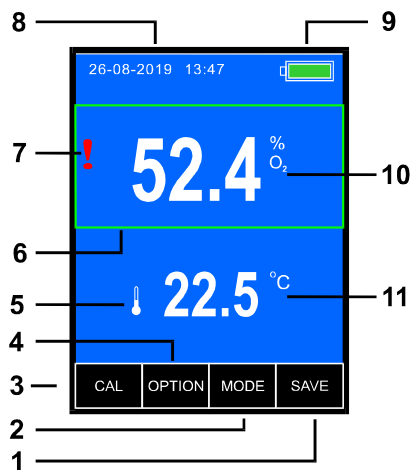


1	Collect results key
2	Parameters selection key
3	Key to enter the calibration mode (hold down for 2s)
4	Key for selecting the measurement function
5	Symbol of the connected temperature sensor
6	Fixed measurement frame
7	Oxygen sensor warning symbol
8	Date and time
9	Battery status
10	Oxygen measurement result with unit
11	Temperature measurement result with unit

Parameter screen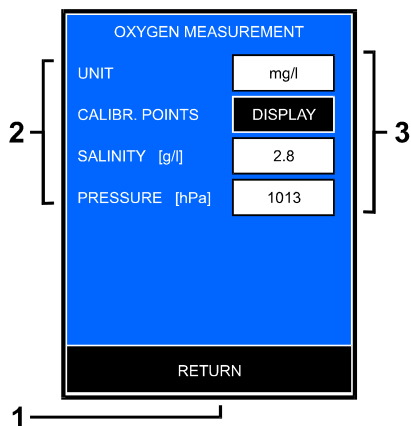

1	Key to return to the previous screen
2	Description of parameters
3	Parameter selection keys

Calibration screen

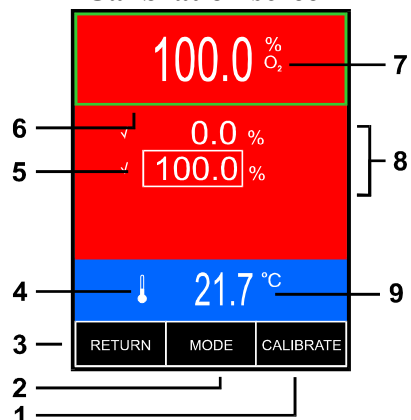1	Calibration result confirmation key
2	Parameter selection key
3	Key to return to the previous screen
4	Symbol of the connected temperature sensor
5	Confirmation of calibration in point
6	Fixed measurement frame
7	Oxygen measurement result in calibration mode
8	Calibration points
9	Temperature measurement result with unit

TROUBLESHOOTING

Problem	Cause	Solution
Meter not switching on / No power	Discharged batteries	Charge batteries
	Incorrect battery polarity	Check the polarity
	Battery failure	Replace batteries with AA (R6) Rechargeable type
	Power adapter failure	Check with other 5V USB power adapter (for mobile)
Unstable reading, lack of reading	Poor connection	Check sensor connection to the unit
	Air bubble in the membrane cap	Check for air bubbles, refill electrolyte or replace the membrane
	DO sensor membrane contaminated / broken	Clean the sensor according to its manual or replace the membrane
	Anode or cathode is contaminated	Clean zinc anode with sandpaper. Clean silver cathode, on top, only with smooth cloth!
	Break in cable, connector or internally in sensor	Check cable and connector, replace the sensor
	DO sensor defective	Replace sensor
Meter is not calibrating	Membrane polluted	Clean or replace the membrane
	Membrane defective - broken	Replace membrane
	Electrolyte used up	Refill electrolyte
	Calibration mode not entered properly	Press and hold the CAL button till displaying Cal symbol
	Intermittent or no connection of DO sensor	Check sensor connection to the unit
	Unable to calibrate to 100%	Clean, refill or replace the membrane
	Stabilisation time not long enough	Wait longer for result stabilisation
	DO sensor defective	Replace sensor
Result different than expected	DO sensor not calibrated	Calibrate the sensor
	Membrane polluted	Clean or replace the membrane
	Electrolyte used up	Refill electrolyte
	Calibration not correct	Make 2 points calibration
	Reading stabilisation time too short	Longer wait for result stabilisation
	0% DO concentration solution used up	Add sodium sulphite or prepare new solution
	Wrong salinity value introduced	Delete or change introduced salinity value
	Wrong air pressure value introduced	Delete or change introduced air pressure value
	Stabilisation time not long enough	Wait longer for result stabilisation
Result drifting down	DO sensor defective	Replace DO sensor
	Air consumption by sensor	Use stirrer or slowly move the sensor in the vessel
Result drifting up	To intensive stirring, sensor close to air source	Slow down stirring, place the sensor in a different place
Temp. value incorrect or  symbol displayed	Temperature sensor not connected or poor connection	Connect the temp. sensor / check the connection
	Breakage in temp sensor cable or connector / defective sensor	Check connection / replace sensor
Information: oxygen sensor calibration readout error!	DO sensor calibration data loss	Recalibrate the DO sensor
	If the situation appears after each switching on	Contact distributor or service agent
Information: Manufacturer's settings readout error!	Internal hardware or software problem / Losing factory settings	Contact distributor or service agent

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