Instruction manual

VWR® Fixed Speed, Analogue, Digital, & Pulsing Vortex Mixer

EU cat. no

Fixed Speed Vortex Mixers

Euro Plug And Swiss Adapter: 444-0994

Uk Plug: 444-0995

Analogue Vortex Mixers

Euro Plug And Swiss Adapter: 444-0996

Uk Plug: 444-0997

Digital Vortex Mixers

Euro Plug And Swiss Adapter: 444-0998

Uk Plug: 444-0999

Pulsing Vortex Mixers

Euro Plug And Swiss Adapter: 444-1041

Uk Plug: 444-1042

EN	- English		1
FR	- Français		16
ES	- Español		32
DE	- Deutsch		48
IT	- Italiano		63
PT	- Português		80
BG	- Български		96
CZ	- Čeština		99
DK	- Dansk		102
EE	- Eesti		105
FI	- Suomi		108
GR	- Ελληνικά		111
HR	- Hrvatska		114
HU	- Magyar		117
IS	- Íslenska		120
LT	- Lietuvių		123
LV	- Latviešu		126
NL	- Nederlands		129
NO	- Norsk		132
PL	- Polski		135
RO	- Română		138
SI	- Slovenščina		141
SK	- Slovenčina		144
SV	- Svenska		147

Legal address of Manufacturer

Europe

VWR International bv,
Researchpark Haasrode 2020
Geldenaaksebaan 464 B-3001
Leuven
+3216385011
be.vwr.com

UK Importer:

VWR International Ltd
Hunter Boulevard, Magna Park
Lutterworth, Leicestershire, LE17 4XN
uk.vwr.com

Country of Origin

United States

Table of Contents

Introduction	4
Safety Information	4
Package Contents	4
Installation	4
Intended Use	5
Maintenance & Servicing	5
Environmental Conditions	5
Product Specifications	
<i>Fixed Speed Vortex Mixer</i>	6
<i>Analogue Vortex Mixer</i>	6
<i>Digital Vortex Mixer</i>	6
<i>Pulsing Vortex Mixer</i>	6
Digital and Pulsing Control Panel	7
Vortex Mixer Operating Instructions	
<i>Fixed Speed Vortex Mixer</i>	8
<i>Analogue Vortex Mixer</i>	8
<i>Digital & Pulsing Vortex Mixer</i>	9
Head/Accessory Installation Instructions	
<i>Installing the cup head, 7,6cm head or other head attachments</i>	10
<i>Installing Single Tube Holder</i>	11
Installing Vessel Harness	10
Troubleshooting	12
Technical Service	12
Warranty	12
Compliance with Local Laws and Regulations	12
Disposal	13
Accessories	14
Distributor Contact Information	15

Introduction

This manual contains installation operation and maintenance instructions. Please read the manual completely before using.

Safety Instructions

Please read the entire instruction manual before operating the Vortex Mixer.

WARNING! DO NOT use the Vortex Mixer in a hazardous atmosphere or with hazardous materials for which the unit was not designed. Also, the user should be aware that the protection provided by the equipment may be impaired if used with accessories not provided or recommended by the manufacturer or used in a manner not specified by the manufacturer.

Always operate unit on a level surface for best performance and maximum safety.

DO NOT lift the Vortex Mixer by the head. All heads, including cup head, are removable. They can pop off if lifted by them.

CAUTION! To avoid electrical shock, completely cut off power to the unit by disconnecting the power cord from the wall outlet. Disconnect unit from the power supply prior to maintenance and servicing.

Spills should be removed promptly. **DO NOT** immerse the unit for cleaning.

DO NOT operate the unit if it shows signs of electrical or mechanical damage.

Earth Ground - Protective Conductor Terminal

Alternating Current

Package Contents

Fixed Speed, Analog, Digital or Pulsing Vortex Mixer

Description	Quantity
183cm power cord with Euro plug and UK plug or Swiss adapter	1
cup head*	1
7,6cm head*	1
7,6cm rubber head cover*	1

* These accessories come with all units except the Pulsing Vortex Mixer which is also supplied with a 1,5 to 2mL Micro-Tube Holder with built-in cup head.

Installation

Upon receiving the VWR Vortex Mixer, check to ensure that no damage has occurred during shipment. It is important that any damage that occurred in transport is detected at the time of unpacking. If you do find such damage the carrier must be notified immediately.

After unpacking, place the Vortex Mixer on a level bench or table, away from explosive vapors. Ensure that the surface on which the unit is placed will withstand typical heat produced by the unit. Always place the unit on a sturdy work surface.

The Vortex Mixer is supplied with a power cord that should be plugged into a properly grounded outlet. If the supplied power cord does not meet your needs, please use an approved power cord that suits local codes and electric supply. Replacement of the plug must be made by a qualified electrician.

Intended Use

The VWR Vortex Mixers are intended for general laboratory use. Safety cannot be guaranteed if used outside of the intended use.

Maintenance & Servicing

The Vortex Mixer is built for long, trouble-free, dependable service. No lubrication or other technical user maintenance is required.

The unit should be given the care normally required for any electrical appliance. Avoid wetting or unnecessary exposure to fumes. Spills should be removed promptly. **DO NOT** use a cleaning agent or solvent on the front panel which is abrasive or harmful to plastics, nor one which is flammable. Always ensure the power is disconnected from the unit prior to any cleaning. If the unit ever requires service, contact your VWR International representative.

Environmental Conditions

Operating Conditions: Indoor use only.

Temperature: 4 to 40°C

Humidity: 20% to 85% relative humidity, non-condensing

Altitude: 0 to 2000 M above sea level

Non-Operating Storage:

Temperature: -20 to 65°C

Humidity: 20% to 85% relative humidity, non-condensing

Installation Category II and Pollution Degree 2 in accordance with IEC 664.

Product Specifications

Fixed Speed Vortex Mixer

Overall dimensions (L x W x H):	17,8 x 11,2 x 13,5 cm
Electrical (50/60 Hz):	0,6 amps, 135 watts
Speed range:	2500rpm
Orbit:	4,9mm
Duty rating:	intermittent duty
Controls:	none
Ship weight:	4,5kg

Digital Vortex Mixer

Overall dimensions (L x W x H):	17,8 x 11,2 x 13,5 cm
Electrical (50/60 Hz):	0,6 amps, 135 watts
Speed range:	500 to 2500rpm
Timer:	1 second to 9999 minutes
Orbit:	4,9mm
Duty rating:	intermittent duty
Controls:	see page 7
Ship weight:	4,5kg

Analog Vortex Mixer


Overall dimensions (L x W x H):	17,8 x 11,2 x 13,5 cm
Electrical (50/60 Hz):	0,6 amps, 135 watts
Speed range*:	300 to 2500rpm
Orbit:	4,9mm
Duty rating:	intermittent duty
Controls:	auto/off/on rocker switch, speed knob: variable 1 to 10 dial marks
Ship weight:	4,5kg

Pulsing Vortex Mixer

Overall dimensions (L x W x H):	17,8 x 11,2 x 13,5 cm
Electrical (50/60 Hz):	0,6 amps, 135 watts
Speed range*:	500 to 2500rpm
Timer:	1 second to 9999 minutes
Orbit:	2,5mm
Duty rating:	intermittent duty
Controls:	see page 7
Ship weight:	4,5kg

** Maximum speed will vary depending on accessory being used.*

Digital & Pulsing Vortex Mixer Control Panel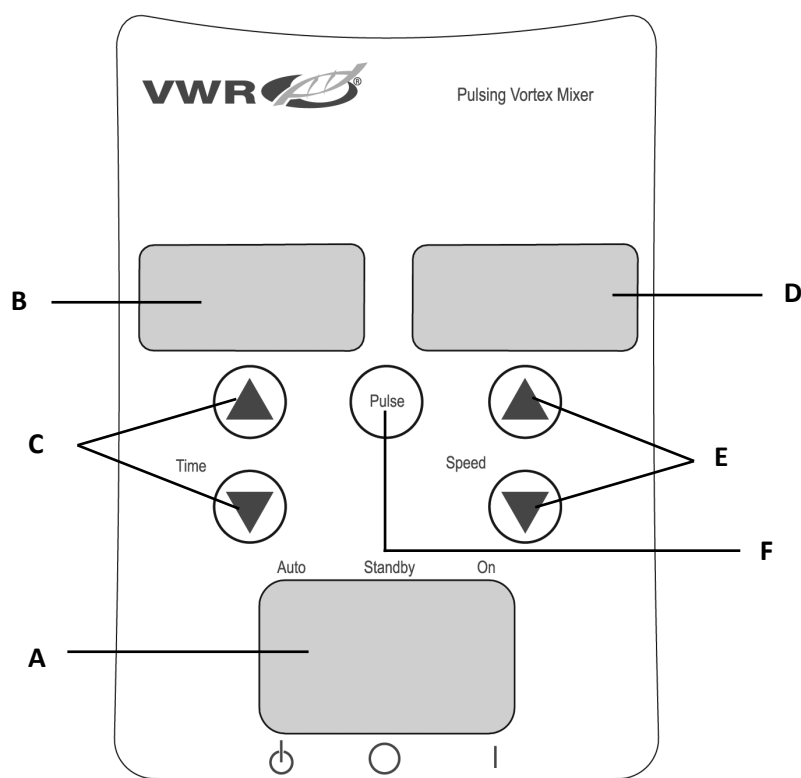

**Same controls for the Digital Vortex Mixer but without the 'pulse' button.*

The front panel of the Digital/Pulsing Vortex Mixer contains all the switches, controls and displays needed to operate the unit.

A. 3-way rocker switch: Auto/standby/on rocker switch starts/stops the vortexing function.

B. Time display: Displays accumulated time (continuous mode) or how much time is remaining (timed mode). The display range is from 0 to 9,999 minutes in one (1) second increments. The display will indicate minutes and seconds until the time reaches 99 minutes and 59 seconds (99:59), then the display will automatically display minutes up to 9,999.

C. Up/down arrows for set-point control.

D. Speed display: Displays the speed of the Vortex Mixer.

E. Up/down arrows for set-point control.

F. Pulse button: Starts the pulsing mode.

Vortex Mixer Introduction

Your Vortex Mixer is ready for most one-handed applications. The Vortex Mixer operates by mixing tubes just prior to testing.

Operating Tips

These units are not intended for continuous use over extended periods of time. A built-in temperature sensor will prevent the unit from overheating by shutting the motor down. If this occurs, allow the motor to cool down and the unit will be able to resume normal operation. Decrease the duration of the time unit is used between rest periods.

See page 10-11 for 'Head/Accessory Installation Instructions'.

Fixed Speed Vortex Mixer Operating Instructions

The Fixed Speed Vortex Mixer runs only at 2500rpm.

1. Plug the power cord into a properly grounded outlet. The unit is now on and in a standby mode.
2. The vessel you are using must be pushed down on the head attachment to achieve motion. To get the desired mix, vary the angle of contact and pressure against the head.
3. The mixer will return to standby mode when the vessel is lifted off the head.
4. The unit can become warm to the touch with constant use. There is nothing wrong with it. At high performance settings the motor (UL recognized component) simply puts out more heat. **IMPORTANT: DO NOT** use large accessories on the Fixed Speed Vortex Mixer. This unit is designed for use with cup head or 7,6cm head with 7,6cm rubber head cover only.
5. To completely cut off power to the unit, disconnect the power cord from the wall outlet.

Analog Vortex Mixer Operating Instructions

1. Make sure the 3-way rocker switch is in the center, off position. Plug the power cord into a properly grounded outlet.
2. For continuous operation, push the rocker switch to the right, on position. For intermittent/touch operation, push the rocker switch to the left, auto position.
3. In either case, turn the speed knob to 1. In the on position, you will see the head in motion immediately. In the auto position, the vessel you are using must be pushed down on the head attachment to achieve motion. To get the desired mix, vary the speed, using the speed knob, and/or vary the angle of contact and pressure against the head.
4. When finished with either continuous or intermittent operation, return rocker switch to the center, off position.
5. The continuous, on mode is for the larger accessories. In continuous on mode, **DO NOT** run the unit above 900rpm (analog speed setting of approximately 5).
6. The intermittent, auto mode is intended for short mixing times of one (1) minute or less at full speed. A rest period of two (2) minutes is recommended to minimize the possibility of overheating the motor. The unit can become warm to the touch with constant use. There is nothing wrong with it. At high performance settings the motor (UL recognized component) simply puts out more heat. **IMPORTANT:** When using larger accessories (Insert Retainer with foam inserts or Ampule Tube Holders), **DO NOT** run the unit above 900rpm, (analog speed setting of approximately 5). **UNIT DAMAGE WILL OCCUR.**
7. To completely cut off power to the unit, disconnect the power cord from the wall outlet.

NOTE: Actual speed is subject to operating conditions and application. When selecting the ideal analog speed setting for your application, please evaluate loading and duty cycle as these may impact the unit's actual motor speed.

Digital & Pulsing Vortex Mixer Operating Instructions

1. Make sure the 3-way rocker switch is in the center, standby position. Plug the power cord into a properly grounded outlet. The unit is now on and in a standby mode.
2. For continuous operation, push the rocker switch to the right, on position. For intermittent/touch operation, push the rocker switch to the left, auto position.
3. In either case, set the speed. In the on position, you will see the head in motion immediately. In the auto position, the vessel you are using must be pushed down on the head attachment to achieve motion. To get the desired mix, vary the speed, using the speed up/down arrows, and/or vary the angle of contact and pressure against the head.

Setting speed:

To set the speed press the up/down arrows below the speed display until you reach the desired speed.

- Speed while running in the on position is 500 to 2500rpm.
- Speed while running in the auto position is 1000 to 2500rpm.
- Speed while running in the pulse mode is 2500rpm, running for 30 seconds, resting 10 seconds.

Setting time:

Time operation works in both auto and on positions.

- To run in timed mode (programmed time), start by pressing the rocker switch to the left, auto position. Press the up/down arrows below the time display until you reach the desired time. While running unit in timed mode, the time display will show time remaining, counting down to zero (0:00).
- To run in continuous mode (accumulated time), reset the timer to zero (0:00) before running. Time display will show the accumulated time and run until you press the rocker switch to the center, standby position.

Pulse operation: (for Pulsing Vortex Mixer only)


- Set the time if desired.
 - Move the rocker switch to the right, on position. Press the pulse button, in the center of the front panel. The speed display will show "PULSE". The unit will run at 2500rpm, cycling a run time of 30 seconds, followed by a resting time of 10 seconds. The unit will continue vortexing until the timer reaches zero (0:00) or you stop the unit by moving the rocker switch to the center, standby position.
4. When finished with either continuous, intermittent or pulse operation, return rocker switch to the center, standby position.
 5. The continuous, on mode is for the larger accessories. In continuous on mode, **DO NOT** run the unit above 900rpm.
 6. The intermittent, auto mode is intended for short mixing times of one (1) minute or less at full speed. A rest period of two (2) minutes is recommended to minimize the possibility of overheating the motor. The unit can become warm to the touch with constant use. There is nothing wrong with it. At high performance settings the motor (UL recognized component) simply puts out more heat. **IMPORTANT:** When using larger accessories (Insert Retainer with foam inserts or Ampule Tube Holders), **DO NOT** run the unit above 900rpm. **UNIT DAMAGE WILL OCCUR.**
 7. To completely cut off power to the unit, disconnect the power cord from the wall outlet.

NOTE: In the event of a power outage or hazard that will interrupt unit operation, the unit will restart at set speed and/or temperature when a power outage or hazard is restored.

Head/Accessory Installation Instructions

Installing the cup head, 7,6cm head or other head attachments:

1. Move the 3-way rocker switch to the center, off/standby position. For the Fixed Speed Vortex Mixer disconnect the power cord from the wall outlet.
2. If using an Insert Retainer with tube holder, insert foam tube holder into Insert Retainer by lining up the recessed notches of the foam with the fingers of the retainer. Push foam to the bottom of the retainer.
3. Place head on post of the unit, aligning the flat "alignment mark" inside of the head mounting hole with the flat "alignment mark" on the post.
4. Press center of head down firmly until it snaps into place. **DO NOT** use a 7,6cm head without the 7,6cm rubber head cover. Add samples to attachments after the head is secured onto the Vortex Mixer. Be sure samples are secure before turning unit on.
5. To remove an installed head, hold the Vortex Mixer housing with one hand and firmly pull the head straight up. Larger heads may require using both hands to force head off post. Remove samples before removing attachments.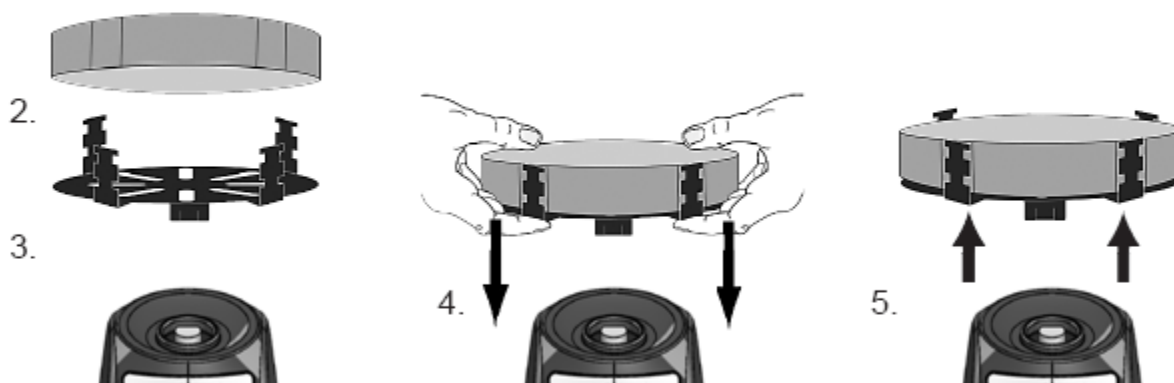


Installing Vessel Harness:

1. Install Insert Retainer without foam tube holder onto Vortex Mixer.
2. Stretch vessel harness elastic bands over opposite fingers of the Insert Retainer. Place bands on either upper or lower notches of the retainer.
3. Place sample in center of the harness bands. Be sure sample is secure before turning unit on.

Head/Accessory Installation Instructions

Installing Single Tube Holder:

1. Move the 3-way rocker switch to the center, off/standby position. Single Tube Holder should **NOT** be used on the Fixed Speed Vortex Mixer.
2. Remove cup head, or any other accessory from the Vortex Mixer.
3. Remove the adhesive backing from adapter plate and line up the opening of the adapter plate over the opening of the housing. Press down firmly to ensure a proper bond between the adaptor and the unit. Align the base plate of the Single Tube Holder with the adaptor plate from the previous step. Keep the center openings of the plates concentric to avoid mechanical interference. The magnets will hold the Single Tube Holder in place. To realign this attachment, pull up from the handle until the magnets release, check the alignment and place again.
4. Align the base plate of the Single Tube Holder over the post of the Vortex Mixer with the two Single Tube Holder alignment posts at the back of the unit. Lower Single Tube Holder onto the adapter plate, the magnets will hold the Single Tube Holder in place. To realign this attachment, pull up from the handle until the magnets release, check the alignment and place again.
5. Now that the Single Tube Holder is in place, attach the cup head on the post of the unit, aligning the orientation arrow towards the rear of the unit.
6. Lift up on the spring-loaded upper portion of the Single Tube Holder and slide a 6,4 to 11,4cm long tube on the cup head so that the tube is aligned in the middle. The 0,5mL Micro-Tube Holder or the 1,5 to 2mL Micro-Tube Holder will also fit into the Single Tube Holder. You are now ready to mix your sample.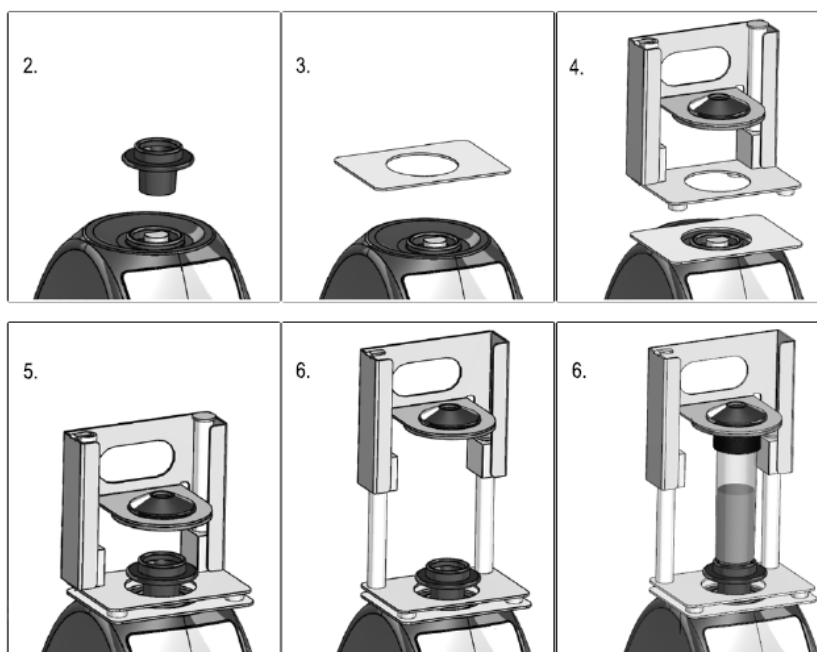

Troubleshooting

Review the information in the table below to troubleshoot operating problems.

Problem	Cause	Solution
Unit will not run	Mechanical obstruction or Motor obstruction	Remove obstruction and press down on cup head several times to confirm no obstruction exists. If problem persists, please contact your VWR International representative for repair.
Unit is excessively noisy	Sensor fan misaligned or Motor misaligned	This error cannot be fixed by the end user. Please contact your VWR International representative for repair.

Technical service

Web Resources

Visit the VWR website at www.vwr.com for:

- Complete technical service contact information
- Access to the VWR Online Catalogue, and information about accessories and related products
- Additional product information and special offers

Contact us For information or technical assistance contact your local VWR representative or visit. www.vwr.com.

Warranty


VWR warrants that this product will be free from defects in material and workmanship for a period of five (5) years from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non-performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

Compliance with local laws and regulations

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorizations necessary to run or use the Product in its local environment. VWR will not be held liable for any related omission or for not obtaining the required approval or authorization unless any refusal is due to a defect of the product.

Equipment disposal

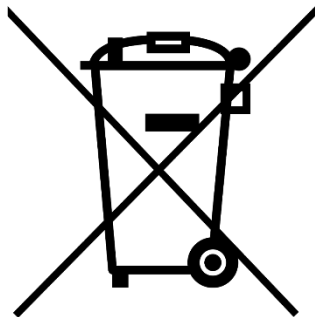This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste.

Instead it's your responsibility to correctly dispose of your equipment at lifecycle -end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

Thank you

Accessories

Micro-Tube Holder - Mixes (48) 0,25 to 2mL microtubes. Requires Insert Retainer.


Description	Part Number
Micro-Tube Holder (2 pk)	444-0218

Insert Retainer	444-7048
-----------------	----------

Micro-Well Plate Holder - Ideal for mixing 96 well plates or deep well blocks. Requires Insert Retainer.

Description	Part Number
Micro-Well Plate Holder (2 pk)	444-0219

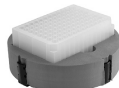Insert Retainer	444-7048
-----------------	----------

9 to 13mm Tube Holder - Ideal for mixing 5mL culture tubes and micro-vials. Requires Insert Retainer.

Description	Part Number
9 to 13mm Tube Holder (2 pk)	444-7051

Insert Retainer	444-7048
-----------------	----------

14 to 19mm Tube Holder - Ideal for mixing up to (8) 15mL centrifuge tubes. Requires Insert Retainer.

Description	Part Number
14 to 19mm Tube Holder (2 pk)	444-7052

Insert Retainer	444-7048
-----------------	----------

Ampule Tube Holder - Mixes up to 4 storage vials and test tubes.

Description	Part Number
15-17mm Ampule Tube Holder	444-0212

10-17mm Ampule Tube Holder	444-0213
----------------------------	----------

Tube Holder Kit

Description	Part Number
Includes: (1) 9-13mm Tube Holder, (1) 14-19mm Tube Holder, (1) 20-25mm Tube Holder, 1 Flat Foam Insert, 2 Vessel Harness, and 1 Insert Retainer	444-7047

Cup Head - Designed for mixing one tube at a time.

Description	Part Number
Cup Head	444-7060

1,5 to 2mL Micro-Tube Holder - Mixes (18) 1,5 to 2mL micro-tubes. For use with Single Tube Holder.

Description	Part Number
--------------------	--------------------

20 to 25mm Tube Holder - Ideal for mixing up to (8) 50mL centrifuge tubes. Requires Insert Retainer.


Description	Part Number
20 to 25mm Tube Holder (2 pk)	444-7053

Insert Retainer	444-7048
-----------------	----------

Flat Foam Insert - Ideal for custom applications. Can be cut or drilled to fit your specifications. Requires Insert Retainer.

Description	Part Number
Flat Foam Insert (2 pk)	444-0211

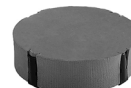Insert Retainer	444-7048
-----------------	----------

Vessel Holder - Mixes Erlenmeyer flasks and media bottles. Requires Insert Retainer.

Description	Part Number
Vessel Harness (2 pk)	444-7050

Insert Retainer	444-7048
-----------------	----------

Stainless Steel Micro-Tube Holder - Mixes up to (12) 1,5 to 2mL micro-tubes. Stainless steel construction.

Description	Part Number
Micro-Tube Holder	444-0214

Single Tube Holder - Single tube, hands free mixing. Easily attaches to the top of any Vortex Mixer with the use of a strong magnetic base. Accepts tubes from 6,4 to 11,4cm in length. Minimum tube diameter of 1,9cm.

Description	Part Number
Single Tube Holder	444-0217

7,6cm Rubber Head Cover & 7,6cm Head - Designed for mixing irregular shaped objects.

Description	Part Number
7,6cm Rubber Head Cover	444-7059

7,6cm Head	444-7058
------------	----------

7,6cm Head with Rubber Cover	444-0208
------------------------------	----------

0,5mL Micro-Tube Holder - Mixes (24) 0,5mL micro-tubes. For use with Single Tube Holder.

Description	Part Number
0,5mL Micro-Tube Holder	444-0215

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India

VWR Lab Products Private Limited
No.139. BDA Industrial Suburb,
6th Main, Tumkur Road, Peenya Post,
Bangalore, India – 560058
Tel.: +91 80 28078400
Email: vwr_india@vwr.com

Ireland

VWR International Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: +353 01 88 22 222
Email: sales.ie@vwr.com

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