European Catalogue Number:
390-0963 INCU-Line Tower

Intended use

VWR INCU-Line Tower is a general purpose convection incubator with two separate chambers, capable of uniformly cooling and heating samples from 10 °C below ambient up to a maximum of 50 °C. The unit has two transparent doors, so samples can be observed without opening the doors, thus preventing temperature changes in the chambers.

Overview

Description of front panel

POWER - power switch: For turning the incubator ON and OFF. It lights when the incubator is switched ON.

▲ - UP key: For increasing the set-temperature.

▼ - DOWN key: For decreasing the set-temperature.

Display: For displaying the actual temperature (dot blinks when you change the set temperature).

Getting Started

Check the data plate for voltage, cycle, wattage and ampere requirements. If matched to your power source, plug the power cord into a grounded outlet. Voltage should not vary more than 10 % from the data plate rating.

In selecting a location, consider all conditions which might affect performance, such as heat from radiators, autoclaves, etc. Avoid direct sun, fast-moving air currents, heating/cooling ducts and high traffic areas. Allow a minimum of 50 mm between the unit and walls or partitions which might obstruct free airflow.

Operation

Do not use this incubator in explosive or flammable environments. Do not heat or incubate flammable, explosive or highly reactive materials in this incubator as serious injury may result.

Place the unit on a level surface and plug the unit into a properly grounded outlet of appropriate voltage.

Install the shelves into both chambers, at the desired height, with the shelf edges up. If you intend to place the samples on the bottom of the chamber, install the shelf so that it sits with the edges down on the chamber bottom, thus creating an air gap between the chamber bottom and your samples.

Place the samples in one or both chambers and close the doors. Make sure that both doors are securely closed.

Turn the power switch to the ON position. The switch should illuminate. On both displays all "888" are displayed at first, then software version "VX.X", and then the current temperatures for both chambers. Incubator starts to cool or to heat individual chambers independently, regarding to the set temperature for each chamber.

Display and keys on the left side (above the wave line) are used to control the upper chamber. Display and keys on the right side (below the wave line) are used to control the lower chamber.

Set desired temperature for each chamber with the corresponding UP or DOWN key. You must press and hold the UP or DOWN key for more than 2 seconds, until the DOT on the DISPLAY starts to blink. When the DOT starts to blink (it blinks for 3 seconds), set the temperature to the desired value, between 10 °C below ambient and up to a maximum of 50 °C. When the desired temperature is set, the display will automatically show the current temperature after 3 seconds (DOT stops to blink).

You can always display the [SET temperature](#) by pressing and holding the relevant UP or DOWN key for more than 2 seconds.

If you want to use only one chamber, you can turn the other chamber OFF by simultaneously pressing (short press) the corresponding UP and DOWN keys. Message "OFF" appears on the display to indicate that the chamber is turned OFF. When you want to use the chamber again, turn it ON by simultaneously pressing (short press) the corresponding UP and DOWN keys. Current temperature appears on display and cooling or heating of the chamber starts.

Temperature readjustment

Perform the following procedure individually for each chamber.

The incubator has software for temperature readjustment. You must first measure the temperature in the middle of the chamber or in the sample for a minimum of 2 hours. You must measure the temperature with a certificated reference temperature measuring instrument with 0.1 °C precision or better. After 2 hours or more read the temperature on a thermometer and compare it with the temperature on the LED display. Difference between thermometer and display is the value that you will enter in the software.

You can insert the thermometer into the upper chamber through the opening on the top of the incubator, and into the lower chamber through the opening on the left side of the incubator.

Example 1: Reference temperature is 37.9 °C, temperature on display is 37 °C.

Difference is 37.9 - 37 = 0.9 Value is 0.9 (value which you enter in the software).

Example 2: Reference temperature is 36.2 °C, temperature on display is 37 °C.

Difference is 36.2 - 37 = -0.8 Value is -0.8 (value which you enter in the software).

Procedure for temperature readjustment:

Press the UP and DOWN keys simultaneously for more than 5 seconds (while current temperature is displayed on display) - message "COR" (correction) appears on display and the DOT starts to blink.

With the relevant UP or DOWN key set the value (while DOT is blinking). Pay attention to the sign of the value (plus or minus). After 3 seconds, when you have set the temperature, display will automatically show the current temperature.

For more details, please consult the full manual.

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