
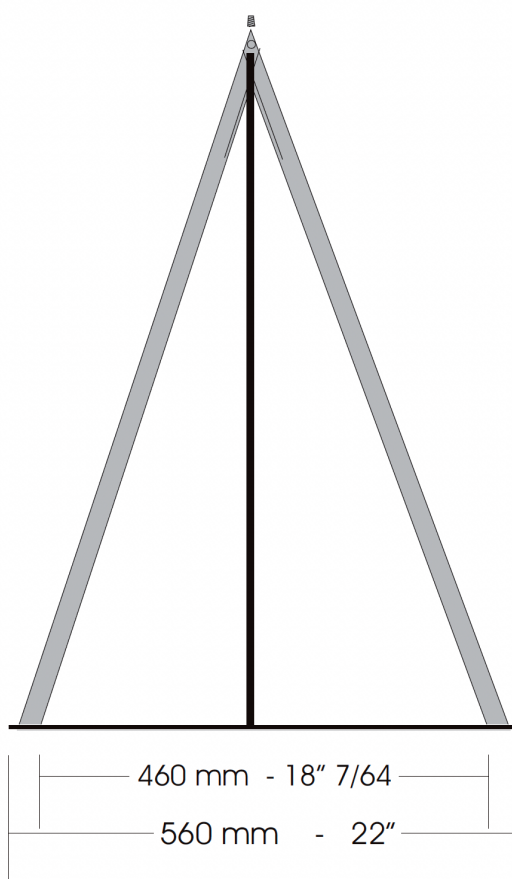
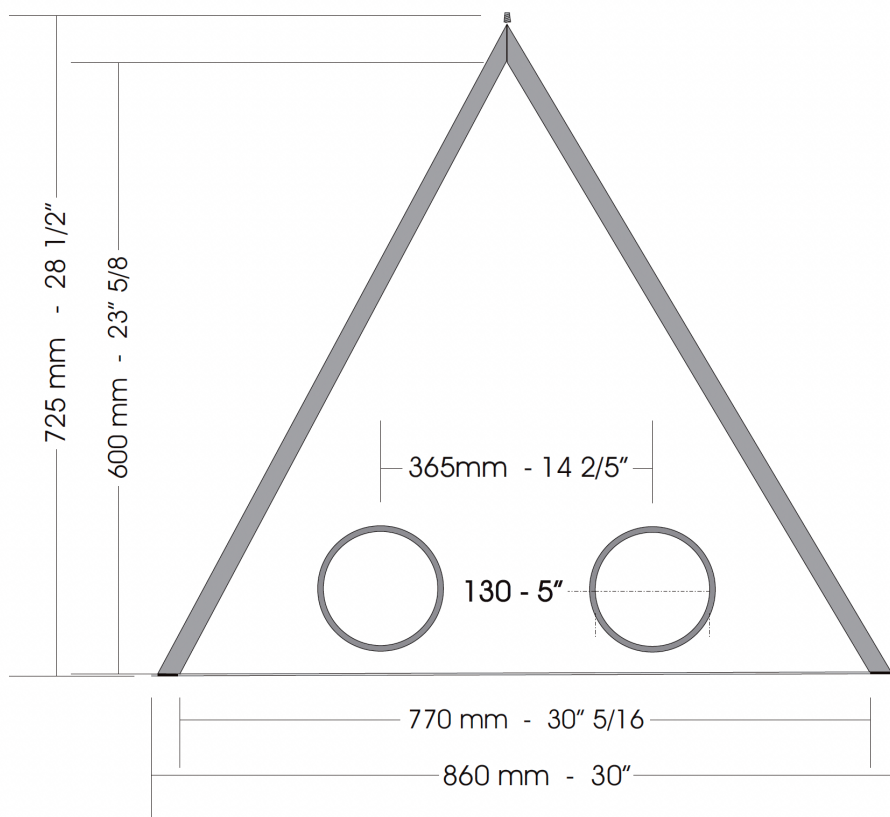
Pyramid

Erlab portable glove bag
La boîte à gants Erlab
Glovebox portatile Erlab
Caja de guantes transportable Erlab
Erlab tragbarer Handschuhkasten

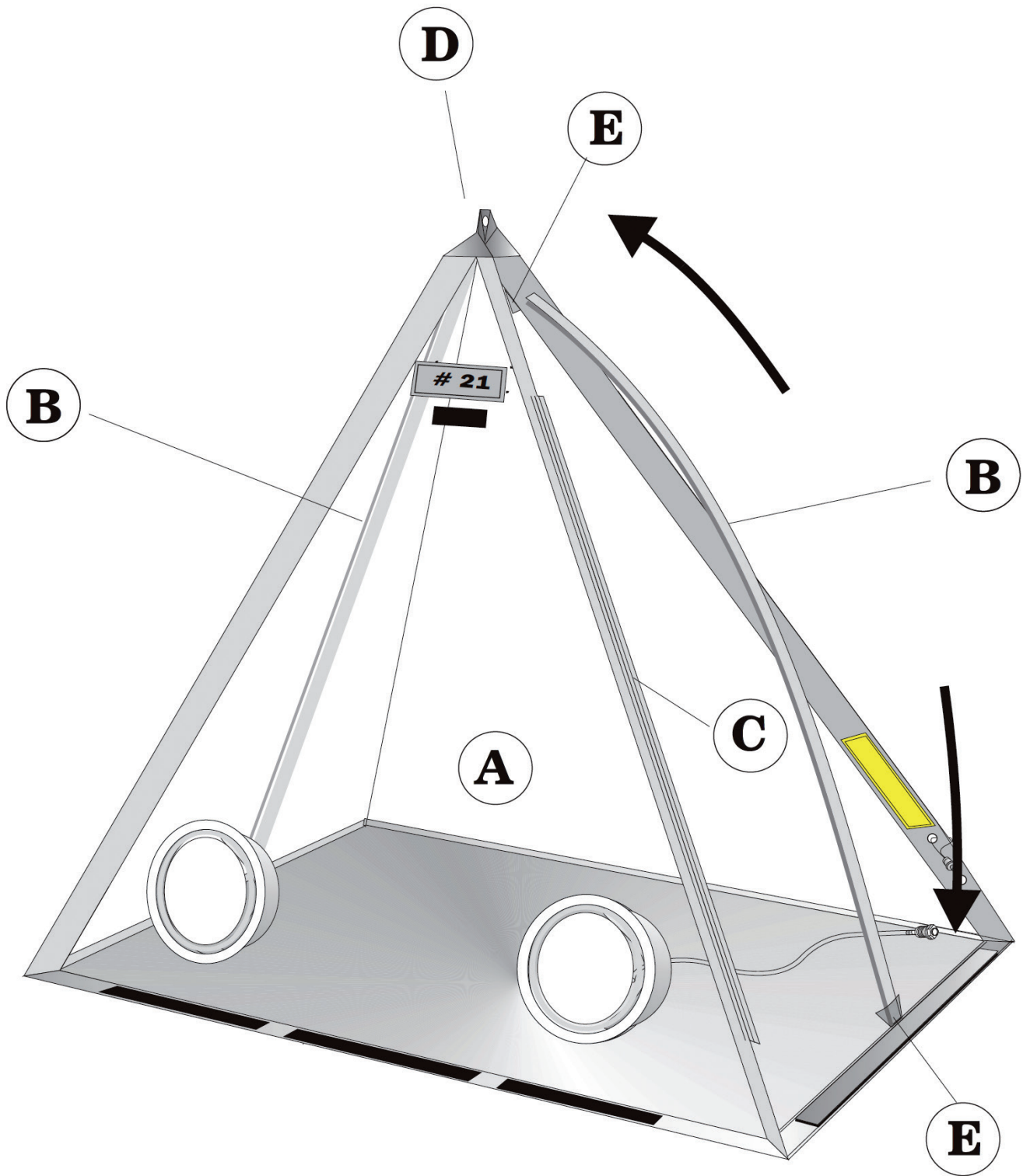Assembly Instructions **Notice de Montage** **Instrucciones de Montaje** **Montage Beschreibung** **Istruzioni per il Montaggio**

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IMPORTANT :

- a)** The Pyramid must always be carried by lifting the enclosure at its top.
- b)** After placing samples within the Pyramid, always check that the zip **C** is properly sealed prior to performing any manipulations. Place the thumb and the index fingers on each side of the Zip **C** and exert a constant pressure by sliding both fingers from top to bottom and from bottom to top. Repeat this procedure as many times as necessary.
- c)** When carried away, the Pyramid must always remain assembled.
- d)** Avoid pressure on the side walls once samples have been introduced so as to preserve their integrity before a full laboratory investigation can be made.

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Erlab's state of the art Research & Development Laboratory relies exclusively on filtration

About Erlab

We provide safety, we protect your health

Erlab invented the ductless fume hood in 1968. With more than 50 years of experience in the field of chemical filtration and protection of laboratory personnel; we know the formula for safety. With Erlab, you will never have to wonder or worry if our products are safe. We build each one of the following 7 ingredients into our products, and without all of them, your health and safety will be compromised.

1 Erlab R&D Laboratory

The engineers and chemists in our state-of-the-art R&D laboratory understand molecular filtration. We are committed to designing products that are safe and of the highest quality, strive to improve our products, and continuously develop new products that provide greater protection in the laboratory.

2 Strict Safety Standards

We hold ourselves to the highest standard and adhere to the strict AFNOR NF X 15-211: 2009 filtration safety standard as endorsed by ANSI Z9.5-2012.

3 A Published Chemical Listing

It all begins here. Without this listing, we are not compliant with AFNOR NFX 15-211. Our in-house laboratory tests, as well as independent testing, to verify the retention capacity of over 700 chemicals for our filters.

4 Independent Testing

Erlab filters have been independently tested multiple times at various concentrations guaranteeing that our safety solutions all adhere to the strict performance criteria of the AFNOR NF X 15-211:2009 standard assuring that the emission concentration at the filter exhaust will always be lower than 1% of the TLV.

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Our laboratory specialists will recommend the appropriate filtration fume hood, type of filter, and personalized advice.

6 Certificate of Validation for the chemicals used in the hood

A certified PhD chemist issues a Certificate of Validation with a list of the chemicals approved for use in the hood.

7 Our Safety Program

We back up our products 100%. This program includes your specialized chemical evaluation, validation of your hood upon installation, and a filtration safety specialist at your service to ensure that your hood is operating to its full potential.