pHenomenal VWR[®] MD 8000 H

European Catalogue Number(s):

PH SET:	665-0494
COND SET:	665-0495
DI.OX SET:	665-0496

Version: 1
Issued: November 28, 2018 8:37 am

Legal Address of Manufacturer:

VWR International bvba
Researchpark Haasrode 2020
Geldenaaksebaan 464
B-3001 Leuven
+32 16 385011
<http://be.vwr.com>

Country of origin: Germany.

Contents

EN English	4
DE Deutsch.....	6
ES Español.....	8
FR Français	10
IT Italiano	12
PT Português	14
CS Český	16
DA Dansk	18
EL Ελληνικά	20
FI Suomi	22
HU Magyar	24
NL Nederlands	26
NO Norsk	28
PL Polski.....	30
RO Română	32
RU Русский	34
SK Slovenský	36
SV Svenska	38
TR Türkçe	40

MD 8000 H

Overview

The MD 8000 H meter enables you to perform measurements (pH, conductivity, D.O.) quickly and reliably.

The MD 8000 H meter provides a maximum degree of operating comfort, reliability and measuring certainty.


Keypad

Keys are indicated by brackets <..>.

The key symbol (e.g. <ENTER>) generally indicates a short keystroke (under 2 sec). A long keystroke (approx. 2 sec) is indicated by the underscore behind the key symbol (e.g. <ENTER__>).


<F1>	Situation dependent function (is shown on the display) e.g.: <F1>/[Menu]: Opens the menu for measurement settings <F1__>/[Menu]: Opens the menu for system settings
<F1__>	
<F2>	
<F2__>	
<On/Off>	Switches the meter on or off
<M>	Selects the measured parameter
<CAL>	Calls up the calibration procedure Displays the calibration data
<CAL__>	
<STO>	Saves a measured value manually Opens the menu for the automatic save function
<STO__>	
<RCL>	Displays the manually stored measured values Displays the automatically stored measured values
<RCL__>	
<▲><▼> <◀><▶>	Menu control, navigation
<ENTER>	Opens the menu for measurement settings / confirms entries Opens the menu for system settings
<ENTER__>	
<AR>	Freezes the measured value (HOLD function) Switches the AutoRead measurement on or off

Display (example: pH measurement)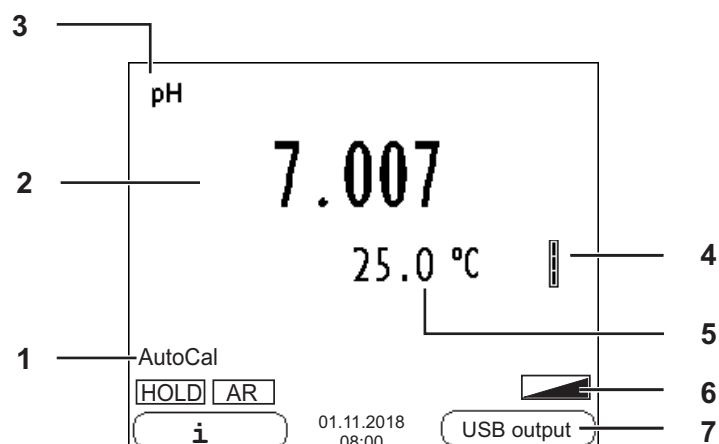

1	Status information
2	Measured value (with unit)
3	Measured parameter
4	Sensor symbol (calibration evaluation)
5	Measured temperature (with unit)
6	Further status information
7	Softkeys and date/time

Jack field /connections

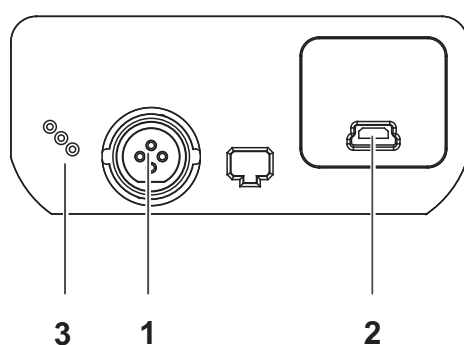1	Digital sensors (IDP) (pH, conductivity, D.O.)
2	USB B interface (device)
3	Service interface

Initial
commissioning

- 1 Insert the supplied batteries.
Make sure that the poles are in the correct position.
- 2 Connect the power pack if necessary.
- 3 Switch on the meter with the <On/Off> key.

Caution

Follow the safety instructions of the used sensors.
The operating manuals of the sensors are available on the supplied CD.

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VWR International AB
VWR International AG
VWR International Ltd