Instrucciones

INSTALACIÓN:

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United States
VWR International, LLC
Radnor Corporate Center, Building
One, Suite 200
100 Matsonford Road
Radnor, PA 19087
800-932-5000
<http://www.vwr.com>

Instructions

4-Channel Electronic Timer and Clock

North American Catalog Number 89203-042

European Catalogue Number 609-0181

INSTALLATION:

1. Open the battery cover.
2. Install battery with correct polarity positioning.
3. Close the battery cover.
4. Peel off display protective sheet.
5. Press and hold [CLOCK] until display flashing.
6. Press [12/24] button on the back to select 12 or 24 hours time format.
7. Press [HR], [MIN] and [SEC] to set the correct time. *
8. Press [CLOCK] to finish clock setting.

OPERATION:

COUNT DOWN

1. Press [TI] to select timer 1.
2. Press [HR], [MIN] and [SEC] to set the desired count down time. *
3. Press [START/STOP/MEMORY] to start TIMER 1's [TI] count down.
4. Press the [START/STOP/MEMORY] button again to pause.
5. Press the [START/STOP/MEMORY] button once more to resume count down.
6. The timer will sound when display value becomes 0:00:00 and will count up to show the elapsed time.
7. Press [START/STOP/MEMORY] to stop TIMER 1.
8. Press the button again to recall the memorized preset value.
9. Press [CLEAR] to clear all values.
10. Press T2, T3 or T4 to select the desired timer.
11. Repeat step 2 to 9 to set and run the selected timer.

COUNT UP

1. Press TI, T2, T3, or T4 to select the desired timer.
2. Press [CLEAR] to clear display values.
3. Press [START/STOP/MEMORY] to start count up.
4. Press the [START/STOP/MEMORY] button again to pause.
5. Press the [START/STOP/MEMORY] button once more to resume count up.

*During setting the values, press and hold the [HR], [MIN] or [SEC] button will increase the values automatically.

Battery Handling/Packaging

- Remove batteries from equipment immediately after they fail to operate the equipment.
- If equipment uses two or more batteries, always replace batteries in complete sets.
- Do not attempt to recharge alkaline batteries, they are not rechargeable.
- Do not handle hot or warm batteries.
- Do not heat, incinerate, crush, puncture or mutilate batteries.
- Wear personal protective equipment if batteries show signs of leakage, bulging, swelling or deformity.

Warning: If potassium hydroxide electrolyte from an alkaline battery comes in contact with the skin, do not try to neutralize the electrolyte with vinegar or any other acidic solutions. Neutralization may trap electrolyte on the skin. Flush the affected skin area with copious amounts of water. If the battery electrolyte gets into your eyes, it can cause severe damage and/or blindness.

Equipment disposal

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste. Instead it's your responsibility to correctly dispose of your equipment at lifecycle -end by handing it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment. For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment. By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health. Thank you

Technical service

Web Resources

Visit the VWR website at www.vwr.com for:

- Complete technical service contact information
- Access to the VWR Online Catalogue, and information about accessories and related products
- Additional product information and special offers

Contact us For information or technical assistance contact your local VWR representative or visit www.vwr.com.

Warranty

VWR warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non performance. Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

Compliance with local laws and regulations

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorizations necessary to run or use the product in its local environment. VWR will not be held liable for any related omission or for not obtaining the required approval or authorization, unless any refusal is due to a defect of the product.

Manufactured for

Europe
VWR International bvba
Researchpark Haasrode 2020
Geldenaaksebaan 464
B-3001 Leuven
+ 32 16 385011
<http://www.vwr.com>

United States
VWR International, LLC
Radnor Corporate Center, Building
One, Suite 200
100 Matsonford Road
Radnor, PA 19087
800-932-5000
<http://www.vwr.com>

Post Consumer Waste
996050004 Rev. 7
Printed in the USA

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