VWR® Thermometer

INSTRUCTION MANUAL

North American Catalog Numbers: 62344-734

European Catalogue Numbers: 620-1673

Version: 1

Issued: 3rd November 2014

Legal Address of Manufacturer

United States

VWR International
Radnor Corporate Center
Building One, Suite 200
100 Matsonford Road, P.O. Box 6660
Radnor, PA 19087
800-932-5000
<http://www.vwr.com>
Country of Origin – Hong Kong

Europe

VWR International bvba
Researchpark Haasrode 2020
Geldenaaksebaan 464
B-3001 Leuven
+ 32 16 385011
<http://be.vwr.com>

Intended use

This item is intended for general laboratory use to monitor time, temperature, and humidity.
Before using this product, open the battery compartment and remove the battery tab.

Product Specifications

Range: 32.0 to 122.0°F (0.0 to 50.0°C) Resolution: 0.1° Accuracy: $\pm 1^{\circ}\text{C}$ RH range: 20 to 90% Resolution: 1% RH
Accuracy: $\pm 5\%$ RH mid-range, otherwise 8% Dimensions: 108 x 57 x 12.7 mm Weight: 71g Description of Buttons and
Switches / Operational keys

Key Functions Location

Clear Clears the memory Front Humidity Memory Displays the Minimum and Maximum Humidity Front TEMP Memory
Displays the Minimum and Maximum Temperature Front HR Sets the Hour Back MIN Sets the Minutes Back Clock
Set/LOCK Switch set to Clock Set allows the user to set the time; LOCK locks the display to read the current time of
day Back 12HR/24HR Switch between 12hr (AM/PM) time and 24hr (military/or European) time Back °C/°F Switch
between Celsius to Fahrenheit Back

Getting Started

Before using this hygrometer, open the battery compartment and remove the battery tab. If probe temperature
display is desired, plug the supplied temperature probe into the socket located on the side of the hygrometer.
Operation

SETTING TIME-OF-DAY

Slide the CLOCK SET switch, located on the back of the unit, to the CLOCK SET position. Press the HR button to set
the hours. Press the MIN button to set the minutes. Press and hold the button to rapidly advance the display. With
the desired time-of-day indicated on the display, slide the CLOCK SET switch to the LOCK position.

12/24 HOUR DISPLAY

Slide the 12HR/24HR switch, located on the back of the unit, to the desired position. In 12 hour format, "AM/PM" will appear on the display.

SELECTING °C OR °F

To display the temperature reading in Fahrenheit or Celsius, slide the °C/°F switch, located on the back of the unit, to the desired position.

ABOUT THE MEMORY

Minimum and maximum memories are NOT programmable. The minimum temperature and humidity recorded into memory are the minimum temperature and humidity achieved since the last time the memory was cleared. The maximum temperature and humidity recorded into memory are the maximum temperature and humidity achieved since the last time the memory was cleared. The minimum and maximum memories are maintained over the period since the memory was cleared.

VIEWING THE HUMIDITY MEMORY

To view the humidity memory, press the HUMIDITY MEMORY button. The maximum humidity reading achieved will be displayed ("MAX" will appear on the display). Press the HUMIDITY MEMORY button a second time, within 10 seconds, to view the minimum humidity achieved ("MIN" will appear on the display). **Note:** Ten (10) seconds after pressing the HUMIDITY MEMORY button, the unit automatically returns to the current temperature/humidity reading. Press the HUMIDITY MEMORY button a third time, within 10 seconds, to return to the current humidity reading.

CLEARING THE HUMIDITY MEMORY

The minimum and maximum humidity memory may be cleared independently. To clear the desired humidity memory (MIN or MAX), press the CLEAR button while that memory value is being displayed (see the "Viewing the Humidity Memory" section). The display will show "--" for approximately 8 seconds and will then display the current humidity reading. **VIEWING THE TEMPERATURE MEMORY** To view the temperature memory, press the TEMP MEMORY button. The maximum temperature reading achieved will be displayed in the top display ("MAX" will appear to the right of the value). The minimum temperature reading achieved will be displayed in the middle display ("MIN" will appear to the right of the value). Press the TEMP MEMORY button a second time, within 5 seconds, to return the current temperature reading and time-of-day display.

Note: Five (5) seconds after pressing the TEMP MEMORY button, the unit automatically returns to the current temperature/time-of-day display.

CLEARING THE TEMPERATURE MEMORY

To clear the temperature memory (MIN and MAX), press the CLEAR button while the temperature memory is being displayed (see the "Viewing the Temperature Memory" section). The display will show "--" for approximately 8 seconds and will then display the current temperature and time-of-day.

Troubleshooting

If this unit does not function properly for any reason, replace the battery with a new high quality battery (see the “Battery Replacement” section). Low battery power can occasionally cause any number of “apparent” operational difficulties. Replacing the battery with a new fresh battery will solve most difficulties.

BATTERY REPLACEMENT

Erratic readings, faint readings, or no display are all indications that the battery must be replaced. Remove the battery cover by sliding it in the direction indicated by the arrow. Remove the exhausted battery and replace it with a new AAA alkaline battery. Make certain to insert the new battery with the proper polarity as indicated in the illustration in the battery compartment. Replace the battery cover. Technical service

Web Resources Visit the VWR’s website at www.vwr.com for:

- Complete technical service contact information
- Access to VWR’s Online Catalogue, and information about accessories and related products
- Additional product information and special offers

Contact us For information or technical assistance contact your local VWR representative or visit. www.vwr.com.

Warranty VWR International warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of purchase. If a defect is present, VWR will, at its option, repair, replace, or refund the purchase price of this product at no charge to you, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. For your protection, items being returned must be insured against possible damage or loss. This warranty shall be limited to the replacement of defective products. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

Equipment disposal

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste. Instead it's your responsibility to correctly dispose of your equipment at lifecycle -end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment. For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment. By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

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Local VWR offices in Europe and Asia Pacific

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India

VWR Lab Products Private Limited
135/12, Brigade Towers, 2nd Floor
Front wing, Brigade Road,
Bengaluru, India – 560 025
Tel: +91-2522-647911/922 (Mumbai)
Tel: +91-80-41117125/26 (Bangalore)
Fax +91-80-41117120
E-mail: vwr_india@vwr.com

Ireland / Northern Ireland

VWR International Ltd / VWR International
(Northern Ireland) Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: 01 88 22 222
Fax: 01 88 22 333
E-mail sales@ie.vwr.com

Italy

VWR International PBI S.r.l.
Via San Giusto 85
20153 Milano (MI)
Tel.: 02-3320311/02-487791
Fax: 02-332031307/02-40090010
E-mail: info@it.vwr.com

The Netherlands

VWR International B.V.
Postbus 8198
1005 AD Amsterdam
Tel.: 020 4808 400
Fax: 020 4808 480
E-mail: info@nl.vwr.com

New Zealand

VWR International LP
241 Bush Road
Albany
0632 New Zealand

Norway

VWR International AS
Haavard Martinsens vei 30
0978 Oslo
Tel.: 02290
Fax: 815 00 940
E-mail: info@no.vwr.com

Poland

VWR International Sp. z o.o.
Limbowa 5
80-175 Gdansk
Tel.: 058 32 38 210
Fax: 058 32 38 205
E-mail: labart@pl.vwr.com

Portugal

VWR International - Material de
Laboratório, Lda
Edifício Neopark
Av. Tomás Ribeiro, 43- 3 D
2790-221 Carnaxide
Tel.: 21 3600 770
Fax: 21 3600 798/9
E-mail: info@pt.vwr.com

Singapore

VWR Singapore Pte Ltd
18 Gul Drive
Singapore 629468
Tel: +65 6505 0760
Fax: +65 6264 3780
E-mail: sales@sg.vwr.com

Spain

VWR International Eurolab S.L.
C/ Tecnología 5-17
A-7 Llinars Park
08450 - Llinars del Vallès
Barcelona
Tel.: 902 222 897
Fax: 902 430 657
E-mail: info@es.vwr.com

Sweden

VWR International AB
Fagerstagatan 18a
163 94 Stockholm
Tel.: 08 621 34 00
Fax: 08 621 34 66
E-mail: info@se.vwr.com

Switzerland

VWR International GmbH
Lerzenstrasse 16/18
8953 Dietikon
Tel.: 044 745 13 13
Fax: 044 745 13 10
E-mail: info@ch.vwr.com

Turkey

Pro Lab Laboratuvar Teknolojileri Ltd.Şti.
a VWR International Company
Orta Mah. Cemal Gürsel Caddesi
Ördekcioglu İşmerkezi No.32/1
34896 Pendik - Istanbul
Tel.: +90216 598 2900
Fax: +90216 598 2907
Email: info@pro-lab.com.tr

UK

VWR International Ltd
Customer Service Centre
Hunter Boulevard
Magna Park
Lutterworth
Leicestershire
LE17 4XN
Tel.: 0800 22 33 44
Fax: 01455 55 85 86
E-mail: uksales@uk.vwr.com