07533-50 with Pump Head 07518-02

OPERATING MANUAL:
PUMP DRIVES

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07533-80

WARNING: PRODUCT USE LIMITATION

These products are not designed for, nor intended for use in, patient-connected applications, including, but not limited to, medical and dental use and, accordingly, have not been submitted for FDA approval.

Variety of Pump Heads Accepted

Mount 1 or 2 MASTERFLEX® L/S® Pump Heads, and all MASTERFLEX-compatible pump heads.

Setup and Drive Operation

1. Mount pump head and load tubing. (See pump head manual.)
2. Connect 16 ga. power input leads to power supply.
 - Positive input: Red
 - Negative input: Black
 - Case Ground: Green

Optional

3. Connect tachometer and/or cooling fan to rear motor shaft.

NOTE:

Shaft dimensions: 1.0 in ± 0.050 in long;
0.375 in ± 0.010 in diameter.

Shaft speed: approx. 33 times pump head shaft speed (low-speed units) approx. 6 times pump head shaft speed (high-speed units).

4. Provide enclosure.

OEM Customers: Contact your dealer for a complete set of mechanical and electrical considerations.

Replacement Parts

- 07553-06, Gear Service Kit, High-speed units
- 07553-08, Gear Service Kit, Low-speed units
- 07553-09, Gear only, High-speed units
- 07533-01, Motor Brushes, set of two
- 07520-03, Brush cap holder

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SPECIFICATIONS
SPÉCIFICATIONS
TECHNISCHE DATEN
ESPECIFICACIONES
SPECIFICHE

WARRANTY

Use only MASTERFLEX precision tubing with MASTERFLEX pumps to ensure optimum performance. Use of other tubing may void applicable warranties.

The Manufacturer warrants this product to be free from significant deviations from published specifications. If repair or adjustment is necessary within the warranty period, the problem will be corrected at no charge if it is not due to misuse or abuse on your part, as determined by the Manufacturer. Repair costs outside the warranty period, or those resulting from product misuse or abuse, may be invoiced to you.

The warranty period for this product is two (2) years from date of purchase.

PRODUCT RETURN

To limit charges and delays, contact the seller or Manufacturer for authorization and shipping instructions before returning the product, either within or outside of the warranty period. When returning the product, please state the reason for the return. For your protection, pack the product carefully and insure it against possible damage or loss. Any damages resulting from improper packaging are your responsibility.

TECHNICAL ASSISTANCE

If you have any questions about the use of this product, contact the Manufacturer or authorized seller.

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Masterflex LLC