

SPECIFICATIONS


Range: 14.0 to 230.0 °F / -10.0 to 110.0 °C

Resolution: 0.1°

Display Update Rate: 10 seconds

OPERATION

1. Attach the probe and secure the connection by tightening. The connector and plug are keyed. Make certain that the plug is inserted correctly, do not insert the plug incorrectly and force it.

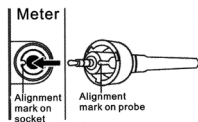

2. Press the  button to turn the unit on.
3. Insert the probe into the material to be measured and read the display.
4. Press the  button to turn the unit off. To conserve battery life, always turn the unit off when not in use.


SELECTING °F OR °C

1. Press the  button to turn the unit off.
2. Press and hold the  button for 8 seconds.
3. Continue to press the  button when the display shows a reading. The display will alternate between showing the temperature in °F and °C.
4. Release the  button once the desired measurement unit is displayed.

PROBE USE

The probe is supplied with a holder and ferrel. The ferrel may be tightened so that while using with the holder the probe may be set at the desired depth.


The probe and cable are waterproof; however, for extended measurement periods, do not leave the probe handle and cable submersed. Utilise the holder and ferrel for extended measurement.

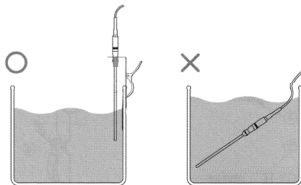

DISPLAY MESSAGES

- E 0 2** Indicates that the temperature being measured is below the measurement range of the thermometer or that the probe and/or cable has been damaged.
- E 0 3** Indicates that the temperature being measured is above the measurement range of the thermometer or that the probe and/or cable has been damaged.
- E 3 1** Indicates that the electronics have been damaged.

WALL MOUNT HANGER

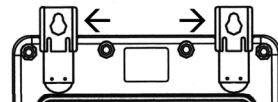

ALL OPERATIONAL DIFFICULTIES

If this thermometer does not function properly for any reason, replace the batteries with new high quality batteries (see the "Battery Replacement" section). Low battery power can occasionally cause any number of "apparent" operational difficulties. Replacing the batteries with new fresh batteries will solve most difficulties.

BATTERY REPLACEMENT

Erratic readings, a faint display, no display or  appearing on the display are all indications that the batteries must be replaced. Remove the battery cover by removing the six battery cover screws. Remove the exhausted batteries and replace them with 4 new AAA Alkaline batteries. Make certain to install the new batteries in the directions indicated in the battery compartment (improper battery placement may cause damage to the electronics). Replace battery cover and tighten the 6 screws securely.

Technical service

Web Resources

Visit the VWR's website at www.vwr.com for:

- Français, Español, Deutsch, Italiano, Português version
- Complete technical service contact information
- Access to VWR's Online Catalogue, and information about accessories and related products
- Additional product information and special offers

Contact us for information or technical assistance contact your local VWR representative or visit: www.vwr.com.

Warranty

VWR International warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of purchase. If a defect is present, VWR will, at its option, repair, replace, or refund the purchase price of this product at no charge to you, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. For your protection, items being returned must be insured against possible damage or loss. This warranty shall be limited to the replacement of defective products. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

**GO TO VWR.COM FOR THE LATEST NEWS,
SPECIAL OFFERS AND DETAILS OF YOUR LOCAL
VWR DISTRIBUTOR**

Cat. No. 89369-136 / 620-1975

©2015 92-6406-60 Rev. 1 261015



**TRACEABLE®
EXTRA-LONG
PROBE
WATERPROOF
THERMOMETER
INSTRUCTIONS**