VWR

Stereomicroscope 

INSTRUCTION MANUAL

Model	European Catalogue Number
STB150	630-2942

Version: 1.4

Issued: 18, 12, 2023

Legal Address of Manufacturer

Europe

VWR International bvba

Researchpark Haasrode 2020

Geldenaaksebaan 464

B-3001 Leuven

+ 32 16 385011

<http://be.vwr.com>

Country of origin: **ITALY**

Table of Contents

1.	Warning	3
2.	Safety Information	3
3.	Package Contents	3
4.	Unpacking	3
5.	Intended use	3
6.	Symbols and conventions	3
7.	Product specifications	4
8.	Instrument description	4
8.1	Assembling procedure	4
9.	Use of the microscope	5
9.1	Adjust interpupillary distance	5
9.2	Focus	5
9.3	Diopter compensation	5
9.4	Magnification	6
9.5	Contrast disc	6
9.6	Inserting the batteries	6
9.7	Use of illumination	7
10.	Troubleshooting	8
11.	Repair and maintenance	9
12.	Technical service	9
13.	Warranty	9
14.	Compliance with local laws and regulations	9
	Disposal	10

1. Warning

This microscope is a scientific precision instrument designed to last for many years with a minimum of maintenance. It is built to high optical and mechanical standards and to withstand daily use.

We remind you that this manual contains important information on safety and maintenance, and that it must therefore be made accessible to the instrument users.

We decline any responsibility deriving from incorrect instrument use that does not comply with this manual.

2. Safety Information

Avoiding Electrical Shock

Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off position.

Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users have full responsibility to use this equipment safely.

Please follow the guidelines below, and read this manual in its entirety to ensure safe operation of the unit.

3. Package Contents

DESCRIPTION	QUANTITY
Stereomicroscope	1
Stereo head -) Binocular, fixed zoom	1
Eyepiece WF10x/20	1 PAIR
Transmitted light illuminator	1
Incident light illuminator	1
Stand with integrated focus system	1
Stage clips	1 PAIR
Power supply output 5V dc multiplug	1
Dust cover	1

4. Unpacking

The microscope is housed in a moulded Styrofoam container. Remove the tape from the edge of the container and lift the top half of the container. Take care to ensure that the optical items (objectives and eyepieces) do not fall out and get damaged. Using both hands (one around the neck and one around the base), lift the microscope from the container and put it on a stable desk.

5. Intended use

For research and teaching use only. Not intended for any animal or human therapeutic or diagnostic use.

6. Symbols and conventions

The following chart is an illustrated glossary of the symbols that are used in this manual.

CAUTION

This symbol indicates a potential risk and alerts you to proceed with caution


ELECTRICAL SHOCK

This symbol indicates a risk of electrical shock.

7. Product specifications

Model	Head	Eyepiece	Objectives	Working distance	Stand	Illuminator
STB150	Binocular, 45° inclined, fixed	WF 10x/20	2x – 4x selectable	57 mm	Fixed with focus and handle	Incident: 1 W LED Transmitted: 1 W LED Touch brightness control Rechargeable batteries

8. Instrument description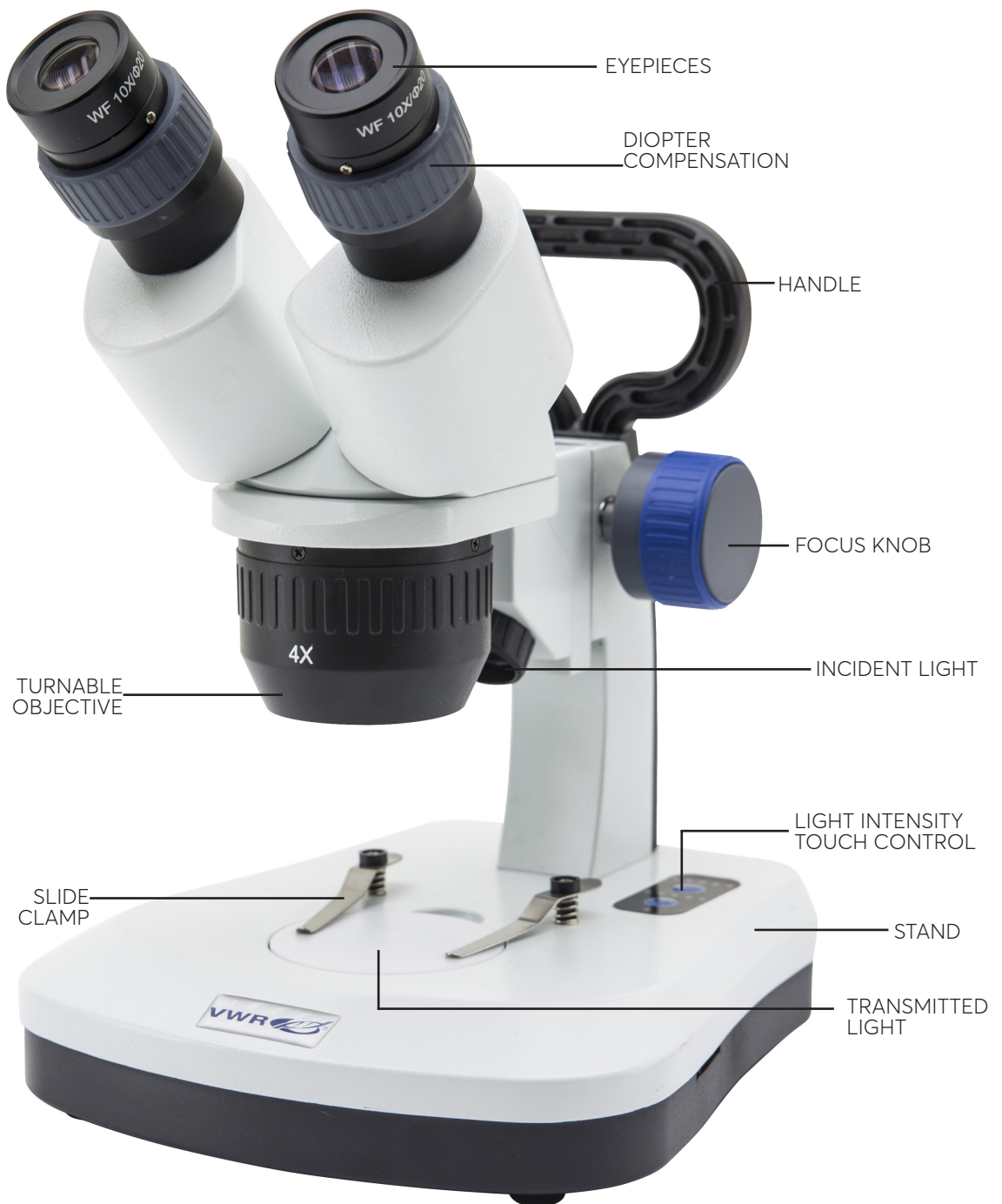

8.1 Assembling procedure


No installation procedure is required for this model.

The only thing you need to do is to connect the power supply to the microscope and the wall socket.

9. Use of the microscope

9.1 Adjust interpupillary distance

1. Hold the right and left parts of the observation head by both hands and adjust the interpupillary distance by turning the two parts (Fig. 1) until one circle of light can be seen.
 - If two circles appear, the interpupillary distance is too big.
 - If two overlapped circles appear, the interpupillary distance is too small.

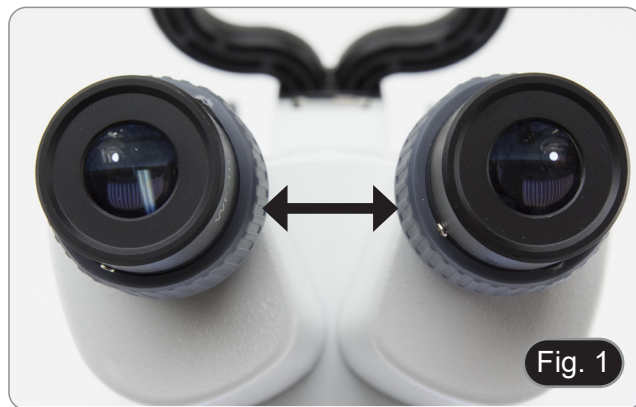9.2 Focus

Put the sample to be observed on the stage plate and focus the sample using the focusing knobs of the stand ①. (Fig. 2)

9.3 Diopter compensation

- This compensation makes possible for people wearing glasses to adjust the microscope to their eyes and use the microscope without glasses.
1. Insert the 2x objective and focus the specimen with the focusing knob. (Fig. 3)
 2. Insert the 4x objective and focus the specimen with the focusing knob.
 3. Return to the lowest magnification: the specimen will be out of focus.
-
4. Adjust the diopter compensation ring of the right eyepiece until the image of the right eyepiece is clear and sharp (Fig. 4). Repeat the procedure for the left eyepiece.
 5. Then, check the focus of the image for the whole zoom range. It should now be perfectly parfocal (focus is always maintained during the change of magnification).

9.4 Magnification

1. Rotate the objective to insert the desired lens in the optical path. (Fig. 5)
- Rotation in clockwise direction will move from 2X to 4X lens.
 - Counterclockwise rotation will change from 4X to 2X.

Fig. 5

9.5 Contrast disc

You can use the black/white disc in order to enhance the contrast of the image when using the incident illumination.

If you want to look at bright opaque objects, place the disc with the black side up. (Fig. 6)

Fig. 6

9.6 Inserting the batteries

- **Rechargeable batteries are not provided; it is necessary to use 3x AA rechargeable NiMh 1.2V.**
 - Before you start using the instrument, you must insert the rechargeable batteries.
1. Open the battery cover (located on the bottom of the instrument) and insert the provided batteries, taking care to respect the polarities. (Fig. 7-8)
 2. When finished, close the battery cover again.

Fig. 7

Fig. 8

9.7 Use of illumination

LED ① indicates the power status of the microscope. (Fig. 9)

- LED ON (green): power supply connected or charged batteries inserted.
 - LED OFF (no color): Power supply not connected or batteries discharged or not inserted.
1. Press the "Inc" key or the "Tra" key to increase the light intensity.
 2. The green LEDs will light up with each press of the key, and the light intensity will increase accordingly.
- **Once the fifth LED is lit, a subsequent press of the key will turn off the LEDs and also the light of the microscope. This is an intended effect and is not a defect.**

10. Troubleshooting

Review the information in the table below to troubleshoot operating problems.

PROBLEM	CAUSE	SOLUTION
I. Optical Section:		
The illumination is ON, but the field of view is dark	The plug is not connected to the illumination	Connect the cable
	The brightness is too low	Adjust to a proper setting
The edge of the field of view is vignetted or the brightness is asymmetric	The incident illuminator is not correctly oriented	Change the angle of the incident illuminator
Dust and stains can be seen in the field of view	There are stains and dust on the specimen	Clean the specimen
	There are stains and dust on the eyepiece	Clean the eyepiece
Poor image quality: • The image is not sharp • The contrast is not high • The details are not clear • Image glares	The lenses (additional lens, objective, eyepieces) are dirt	Thoroughly clean all the optical system
One side of the image is out of focus	The specimen is out of place (tilted)	Place the specimen flat on the stage
II. Mechanical Section:		
The focus knob is hard to turn	The tension adjustment collar is too tight	Loosen the tension adjustment collar
The focus is unstable	The tension adjustment collar is too loose	Tighten the tension adjustment collar
III. Electric Section:		
The LED doesn't turn on	No power supply	Check the power cord connection
The brightness is not enough	The brightness adjustment is low	Adjust the brightness
The light blinks	The power cord is poorly connected	Check the power cord
IV. Viewing tube assembly:		
The field of view of the two eyes is different	The interpupillary distance is not correct	Adjust the interpupillary distance
	The diopter correction is not right	Adjust the diopter correction
	The viewing technique is not correct, and the operator is straining the eyesight	When look into the eyepieces, do not stare at the specimen but look at the whole field of view. Periodically, move the eyes away to look at a distant object, then back into the eyepieces

11. Repair and maintenance

Microscopy environment

This microscope is recommended for use in a clean, dry and shock free environment with a temperature of 0-40°C and a maximum relative humidity of 85 % (non condensing). Use a dehumidifier if needed.

To think about when and after using the microscope

The microscope should always be kept vertical when moving it so that no moving parts, such as the eyepieces, fall out. Never mishandle or impose unnecessary force on the microscope.

Never attempt to service the microscope yourself.

After use, turn down the illumination intensity control and turn the light off. Cover the microscope with included the dust cover, and keep it in a dry and clean place.

Electrical safety precautions

Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off-position.

Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users do have full responsibility to use this equipment safely.

Cleaning the optics

If the optical parts require cleaning first use compressed air.

If that is not sufficient use a soft lint-free cloth with water and a mild detergent.

And as a final option use the piece of cloth moistened with a 3:7 mixture of ethanol and ether.

- **Note: ethanol and ether are highly flammable liquids. Do not use them near a heat source, near sparks or near electric equipment. Use these chemicals in a well ventilated room.**

Remember to never wipe the surface of any optical items with your hands. Fingerprints can damage the optics. Do not disassemble objectives or eyepieces attempting to clean them.

For the best results, use the VWR cleaning kit.

If you need to send the microscope to manufacturer for maintenance, please use the original packaging.

12. Technical service

Web Resource

Visit the VWR's website at **www.vwr.com** for:

- Complete technical service contact information
- Access to VWR's Online Catalogue, and information about accessories and related products
- Additional product information and special offers

Contact us For information or technical assistance contact your local VWR representative or visit **www.vwr.com**.

13. Warranty


VWR International warrants that this product will be free from defects in material and workmanship for a period of five (5) years from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non-performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

14. Compliance with local laws and regulations

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorizations necessary to run or use the Product in its local environment. VWR will not be held liable for any related omission or for not obtaining the required approval or authorization, unless any refusal is due to a defect of the product.

Disposal

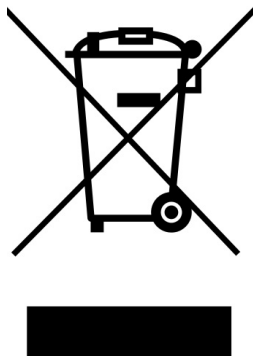This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste.

Instead it is your responsibility to correctly dispose of your equipment the end of its life cycle by handling it over to an authorized facility for separate collection and recycling. It is also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

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Ireland / Northern Ireland

VWR International Ltd /
VWR International (Northern Ireland) Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: 01 88 22 222
Fax: 01 88 22 333
E-mail: sales@ie.vwr.com

Italy

VWR International PBI S.r.l.
Via San Giusto 85
20153 Milano (MI)
Tel.: 02-3320311/02-487791
Fax: 800 152999/02-40090010
E-mail: info@it.vwr.com

The Netherlands

VWR International B.V.
Postbus 8198
1005 AD Amsterdam
Tel.: 020 4808 400
Fax: 020 4808 480
E-mail: info@nl.vwr.com

New Zealand

Global Science - A VWR Company
241 Bush Road
Albany 0632, Auckland
Tel.: 0800 734 100
Fax: 0800 999 002
E-mail: sales@globalscience.co.nz

Norway

VWR International AS
Haavard Martinsens vei 30
0978 Oslo
Tel.: 02290
Fax: 815 00 940
E-mail: info@no.vwr.com

Poland

VWR International Sp. z o.o.
Limbowa 5
80-175 Gdansk
Tel.: 058 32 38 200 do 204
Fax: 058 32 38 205
E-mail: info@pl.vwr.com

Portugal

VWR International -
Material de Laboratório, Lda
Edifício Neopark
Av. Tomás Ribeiro, 43- 3 D
2790-221 Carnaxide
Tel.: 21 3600 770
Fax: 21 3600 798/9
E-mail: info@pt.vwr.com

Singapore

VWR Singapore Pte Ltd
18 Gul Drive
Singapore 629468
Tel.: +65 6505 0760
Fax: +65 6264 3780
E-mail: sales@sg.vwr.com

Spain

VWR International Eurolab S.L.
C/ Tecnología 5-17
A-7 Llinars Park
08450 - Llinars del Vallès
Barcelona
Tel.: 902 222 897
Fax: 902 430 657
E-mail: info@es.vwr.com

Sweden

VWR International AB
Fagerstagatan 18a
163 94 Stockholm
Tel.: 08 621 34 00
Fax: 08 621 34 66
E-mail: kundservice@se.vwr.com

Switzerland

VWR International GmbH
Lerzenstrasse 16/18
8953 Dietikon
Tel.: 044 745 13 13
Fax: 044 745 13 10
E-mail: info@ch.vwr.com

Turkey

Pro Lab Laboratuvar Teknolojileri Ltd.Şti.
a VWR International Company
Orta Mah. Cemal Gürsel Caddesi
Ördekcioglu İşmerkezi No.32/1
34896 Pendik - Istanbul
Tel.: +90216 598 2900
Fax: +90216 598 2907
Email: info@pro-lab.com.tr

UK

VWR International Ltd
Customer Service Centre
Hunter Boulevard - Magna Park
Lutterworth
Leicestershire
LE17 4XN
Tel.: 0800 22 33 44
Fax: 01455 55 85 86
E-mail: uksales@uk.vwr.com

**GO TO VWR.COM FOR THE
LATEST NEWS, SPECIAL OFFERS
AND DETAILS OF YOUR LOCAL
VWR DISTRIBUTOR**

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