Short user's manual

VWR Refrigerator +5C 118L sparkfree

VWR Refrigerator +5C 118L

VWR Refrigerator +5C 310L sparkfree

VWR Refrigerator +5C 310L

VWR Freezer -25C 92L sparkfree

VWR Freezer -25C 92L

EU cat. No. 471-1307 / 471-1308 / 471-1309 / 471-1310 / 471-1311/ 471-1312

Manufactured in China

EN:

Temperature setting

The setpoint is pre-set at 5°C .

- ① Press "SET" for 5 seconds, "TS" displays.
- ② Press "SET" again, current setpoint flashes.
- ③ Use "+" or "-" to change the setpoint, then press "SET" to save.

Check max / min temperature

Press both "Alarm" and "+" to check max temperature.

Press both "Alarm" and "-" to check min temperature.

Precautions

- This product shall be used under specified ambient temperature. In an environment of high humidity and temperature, there may be condensation on the glass door, which is normal without any influence on storage temperature inside the refrigerator. In case of condensation, please wipe with dry cloth to reduce the condensation, please improve the ventilation and drop the ambient temperature if possible.
- This product is required to be placed in a location that is dry, ventilated and sheltered from sunlight, a minimum of 10cm space must be kept around the refrigerator for ventilation and heat dissipation and the top space shall not be less than 30cm.
- Do not immediately turn on the refrigerator after it is installed in place. Let it sit for 24 hours before power up to ensure a normal operation.
- At the time of the initial use, please run the equipment without load for 5 hours before putting in articles. When the temperature has not reached specified value, it is normal for the appearance of high-temperature alarm.
- Please refer to location of "load line" inside the refrigerator at the time of article storage in order to not block the interior air flow.
- Before loading the refrigerators, make sure the unit is at set temperature. Do not load in more than 1/3 of the storage volume to avoid thermally overloading the unit.
- If the door is opened, the temperature of the refrigerator will warm up. That is normal.
- The temperature will recover to a stable condition in a short time. To reduce the possibility of temperature fluctuation in the refrigerator, please attempt to cut down the time to open the door to remove and load products.
- Clean the condenser regularly (at least each 90 days) to ensure that the condenser functions properly.
- In case of alarm or other failures, please check the manual for troubleshooting or contact service technicians. Please do not disassemble by yourself.
- Do not store corrosive items such as acids and alkalis in the refrigerator and freezer. Failure to do so may damage the refrigerator-freezer internal components or electrical parts.
- Electric cabinet can only be opened by service technicians.

For more information, please consult the full manual.

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