

Protein Extraction and Lysis Buffers (PE LB™)

ABSTRACT

Use of PE LB™ Buffers for Extraction of Carbonic Anhydrase and Alkaline Phosphatase from *E. coli*, yeast, mouse pancreases and cultured human cells is evaluated.

Lysis and extraction of protein from cellular and tissue samples is the first critical step for biochemical analysis. Selection of lysis and extraction buffers requires considerations such as recovery of the protein of interest as well as stability of biological activity (1,2). G-Biosciences has developed a series of Protein Extraction & Lysis Buffers (PE LB™), which ensure good protein recovery while maintaining biological activity.

The following PE LB™ series offers a wide selection of buffers for lysis and extraction of protein from bacteria, yeast and animal cells and tissues. PE LB™ buffer system is based on a proprietary combination of organic buffering agents, mild non-ionic detergents, and a combination of various salts and agents to enhance extraction of protein and maintain stability of biological activities of the proteins. Depending on application, additional agents such as chelating agents, reducing agents and protease inhibitors may be added in to the PE LB™ buffer system.

The PE LB™ series of buffers is compatible with most downstream applications including running various chromatography, gel electrophoresis applications, and protein folding procedures. It is also compatible for protein estimation with NI™ (Non-Interfering) Protein Assay (Cat. No. 82022-356).

In this study, PE LB™ buffers are evaluated for the extraction of carbonic anhydrase and alkaline phosphatase from *E. coli*, yeast, mouse pancreases, and culture human cells. Protocols for the individual PE LB™ buffers were followed.

Bacterial PE LB™

Bacterial-PE LB™ (Cat. No. 82021-546) has been specifically developed for the extraction of total soluble proteins and inclusion bodies from bacterial cells. Bacterial-PE LB™ is a proprietary improvement on the lysozyme based lysis method.

Method: Freshly grown *E. coli* from a 1.5ml culture was used. The bacterial pellet was suspended in PE LB™ and an appropriate amount of LongLife™ Lysozyme (Cat. No. 82021-516), supplied with the kit, was added to the suspension. Protease cocktail ProteaseArrest™ (Cat. No.

82022-478) was added to prevent proteolysis damage to the proteins. The suspension was incubated in an ice bucket for 30 minutes. After 30 minutes, the suspension was vigorously vortexed and centrifuged to collect a clear lysate for further analysis.

Yeast PE LB™

Yeast-PE LB™ (Cat. No. 82023-266) is a proprietary improvement on the Zymolyase® based spheroplast preparation and extraction of soluble proteins from yeast cells.

Method: Freshly grown yeast from a 1.5ml culture was used. The yeast pellet was suspended in PE LB™ buffer and an appropriate amount of LongLife™ Zymolyase® (Cat. No.82021-514), provided with each kit, was added. Protease cocktail ProteaseArrest™ (Cat. No. 82022-478) was added to prevent proteolysis damage to the proteins. The suspension was incubated in an ice bucket for 30 minutes. After 30 minutes, the suspension was vigorously vortexed and centrifuged to collect a clear lysate for further analysis

Mammalian Cell PE LB™

Mammalian Cell-PE LB™ (Cat. No.82022-596) is useful for extraction of total soluble proteins from mammalian cultured cells.

Method: Freshly grown 5 million Jurkat cells were used in this study. The cell pellet was suspended in 0.5ml PE LB™. A protease cocktail, ProteaseArrest™ (Cat. No. 82022-478), was added into PE LB™ to prevent proteolysis damage to the proteins. The cell suspension was vortexed and pulled up and down the pipetor a few times. The suspension was centrifuged and a clear lysate was collected for analysis.

Tissue PE LB™

Tissue PE LB™ (Cat. No. 82022-598) has been specifically developed for the extraction of total soluble protein from animal tissues.

Method: 50mg mouse pancreas tissue was used in this study. The tissue sample was homogenized in 0.5ml PE LB™. Protease cocktail ProteaseArrest™ (Cat. No. 82022-478) was added into PE LB™ to prevent proteolysis damage to the proteins. The homogenate was centrifuged and a clear lysate was collected for analysis.



think proteins! think G-Biosciences!



RESULTS AND DISCUSSION

The results of carbonic anhydrase and alkaline phosphatase activities are shown in Figure 1. Bacterial PE LB™ eliminates the need for laborious mechanical lysis of bacterial cells and removal of nucleic acids. Bacterial PE LB™ allows extraction of soluble proteins and concurrent removal of nucleic acids released during cell lysis. The lysis by Bacterial PE LB™ eliminates viscosity build up.

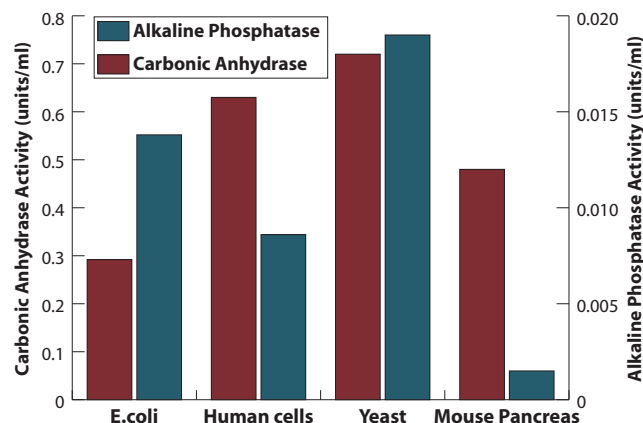


Figure 1: PE LB™ System maintains the biological activity of proteins. Extraction of carbonic anhydrase or alkaline phosphatase from E.coli, human cells, yeast and mouse pancreas with Bacterial, Mammalian Cell, Yeast and Tissue PE LB™ respectively. The resulting lysates were submitted to enzyme assays and both enzymes retain their biological activity.

Yeast PE LB™ buffer kit provides a simple and versatile method of yeast protein extraction and eliminates the need of glass beads for lysis of yeast cells. Yeast PE LB™ buffer eliminates viscosity build up due to the release of genomic nucleic acids, allowing effective clarification of the lysate.

Mammalian Cell PE LB™ and Tissue PE LB™ are simple to use, quickly extract proteins, and release carbonic anhydrase and alkaline phosphatase activities.

Insect PE LB™ has been developed to provide a simple and versatile method of extraction of total biologically active, soluble proteins from adherent or suspension cultured insect cells, including Sf9 and Sf21.

The protein extract prepared by using PE LB™ buffers are suitable for enzyme studies and other downstream applications such as protein purification, electrophoresis, immuno-assays, and so forth.

CONCLUSION

PE LB™ extraction and lysis buffers ensure good recovery of carbonic anhydrase and alkaline phosphatase activities while maintaining biological activity.

ORDERING INFORMATION

VWR Cat. No.	Description	Size
82021-546	Bacterial PE LB™ Kit with Lysozyme	100 x 50µl cell pellets
82022-592	Bacterial PE LB™ Buffer	500 mL
82023-266	Yeast PE LB™ Kit with Zymolyase®	100 x 50µl cell pellets
82022-594	Yeast PE LB™ Buffer	500 mL
82022-596	Mammalian PE LB™ Buffer	500 mL
82022-598	Tissue PE LB™ Buffer	500 mL
95057-584	Insect PE LB™	250 mL

REFERENCES

Tissue PE LB™

1. Kavanagh, K. et al (2012) J Gerontol A Biol Sci Med Sci. 10:1093
2. Kavanagh, K. et al (2011) Am J Physiol Endocrinol Metab 300:E894
3. Vukelic, S. et al (2011) J Biol Chem 286:10265
4. Ray, S. et al (2008) Mol Endocrinol 22:1125
5. Stein, D. et al (2008) J Antimicrob Chemother 62:555
6. Wang, Z. et al (2007) J Neurosci 27:3686
7. Shariat-Madar, Z. et al (2006) Blood. 108:192
8. Yoshino, O. et al (2006) PNAS 103:10678
9. Yao, L. et al (2005) Blood. 106:4093
10. Mangino, M. et al (2004) Am J Physiol Renal Physiol 286:F838

Mammalian Cell PE LB™

1. Zou, X. et al (2011) J. Biol. Chem. 286:1301
2. Yu, J. et al (2008) PNAS. 105:19300
3. Zhang, L. et al (2006) Cancer Gene Ther. 13:74
4. Valverde, P. et al (2004) Exp Eye Res. 78:27
5. Qin, M. et al (2003) Clin. Cancer Res. 9:4992

Yeast PE LB™

1. Saribas, A. et al (2004) Glycobiology 14:1217



1.800.932.5000
vwr.com



Prices and product details are current when published; subject to change without notice. | Certain products may be limited by federal, state, provincial, or local regulations. | VWR makes no claims or warranties concerning sustainable/green products. Any claims concerning sustainable/green products are the sole claims of the manufacturer and not those of VWR International, LLC. All prices are in US dollars unless otherwise noted. Offers valid in US and Canada, void where prohibited by law or company policy, while supplies last. | VWR, the VWR logo and variations on the foregoing are registered (®) or unregistered trademarks and service marks of VWR International, LLC and its related companies. All other marks referenced are registered by their respective owner(s). | Visit vwr.com to view our privacy policy, trademark owners and additional disclaimers. © 2012 VWR International, LLC. All rights reserved.