

Short user's manual

VWR® MINI VORTEX MIXER

EU cat. No 444-1337

Manufactured in China

Specifications

Speed	3000rpm
Orbital Diameter	4.5mm
Capacity	Max. 80ml
Tube Diameter	Max. 30mm



Operation Instructions

1. Please connect the device to the power supply.
2. Place the tube on the top socket.
3. Press the tube slightly to active shaking.
4. Lift the tube, and the device will stop shaking.

Cautions for Usage

1. Use only approved accessories.
2. Do not modify the power cord provided with the device.
3. Please only use the original power supply adaptor in the device.
4. Place the device on a non-slip and sturdy surface.
5. Be aware of the risk of squirting liquids while stirring.
6. Please avoid using more than 0.8 kg of pressure to activate the device.
7. Do not mix the liquid will react dangerously or use any flammable or any hazardous liquids.

For more information, please consult the full manual