

VWR[®] P Water System

INSTRUCTION MANUAL

European Catalogue Numbers:

171-1247	VWR[®] P Water System
171-1248	VWR[®] P Water System with TOC



Legal Address of Manufacturer

Europe

VWR International bv
Researchpark Haasrode 2020
Geldenaaksebaan 464
B-3001 Leuven
+ 32 16 385011
be.vwr.com

UK Importer:

VWR International Ltd
Hunter Boulevard, Magna Park
Lutterworth, Leicestershire, LE17 4XN
uk.vwr.com

Country of origin China

Table of Contents

1 INTRODUCTION

- 1.1 System Exterior
- 1.2 Product Features
- 1.3 Main Applications
- 1.4 Specifications
- 1.5 Operation Principles
- 1.6 Technical Specifications

2 INSTALLATION

- 2.1 Installation Site Requirements
- 2.2 Feed Water Requirements
- 2.3 Feed Water Quality Measurements
- 2.4 Installation

3 SYSTEM STARTUP AND OPERATION

- 3.1 Check List Prior to System Startup
- 3.2 System Initial Startup
- 3.3 System Setting-up
- 3.4 Information Menu
- 3.5 Maintenance Menu

4 DISPENSING WATER

- 4.1 Dispensing Water
- 4.2 Adjusting Dispense Rate
- 4.3 Volumetric Dispensing

5 MAINTENANCE

- 5.1 Basic Alerts
- 5.2 Basic Troubleshooting
- 5.3 Replacing Consumables
- 5.4 History Inquiry and Export

6 PARTS AND ORDER INFORMATION

7 TECHNICAL SERVICE

8 WARRANTY

9 DISPOSAL

1 INTRODUCTION

The VWR P water system is designed to produce ultrapure water (Type I) from pure water. Water quality produced meets or exceeds the corresponding ASTM, CAP, CLSI and ISO 3696 / BS 3997 water standards.

This manual describes in detail about system performance characteristics, installation, operation, and routine maintenance. Please read this manual thoroughly for its instructions on installation, use and maintenance. Proper installation, use and maintenance guarantee the continuous flow of high quality ultrapure water.

Please contact us or your local distributor if you encounter any issues during installation and use. Professional engineers are fully trained to support you.


Safety Information



WARNING!

- 1) Always power down the system before plug or unplug any hardware. Never hot swap (hot plug) any hardware, including dispenser, monitor as may damage control boards.
- 2) Disconnect the unit from the power supply prior to perform service work.
- 3) Refer service to qualified personnel.

1.1 System Exterior

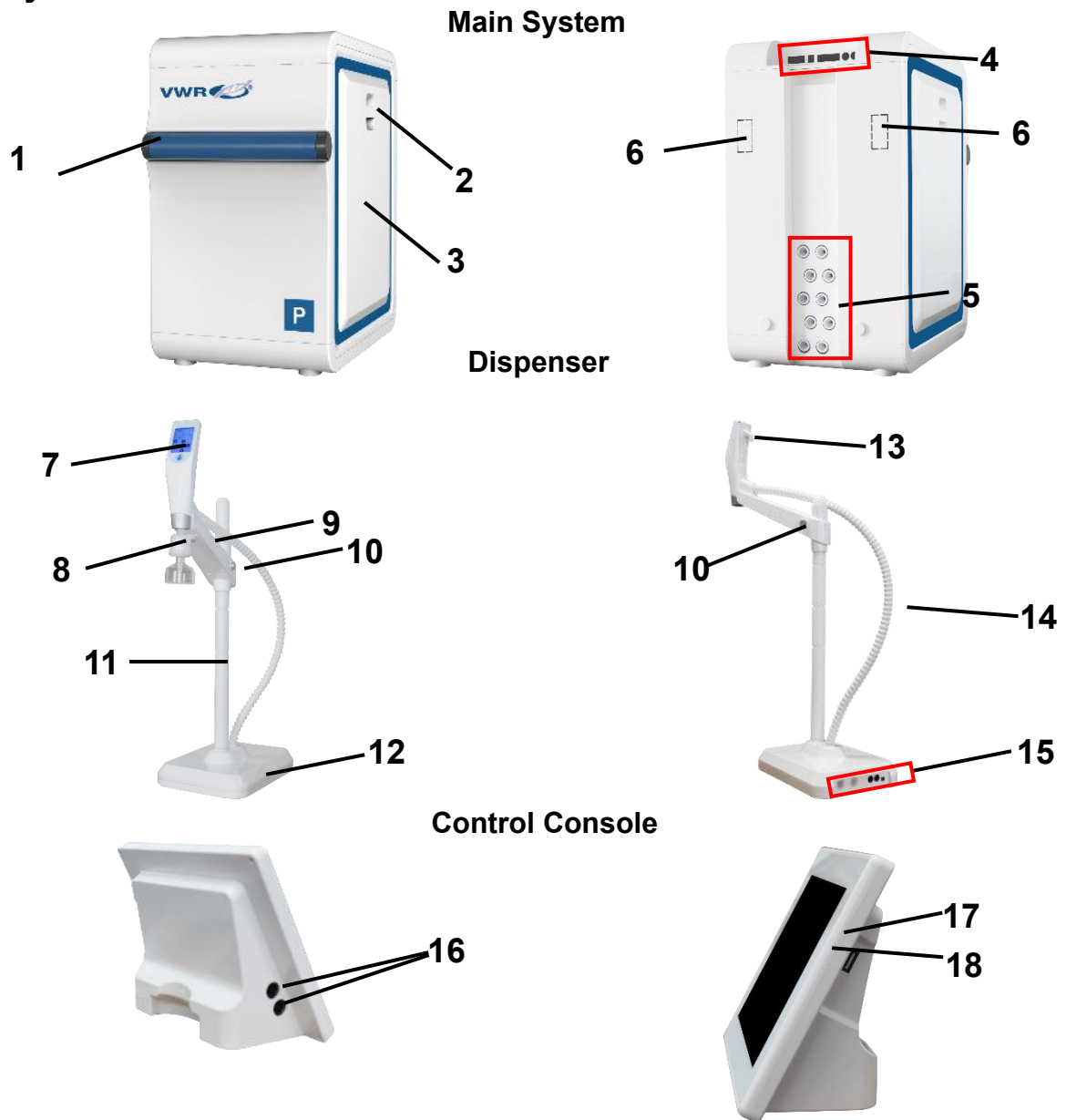

1	Cross bar and dispenser rest on with optional accessories	2	Side panel handle
3	Side panel	4	Electrical connections
5	Water outlets	6	Wall mount ports
7	Touchscreen	8	Final filter
9	Crossbar	10	Crossbar release button
11	Center post	12	Dispenser base
13	Finger rest hook	14	Corrugated tubing
15	Cable and water outlets	16	Outlets for CAN cable or power adapter
17	USB Port	18	SD card port

Description of the system

The VWR P manages the production and distribution of ultrapure water from purified water. It consists of three main units:

- ◆ Main System: manages the production of ultrapure water.
- ◆ Dispenser: integrates the touchscreen and manages the dispensing.
- ◆ Control Console: controls and monitors the water system and other components on the 8-inch touchscreen.

1.2 Product Features

The VWR P ultrapure water system provides an integrated solution for lab water supply. This system is easy to install, easy to operate and easy to maintain.

This system has the following features:

- ◆ Type I ultrapure water is produced from pure water.
- ◆ The system can be link to multiple dispensers via cable or wireless.
- ◆ Ultra purification cartridge is filled with media to remove trace ions and organics.
- ◆ Built-in 185/254 nm dual wavelength lamp reduces TOC level in ultrapure water.
- ◆ On-line TOC monitor is based on complete oxidation method to obtain more accurate measurement of the TOC.
- ◆ Various final filters (optional) can ensure ultrapure water without particles, bacteria or pyrogen.
- ◆ RFID tags ensure perfect placement of consumables and trace their performance.
- ◆ Manual and volumetric dispensing, adjustable dispensing rate, and water quality monitoring are built into the dispenser screen.
- ◆ The control console is an 8-inch touchscreen. The console controls system and peripheral devices (such as dispensers and tank sanitization module). All operations can be done on the console.
- ◆ Water quality, operation parameters, the status of the system, dispensers, components, and peripheral devices are stored and displayed on the large color touchscreen monitor.
- ◆ The console screen and dispenser screens are water-proof. You may operate the console and dispenser with wet latex gloves on.
- ◆ Signature verification is required for maintenance and service.
- ◆ VWR provides full documents support to meet user's GMP, GSP, GAP, GCP, GLP requirements.

1.3 Main Applications

Ultrapure water can be used in many areas. Here are some typical applications.


- HPLC (High Performance Liquid Chromatography) mobile phase preparation
- Preparation of reagent blank solution
- As sample diluents for GC, HPLC, AA, ICP-MS and other analytical techniques
- Preparation of buffer and culture media for mammalian cell culture
- Nano-material science
- Preparation of molecular biology reagents, etc.

1.4 Specifications

Ultrapure Water (Type I)	
Dispensing Rate	Variable speed dispensing up to 2 L/min
Resistivity	18.2 MΩ.cm (at 25 °C)
TOC	< 5 ppb
Particles (>0.2 µm)	<1 /mL (with a 0.2 µm final filter)
Microorganism	< 0.01 cfu/mL (with a 0.2 µm final filter or Bio-filter)
Pyrogen Content	< 0.001 EU/mL (with Bio-filter)
RNAses	< 0.5 pg/mL (with Bio-filter)
DNAses	< 10 pg/mL (with Bio-filter)
Electrical Connections and Specifications	
Operating Voltage	24 VDC
Input Voltage	100 – 240 VAC
Main System Power	< 200 W
Dimensions and Weights	
Main System Width × depth × height	32 × 44 × 54 cm (12.6 × 17.3 × 21.3 in)
System Weight	17 kg
Control Console Length × width	22 × 17 cm (8.7 × 6.7 in)
Control Console Weight	0.75 kg
Dispenser Width × depth × height	21 × 29 × 61 cm (8.3 × 11.4 × 24.0 in)
Dispenser Weight	5 kg

1.5 Operation Principles

VWR P ultrapure water system produces ultrapure water from purified water. Pure water flows through H Pack for further deionization, through a UV lamp chamber with a dual wavelength UV lamp to destroy trace organic pollutants in water, then through a polishing cartridge to remove the last trace of ions, then a 0.2 µm final filter to the outlet.

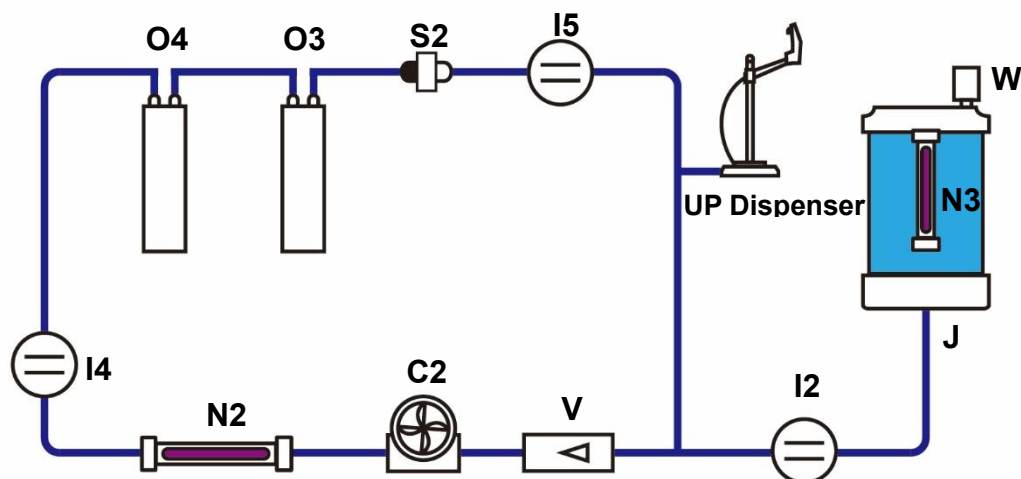

No.	Description	No.	Description
C2	Distribution Pump	I2	Feed Water Conductivity Sensor
I4	TOC Sensor	I5	UP Resistivity Sensor
J	Pure Water Tank (optional)	N2	185 nm UV Lamp
N3	Tank UV Lamp	O4	H Pack
O3	U Pack	S2	Inline Filter
V	Flow Meter	W	Tank Vent Filter

1.6 Technical Specifications

Measurement Range	Channel A (Feed Water): 0 – 200 µS/cm Channel B (UP): 1 – 18.2 MΩ·cm
Pressure Range	0 – 1.6 MPa
Flow Meter Range	0.25 – 3 L/min
Temperature Range	0 – 100 °C
Tank Level	0 – 2 m Continuous display, or 0 – 3 m submersible sensor
Temperature Compensation Range	Automatic temperature compensation of readings Temperature compensation range: 5 – 45 °C to 25 °C

2 INSTALLATION

2.1 Installation Site Requirements

Item	Requirements
Installation space	≥ 50 cm × 50 cm total
Feed water	Purified water: EDI, RO, DI or distilled water
Working Temperature	5 – 45 °C
Power	100 – 240 VAC ± 10%, 1 KW (5 A)
Humidity	20 – 80%

2.2 Feed Water Requirements

Parameter	Value or Range
Pressure	0 – 15 psi (0 - 1 bar)
Conductivity at 25 °C	< 100 µS/cm
Water temperature	5 – 45 °C
Total Organic Carbon (TOC)	< 50 ppb

2.3 Feed Water Quality Measurements

Parameter	Test Method
Pressure	Use a pressure gauge
Conductivity at 25 °C	Use a conductivity meter
Water temperature	Use a thermometer
Total Organic Carbon (TOC)	Use a TOC analyzer

2.4 Installation

2.4.1 Preparation for Installation

2.4.1.1 Power Supply

Power supply must be properly grounded

2.4.1.2 Feed Water

Twenty liters of purified water

2.4.2 Items Included

Package			
No.	Name	Quantity	Note
1	VWR P Main System	1	
2	Control Console	1	Including a cable connected to main system
3	Accessories Bag	1	Including 8 mm PE tubing (8 meter), one roll of Teflon tape, a male connector, and a power adapter (with power cords)
4	Instruction Manual	1	On USB
5	Quick Installation Guide	1	

2.4.3 System Structure and Flow Diagrams

2.4.3.1 Main System Back Side View


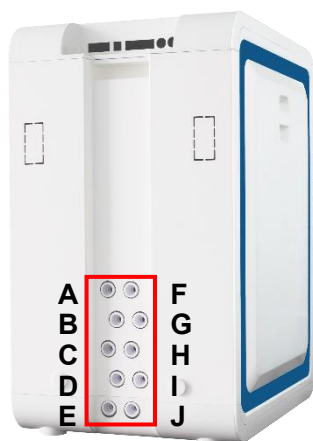


Illustration of water inlets and outlets


No.	Label	Description	Tubing Diameter	Connect to
A	HP Dispenser A	Spare	8 mm	Not used
B	HP Dispenser B	Spare	8 mm	Not used
C	EDI Drain	Spare	8 mm	Not used
D	Pure Water In	Pure water inlet	8 mm	Pure water tank
E	Product To Tank	Spare	8 mm	Not used
F	UP Dispenser A	For ultrapure water	8 mm	Dispenser
G	UP Dispenser B	For ultrapure water	8 mm	Dispenser
H	RO Drain	Spare	8 mm	Not used
I	Tap Water In	Spare	8 mm	Not used
J	Spare	Spare	8 mm	

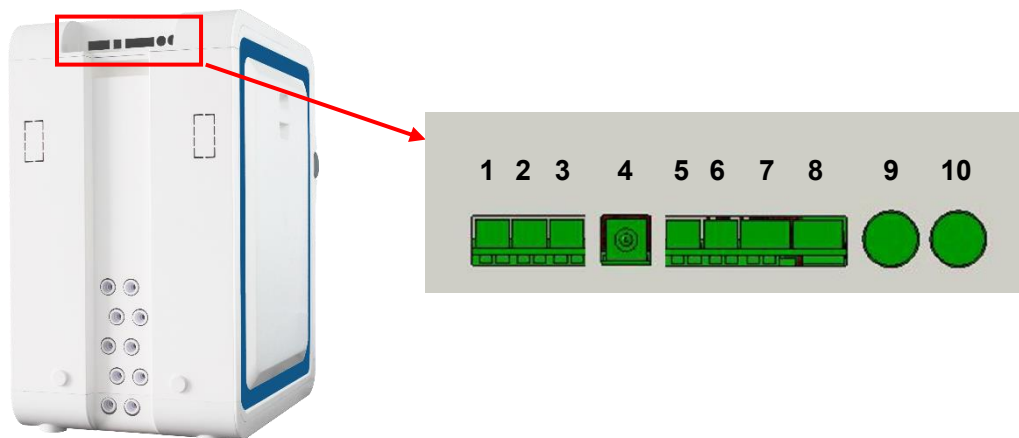

Description of electrical outlets

No.	Label	Description	Connect to
1	Feed Valve	For the feed solenoid valve control line	Feed solenoid valve (Optional)
2	Tank UV	For the sanitization module control line	Not in use
3	Feed Pump	For the booster pump control line	Not in use
4	Power	Power cord port	Power adapter
5	Prefilter Sensor	For backwash signal line	Not in use
6	Overflow Sensor	For overflow sensor signal line	Not in use
7	Leak Sensor	For leak sensor signal line	Leak sensor (Optional)
8	Level Sensor	For the pure water tank level sensor control line	Pure water tank level sensor (Optional)
9	CAN A	Data cable port	Dispenser or control console
10	CAN B	Data cable port	Dispenser or control console

2.4.3.2 VWR P External Connection Diagram

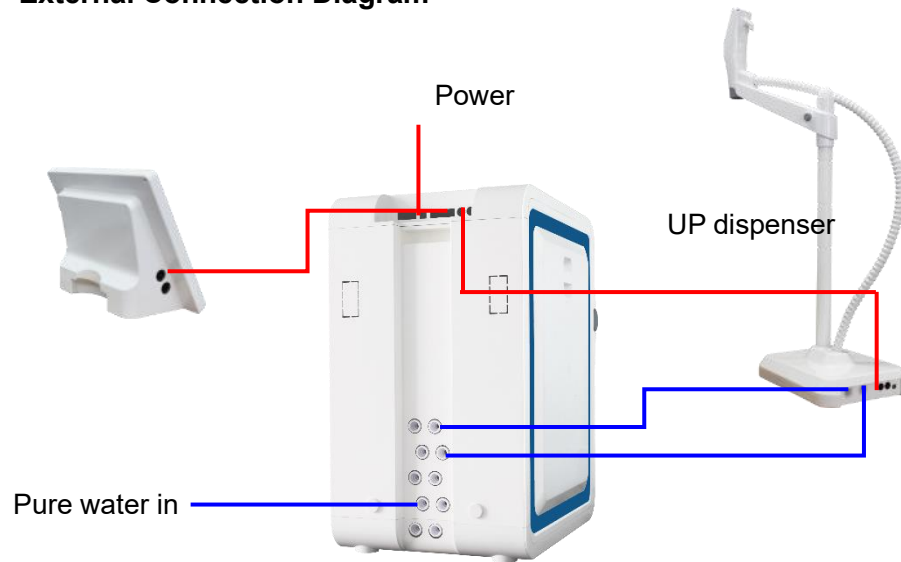

Red Line: Electrical loop


Blue Line: Water loop

2.4.4 Installing the Leak Protector (Optional)


2.4.4.1 Leak Protector Specifications

Catalog No.	171-1476
Operating voltage	24 VDC
Cable length	3 m to solenoid valve
Fitting	1/2 " female BSP (Inlet)-8 mm quick connector (Outlet)
Method of water leak detection	Conductivity
Operating temperature	5 – 45 °C
Maximum inlet pressure	6 bar (90 psi)

2.4.4.2 Illustration of Installation

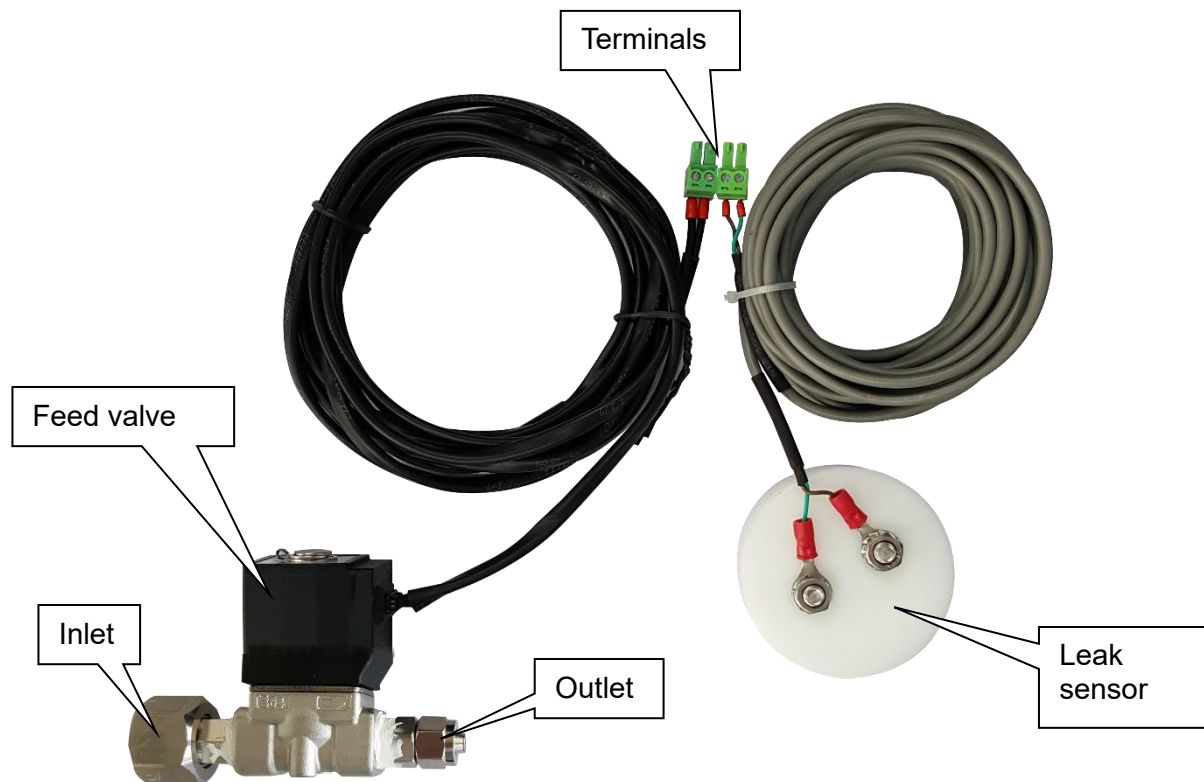

2.4.4.3 Installation

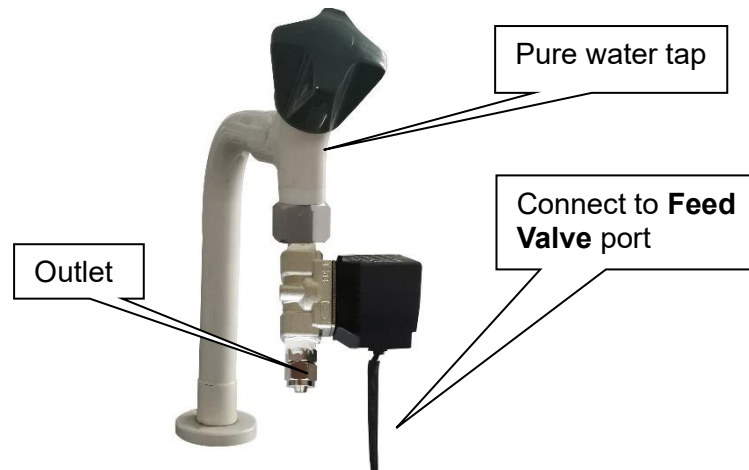

Follow the steps below if the diameter of the faucet is 1/2 inch male thread.

- As the picture shown above, connect the inlet of the feed valve to the faucet.
- Cut an appropriate length of 8 mm PE tubing. Insert one end of the tubing into the port **Pure Water In** on the main system and the other end into the outlet of the feed valve.
- Place the sensor metal probes side down at a lower point on the floor or near the water system.
- Plug the terminal of feed valve into the port **Feed Valve** on the back of the main system, and plug the terminal of leak sensor into the port **Leak Sensor**.

2.4.4.4 Instruction for use


The leak protector is activated automatically when the system powers up.

When it detects water, the system will shut the **Feed Valve** down, and switch to Standby mode automatically. A red icon  will appear on the screen and an alarm message box will pop up. The message "**Leak or Tank Overflow, all valves are closed, system could run after troubleshooting**" will disappear by tapping **Confirm** when the leak is resolved. Switch system from Standby to Ready mode manually so system will work again.

2.4.5 Installing the Dispenser

This water system offers two different dispensers: Dispenser 1 has an anchor base and stand, while Dispenser 3 is anchored on the system itself with no dispenser base.

2.4.5.1 Installing the Dispenser 1

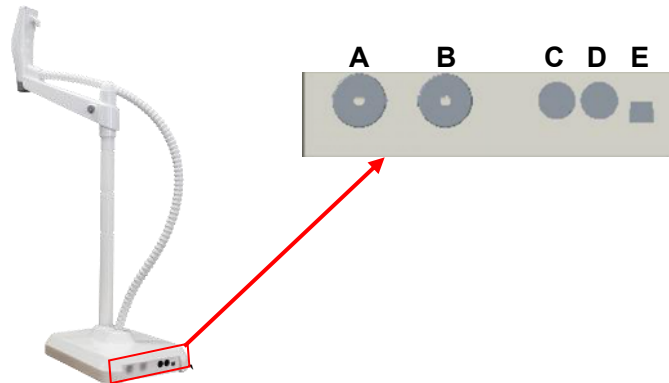

No.	Description	Connect to
A	8 mm tubing port	Main system
B	8 mm tubing port	Main system
C	Data cable port	Main system or power
D	Data cable port	Main system or power
E	Foot switch port	Foot switch (Optional)

2.4.5.1.2 Items included


Package			
No.	Name	Quantity	Note
1	Dispenser base	1	The handle is connected to the base with corrugated pipe.
2	Center post	1	With Screw
3	Crossbar	1	
4	Accessories bag	1	Including one roll of PVC tubing, a male barb connector and an Allen wrench.

2.4.5.1.3 Installing Dispenser Kit (171-1273)


- a) Open the package and take out the dispenser from the package. Set the dispenser base on a stable surface. This kit comes with a 3-meter link bundle.



- b) Install the dispenser center post to the base, then insert the crossbar on the center post, place the dispenser handle on the magnetic base of the dispenser arm.
- c) Take the tubing bundle out from the package, including two pieces of tubing and a cable. Install the tubing and cable as shown in the pictures below.

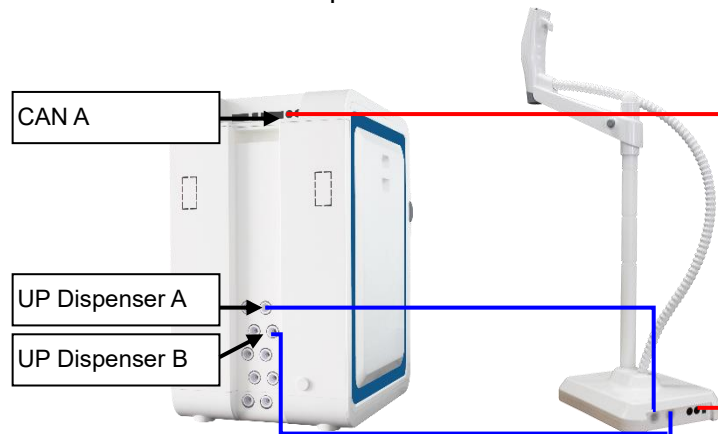

Connection of one dispenser for UP water

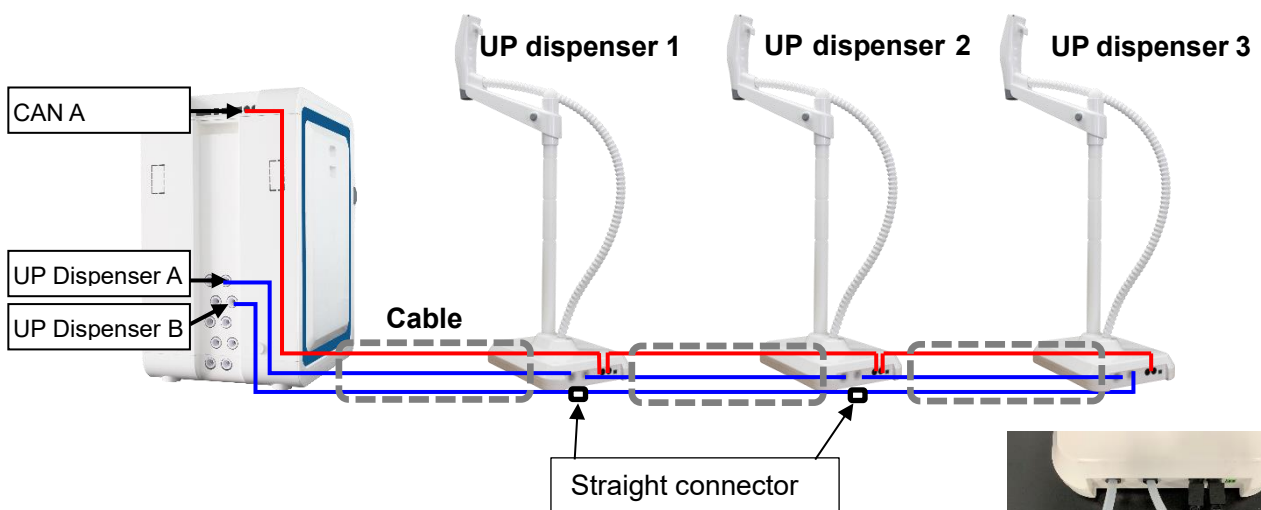

Connection of multiple dispensers for UP water



2.4.5.2 Dispenser 3 (Optional)



No.	Description	Connect to
1	Touch screen	
2	Dispense button	
3	Dispenser tubing connector	Main system
4	Dispenser cable connector	Main system

Package			
No.	Name	Quantity	Note
1	Dispenser	1	
2	Base kit	1	Including screws
3	Accessories bag	1	One dispenser adapter, two 8 mm tee connectors, two 8 mm quick connectors, eight 8 mm lock-clamps, two 6 mm lock-clamps, 1.5 m CAN cable

Installing the dispenser 3 (171-1275)


- a) Take out and assemble the base kit. The base is assembled differently on the left and right sides of the crossbar. Refer to the pictures below to install the base.



Install the base on the left side of the unit



Base kit

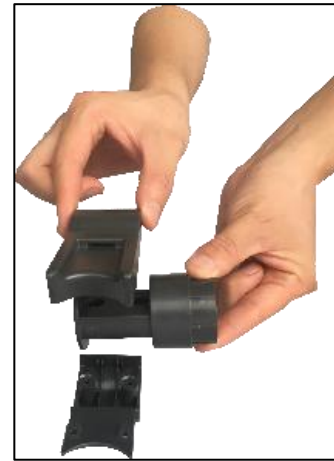

Install the base on the right side of the unit

- b) Open the side panels of the main system on both sides. Remove the screws as shown below.




The left side of the VWR G




The right side of the VWR G

- c) Remove the corresponding caps from the crossbar. Remove the fixed block from the cap.

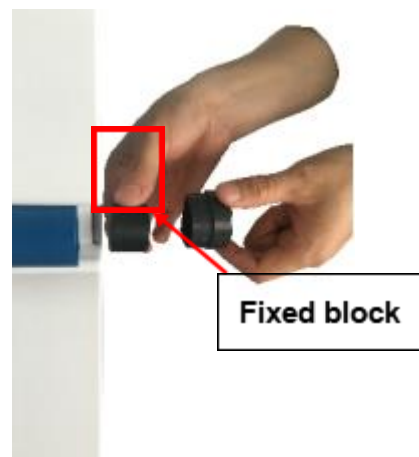


- d) Insert the holding block into the base and then insert the base into the crossbar. Fix it with the screw.

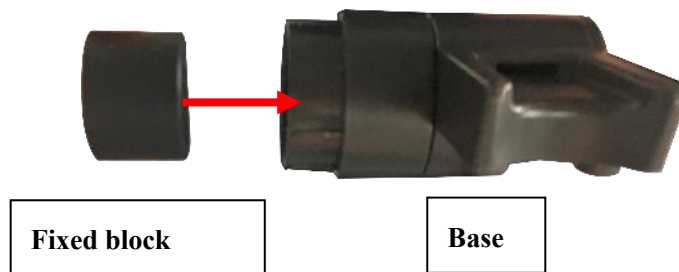


e) Install the tubing and cable.

Ultrapure dispenser: Insert the dispenser tubing connectors into the **UP Dispenser A** port and **UP Dispenser B** port on the back of the main system respectively. Then insert the dispenser cable connector into the **CAN A** or **CAN B** port on the back of the system.

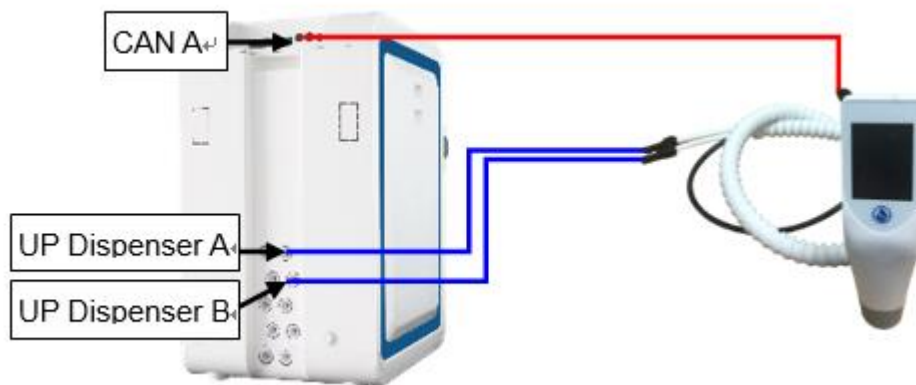


Pure dispenser: Insert the dispenser tubing connectors into the **HP Dispenser A** port and **HP Dispenser B** port on the back of the main system respectively. Then insert the dispenser cable connector into the **CAN A** or **CAN B** port on the back of the system.

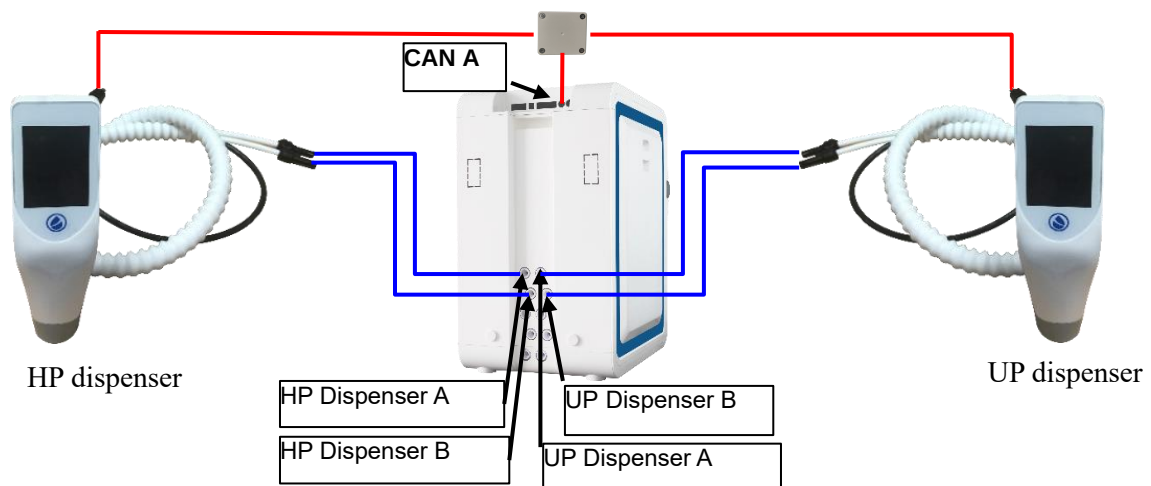


Pure and Ultrapure dispenser: Follow the picture above to connect the pure dispenser and the ultrapure dispenser using the dispenser adapter.

2.4.5.3 Installing the Wireless Dispenser Upgrade Kit

- a) For dispenser connected using wireless, need to use the upgrade kit (**171-1274**) to install the dispenser. The kit comes with a 10-meter link bundle and a power adapter
- b) Assemble the dispenser as that described above.
- a) Connect the tubing and power adapter as shown in the pictures below. Take the tubing out from the package, cut two pieces in proper length and insert them to the system and dispenser base as shown below.

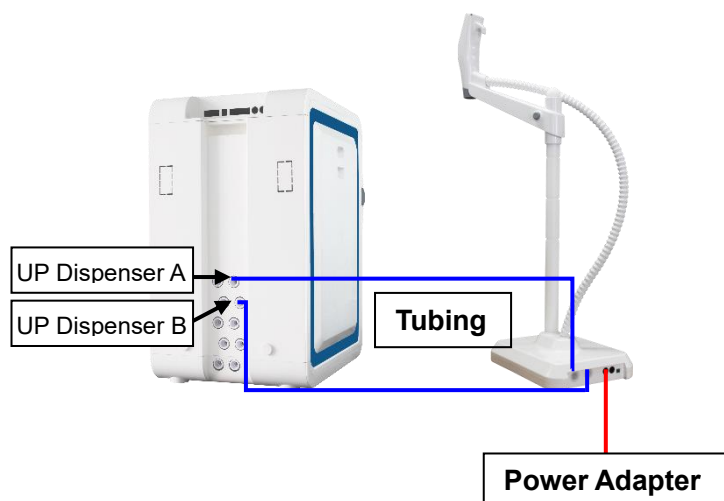

Connection of one dispenser for UP water

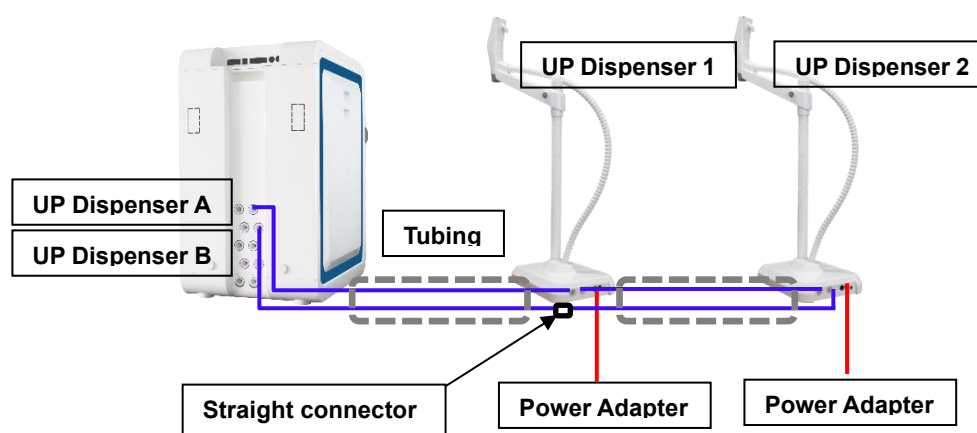


Connection of multiple dispensers for UP water



2.4.5.4 Installing the Dispenser Kit (RG0P0U001) and the Wireless Dispenser Upgrade Kit

- For one dispenser connected via wire and another one dispenser connected using wireless, need to use the upgrade kit (**RG0P0U012**) to install the dispenser. The kit comes with a 10-meter link bundle and a power adapter
- Assemble the dispenser as that described above.
- Connect the tubing and power adapter as shown in the pictures below. Take the tubing out from the package, cut three pieces in proper length and insert them to the system and dispenser base as shown below.

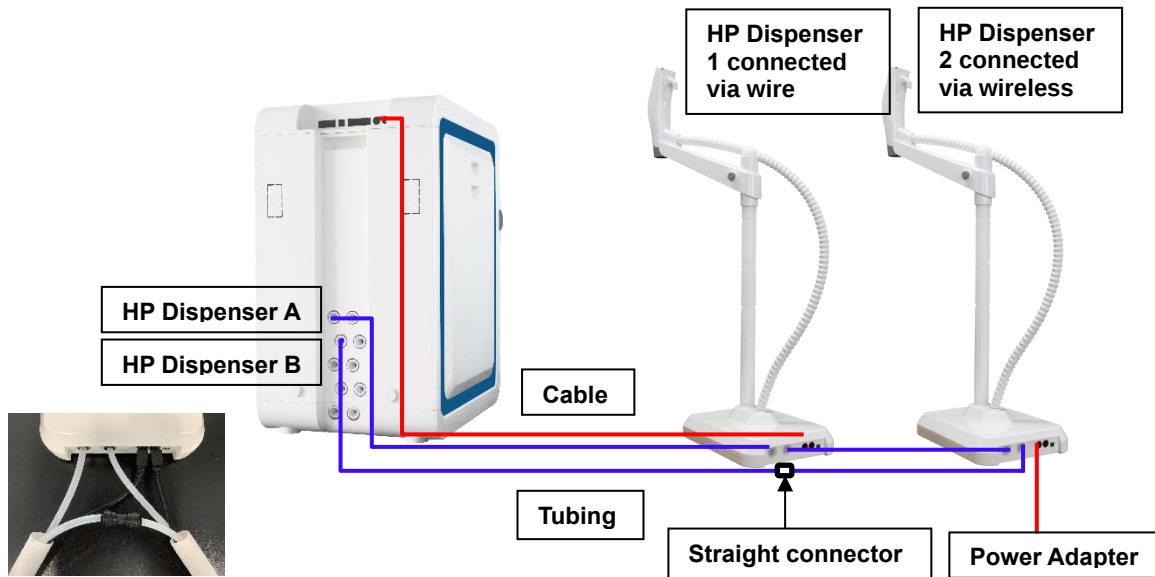


2.4.6 Installing the Water Tubing



2.4.6.1 Connecting the Inlet Feed Tubing

- Remove the stopper on the pure water inlet on the system. Take the 8 mm PE tubing from the accessories bag and cut a piece of the tubing at proper length for feed.
- Insert one end of the 8 mm PE tubing into the water inlet on the system (Mark as **Pure Water In**). Gently pull the tubing outwards to check if the connection is secure.
- Connect the other end of the tubing to a feed water tank; ensure the tubing inlet is immersed to the bottom of the tank.

2.4.7 Installing External Cable Connection

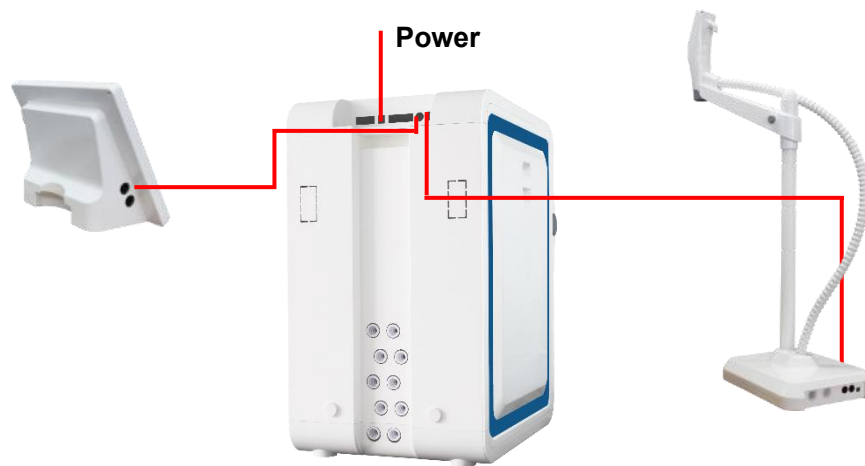

Make sure all power cords and signal cables are connected to their corresponding ports.

3 SYSTEM STARTUP AND OPERATION

3.1 Check List Prior to System Startup

Prior to system startup, use the table below to make sure all parts have been installed and connected, and quality of pure water meets minimal requirements.

Check List	
Feed water quality Pure water: Conductivity < 100 µS/cm (at 25 °C), Total Organic Carbon (TOC) < 50 ppb, Water temperature: 5 – 45 °C, Pressure: 0 – 15 psi (0 – 1 bar)	
Feed water connected	
Control console connected	
Dispenser installed and connected	
Power adapter plugged in	

Note: Cartridges will be installed AFTER the system power-up. DO NOT install them now.




3.2 System Initial Startup

Connect the power adapter to the system to power on the system, set-up menus will show up in order at the initial power-up in a few seconds.

3.2.1 Language Setup

	The language setup menu will show up in the first page. English is default. You may choose the language desired and then tap Next to set up Time & Date.
---	---


3.2.2 Time & Date Setup

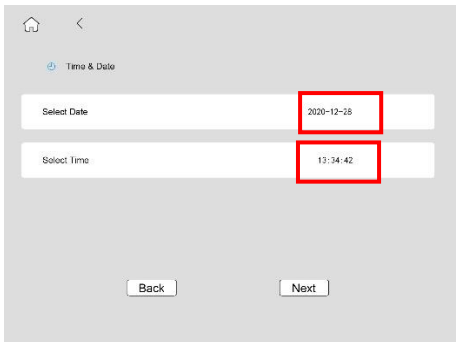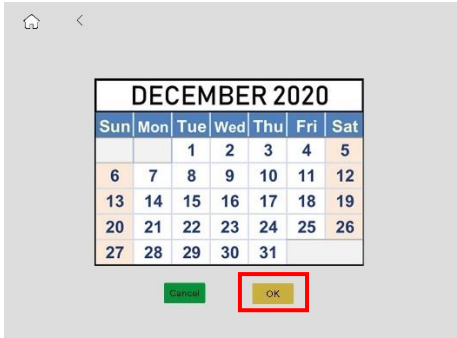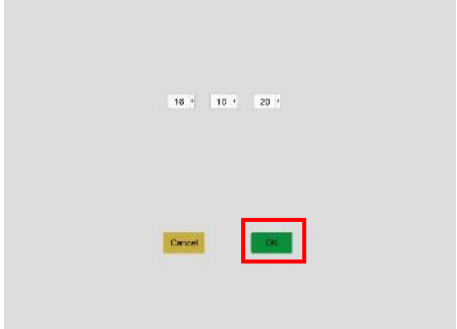	Tap red box to enter set-up menu.
	Select locate date and then tap OK to confirm setting.
	Select locate time and then tap OK to confirm setting. Tap "Next" to enter Set Tank menu.



Note: If the power outage exceeds 3 days, a pop-up box will appear to reset the system Time & Date.

3.2.3 Tank Setup

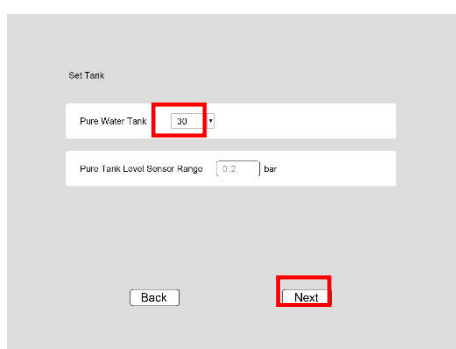


Choose the actual capacity of pure water tank if a tank is installed. Tap **Next** to enter **Other Devices** setup menu.



If the tank is not from VWR, choose **Custom**. Setup the height and volume of the feed water tank. Then enter the Tank Level Sensor Range as pressure. In Bar. Default sensor from VWR is 0.2 bar (Submarine level sensor is 0.3 bar).


3.2.4 Other Devices Setup

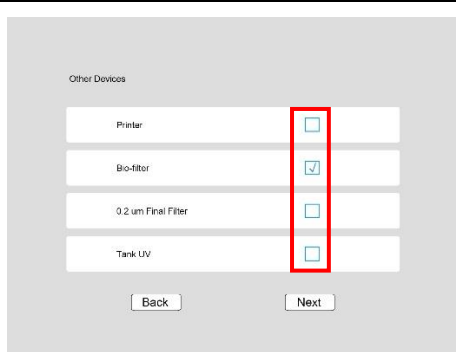

Put “√” in the box of **Printer** if a printer is used.
Put “√” in the box of **Bio-filter** if a Bio-filter is installed.
Put “√” in the box of **0.2 μm Final Filter** if a 0.2 μm final filter is installed.
Put “√” in the box of **Tank UV** if a tank sanitization module is installed.






	Printers supported: Epson Dot matrix VWR recommends Epson's TM-U330D dot matrix printer. Other USB based dot matrix printers may not be supported. After each dispense, the printer will automatically print out dispense time, water quality and other displayed parameters.
---	--

3.2.5 Connectivity Setup

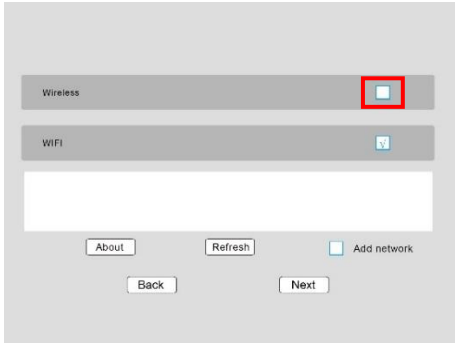	Select Wireless if wireless dispenser is opted. Check “√” in the red box. Tap “Next” to enter consumables installation interface.
---	---


3.2.6 Install Consumables

3.2.6.1 Register Consumables (Not including cartridges)

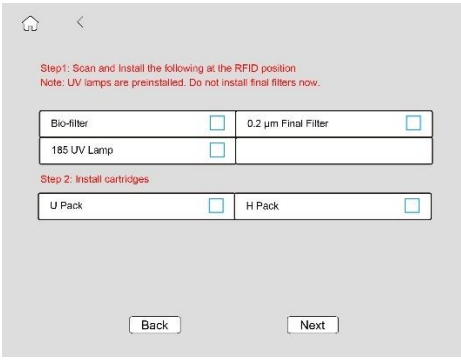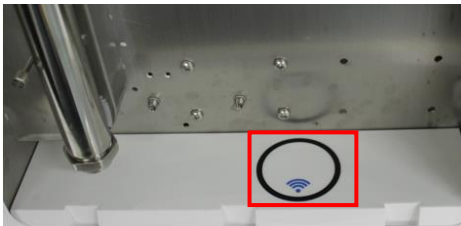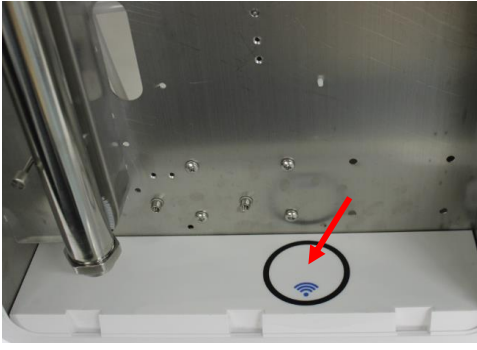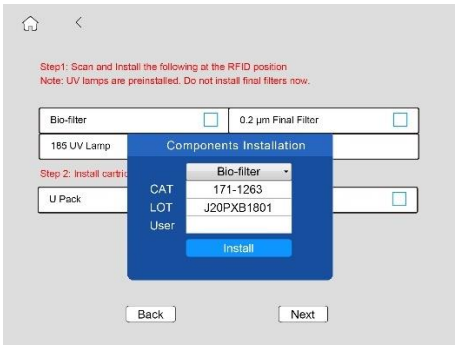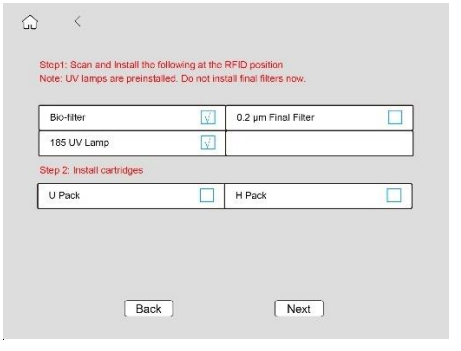 	Follow instructions on the screen, scan RFID tags for the 185 nm UV lamp and final filter at the AC pack position with an antenna sign () on the system. Details are shown below Note: 1. RFID tags are either inside consumable packages, or for preinstalled UV lamps etc, inside USB key package. 2. For traceability, your User Name is required to install/register consumables. For a new installation, enter system manager's User Name.
	Remove the right side panel to find the RFID reader position () as shown left.
	A pop-up dialog will be shown as left after every scan, enter manager's User Name and tap Install to register the consumables
	Follow steps below to install cartridges after finishing the scan.

3.2.6.2 Install and Register Cartridges


VWR P system has two packs, 1 H Pack and 1 U Pack. Each cartridge's position is labelled on the system. Install a cartridge to its corresponding position.

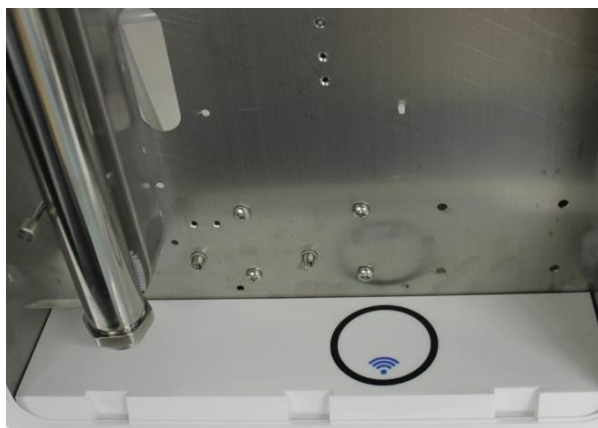

Left Side




Right Side

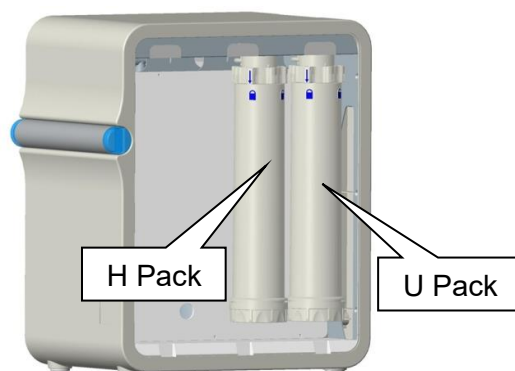

- a) Remove the right side panel from the system as shown below.




- b) Take out the cartridges from their packaging, respectively. Remove caps on the cartridge's inlet and outlet on the system.
- c) Follow the instruction below to install cartridges to their corresponding positions.

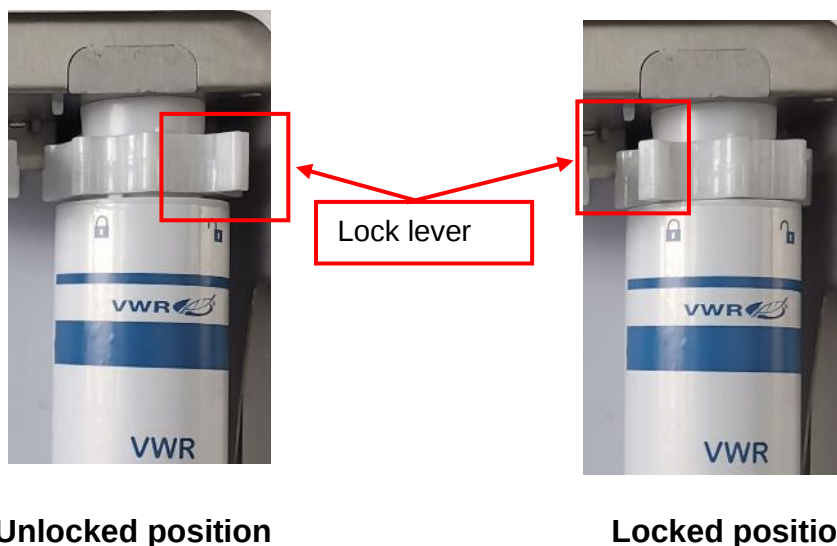

d) Wet the O-rings on the H Pack cartridge with pure water.



- e) Turn the lever on the cartridge adapter to the right side at the unlock position  as shown in the photo
- f) Align the cartridge label facing out. Insert the up end first into the opening on the cartridge adapter, then turn the lever handle to the left  to lock the cartridge.
- g) A pop-up dialog will be shown as below, tap **Install** to complete the installation.

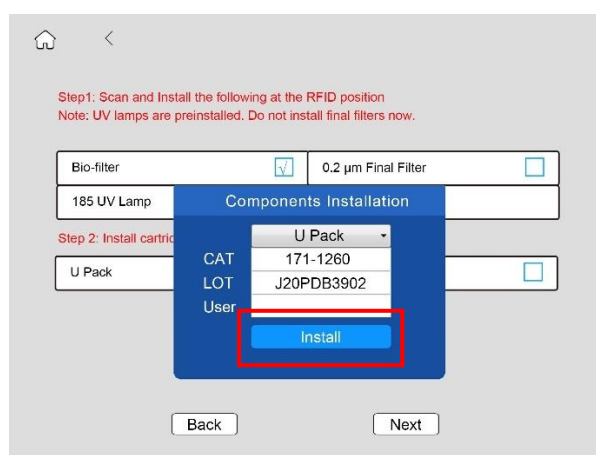


Follow the same procedure to install other cartridges.



Note!

The control console will display “RFID Error!”, and the alarm icon  if a pack is installed incorrectly. Reinstall it to the correct position.

- h) Make sure all consumables have been registered and tap **Next** to enter Dispenser setup menu.

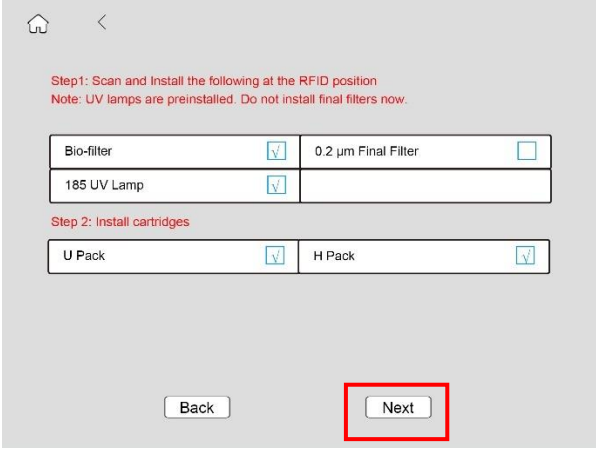

Step 1: Scan and Install the following at the RFID position
Note: UV lamps are preinstalled. Do not install final filters now.


Bio-filter	☒	0.2 µm Final Filter	☐
185 UV Lamp	☒		

Step 2: Install cartridges

U Pack	☒	H Pack	☒
--------	-------------------------------------	--------	-------------------------------------

Back Next

3.2.7 Dispenser Setup

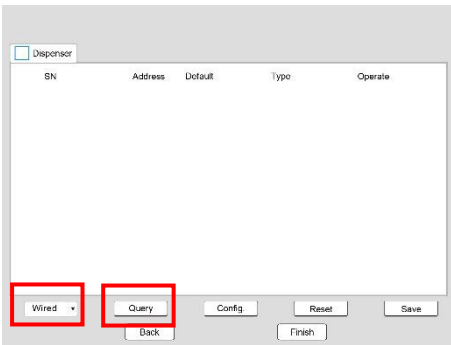

Dispenser

SN	Address	Default	Type	Operate
----	---------	---------	------	---------

Wired Query Config Reset Save Back Finish


Select connection mode for the dispenser from the pull-down menu at the left corner:

Select **Wireless** for wireless connection.

Select **Wired** for the wired connection.

Tap **Query**, the system will automatically search for the dispensers that have been installed.

Note: The dispensers may have prior registered to other systems, causing problems in installation if record in the dispenser is not cleared. Press Query will clear such record.

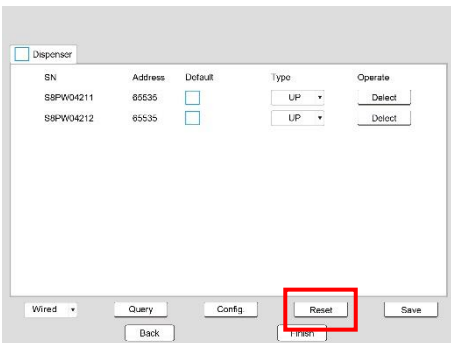


Dispenser

SN	Address	Default	Type	Operate
88PW04211	85535	☒	UP	Detect
88PW04212	85535	☒	UP	Detect


Wired Query Config Reset Save Back Finish

Tap **Reset**, system will automatically reset the address.

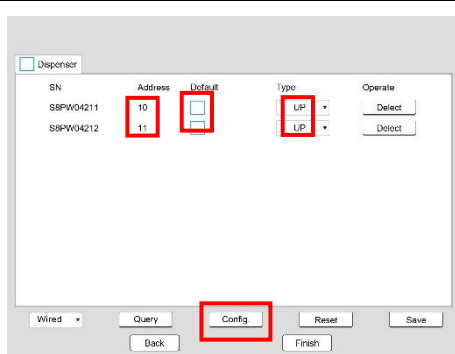

Tap **Config**, system will automatically configure the dispensers. You need to choose a dispenser as default for each type of water. Water will come out from the default dispenser when you tap dispense button on the control console. Choose the Type of the dispenser.

Note: An Address shown between 10 and 20 means the dispenser setup is successful.

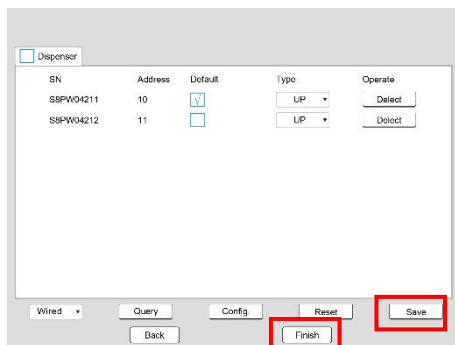

Tap **Save** to save the configuration.

Confirm the configuration by tap **Finish**.




System will enter the home screen automatically.

Now the system is ready to operate

3.3 System Setup

Swipe a finger down the top section of the home screen to see the drop-down bar. Make sure the system is in Ready mode. If not, toggle between Standby  and Ready  modes using the drop-down bar to switch the system from Standby to Ready.

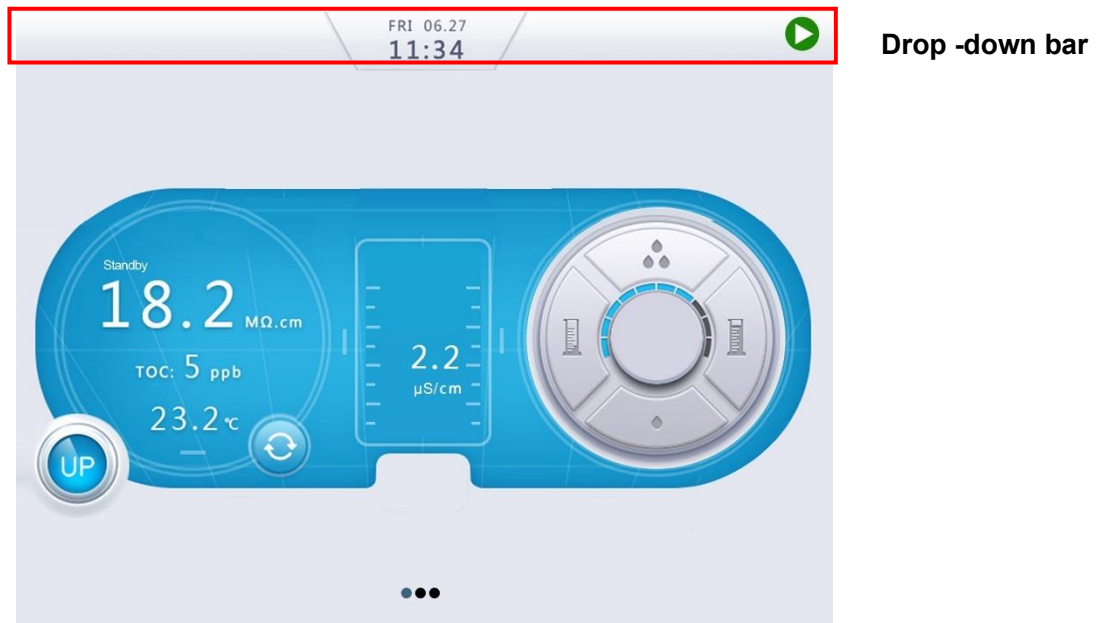

Note!

Alarms  icon and / or alerts  icon may present in the middle of the home screen when setting-up the system. It's normal. Just follow this user manual to finish the rest of operations.

3.3.1 Flushing the H Pack and U Pack.

- Remove stopper from the UP dispenser outlet.
- Screw on the barbed adapter (in the accessory bag) to the dispenser outlet. Connect the PVC tubing inside the accessory bag; insert one end to the adapter and put the other end into the drain.
- Press the dispense button to dispense water. Gas inside the system will be released out of the system.
- Wait till no more gas bubbles coming out of the tubing. Press Dispense button again to stop degassing. Screw off the adapter.


3.3.2 Installing the Final Filter

- Open the attached final filter package.
- Screw the final filter onto the remote dispenser until finger tight (no leak at dispensing). Do not over tighten it as it may damage threads.
- Press the dispense button to flush out gas in the filter.
- Press the dispense button again to stop dispensing.

This completes system hardware installation.

3.4 Information Menu

Control Console Main Page

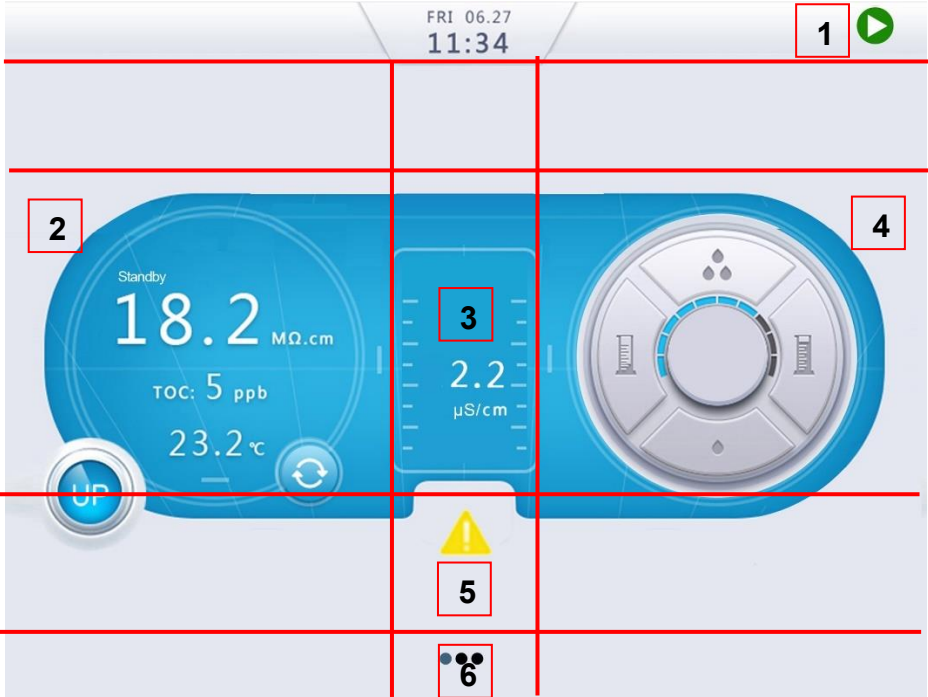

The screenshot shows the Control Console Main Page interface. At the top, it displays 'FRI 06.27' and '11:34'. A green play button icon is in the top right corner. The main display area is divided into several sections: a large blue circular gauge on the left showing 'Standby' and '18.2 MΩ.cm', with 'TOC: 5 ppb' and '23.2 °C' below it; a central section showing '2.2 μS/cm' with a yellow warning triangle icon below it; and a large blue circular gauge on the right. A bar graph is located below the central section. At the bottom, there are three small circular icons: a blue one with '12.5 L', a grey one with '0.0 L', and a grey one with '0.0 L'.

Section 1: Drop-down menu, including the located time and system operating status.

-  Standby
-  Ready

Section 2: UP water quality, TOC, temperature, operation mode and UP water recirculation icon.

Section 3: Feed water resistivity.


Section 4: UP water dispensing set up

-  Volumetric dispensing button, decrease the volume, down to 0.1 L.
-  Volumetric dispensing button, increase the volume, up to 99.9 L.
-  Dispense rate down button, decrease the rate.
-  Dispense rate up button, increase the rate up to 2 L/min.
-  Bar graph shows the relative dispensing rate.
-  Dispensing button, press the button to dispense water after setting up the volume. If it displays 0.0 L, it's non-volumetric dispensing.

Section 5: Alarms  icon and alerts  icon present in the middle to display current alarms or/and alerts.

Section 6:  Indicates multiple screens. Swipe left & right anywhere in blank area on screen to switch pages

Dispenser

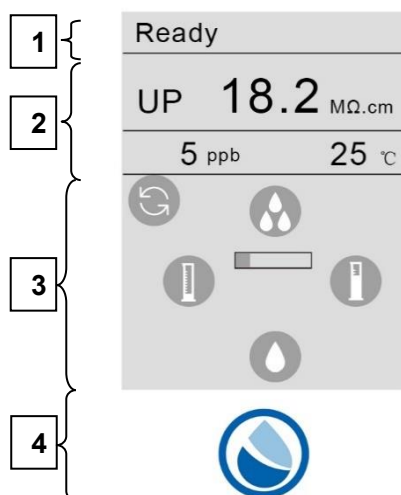

Section 1: System operating status (including Ready, Dispense, Standby, Pressure Relief and Recirculation) and tank level (if a level sensor is installed).

Section 2: Water quality parameters: water resistivity, TOC and temperature.

Section 3: Dispensing set up



Recirculation button



Volumetric dispensing button, decrease the volume, down to 0.1 L.



Volumetric dispensing button, increase the volume, up to 99.9 L.



Dispense rate down button, decrease the rate.




Dispense rate up button, increase the rate up to 2 L/min.



Bar graph shows the relative dispensing rate.


Section 4:  Dispensing button

Control Console - More Pages

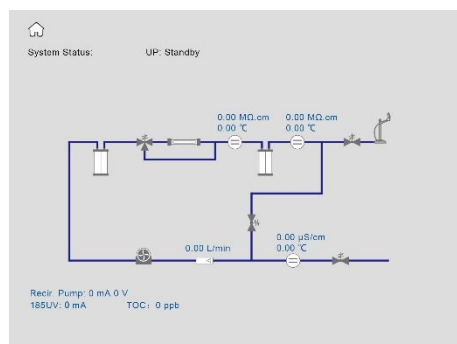

Information Menu I

Swipe the screen to the right. This page shows system operation status and parameters in a flowchart format.

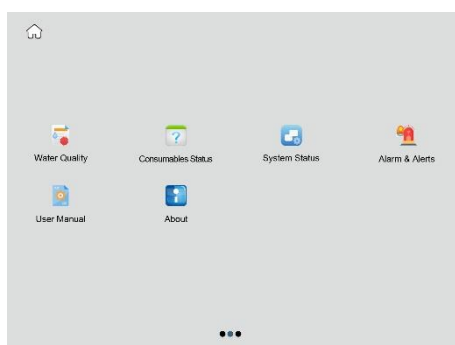

Information Menu II

Swipe the screen to page 2. Page 2 is an Information Menu which contains information about system status and history.

Puro Tank Level	0%	0 L	Feed	2.2 µS/cm	23.2 °C
UP	18.2 MΩ.cm	23.2 °C	TOC	5 ppb	
UP Disp. Rate	0.0 L/min	/			

Water Quality displays measured values about water.

- **UP, TOC and Temp.:** UP resistivity, TOC and temperature are monitored.
- **Feed and Disp. Rate:** Feed water conductivity and UP water dispense rate are monitored.

H Pack	DI	Installation Date 2020-12-28	CAT NO.: 171-1267	
		Replace in 360 Days	LOT NO.: J20PDB3902	
U Pack	DL	Installation Date 2020-12-28	CAT NO.: 171-1260	
		Replace in 360 Days	LOT NO.: J20PDB3902	
185 UV	OH	Installation Date 2020-12-28	CAT NO.: 171-1263	
		Replace in 720 Days	LOT NO.: UV200101B	
Puro filter		Installation Date 2020-12-28	CAT NO.: 171-1265	
		Replace in 3 Days	LOT NO.: J20PXB1801	

Consumables Status

Displays details of consumables installed, including processed volume/time, installation date, remaining life, catalog number and lot number.


You can order consumables by clicking  on this page.



Means the component working normally.



Means the component needing to be replaced.

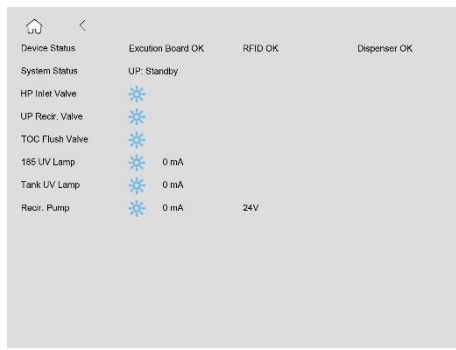

System Status provides information about the components operation status.



Means the component is functional. The screen displays operating current or voltage for UV lamp and pump.



Means the component is off.



Alarms & Alerts:

Alarms sign appears when there is a technical issue. Some **red alarms** stop the system from operating to protect the system.

Alerts appear when a consumable needs to be replaced or a non-critical event occurs.





About






Including: catalog number, serial number, production date, installation date and software version.






You may need to provide some or all of the information when requesting service.




3.5 Maintenance Menu

Notes: Maintenance menu is password protected for system administrators. The log-in information is in the USB stick shipped with the system.

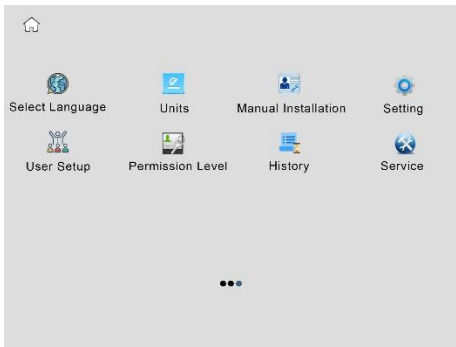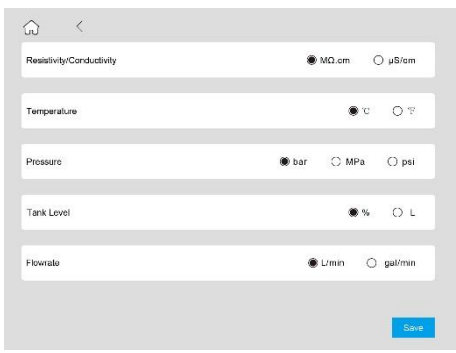	Maintenance Menu Swipe left on the Screen to page 3 in Standby mode. Page 3 is Maintenance Menu which contains parameter setup, RO cleaning, etc. Maintenance Menu is a manager authorized operation interface. When clicking the above options, a password login popup will appear. Enter the user name and password in the popup and click "login" to do the operation. The initial manager user name is "manager", please refer to the "password.txt" file in the USB flash drive for the password.
	Select Language Choose your working language desired.
	Units Change unit of display, including water quality, temperature, pressure and dispense rate.
	Manual installation Users can manually fill in the CAT number and LOT number of consumables in the "manual installation" interface. Click "install" in the next circled box to manually enter information without scanning the RFID. Pay attention to letter case when entering.

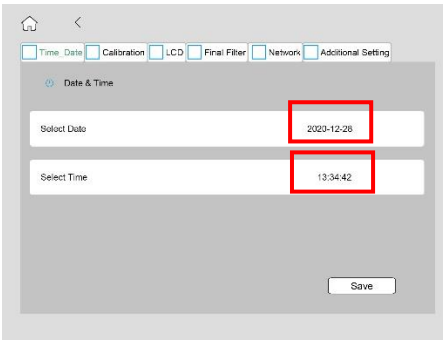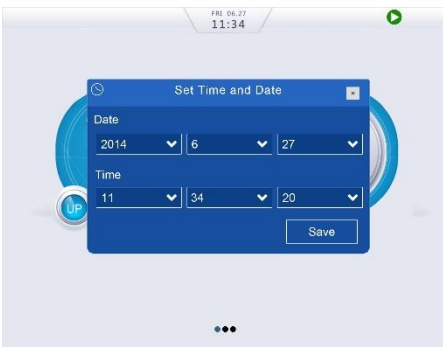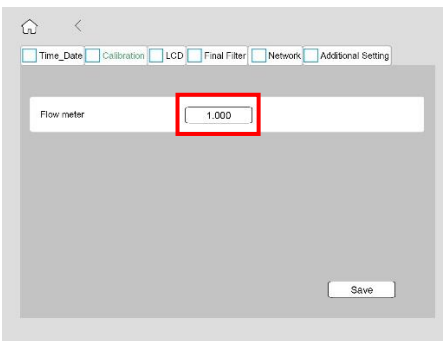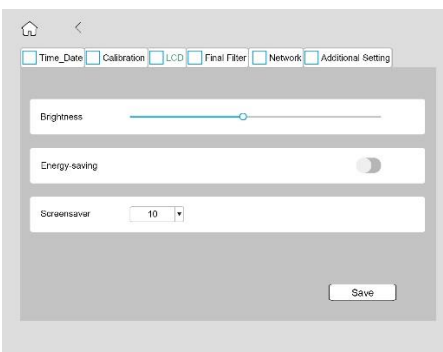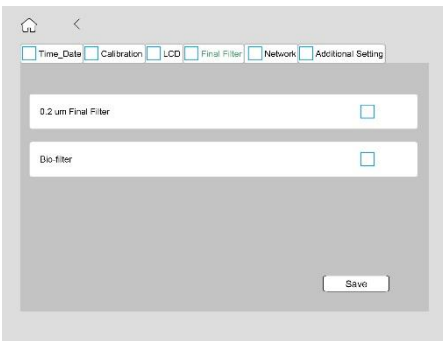	Setting Including time & date, calibration, audio, LCD and additional setting. Time & Date Choose the locate time and date.
	Note: If the power outage exceeds 3 days, a pop-up box will appear to reset the system Time & Date.
	Calibration The calibration menu is used to calibrate water dispense rate from the dispenser.
	LCD Change screen brightness and choose energy-saving if needed. Set the time for screensaver.
	Final Filter Put “√” in the box of 0.2 µm Final Filter if a 0.2 µm final filter will be installed. Put “√” in the box of Bio-filter if a Bio-filter will be installed. Note: In this interface, restarting the system is required after tapping Save, or the setting will not take effect prior to a Restart.

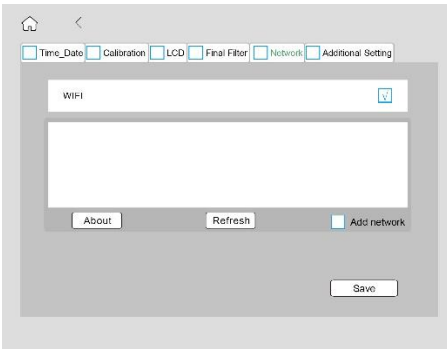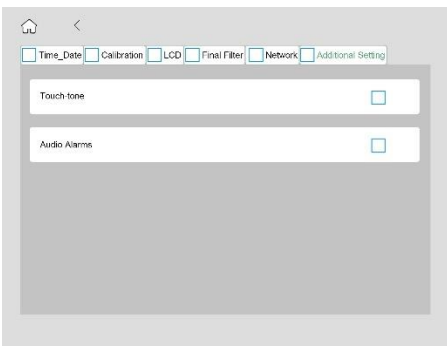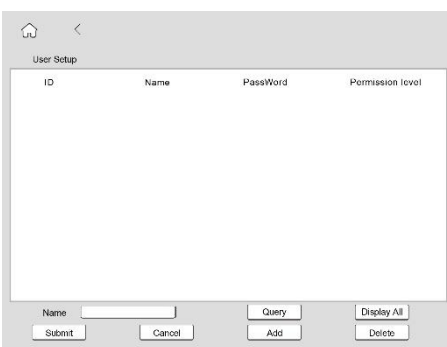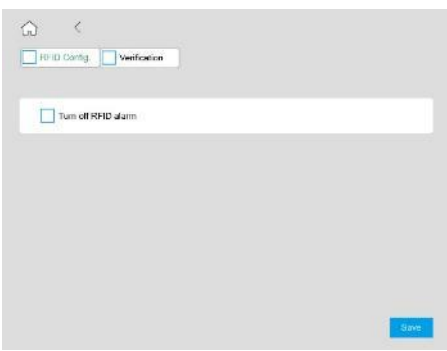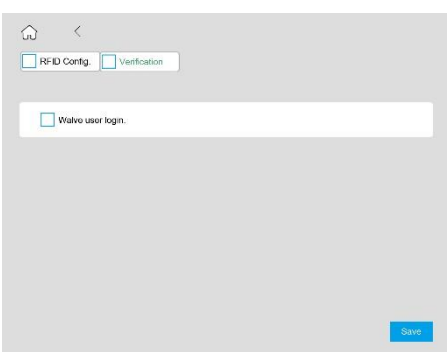	Network In this interface, you can connect the system to Wi-Fi network, make sure there is a “√” in the box, and then tap Refresh to search for your accessible network name (SSID), enter its password. Tap Save to complete the setting. System restart is required after tapping Save. A clip icon  after network name (SSID) means the system is connected to this network at present. Tap About to view IP address for the network if needed.
	Additional Setting Activate Touch-tone, the control console will make a sound when you touch the screen. Also, if activate Audio Alarms, the control console will make beeping sounds when alarm(s) appear(s).
	User Setup Manager can add new account to operate the system. Tap Add, then enter a new username and password under Name and Password to add a new account. Select the user level under Permission. User is for general user, dispense water only. Manager means a user has the right to add new accounts. Then tap Submit to complete the setup.
	Permission Level Including RFID Config. and Verification. RFID Config. Turn on and off RFID alarm. System wouldn't show alarms about consumable status and type.
	Verification Turn on and off user verification. Users don't need to enter a name when replacing consumables if it is checked. It's off by default.

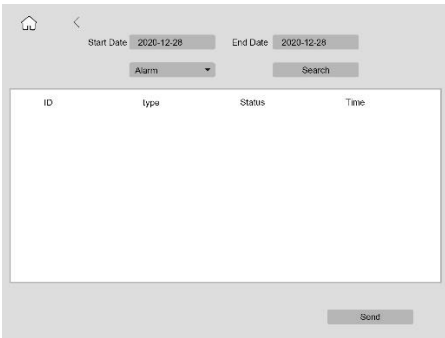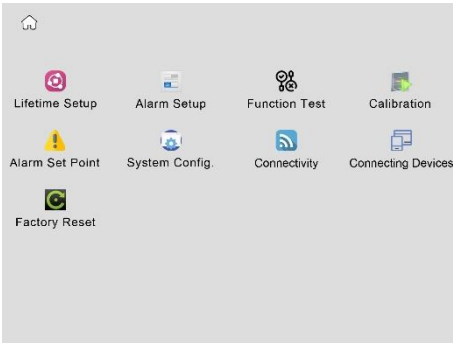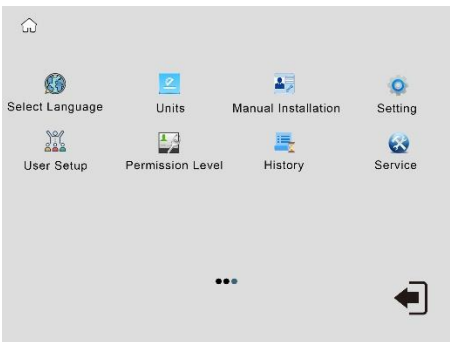	History Provide historical information related to your system. The system data can be filtered by date and the selected data can be exported.
	Service Service menu is an engineer authorized operation interface.
	After the maintenance operation, click the icon  in the lower right corner to exit the login.

4 DISPENSING WATER

Water can be dispensed from the dispenser handle either by pressing the dispense button on the handle or from the main control console.

4.1 Dispensing Water

Tap recirculation icon  on the main console or the dispenser to start recirculation. When the resistivity reaches 18.2 MΩ.cm, press the dispense icon  on the monitor or  on the dispenser to dispense UP water. Press the dispense icon again to stop dispensing.

Note: Dispensers are interlocked. Only one dispenser can be used at a time. Other dispensers will display “In Use” on the dispenser.

4.2 Adjusting Dispense Rate


To change the dispense rate, tap the icon  or  on home screen or dispenser monitor.

4.3 Volumetric Dispensing

	Setting up is the same on the dispenser or the control console. Tap recirculation icon  on the main page to enter Recirculation mode until the resistivity reaches 18.2 MΩ.cm.
	Tap the right cylinder icon to increase or the left cylinder icon to decrease the volume. You can set the volume between 0.1 and 99.9 liters.
	Tap the middle dispensing icon  to start dispensing. Volume dispensed is displayed at the start button position. Volumetric dispensing will stop automatically after completion.

5 MAINTENANCE

5.1 Basic Alerts

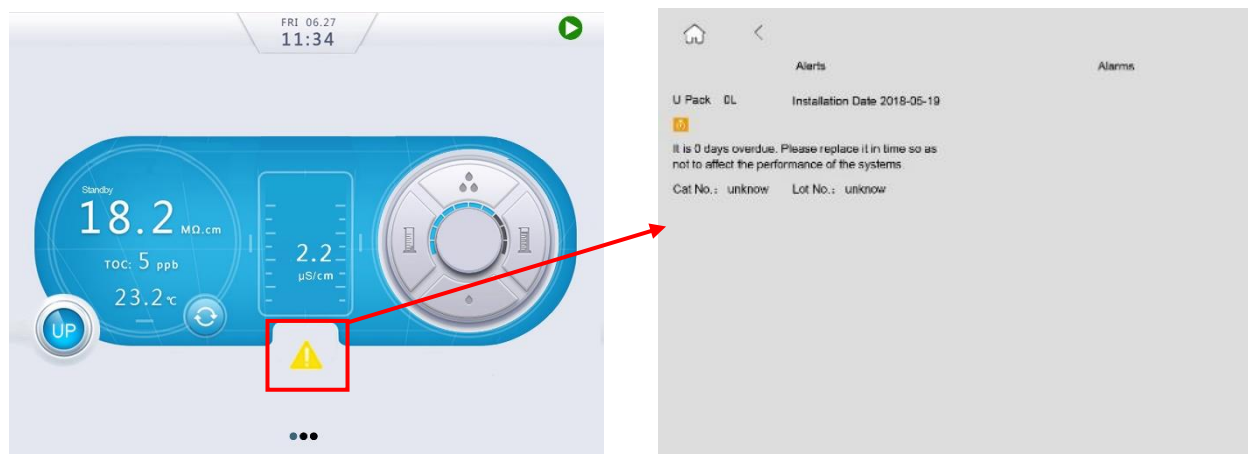

Message	Solution
H-Pack 3000 L Installation Date: XXXX-XX-XX It is XX days overdue. Please replace it in time so as not to affect the performance of the system. CAT NO: 171-1257 LOT NO: S6PDC0101	Exchange the H Pack and reset.
U-Pack 3000 L Installation Date: XXXX-XX-XX It is XX days overdue. Please replace it in time so as not to affect the performance of the system. CAT NO: 171-1258 LOT NO: S6PDC0101	Exchange the U Pack and reset.
185 UV 500 H Installation Date: XXXX-XX-XX It is XX days overdue. Please replace it in time so as not to affect the performance of the system. CAT NO: 171-1283 LOT NO: S6PDC0101	Exchange the 185 nm UV lamp and reset.
Final Filter A 360 D Installation Date: XXXX-XX-XX It is XX days overdue. Please replace it in time so as not to affect the performance of the system. CAT NO: 171-1262 LOT NO: S6PDC0101	Exchange the final filter and reset.

5.2 Basic Troubleshooting

If you encounter any electronic error, **restart the system first before doing any troubleshooting.**

Alarm / Phenomenon	Solution
Cannot dispense water from dispenser. It shows "In Use"	• Another dispenser connected to the same systems is being used. System allows one dispenser working at a time. Wait till the other dispenser stopped dispensing.
Feed Conductivity > upper limit	• Feed water quality is too poor • Contact a VWR professional engineer.
UP product resistivity < SP	• UP product resistivity is below preset value. • Exchange both H Pack and U Pack cartridges or contact a VWR professional engineer.
Low Pure Water Tank Level	• Pure water tank level is below preset value (10%). • Stop dispensing until water tank is refilled.
Check 185 UV	• The 185 nm UV lamp may be burnt. • Exchange the 185 nm UV lamp or contact a VWR professional engineer.
Check Tank UV	• The tank UV lamp may be burnt. • Exchange the tank UV lamp or contact a VWR professional engineer.
H-Pack Error	• Reinstall the H Pack. • Check the RFID chip. • Contact a VWR professional engineer.
U-Pack Error	• Reinstall the U Pack. • Check the RFID chip. • Contact a VWR professional engineer.
H-Pack Not Detected	• Check cable connections. • Reinstall the H Pack. • Contact a VWR professional engineer.
U-Pack Not Detected	• Check cable connections. • Reinstall the U Pack. • Contact a VWR professional engineer.
UP Product Temperature>45°C	• Contact a VWR professional engineer.
UP Product Temperature<5°C	• Contact a VWR professional engineer.
TOC Sensor Temperature>45°C	• Contact a VWR professional engineer.
TOC Sensor Temperature<5°C	• Contact a VWR professional engineer.

TOC Feed Resistivity<15MΩ.cm	• Contact a VWR professional engineer.
Cannot dispense water from dispenser. It shows “In Use”	• Another dispenser connected to the same systems is being used. System allows one dispenser working at a time. Wait till the other dispenser stopped dispensing.
System leakage	• The system detects a leakage. • Shut off power and water supply. • Remove system side panels, turn on water supply and check leaking points. • Reconnect or replace leaking parts.

5.3 Replacing Consumables

Consumable life, in particular cartridges, is heavily dependent upon feed water quality and amount of water used. Below is a general guideline of frequency to replace consumables.

Consumables		Replacement Frequency	Performance Indicator
Dual UV lamp		24 months	Increase in TOC
Final Filter	Remove bacteria and particles	12 months	Reduce in flow rate
		When needed	Flow rate is less than 1 liter per minute
H Pack & U Pack	Polishing pure water	When needed or 12 months	Reduce in resistivity of ultrapure water
Bio-filter	Removes pyrogen, RNAses and DNAses	3 months	Increased pyrogen, RNAses and DNAses levels

System administrator can set Permission for cartridge replacement in Maintenance Menu. System default is OFF. If the **Permission** is **On**, User need to enter a name to verify the installation. Only authorized Users registered in the system are allowed to install or replace consumable.

The following consumables need to be registered to the system on the RFID tag reader at installation.

Scan the RFID chip of the installed 185 nm UV lamp and final filter in their corresponding box to register the consumables. Tap **Install** when a pop-up dialog is appearing if the **Verification** is **Off**.



5.3.1 Replacing the VWR H Pack and VWR U Pack cartridges

- Switch the system to Standby mode, then remove the system right side panel.
- Turn the lever on the cartridge adapter to right to the unlock position, pull the cartridge downwards to disengage the cartridge from its adapter. Install the new one as described in 2.4.4.
- After installation, do not install the system side panel. Power up and degas the system, and check any leakage. If leaks, reseal the cartridge, then test again.
- If there is no leakage, install back the right side panel.
- Remove the final filter before replacing VWR U Pack. Release pressure by dispensing some water. Install the final filter after installing the cartridge.
- Tap **Install** on the pop-up dialog to reset the cartridge life.

5.3.2 Replacing the UP UV Lamp (185 nm)




Warning!

Ultraviolet (UV) radiation is harmful to the eyes and skin. Do not observe the lamp directly when it is illuminated. This system is equipped with a lamp cover to prevent UV light leakage. This cover must be on ALL TIMES when a UV lamp is installed.



Caution!

Keep the UV lamp straight in and out of the stainless-steel chamber during its installation to avoid any action that could cause the lamp to break.

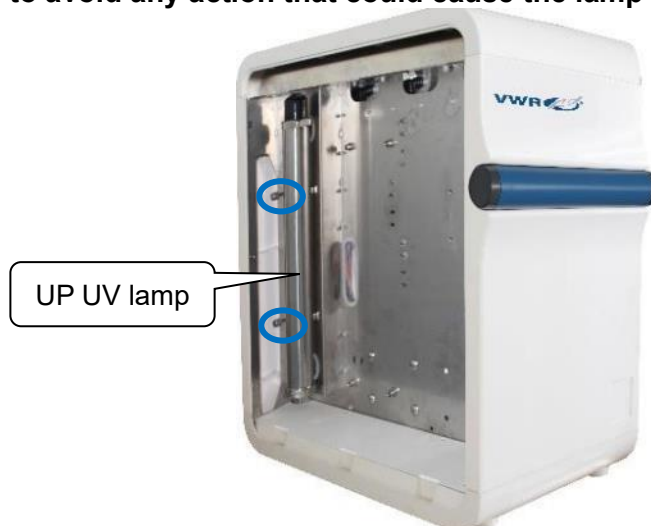

- Turn off the power.
- Remove the left side panel to find the position of 185 nm UV lamp.
- Loosen the UV lamp chamber fixing screws. As shown above, the blue circles indicate the UP water UV lamp chamber fixing screws.
- Tilt the upper end of the UV lamp chamber out.



- e) Wear gloves included in the new UV lamp package. Avoid direct skin contact with the quartz glass of the UV lamp. Remove the UV lamp cover to expose the UV lamp. Unplug the UV lamp from its power cord, and then carefully remove the old UV lamp from the chamber.



- f) Carefully insert the new UV lamp into its chamber. When about 2/3 of the lamp is inserted, hold the UV lamp and connect it to the ballast cable connector (4-pin connector) as shown in the picture, and then gently insert the new UV lamp completely into the chamber.



- g) Cap the UV chamber with the black mask.
h) Install the UV lamp chamber back to the original position.
i) Turn on the system power. Scan the RFID chip of the UV lamp on the system. Tap **Install** on the pop-up dialog to reset the corresponding UV lamp.

5.3.3 Replacing the Final Filter





- a) Remove the final filter
- b) Install the filter into UP outlet firmly, and make sure no leaking. Dispense UP water to degas.
Stop UP dispensing, installation finished.

If a final ultrafiltration cartridge is needed, install the device the same way as the final filter.



Note: Any change of type of final filter, need to reconfigure the final filter types in the setting menu(Maintenance menu→ Setting→ Final Filter)!




5.4 History Inquiry and Export

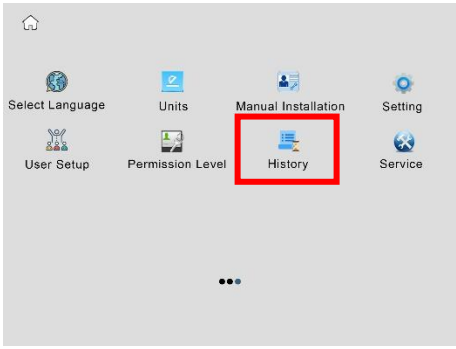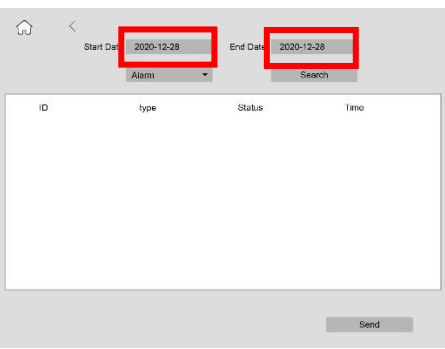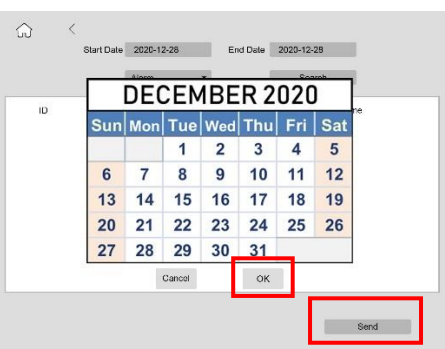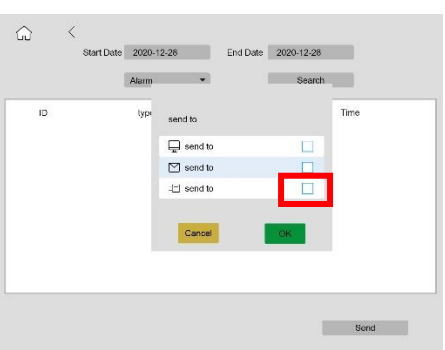
	Insert a USB stick into the port on the control console as shown.
	Swipe left on the Screen to Maintenance Menu in Standby mode. Tap History on Maintenance Menu.
	Tap Start Date and End Date to choose the date that need to check.
	Tap Search to check the history data. The history data include water quality, alarms information, dispensing information, product information and daily record.
	Choose the way to send the data. Check the red box and tap OK, then tap Send.

Read and Print the Exported History Data from a PC Computer:


Exported data are encrypted. The files exported can be read as pdf files or printed using VWR's VWR Data Converter. The converter SW is on the USB that comes with the system.

- Insert a USB stick to the right side of the control console to export the history data as shown above.
- Insert the USB stick to a computer.
- Follow the steps below to read, save or print the data for the first time, then tap the data in the .dcj file in Data to open it directly.

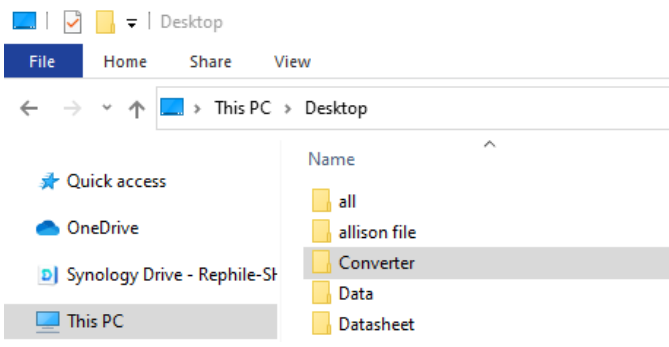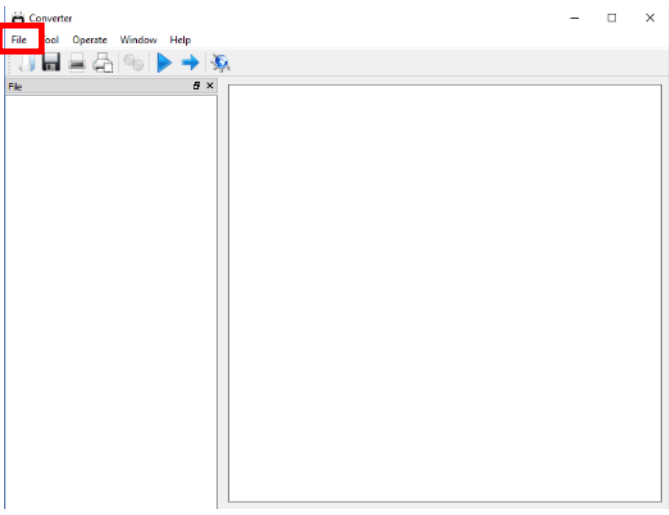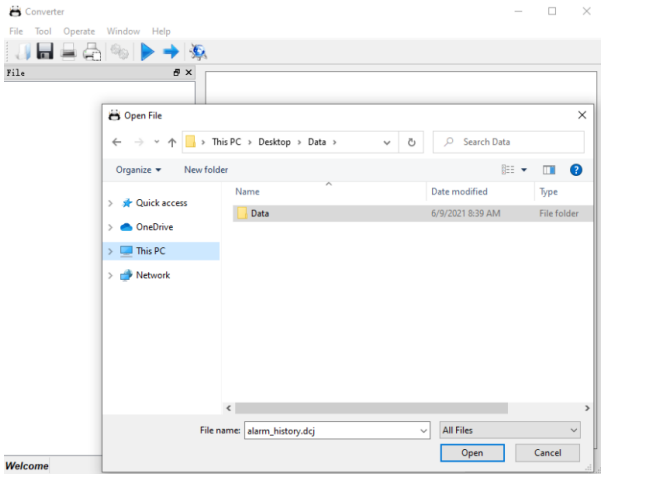
	Copy the VWR Data Converter file to your computer from the SD card. Tap the file "Converter.exe" as an administrator to complete the registration.
	Tap the file "Converter.exe" to open it.
	Tap "File", choose "Open".
	Choose the file folder "Data" from the USB stick to open it.

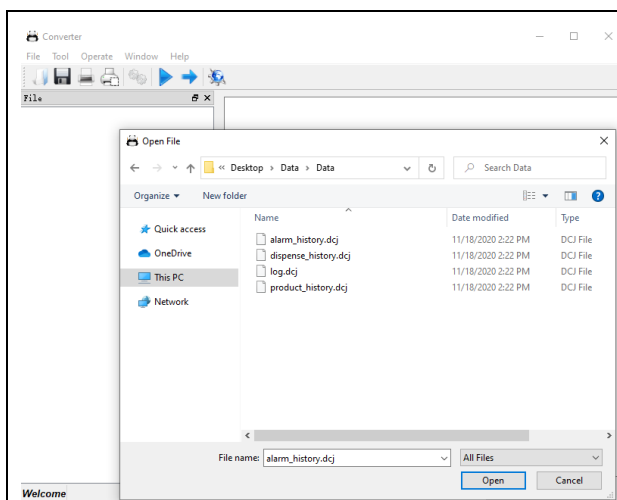

Choose the data file you wish to check and then open it.

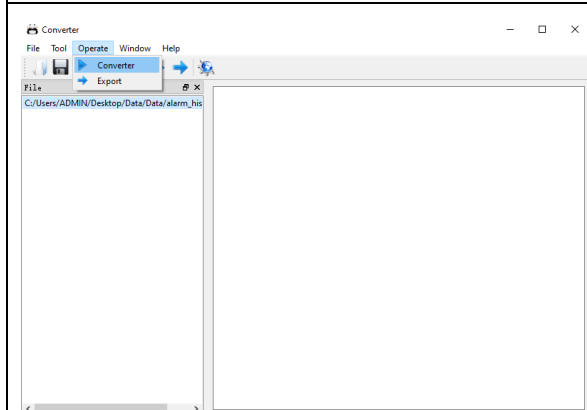

Tap **“Operate”**, then choose **“Convert”**.

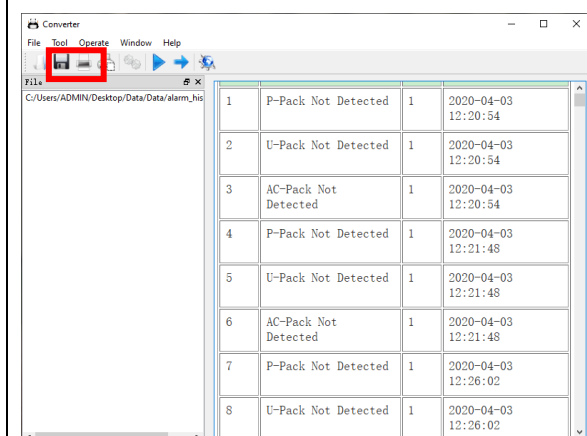

The data are as shown left.

You can **Save** the data in pdf format or **Print** it by tapping the icons as shown in the red boxes.

6 PARTS AND ORDER INFORMATION

VWR P Systems

Catalog Number	VWR P System
171-1247	VWR P Ultrapure water system
171-1248	VWR P Ultrapure water system, with TOC

Commonly Used Consumables

Catalog No.	Product Name	Description	Unit
171-1273	Dispenser with support	With a 3 m corrugated pipe, including data cables and water tubing	Set
171-1274	Upgrade kit for dispenser	With power adapter, three power cords, 20 m water tubing and 10 m corrugated pipe	Set
171-1284	Control console		Each
171-1476	Leak protector		Each
171-1258	VWR U pack		Pack
171-1260	VWR U pack (low TOC)		Pack
171-1257	VWR H pack		Pack
171-1283	185 nm UV lamp		Each
171-1262	Final Filter	0.2 µm PES final capsule filter, 1/pack	Each
171-1263	Bio-filter		Each
171-1280	UP recirculation pump	24 VDC, 1/pack	Each

7 Technical service

Web Resources

Visit the VWR website at www.vwr.com for:

- Complete technical service contact information
- Access to the VWR Online Catalogue, and information about accessories and related products
- Additional product information and special offers

Contact us

For information or technical assistance contact your local VWR representative or visit.

www.vwr.com.

8 Warranty

VWR warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non-performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

Compliance with local laws and regulations

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorisations necessary to run or use the Product in its local environment. VWR will not be held liable for any related omission or for not obtaining the required approval or authorisation, unless any refusal is due to a defect of the product.

9 Equipment disposal



This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste.

Instead it's your responsibility to correctly dispose of your equipment at lifecycle -end by handling it over to an authorized facility for separate and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

Thank you

Local VWR offices in Europe and Asia Pacific

Austria

VWR International GmbH
Graumannsgasse 7
1150 Vienna
Tel.: +43 1 97 002 0
Email: info.at@vwr.com

Belgium

VWR International BV
Researchpark Haasrode 2020
Geldenaaksebaan 464
3001 Leuven
Tel.: +32 016 385 011
Email: vwr.be@vwr.com

China

VWR (Shanghai) Co., Ltd
Bld.No.1, No.3728 Jinke Rd,
Pudong New District
Shanghai, 201203- China
Tel.: +400 821 8006
Email: info_china@vwr.com

Czech Republic

VWR International s. r. o.
Veetee Business Park
Pražská 442
CZ - 281 67 Strábrná Skalice
Tel.: +420 321 570 321
Email: info.cz@vwr.com

Denmark

VWR International A/S
Tobaksvejen 21
2860 Søborg
Tel.: +45 43 86 87 88
Email: info.dk@vwr.com

Finland

VWR International Oy
Valimotie 17-19
00380 Helsinki
Tel.: +358 9 80 45 51
Email: info.fi@vwr.com

France

VWR International S.A.S.
Le Périgares – Bâtiment B
201, rue Carnot
94126 Fontenay-sous-Bois cedex
Tel.: 0 825 02 30 30* (national)
Tel.: +33 (0) 1 45 14 85 00
(international)
Email: info.fr@vwr.com
* 0,18 € TTC/min

Germany

VWR International GmbH
Hilpertstrasse 20a
D - 64295 Darmstadt
Freecall: 0800 702 00 07
Tel.: +49 (0) 6151 3972 0
(international)
Email: info.de@vwr.com

Hungary

VWR International Kft.
Simon László u. 4.
4034 Debrecen
Tel.: +36 (52) 521-130
Email: info.hu@vwr.com

India

VWR Lab Products Private Limited
No.139. BDA Industrial Suburb,
6th Main, Tumkur Road, Peenya
Post,
Bangalore, India – 560058
Tel.: +91-80-28078400
Email: vwr_india@vwr.com

Ireland / Northern Ireland

VWR International Ltd / VWR
International (Northern Ireland) Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: +353 01 88 22 222
Email sales.ie@vwr.com

Italy

VWR International S.r.l.
Via San Giusto 85
20153 Milano (MI)
Tel.: +39 02-3320311/02-487791
Email: info.it@vwr.com

Korea

VWR International ~
17, Daehak 4-ro, Yeongtong-gu,
Suwon-si, Gyeonggi-do
Tel.: +82 31 645 7256
saleskorea@avantorsciences.com

Middle East & Africa

VWR International FZ-LLC
DSP Laboratory Complex
125, Floor 01
Dubai, United Arab Emirates
Tel.: +971 4 5573271
info.mea@vwr.com

The Netherlands

VWR International B.V.
Postbus 8198
1005 AD Amsterdam
Tel.: +31 020 4808 400
Email: info.nl@vwr.com

Norway

VWR International AS
Brynsalleen 4
0667 Oslo
Tel.: +47 22 90 00 00
Email: info.no@vwr.com

Poland

VWR International Sp. z o.o.
Limbowa 5
80-175 Gdansk
Tel.: +48 058 32 38 210
Email: info.pl@vwr.com

Portugal

VWR International - Material de
Laboratório, Lda
Centro Empresarial de Alfragide
Rua da Indústria, nº 6
2610-088 Alfragide
Tel.: +351 21 3600 770
Email: info.pt@vwr.com

Singapore

VWR Singapore Pte Ltd
18 Gul Drive
Singapore 629468
Tel: +65 6505 0760
Email: sales.sg@vwr.com

Spain

VWR International Eurolab S.L.
C/ Tecnología 5-17
A-7 Llinars Park
08450 - Llinars del Vallès
Barcelona
Tel.: +34 902 222 897
Email: info.es@vwr.com

Sweden

VWR International AB
Fagerstagatan 18a
163 94 Stockholm
Tel.: +46 08 621 34 00
Email: info.se@vwr.com

Switzerland

VWR International GmbH
Lerzenstrasse 16/18
8953 Dietikon
Tel.: +41 044 745 13 13
Email: info.ch@vwr.com

UK

VWR International Ltd
Customer Service Centre
Hunter Boulevard
Magna Park
Lutterworth
Leicestershire
LE17 4XN
Tel.: +44 (0) 800 22 33 44
Email: uksales@vwr.com