VWR® Thermometer

INSTRUCTION MANUAL

North American Catalog Number 61161-324

European Catalogue Number 620-2042

Version: 1
Issued: 15th May 2014

Legal Address of Manufacturer

United States

VWR International
Radnor Corporate Center
Building One, Suite 200
100 Matsonford Road, P.O. Box 6660
Radnor, PA 19087
800-932-5000
<http://www.vwr.com>

Europe

VWR International bvba
Researchpark Haasrode 2020
Geldenaaksebaan 464
B-3001 Leuven
+ 32 16 385011
<http://be.vwr.com>

Country of origin - China

INSTALLATION - English

Before using this Thermometer, place your thumb on the ribbed area on the back of the unit and slide the battery cover off. Remove the red tab so that the battery will make proper contact at the terminals. Replace the battery cover.

INTENDED USE

This item is intended for general laboratory use to monitor temperature and humidity.

PRODUCT SPECIFICATIONS

Display: 1½ inch (2.86 cm) LCD
Range Ambient: 23.0 to 122.0 °F (–5.0 to 50.0 °C)
Probe: –58.0 to 158.0 °F (–50.0 to 70.0 °C)
Resolution: 0.1°
Accuracy: ±1°
Sampling Time: 10 seconds
Power: 1 AAA Alkaline Battery
Case: ABS Plastic
Size: 2¾ x 4¼ x ⅝ inches
Accessories: sensor bracket, Velcro™, stand, battery, Traceable® Certificate, instructions.

DESCRIPTION OF BUTTONS AND SWITCHES /OPERATION KEYS

Key	Functions	Location
RESET	Resets the minimum/maximum memory	Front
MIN TEMP	Displays the minimum temperature	Front
MAX TEMP	Displays the maximum temperature	Front

OPERATION

1. Turn the thermometer on by inserting the battery. To insert the battery, slide the Battery door located on the back in the direction of the arrow. Insert the battery as directed by the polarity symbols (+ and -). Replace the Battery Door.
2. The external probe may be used to monitor temperatures in air/gas, liquids, and semi-solids. Both the probe and cable may be totally immersed in water.
3. Select Fahrenheit (°F) or Celsius (°C) by sliding the switch on the back of the unit to the desired setting.
4. Press the RESET button. All segments will be displayed for 3 seconds then the unit will display the current °F or °C temperature.

MEMORY

1. Press the MIN or MAX button once to recall both the ambient and probe minimum/maximum temperature. "MIN" or "MAX" will appear on the display.
2. Press the MIN or MAX button once again to return to the current temperature display. "MIN" or "MAX" will no longer appear on the display.
3. With "MIN" or "MAX" on the display, press the RESET button to reset either the minimum and maximum temperatures. All segments will be displayed for 3 seconds and the minimum or maximum temperatures will be set to the current temperature.

DISPLAY MESSAGES

HHH (Flashing)— Indicates that the temperature being measured is above the range of the unit.

LLL (Flashing)— Indicates that the temperature being measured is below the range of the unit.

MOUNTING

A wire holder is provided with the external probe for permanent mounting. To mount the probe, place a screw or nail through the hole of the wire holder and secure the holder to the wall. The unit is provided with a receptacle for wall mounting. Set a screw into the wall at the location desired. The head of the screw will need to slip into the receptacle on the back of the unit. Do not set the screw flush to the wall. Once the screw has been set properly, hang the unit in place by sliding the receptacle on the back of the unit over the head of the screw. The unit may also be mounted by using the Velcro™ supplied.

BENCH STAND

The unit is supplied with a bench stand that is a part of the case. To use the bench stand, locate the small rectangular opening on either side of the battery cover. Place your finger nail into the opening and flip the stand up. To close the stand, simply snap it shut.

TRACEABLE® CERTIFICATE

To assure accuracy, the unit is supplied with an individually serial-numbered Traceable® Certificate provided from an ISO/IEC 17025 calibration laboratory accredited by A2LA. The recommended recalibration date is noted on the certificate that is included with the product. Annual recalibrations should occur thereafter.

BATTERY REPLACEMENT

Erratic readings, a faint display, or no display are all indicators that the battery is low and must be replaced.

To replace the battery, slide Battery Door located on the back in the direction of the arrow. Insert a new 1.5V AAA alkaline battery, NOT a regular or heavy duty battery, as directed by the polarity symbols (+ and -). Incorrectly installed batteries may damage electronics. Replace the Battery Door.

ALL OPERATIONAL DIFFICULTIES

If this thermometer does not function properly for any reason, please replace the battery with a new high quality battery (see “Battery Replacement” section). Low battery power can occasionally cause any number of “apparent” operational difficulties. Replacing the battery with a new fresh battery will solve most difficulties.

TECHNICAL SERVICE

Web Resources

Visit the VWR's website at www.vwr.com for:

- Complete technical service contact information
- Access to VWR's Online Catalogue, and information about accessories and related products
- Additional product information and special offers

Contact us For information or technical assistance contact your local VWR representative or visit. www.vwr.com.

WARRANTY

VWR International warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non-performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

Equipment disposal

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste. Instead it's your responsibility to correctly dispose of your equipment at lifecycle -end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment. For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

Compliance with local laws and regulations

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorizations necessary to run or use the Product in its local environment. VWR will not be held liable for any related omission or for not obtaining the required approval or authorization, unless any refusal is due to a defect of the product.

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Local VWR offices in Europe and Asia Pacific

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India

VWR Lab Products Private Limited
135/12, Brigade Towers, 2nd Floor
Front wing, Brigade Road,
Bengaluru, India - 560 025
Tel: +91-2522-647911/922 (Mumbai)
Tel: +91-80-41117125/26 (Bangalore)
Fax +91-80-41117120
E-mail: vwr_india@vwr.com

Ireland / Northern Ireland

VWR International Ltd / VWR International
(Northern Ireland) Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: 01 88 22 222
Fax: 01 88 22 333
E-mail sales@ie.vwr.com

Italy

VWR International PBI S.r.l.
Via San Giusto 85
20153 Milano (MI)
Tel.: 02-3320311/02-487791
Fax: 02-332031307/02-40090010
E-mail: info@it.vwr.com

The Netherlands

VWR International B.V.
Postbus 8198
1005 AD Amsterdam
Tel.: 020 4808 400
Fax: 020 4808 480
E-mail: info@nl.vwr.com

New Zealand

VWR International LP
241 Bush Road
Albany
0632 New Zealand

Norway

VWR International AS
Haavard Martinsens vei 30
0978 Oslo
Tel.: 02290
Fax: 815 00 940
E-mail: info@no.vwr.com

Poland

VWR International Sp. z o.o.
Limbowa 5
80-175 Gdansk
Tel.: 058 32 38 210
Fax: 058 32 38 205
E-mail: labart@pl.vwr.com

Portugal

VWR International - Material de
Laboratório, Lda
Edifício Neopark
Av. Tomás Ribeiro, 43-3 D
2790-221 Carnaxide
Tel.: 21 3600 770
Fax: 21 3600 798/9
E-mail: info@pt.vwr.com

Singapore

VWR Singapore Pte Ltd
18 Gul Drive
Singapore 629468
Tel: +65 6505 0760
Fax: +65 6264 3780
E-mail: sales@sg.vwr.com

Spain

VWR International Eurolab S.L.
C/ Tecnología 5-17
A-7 Llinars Park
08450 - Llinars del Vallès
Barcelona
Tel.: 902 222 897
Fax: 902 430 657
E-mail: info@es.vwr.com

Sweden

VWR International AB
Fagerstagatan 18a
163 94 Stockholm
Tel.: 08 621 34 00
Fax: 08 621 34 66
E-mail: info@se.vwr.com

Switzerland

VWR International GmbH
Lerzenstrasse 16/18
8953 Dietikon
Tel.: 044 745 13 13
Fax: 044 745 13 10
E-mail: info@ch.vwr.com

Turkey

Pro Lab Laboratuvar Teknolojileri Ltd.Şti.
a VWR International Company
Orta Mah. Cemal Gürsel Caddesi
Ördekcioglu İşmerkezi No.32/1
34896 Pendik - Istanbul
Tel.: +90216 598 2900
Fax: +90216 598 2907
Email: info@pro-lab.com.tr

UK

VWR International Ltd
Customer Service Centre
Hunter Boulevard
Magna Park
Lutterworth
Leicestershire
LE17 4XN
Tel.: 0800 22 33 44
Fax: 01455 55 85 86
E-mail: uksales@uk.vwr.com