

LAL REAGENT PRODUCTS

For the Detection of Bacterial Endotoxin



TABLE OF CONTENTS

03	FUJIFILM WAKO CHEMICALS U.S.A. CORPORATION	04	THE BACTERIAL ENDOTOXIN TEST
		05	UNDERSTANDING BACTERIAL ENDOTOXIN TESTING
		06	LAL BIOCHEMICAL CASCADE
		07	LAL TECHNIQUES
08	ENDOTOXIN-SPECIFIC LAL REAGENTS	14	SYNTHETIC ALTERNATIVES
09	PYROSTAR™ ES-F SERIES	15	PYROSTAR™ NEO
12	LIMULUS COLOR KY SERIES		
13	LIMULUS PS SINGLE TEST	18	ENDOTOXIN ANALYSIS SYSTEMS
16	PEPTIDOGLYCAN AND BETA GLUCAN DETECTION	19	αBET™ RAPID ENDOTOXIN DETECTION SYSTEM
17	SLP REAGENT	20	TOXIMASTER® QC8
23	ENDOTOXIN TEST-RELATED ACCESSORIES	21	TOXINOMETER® ET-7000
24	BIOCLEAN SERIES	22	BIOTEK MICROPLATE READERS
24	LAL REAGENT WATER	27	ENDOTOXIN TEST-RELATED PRODUCTS
25	GEL CLOT REACTION TUBES	28	CONTROL STANDARD ENDOTOXIN
25	LIMULUS TEST TUBE-S	28	LPS FOR ENDOTOXIN INDICATOR
26	DEPYROGENATED DILUTION TUBES	29	ES BUFFER
26	ALUMINUM CAPS	29	ENDOTOXIN EXTRACTING SOLUTION
30	THE EXPERTS BEHIND THE PRODUCTS	31	FUJIFILM WAKO LAL DISTRIBUTORS

WELCOME TO FUJIFILM WAKO CHEMICALS U.S.A. CORPORATION

FUJIFILM Wako Chemicals U.S.A. Corporation (FUJIFILM Wako) is recognized around the world as a trusted supplier of pure chemicals and reagents. After establishing a sales office in Dallas, Texas, in 1981, we grew significantly in the U.S. and in 1989, relocated our corporate headquarters and manufacturing facility to Richmond, Virginia.

In 2012, we culminated three decades of research and development by introducing our PYROSTAR™ ES-F series to the market. The PYROSTAR™ ES-F series offers a simple, accurate platform for the detection of bacterial endotoxin. This endotoxin-specific LAL reagent is accompanied by a suite of complementary products

to offer a complete solution for your endotoxin testing needs.

Like all FUJIFILM Wako businesses, our LAL division is committed to manufacturing the highest-quality chemical reagents and supplies. Our commitment to delivering exceptional customer service is just as strong. When partnering with our LAL division, you receive personal attention from a team of specialists dedicated to your success. We look forward to providing you with superior service and supporting your laboratory in maintaining the utmost level of quality control.



HORSESHOE CRAB CONSERVATION

FUJIFILM Wako is committed to the sustainability of the Atlantic horseshoe crab (*Limulus polyphemus*) population. We voluntarily follow a set of best practices that involve careful handling and selection of healthy, viable animals. Our procedure is conducted with the utmost diligence to prevent any injury to the donor crabs, and our fishermen always return them to the same waters from which they were collected. FUJIFILM Wako cooperates with the U.S. Fish and Wildlife Service to tag and monitor horseshoe crabs as part of an ongoing conservation effort.

OUR PROMISE

As an FDA-licensed facility, FUJIFILM Wako is committed to ensuring that our production site and reagents comply with the rules, regulations and quality standards set forth by the FDA. We guarantee that the processes used to prepare our products adhere strictly to current Good Manufacturing Practices (cGMPs).



THE BACTERIAL ENDOTOXIN TEST

The Bacterial Endotoxin Test (BET) is used to detect the presence of bacterial endotoxin, which is a membrane component of gram-negative bacteria — a group of bacteria with a number of pathogenic species. Exposure to endotoxin causes the innate immune response of mammals to mount a vigorous defense. In fact, scientists classify endotoxin as pyrogenic because, even when present in small amounts, it induces biological reactions that can escalate rapidly and ultimately result in death. This makes the BET a vital assay in environments where injectables or medical devices are manufactured.

UNDERSTANDING BACTERIAL ENDOTOXIN TESTING

ENDOTOXIN

Endotoxin is a natural component of the outer cell wall of gram-negative bacteria. It is a lipopolysaccharide (LPS), which includes a polysaccharide portion, responsible for the immune response, and a lipid portion, responsible for the biological response. Endotoxin are highly heat resistant and must be exposed to temperatures of 250°C or greater for at least thirty minutes to achieve inactivation. It should be noted that endotoxin can trigger an immune response both when the bacterial cell wall is intact and when fragments are released upon lysis.

ENDOTOXIN CONTAMINATIONS

The BET is required in a number of industries to minimize the risk of human exposure to endotoxin. The goal of testing is to ensure that manufacturing operations prevent endotoxin contamination, which can cause a febrile response to end users, in the final product. This is very important in the pharmaceutical

industry, which produces a variety of drugs, biologics and devices that could expose humans to endotoxin. Though the BET is for end-product testing, it is typically performed throughout the manufacturing process. Endotoxin testing helps pharmaceutical companies avoid releasing defective products to the market. More importantly, it protects vulnerable patients with a low infection threshold.

HISTORY OF THE BET

The BET gets its secondary name — limulus amebocyte lysate (LAL) test — from its reliance on amebocytes derived from the Atlantic horseshoe crab (*Limulus polyphemus*). The blood of this living fossil, which has existed for millions of years, possesses an interesting defense mechanism against gram-negative bacteria. When the animal becomes injured, its blood will immediately bind and form a clot around the invading foreign organisms as an immune response to prevent any additional

infection to the injured area.

In 1953, Frederik B. Bang described the effects of injecting a marine bacterium into *Limulus polyphemus*. When he injected gram-negative bacteria, he observed intravascular clotting; when he injected gram-positive bacteria, no clotting occurred. Bang also observed that the clotting did not require living bacteria; dead cells also triggered the reaction. His research ultimately led to the discovery of the cells responsible for the clotting reaction and how to isolate the active material — an aqueous extract of horseshoe crab blood cells known as limulus amebocyte lysate, or LAL. In the late 1970s, the United States Food and Drug Administration (FDA) approved the LAL test as an acceptable substitution to the rabbit pyrogen test to determine the presence of endotoxin in pharmaceutical drugs, biologics and medical devices.



LAL BIOCHEMICAL CASCADE

There is a biochemical cascade associated with LAL that makes it possible to observe either qualitative or quantitative effects when endotoxin is present. The test itself can be conducted using one of three techniques: gel clot, turbidimetric or chromogenic.


As shown below, the endotoxin interaction initiates the first serine protease precursor (Factor C) from its inactivated form; this in turn activates the second serine protease precursor (Factor B). The activated Factor B stimulates the clotting enzyme, converting it from a proclotting enzyme. The clotting

enzyme cleaves peptide bonds within coagulogen to yield coagulin, the insoluble gel-forming protein produced in the gel clot assay.

The cascade mechanism is identical in the turbidimetric assay, although it is measured differently. In the turbidimetric assay, the rate at which turbidity increases relative to the concentration of endotoxin, is the critical measurement. Although the chromogenic assay has the same protease precursors activated as in both the gel clot and turbidimetric assays, a chromogenic synthetic substrate is cleaved by the clotting enzyme to convert the chromogenic

peptide into a synthetic substrate and a para nitroaniline (pNA). This releases the chromophore, yielding a yellow color whose intensity is directly proportional to the endotoxin concentration.

It should also be noted that LAL can yield a biochemical process in the presence of (1,3)-β-D-glucan by way of an additional factor (Factor G). Because of this, most LAL reagents require buffers to block the (1,3)-β-D-glucan from detection in order to determine the true concentration of endotoxin.

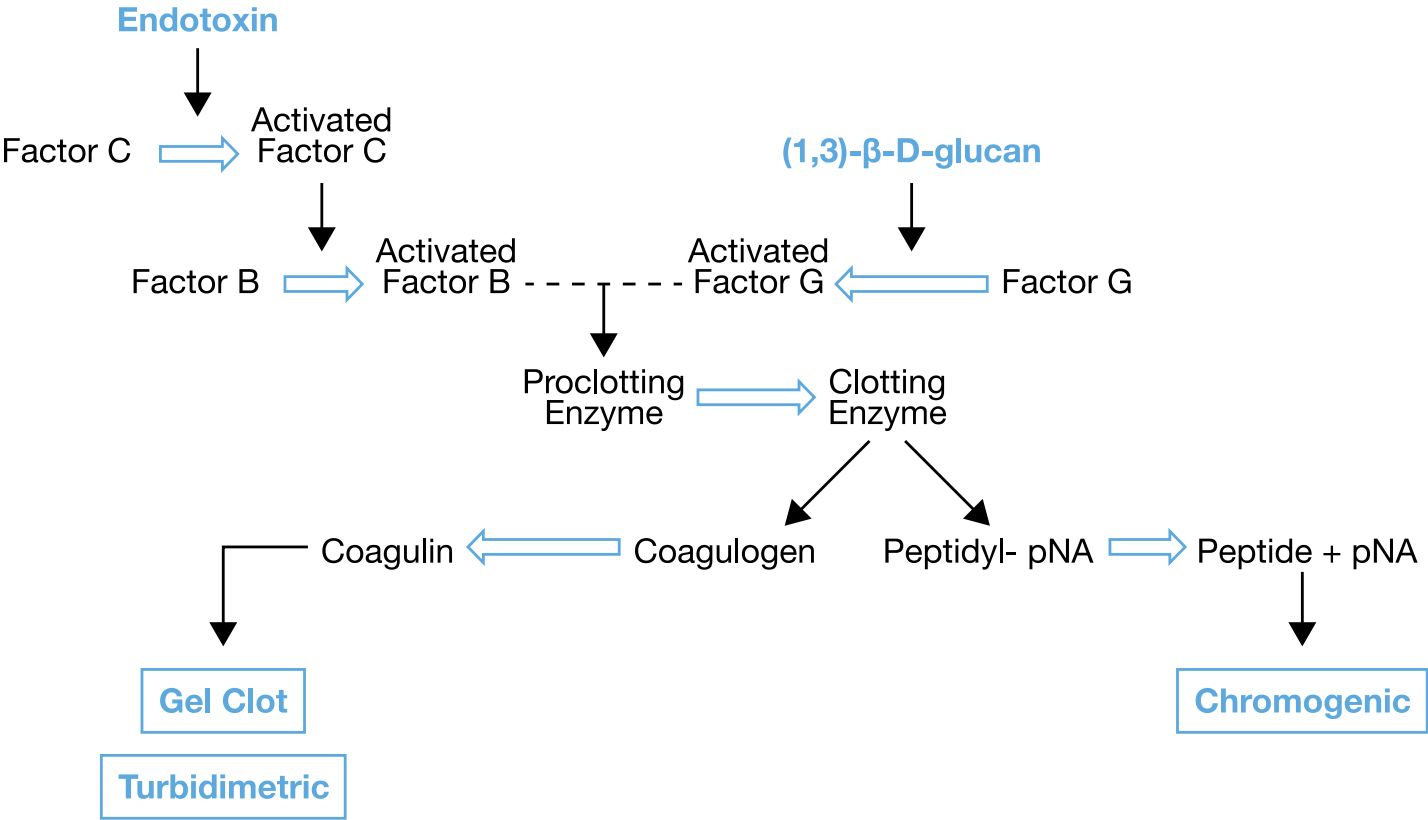

LAL TECHNIQUES

GEL CLOT

Gel clot assays function both qualitatively and semi-quantitatively. Gel clot assays are incubated at 37°C ± 1°C for 60 ± 2 minutes using a water bath, heat block or tube reader. After the incubation time has elapsed, the tube is removed and slowly inverted 180° to determine if there is a firm gel formed at the bottom of the tube. If the integrity of the gel within the tube is intact with no deformation, the result is positive. If the gel collapses, the result is negative.




TURBIDIMETRIC

Turbidimetric assays rely on turbidity change to determine the presence of endotoxin. The endpoint turbidimetric reaction is measured at a determined time after the reagent is mixed with the sample. The kinetic turbidimetric reaction measures the time it takes to reach a predetermined optical density or threshold value. Both methods can be measured around 405 nm incubated at 37°C ± 1°C in a microplate or tube reader, and a linear regression of the standards can be generated.



CHROMOGENIC

Chromogenic assays rely on color change to indicate the presence of endotoxin. The endpoint chromogenic reaction is stopped after a specific incubation period by the addition of acid; the kinetic chromogenic reaction is determined by the time it takes to reach a predetermined optical density or threshold. Regardless of the chromogenic method used, a linear regression of the standards can be generated and the yellow coloration produced can be measured around 405 nm incubated at 37°C ± 1°C in a microplate or tube reader.

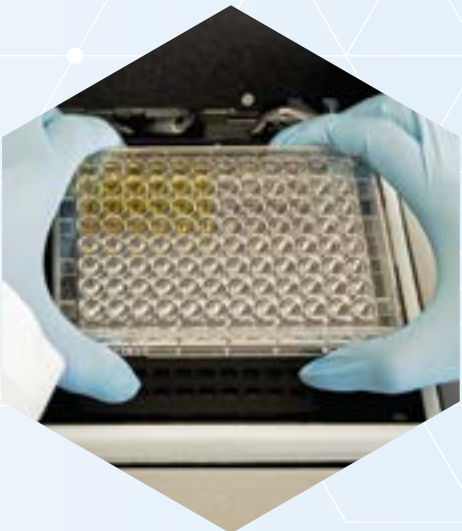

It is crucial that the user has the necessary endotoxin-free accessories to obtain reliable, reproducible and accurate results during LAL testing. Water-for-injection (WFI) or LAL Reagent Water is ideal for endotoxin testing. Reference Standard Endotoxin (RSE) should be used as a standard when testing for endotoxin, though the Control Standard Endotoxin (CSE) is an economic alternative. CSE are standardized against the RSE by each LAL manufacturer, allowing CSE to be used in place of RSE. For more detailed information, refer to regulatory guidance documents for bacterial endotoxin testing.

ENDOTOXIN-SPECIFIC LAL REAGENTS

FUJIFILM Wako provides a line of reagents formulated to perform each of the three endotoxin detection assays: the gel clot assay, the kinetic turbidimetric assay (KTA) and the kinetic chromogenic assay (KCA).

Traditional LAL reagents are untreated, which means that they react not only with endotoxin, but also with (1,3)- β -D-glucan, a fungal cell wall component that initiates the clotting cascade by activating the Factor G pathway (see page 6). The activation of (1,3)- β -D-glucan will generate a false-positive result for endotoxin because it is difficult to determine which pathogen is being measured.

The activation of LAL by (1,3)- β -D-glucan in a sample can be prevented by adding a large amount of carboxymethylated curdlan (CMC), which does not interfere with the quantitation of endotoxin. FUJIFILM Wako first made use of these findings by developing an endotoxin-specific buffer with high concentrations of CMC. Each reagent also contains buffering components that help bring most test mixtures within the pH range needed for accurate and reliable results (6.0–8.0).

PYROSTAR™ | ES-F Series



The PYROSTAR™ ES-F series of reagents are endotoxin-specific and specially formulated to be unreactive to (1,3)- β -D-glucan. In addition, these products are dual-purpose, meaning they are formulated to be used as either a gel clot or kinetic turbidimetric assay. They are available in a single-test or multi-test configuration.

These reagents require only one Certificate of Analysis (COA), regardless of whether they are used as a gel clot or kinetic turbidimetric assay. PYROSTAR™ ES-F/Plate requires a different COA. All reagent kits are matched with a specific Control Standard Endotoxin (CSE). These reagents work well with samples that have color or with samples that present difficulties in quantification by the KCA.



PYROSTAR™ ES-F SINGLE TEST

Single-test vials come with pre-dispensed LAL reagent for a single measurement. This configuration is used by adding 0.2 mL of the sample directly to the reaction vial containing the lyophilized reagent. This single-test vial is designed to be used with our Toxinometer® Measurement System (Toxinometer®) to assist labs transitioning from a gel clot to a kinetic turbidimetric methodology. The single-test configuration is ideal to assay a small number of samples or for a less-experienced user.

Key Features

- Endotoxin-specific reagent avoids false positive results from glucans
- Gel clot sensitivities include 0.015 EU/mL, 0.03 EU/mL, 0.06 EU/mL, 0.125 EU/mL and 0.25 EU/mL
- KTA sensitivities range from 0.001 to 10 EU/mL
- Can be used with the traditional water bath or heat block
- Can be used with the Toxinometer® to facilitate easier transition from gel clot to KTA

PYROSTAR™ ES-F SINGLE TEST KIT 25 single-test vials + 1 vial CSE (500 ng/vial)		
Catalog Number	Gel Clot Sensitivity (EU/mL)	Kinetic Turbidimetric Quantitative Range (EU/mL)
WPESK-0015	0.015	0.001 to 10
WPESK-0003	0.03	0.01 to 10
WPESK-0006	0.06	0.01 to 10
WPESK-0125	0.125	0.01 to 10
WPESK-0025	0.25	0.01 to 10

PYROSTAR™ ES-F MULTI-TEST

This multi-test vial is designed to be used with a water bath, heat block or the Toxinometer®. With this configuration, the user dispenses 0.1 mL of dissolved LAL reagent into the appropriate reaction tube, then adds 0.1 mL of the sample. This multi-test vial is ideal for a user testing a larger number of samples.

PYROSTAR™ ES-F 80 TEST KIT 4 multi-test vials (2.0 mL) + 1 vial CSE (500 ng/vial)		
Catalog Number	Gel Clot Sensitivity (EU/mL)	Kinetic Turbidimetric Quantitative Range (EU/mL)
WPEK4-20015	0.015	0.001 to 10
WPEK4-20003	0.03	0.01 to 10
WPEK4-20006	0.06	0.01 to 10
WPEK4-20125	0.125	0.01 to 10
WPEK4-20025	0.25	0.01 to 10

PYROSTAR™ ES-F 200 TEST KIT 4 multi-test vials (5.2 mL) + 1 vial CSE (500 ng/vial)		
Catalog Number	Gel Clot Sensitivity (EU/mL)	Kinetic Turbidimetric Quantitative Range (EU/mL)
WPEK4-50015	0.015	0.001 to 10
WPEK4-50003	0.03	0.01 to 10
WPEK4-50006	0.06	0.01 to 10
WPEK4-50125	0.125	0.01 to 10
WPEK4-50025	0.25	0.01 to 10

PYROSTAR™ ES-F 2.0 ML BULK KIT 100 multi-test vials (2.0 mL)		
Catalog Number	Gel Clot Sensitivity (EU/mL)	Kinetic Turbidimetric Quantitative Range (EU/mL)
WPEM-20015	0.015	0.001 to 10
WPEM-20003	0.03	0.01 to 10
WPEM-20006	0.06	0.01 to 10
WPEM-20125	0.125	0.01 to 10
WPEM-20025	0.25	0.01 to 10

PYROSTAR™ ES-F 5.2 ML BULK KIT 100 multi-test vials (5.2 mL)		
Catalog Number	Gel Clot Sensitivity (EU/mL)	Kinetic Turbidimetric Quantitative Range (EU/mL)
WPEM-50015	0.015	0.001 to 10
WPEM-50003	0.03	0.01 to 10
WPEM-50006	0.06	0.01 to 10
WPEM-50125	0.125	0.01 to 10
WPEM-50025	0.25	0.01 to 10

Key Features

- Endotoxin-specific reagent avoids false positive results from glucans
- Dual-purpose reagent can perform either a gel clot or kinetic turbidimetric assay (KTA)
- Gel clot sensitivities range from 0.015 to 0.25 EU/mL
- Kinetic turbidimetric sensitivities range from 0.001 to 10 EU/mL
- Due to its high sensitivity, chances of interference are reduced
- Facilitates easier transition from gel clot to KTA



PYROSTAR™ ES-F/PLATE

This multi-test vial is designed to be used with the microplate reader. To use this configuration, the user dispenses 0.05 mL of the sample into the microplate followed by 0.05 mL of dissolved LAL reagent. This assay is ideal for testing a larger number of samples in one run.

PYROSTAR™ ES-F/PLATE TEST KIT 4 multi-test vials + 1 vial CSE (500 ng/vial)		
Catalog Number	Volume (mL)	Kinetic Turbidimetric Quantitative Range (EU/mL)
WPEPK4-20015	2.0	0.01 to 10
WPEPK4-50015	5.2	0.01 to 10

PYROSTAR™ ES-F/PLATE BULK KIT 100 multi-test vials		
Catalog Number	Volume (mL)	Kinetic Turbidimetric Quantitative Range (EU/mL)
WPEPM-20015	2.0	0.01 to 10
WPEPM-50015	5.2	0.01 to 10

Key Features

- Endotoxin-specific reagent avoids false positive results from glucans
- 2.0 mL volume equates to approximately 40 tests per vial
- 5.2 mL volume equates to approximately 100 tests per vial
- The sensitivity ranges from 0.01 to 10 EU/mL

LIMULUS COLOR KY SERIES

The Limulus Color KY Series are quantitative kinetic chromogenic assays (KCA) that are endotoxin-specific and unreactive to (1,3)-β-D-glucan. This series includes both a multi-test kit and a single-test kit, which utilize a synthetic substrate that produces a yellow color to detect endotoxin with high sensitivity. All reagent kits are matched with a Control Standard Endotoxin (CSE). These reagents work well with samples that are very turbid or show difficulty in quantification by the kinetic turbidimetric assay (KTA).



LIMULUS COLOR KY SINGLE TEST

Limulus Color KY single-test vials come with pre-dispensed LAL reagent for a single measurement. This configuration is used by adding 0.2 mL of the sample directly to the reaction vial containing the lyophilized reagent. This single-test vial is ideal for a user testing a smaller number of samples or one who is less experienced.

LIMULUS COLOR KY TEST KIT 25 single-test vials + 1 vial CSE (500 ng/vial)		
Catalog Number	Kinetic Chromogenic Quantitative Range (EU/mL)	Number of Tests
291-53601	0.0002 to 5	25 tests
Key Features		
• KCA quantitative range detection limit of 0.0002 to 5 EU/mL • Reagent reacts with endotoxin over a wide range of concentrations • High sensitivity to reduce the chance of interfering factors • Can be used with the Toxinometer® • Reduces the chance of contamination and no wasted reagent during testing		

LIMULUS COLOR KY MULTI TEST

This multi-test vial is designed to be used with the Toxinometer® or a microplate reader. To use this configuration with a microplate reader, the user dispenses 0.05 mL of the sample into a microplate containing 0.05 mL of dissolved LAL reagent. When utilizing the Toxinometer®, the user dispenses 0.1 mL of dissolved LAL reagent into the appropriate reaction tube followed by 0.1 mL of the sample. The multi-test vial is ideal for a user with a larger number of samples that are turbid.

LIMULUS COLOR KY TEST KIT 3 multi-test vials (2.0 mL) + 1 vial CSE (500 ng/vial)	
Catalog Number	Kinetic Chromogenic Quantitative Range (EU/mL)
291-53101	0.0005 to 5

Key Features

- Reagent reacts with endotoxin over a wide range of concentrations
- Reduces the chance of interfering factors due to high sensitivity
- Can be used with the Toxinometer® or microplate reader
- KCA quantitative range detection limit of 0.0005 to 5 EU/mL
- The sample size is 0.1 mL when used with tube reader; 0.05 mL sample size when used with microplate reader
- The kit will perform 60 tests on our Toxinometer or 120 tests on a microplate reader



LIMULUS PS SINGLE TEST

The Limulus PS Single Test contains endotoxin-specific single-test vials and an affinity resin suspension called PyroSep™, which is designed to overcome any product interference by adsorbing potential endotoxin in samples while washing away the inhibitory components.

Case Studies: Using the Limulus PS Single Test

Endotoxin Detection in Peritoneal Dialysate/Replacement Fluid

5 mL samples of commercially available dialysate (Dialysate X) and two kinds of replacement fluids for artificial kidneys (Fluids Y and Z) were spiked with two different concentrations of Reference Standard Endotoxin: 0.1 EU/mL and 1.0 EU/mL. The samples were analyzed using the PyroSep™ method. All endotoxin spikes showed recovery within the acceptable range of 50% to 200%.

Endotoxin Recovery from Peritoneal Dialysate/Replacement Fluid Using the PyroSep™ Method

SAMPLE	ENDOTOXIN SPIKE (EU/mL)	ENDOTOXIN RECOVERY (%)
Dialysate X	0.1	80
	1.0	72
Fluid Y	0.1	144
	1.0	102
Fluid Z	0.1	108
	1.0	118



LIMULUS PS SINGLE TEST		
Catalog Number	Contents	
299-54501	- 20 LAL ES single-test vials - PyroSep™ resin suspension - LAL reconstitution solution - Wash solution - Sample diluent - 20 glass capillary columns	

PS ACCESSORY KIT (OPTIONAL)		
Catalog Number	Contents	
294-33311	- 12 syringes (20 mL) - 12 T-shape stopcocks R Type - 20 Dedicated adapters - 12 Bulldog clips - 25 Polystyrene tubes (5 mL) - 25 Polystyrene tubes (14 mL) - Test tube rack	

Key Features

- Endotoxin-specific reagent avoids false positive results from glucans
- This kit is in single-test configuration to avoid contamination
- KTA quantitation for this product is performed on the Toxinometer®
- Allows endotoxin-specific measurement in samples that typically have inhibitory components: ethanol, silicone oil, soybean oil, olive oil, stearic acid, egg yolk lecithin, oil-adjuvanted vaccine, albumin products, coagulation factor products (factor 8), interferon formulation, antithrombin III, immunoglobulin product and other ethanol-soluble samples

SYNTHETIC ALTERNATIVES

PYROSTAR™ Neo is based on a genetically engineered approach to produce reactive factors required for endotoxin detection at pharmaceutical and medical equipment manufacturing sites. It also facilitates the same sensitivity in endotoxin detection as our limulus amebocyte lysate (LAL) reagent, PYROSTAR™ ES-F.

This reagent utilizes the Kinetic Chromogenic Assay (KCA) to measure endotoxin content. It is composed of recombinant reactive factors equivalent to Factor C, Factor B and a proclotting enzyme, which constitute the horseshoe crab's natural endotoxin reaction system, allowing for endotoxin detection that offers the same sensitivity as the LAL reagent. In addition, the absence of factor G eliminates interference due to reaction with β -glucan, enabling improved specificity.



SYNTHETIC ALTERNATIVES


SYNTHETIC ALTERNATIVES

PYROSTAR™ Neo is a new line of endotoxin detection reagents developed by recombinant technology. It mimics traditional LAL reagents with three recombinant factors (Factor C, Factor B, and proclotting enzyme) and the pNA chromogenic group, offering an assay with high sensitivity and endotoxin specificity.

Key Features

- Colorimetric method, can be used with an absorbance plate reader
- 3-Factor system mimics the same cascade reaction as traditional LAL
- Endotoxin-specific reagent eliminates the risk of false positives from (1,3)- β -D-glucan
- 100% animal free
- Quantitative range: 0.001 to 50EU/mL
- High sensitivity with less lot-to-lot variation
- Stable storage after dissolution (4hrs. at 2-8°C and 2 weeks at 30°C)
- Validation package provided by the manufacturer

Catalog Number	Kinetic Chromogenic Quantitative Range (EU/mL)	Number of Tests
294-36731	0.001 - 50EU/mL	50 tests

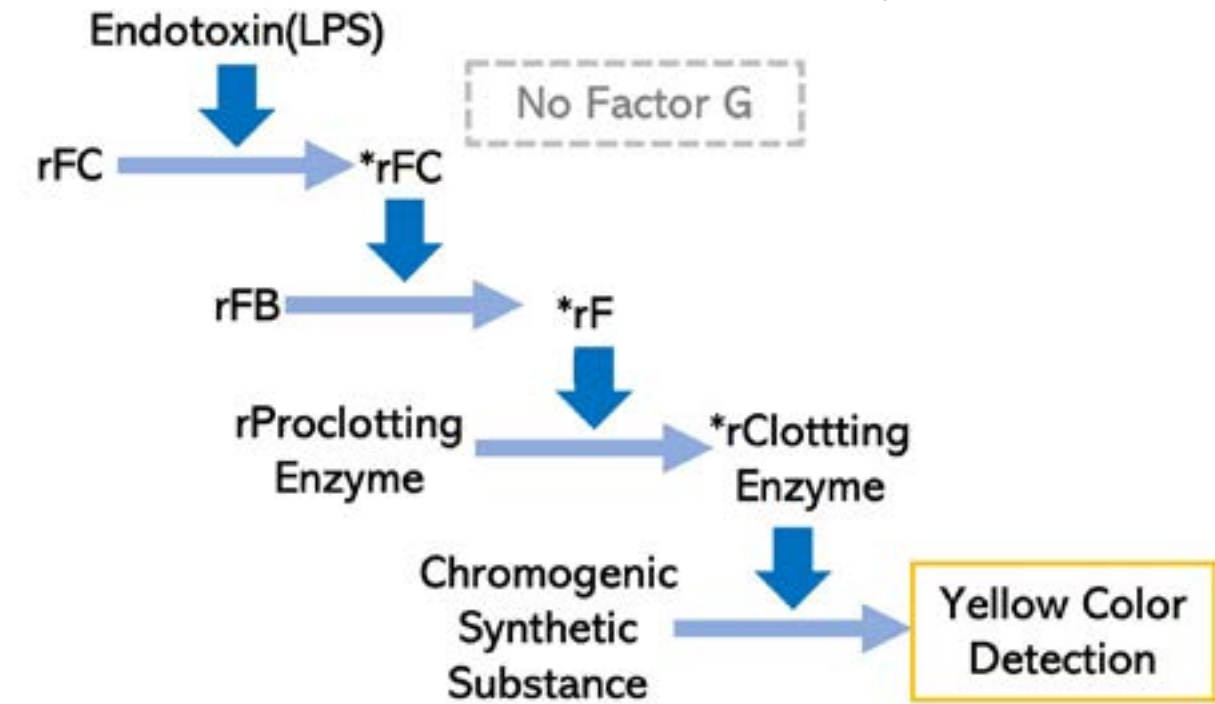

PEPTIDOGLYCAN AND BETA GLUCAN DETECTION

Quantitative determination of peptidoglycan (PG) and (1,3)-β-D-glucan (BDG) in parenteral solutions and active pharmaceuticals.

The SLP-HS Single Reagent Set II is able to detect both peptidoglycan (PG) and (1,3)-β-D-glucan (BDG). The product utilizes a self-defense mechanism found within the hemolymph of the silkworm Bombyx mori, which results in a color change by generating a black melanin pigment in the presence of microbial contamination. Since PG is found in all bacterial cell walls and BDG in most fungi, the SLP (silkworm larvae plasma) reagent constitutes a highly sensitive and broadband method for detection of microbial contamination.

PEPTIDOGLYCAN AND BETA GLUCAN DETECTION

SLP-HS SINGLE REAGENT SET II

Quantitative determination of peptidoglycan (PG) and (1,3)-β-D-glucan (BDG) in parental solutions and active pharmaceutical.

The SLP-HS Single Reagent Set II is able to detect both peptidoglycan (PG) and (1,3)-β-D-glucan (BDG). The product uses a self-defense mechanism of silkworm Bombyx mori which leads to a color change by building a black melanin pigment in the presence of microbial contamination. Since PG is found in all bacterial cell walls and BDG in most fungi, the SLP (silkworm larvae plasma)-reagent constitutes a highly sensitive and broadband method for detection of microbial contamination.

Key Features

- SLP (Silkworm Larvae Plasma reagent) based test principle
- Kinetic Colormetric Assay
- PG isolated from S. aureus used as standard
- BDG Detection Limit: 1pg/mL
- PG Detection Limit: 10pg/mL
- Analysis automated through Toximaster® software
- Measurement time: 120 minutes at 30 °C
- Testing for microbial contamination in dialysate and its raw materials
- Detection of fungal compounds in pharmaceuticals, medical devices, biologics, and genetically engineered products
- Investigation of bacterial pollutants in commercial water

Catalog Number	Product Name	Number of Tests
296-81001	SLP-HS Single Reagent Set II	20 tests



ENDOTOXIN ANALYSIS SYSTEMS

FUJIFILM Wako provides analysis systems with software that is user friendly and capable of performing all methodologies for LAL testing.

ENDOTOXIN ANALYSIS SYSTEMS



αBET™ RAPID ENDOTOXIN DETECTION SYSTEM

The CMD αBET™ system is a fully integrated testing system that provides a rapid and sustainable solution to endotoxin testing without compromising on sensitivity or performance. Using our FDA-licensed turbidimetric Limulus amoebocyte lysate (LAL) reagent, quantitative results are provided in under half the time of other testing formats, with less reagent consumption and minimal user input.

Catalog Number	Product Name	Contents
CMD-A-0007-EN	CMD αBET™ Instrument	Instrument, power cord, universal adaptor, user's manual, barcoded key cards
CMD-A-0006-C	CMD αBET™ Consumable Pack	25 x Cartridge + 5 x CMD Reagent

Key Features

- Portable and fully integrated bench top system with FDA 21 CFR Part 11 compliant software
- Data back up to PC by binary code, preventing the ability of data modification
- Backups made in two formats-native firmware and HTML for ease of viewing records
- 7-inch Touchscreen
- QR Code access cards included (1 administrator card, 10 user cards)
- Run times of no more than 30 minutes per assay can achieve 0.001EU/mL
- Uses less reagent than traditional methods
- Minimal user input needed beyond addition of sample
- Guided steps for performing various assays
- Compatible with color and opaque samples
- For use with PYROSTAR™ ES-F 0.015EU/mL reagent.

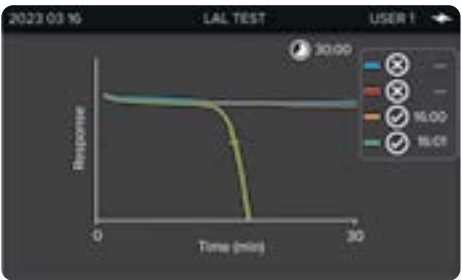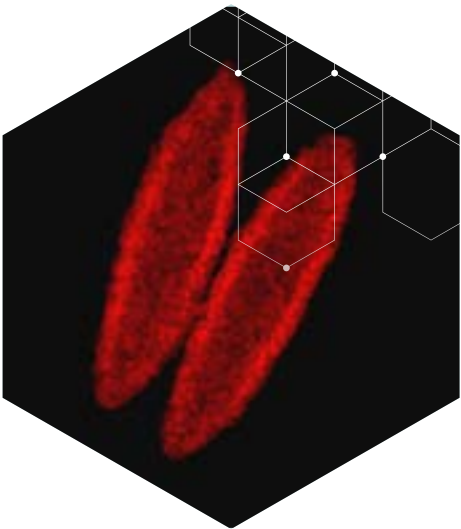

Calibration ID	Calibration Date	Test Date / Time	Instrument ID
0001	2023/03/16	2023/03/16	0001
0002	2023/03/16	2023/03/16	0002
0003	2023/03/16	2023/03/16	0003
0004	2023/03/16	2023/03/16	0004
0005	2023/03/16	2023/03/16	0005
0006	2023/03/16	2023/03/16	0006
0007	2023/03/16	2023/03/16	0007
0008	2023/03/16	2023/03/16	0008
0009	2023/03/16	2023/03/16	0009
0010	2023/03/16	2023/03/16	0010

TOXIMASTER® QC8 SOFTWARE

The Toximaster® QC8 software supports data processing based on the protocols complying with three types of pharmacopeia (USP/EP/JP) for bacterial endotoxin testing. This software is compliant with FDA 21 CFR Part 11 ERES (electronic records/electronic signature), which requires the proper information from all parties involved in the testing of the sample. Toximaster® QC8 offers excellent audit trail capabilities, can provide statistical processing of means and standard deviations, and has robust data processing functions. These data processing functions can be performed by defining sample types such as standards, controls and test samples.

TOXINOMETER® ET-7000 – 21 CFR PART 11 (ERES) COMPLIANT SOFTWARE


Catalog Number	Model	Contents
295-36141	Part 11 software only	Toximaster® QC8 Software

BIOTEK ELX808IU MICROPLATE READER – 21 CFR PART 11 (ERES) COMPLIANT SOFTWARE

Catalog Number	Model	Contents
294-36211	Part 11 software only	Toximaster® QC8 Software

Key Features

- FDA 21 CFR Part 11 ERES compliant
- Endotoxin determination in compliance with pharmacopoeias (USP/EP/JP) for BET
- Capable of creating three types of curves: internal standard curve, manual input curve and a measured curve
- Product endotoxin limit and MVD generation
- Allows trending of product results over time and early detection of potential product failures
- Generates hard copy printouts with all pertinent information for routine audits through three unique audit trails
- Creation of a standard workflow to be reviewed, confirmed and approved prior to operating
- Security functions to lock the application, disable an account and lock out the system
- Backup of the user management database and automatic backup of the system information database
- Improved precision and accuracy over the traditional gel clot method
- Ideal software for comparative testing and validating most LAL

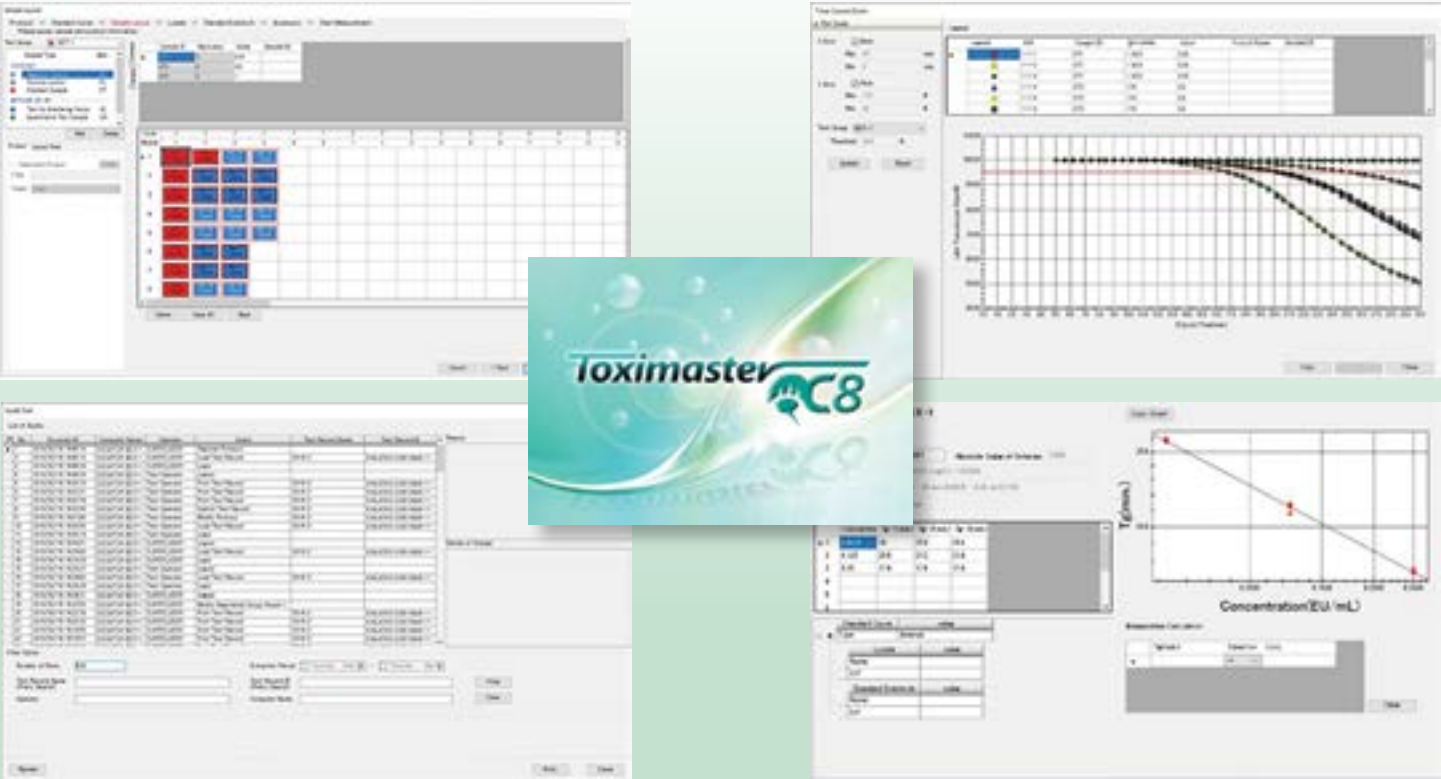

Key Features

- Single assay module can simultaneously measure up to 16 samples
- Expansion modules are available (up to 8) to extend the number of samples in multiples of 16
- Single-test configuration avoids the “hot wells” phenomenon associated with microplate readers by evenly distributing the temperature in all single wells
- Reduces the chance of contamination due to the single-well formats
- Ideal for users who are converting from gel clot method to kinetic turbidimetric assays

TOXINOMETER® ET-7000 ENDOTOXIN MEASUREMENT SYSTEM

The Toxinometer® ET-7000 is a computer-operated kinetic incubating tube reader, designed to be exceptionally user-friendly. Our state-of-the-art expansion modules can be connected to allow for endotoxin testing in a wide range of fields and sample quantities.

The Toxinometer® ET-7000 can perform all LAL testing methodologies (gel clot, kinetic turbidimetric and kinetic chromogenic assays). The instrument has 16 wells for single-test configuration in order to easily detect erroneous readings or contamination associated with each well. The kinetic incubating tube reader has uniformity temperature settings of 30°C and 37°C.

TOXINOMETER® ET-7000 + 21 CFR PART 11 (ERES) COMPLIANT SET

Catalog Number	Model	Power Source	Contents
294-35871	Toxinometer® ET-7000/U Part 11 Set	100–120 ± 10% VAC (USA)	- 1 Toxinometer® ET-7000 - Toximaster® QC8 Software
299-35821	Toxinometer® ET-7000/E Part 11 Set	220–240 ± 10% VAC (Europe)	- System Validation Doc.

TOXINOMETER® ET-7000 ANALYSIS MODULE

Catalog Number	Model	Contents
293-36061	Toxinometer® ET-7000 Expansion Module	16-well expansion unit



BIOTEK MICROPLATE READERS

The BioTek microplate readers provide the flexibility to manipulate and analyze your data. This multi-channel reader continues BioTek’s tradition of offering the strictest specifications of any microplate reader that ensure both accurate and repeatable results. The microplate readers offer a superior 4-zone™ incubator, providing excellent stability for temperature-sensitive assays such as endotoxin analysis and long-term bacterial and yeast growth studies. FUJIFILM Wako now offers this highly reliable instrument with a convenient laptop and software package specifically dedicated to LAL testing.

- Key Features
- 96-well configuration, ideal for high throughput of samples
 - Superior 4-Zone™ incubation
 - Fast kinetics, endpoint and linear well scanning
 - Optical design eliminates crosstalk
 - Extensive endotoxin analysis with Toximaster® QC8 software available for the Biotek Elx808 model

BIOTEK ELX808 MICROPLATE READER		
Catalog Number	Model	Contents
BIOT-808IU	ELx808	- 1 BioTek ELx808 unit
BIOT-35931	ELx808	- 1 BioTek ELx808 unit - Toximaster® QC8 Software



BIOTEK EPOCH 2 MICROPLATE READER		
Catalog Number	Model	Contents
BIOT-EP2GS	Epoch 2	- 1 BioTek Epoch 2 unit



BIOTEK 800TS MICROPLATE READER		
Catalog Number	Model	Contents
BIOT-800GS	800TS	- 1 BioTek 800TS unit



ENDOTOXIN TEST-RELATED ACCESSORIES

FUJIFILM Wako offers a premiere collection of endotoxin-free accessories including pipette tips, microplates, test tubes, caps and LAL Reagent Water designed for worry-free endotoxin testing.

ENDOTOXIN TEST-RELATED ACCESSORIES

BIOCLEAN SERIES

BioClean Series pipette tips and microplates are endotoxin-free. These micropipette tips and microplates are individually wrapped to reduce the chance of contamination and are for all types of LAL testing. BioCleanTip Wako® are available in both 200 µL and 1000 µL volumes, with the 200 µL volume tips being offered in standard and extended sizes. The Extend S tips are used for aliquoting multi-test reagent into Gel Clot Reaction Tubes, Limulus Test Tube-S or BioCleanPlate Wako™. This pipette tip can deliver the reagent to the bottom of the tubes/wells without introducing contamination.

Catalog Number	Product Name	Volume Size	Quantity
291-35021	BioCleanTip Wako® 200 II	200 µL/tip	100 tips
298-35031	BioCleanTip Wako® 1000 II	1000 µL/tip	100 tips
294-35011	BioCleanTip Wako® Extend S	200 µL/tip	100 tips
293-35221	BioCleanPlate Wako™	96 wells/plate	50 plates

Key Features

- Endotoxin-free pipette tip
- Endotoxin-free 96-well microplate
- Individually packaged, ideal for use in clean rooms
- Packaging film features dust-free protection
- Pipette tips fit most micropipettes
- Plates are designed for use with microplate readers

LAL REAGENT WATER

LAL Reagent Water can be used to fulfill all aspects of endotoxin testing. From reconstitution of both the CSE and LAL for creation of standard curves to the making of product dilutions, LAL Reagent Water provides reliable and reproducible data when paired with our endotoxin-specific LAL reagents.

Catalog Number	Product Name	Quantity
LRW-2030	LAL Reagent Water, 30 mL	20 x 30 mL glass bottles
LRW-12125	LAL Reagent Water, 125 mL	12 x 125 mL plastic bottles

Key Features

- Endotoxin-free water
- Steam-sterilized by USP standards
- Non-LAL reactive
- Great for reconstitution of LAL, RSE and CSE
- Used to perform dilution series of RSE and CSE
- Used for sample dilutions
- 30mL lysate reagent water is suitable for BET
- 125mL lysate reagent water is suitable for BET, Peptidoglycan, and β-glucan testing



GEL CLOT REACTION TUBES

Our Gel Clot Reaction Tubes are made of high-quality borosilicate glass, depyrogenated at 250°C. This results in an endotoxin-free tube that is designed to perform successful LAL gel clot assays with heat blocks and water baths.

Catalog Number	Product Name	Quantity
CT-1075	10 x 75 mm Borosilicate Test Tubes	200 tubes/pkg
CT-1075B	10 x 75 mm Borosilicate Test Tubes, Bulk	1250 tubes/pkg

Key Features

- Endotoxin-free test tubes
- Endotoxin-free tubes depyrogenated at 250°C
- Made with borosilicate glass and designed for gel clot testing



LIMULUS TEST TUBE-S

Our specialized Limulus Test Tube-S are made of high-quality borosilicate glass depyrogenated at 250°C. These tubes have a special curvature that allows the appropriate amount of light to transmit through the sample to measure the endotoxin concentration based on the method used. This glass tube is designed for the Toxinometer® Measurement System.

Catalog Number	Product Name	Dimensions	Quantity
292-32751	Limulus Test Tube-S with Aluminum Cap	12 x 75 mm	80 caps/tubes
293-26551	Limulus Test Tube-S	12 x 75 mm	100 tubes

Key Features

- Endotoxin-free test tubes
- Designed for gel clot, KTA and KCA testing on the Toxinometer®
- Made with borosilicate glass

DEPYROGENATED DILUTION TUBES

Our Depyrogenated Dilution Tubes are made of high-quality borosilicate glass depyrogenated at 250°C. This results in endotoxin-free tubes that can perform successful LAL dilution series, from creating the standard curve to sample dilutions, with minimal chance of potential contamination.

Catalog Number	Product Name	Quantity
DL-13100	13 x 100 mm Borosilicate Test Tubes	50 tubes

Key Features


- Endotoxin-free test tubes
- Made with borosilicate glass



ALUMINUM CAPS

Our aluminum caps assist in reducing the chance of contamination, preventing any additional substances or material from entering the testing tubes.

Catalog Number	Product Name	Dimensions	Quantity
CP-1075	Aluminum Caps for 10 x 75mm	11.2 x 11.5 mm	100 caps/pkg
CP-13100	Aluminum Caps for 13 x 100mm	14.5 x 18 mm	50 caps/pkg
293-28251	Aluminum Cap-S	14.7 x 18 mm	100 caps/pkg

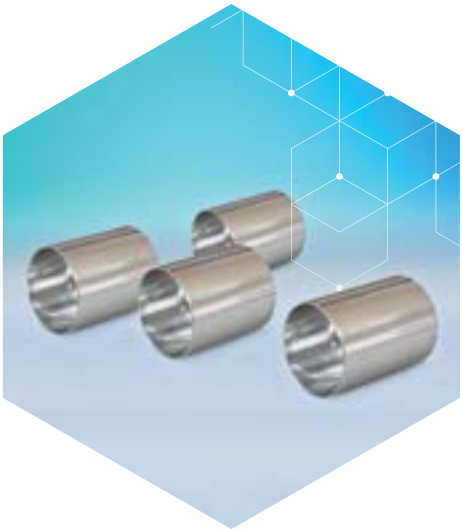

ENDOTOXIN TEST-RELATED PRODUCTS

Our Endotoxin Test-Related Products are additional specialty kits that assist with the successful completion of bacterial endotoxin tests.



ENDOTOXIN TEST-RELATED PRODUCTS

CONTROL STANDARD ENDOTOXIN

Our Control Standard Endotoxin (CSE) is a lyophilized product composed of a 500 ng vial of endotoxin, purified from *E. coli* UKT-B strain that serves as an accurate and dependable standard for endotoxin testing.

Catalog Number	Contents
CSE4037-5006	6 vials (500 ng/vial)

Key Features

- Endotoxin derived from *E. coli* UKT-B
- Can be used to prepare controls and standard curves
- Potency of CSE supplied to match each lot of LAL
- Reconstituted CSE can be stored at 2°C–10°C for one month

LPS FOR ENDOTOXIN INDICATOR

The LPS for Endotoxin Indicator is composed of *E. coli* O55:B5 (>100,000 EU/vial) and serves as a stock solution to prepare multiple endotoxin indicators for depyrogenation validation studies. The prepared endotoxin indicators may be tested using reagents from our PYROSTAR™ ES-F series.

Catalog Number	Contents
WLPS-0100K	6 vials (>100,000 EU/vial)

Key Features

- *E. coli* O55:B5 (>100,000 EU/vial)
- Can make multiple endotoxin indicators from one vial
- Used for validation studies



ES BUFFER

The ES Buffer is an endotoxin-specific buffer that is ideal for untreated LAL. Though all LAL reagents produced by FUJIFILM Wako are endotoxin-specific, other commercially available LAL reagents are not. Without this buffer, tests can activate (1,3)-β-D-glucan when trying to determine the endotoxin concentration in the sample.

Catalog Number	Contents
ESB-0006	6 vials x 6.0 mL

Key Features

- Each vial has solution for use in reconstituting a maximum of 5.2 mL of reagent
- Ideal for LAL that has the potential to react with glucans, thus making an endotoxin-specific reagent

ENDOTOXIN EXTRACTING SOLUTION

Traditionally, water or saline solution has been used to extract endotoxin in tests involving medical devices and equipment; however, the efficacy of this extraction method is debated. In order to provide our customers with a more reliable method, FUJIFILM Wako has developed a solution containing human serum albumin (HSA) that is capable of extracting endotoxins that cannot be extracted in water or saline solution.

Catalog Number	Contents
293-51601	4 vials x 10 mL

Key Features

- Capable of extracting endotoxins that are unable to be extracted in water or saline
- Recommended for use in endotoxin testing on equipment and devices that may come into contact with fluids containing blood or protein

THE EXPERTS BEHIND
THE PRODUCTS

Our team at FUJIFILM Wako would like to thank you for your interest, and we look forward to working with you!



CONTACT OUR DISTRIBUTORS

Looking for a distributor in your region?

Please see our list of distributors around the globe:

ALGERIA

Pharmascience

Cooperative Computer, Villa N°37
Douaouda, Tipaza, 42 445-Algeria
Contact: Dr. Ali Melzi and Dr. Salim Ould-Larbi
Mobile: +213 667 899 062
Phone: +213 (0) 24 216 172 / +213 (0) 24 216 173
Email: contact@pharmascience-dz.com
salim.ouldlarbi@pharmascience-dz.com
Website: http://pharmascience-dz.com/

ARGENTINA

Endolab S.A.

Emilio Lamarca 381/397
Buenos Aires, Argentina
Contact: Osvaldo Fernandez Poblet
Phone: +54 011 4733 4047
+54 011 4792 7820
Email: ventas@endolabsa.com.ar
Website: www.endolabsa.com.ar

2MS S.R.L.

Llerena 3192
CABA
CP 1427.
Argentina
Contact: Dr. Marcelo Madeo
Phone: +5491145247878/7979
Email: 2ms@2ms.com.ar
Website: www.2ms.com.ar

BANGLADESH

Bioaid

1/B/1 (Flat A1) Road#01,
Shyamoli, Dhaka-1207, Bangladesh
Contact: Md. Abdul Latif
Phone: +880 29134673
Email: bioaidbd@gmail.com

M/S RIFA Enterprise

1 New Eskaton Rd.
Razzak Plaza, Floor 10, Dhaka-1000
Contact: AKM Faizur Rahman
Phone: +880 1911 367920
+880 2 934250
Email: faizur7@dhaka.net

BELARUS

DP Unilab

220094 Minsk, Gorovtza str.24, Apartment 3, Office 4
Republic of Belarus
Contact: Eugene Kostoglod
Phone: +375 29 395 25 90
Fax: +375 17 319 57 91
Email: info@unilab.by
Website: www.unilab.by
www.dpunilab.by

BRAZIL

Brazil Nova Analítica

Av. Fábio Eduardo Ramos Esquivel, 2100 Centro,
Diadema - SP, 09941-202, Brazil
Phone: +55 11 2162 8080/ (11) 970277918
Email: reagentes@novanalitica.com.br
Website: www.analiticaweb.com.br

CANADA

Fisher Scientific

Phone: +1 800 234 7437
Website: www.fishersci.ca

COLOMBIA

ADN Internacional

Carrera 58 #127-59 of 378
Colombia
Contact: Ximena Moreno / Sandra Cortes
Phone: +57 601 9261278 / +57 3173698044
Email: farma@adninternacional.co / gerencia@adninternacional.co
Website: www.adninternacional.co

COSTA RICA

Makol OCR S.A.

Heredia, Costa Rica
Contact: Angela Rauff
Phone: +506 22939694
Fax: +506 22930562
Email: angelar@makolcr.com
Website: www.makolcr.com

DOMINICAN REPUBLIC, PUERTO RICO, JAMAICA,
TRINIDAD AND TOBAGO, AND CURACAO

BDC Serralles, srl

Ave. Gregorio Luperon 51
Santo Domingo, DN 10137
Dominican Republic
Contact: Manuel Bros
Phone: +1 809 338 8888
Fax: +1 809 338 8889
Email: bdcint@bdcint.com.do
Website: www.bdcint.com.do

Commercial Manager

Contact: Tomas Garcia
Phone: +1 809 338 8888 ext. 100
Email: tomas@bdcint.com.do

Logistics

Contact: Samuel Fersobe
Phone: +1 809 338 8888 ext. 300
Email: purchasing1@bdcint.com.do

EGYPT

BioScience Group B.S.G

Al-Mansouria, Monshaat Al-Qanatar, Giza, Egypt
Contact: Mohamed Lamie
Phone: +20 1021011177
Fax: +20 38728772
Email: inquiries@bioScience-bsg.com
Website: www.bioscience-bsg.com/

Femto Trade Co.

21 Dr Michel Bakhoum Street,
Dokki, Giza, Egypt
Contact: Dr. Essam Abdallah
Phone: +20 1005126629
Fax: +20 233928424
Email: ahmed@femtotrade.com / essamfemto@yahoo.com
Website: www.femtotrade.com

ETHIOPIA

Binocular Trading

Bole Sub City, Woreda 5, Mulugeta Zeleke,
Building 108
Addis Ababa, Ethiopia
Contact: Zenaw Workneh
Phone: +251 911 733801
Email: zenaww@gmail.com

EUROPE

FUJIFILM Wako Chemicals
Europe GmbH

Fuggerstr. 12, 41468 Neuss, Germany
Contact: Yannic Werner
Phone: +49 2131 311 101
Fax: +49 2131 311 110
Email: labchem_wkeu@fujifilm.com
Website: www.labchem-wako.fujifilm.com/europe/index.html

INDIA

BioSolutions

102, Ashok Heights, Dr. S. Radhakrishnan Marg
Andheri (East), Mumbai-400069, India
Contact: Mr. Prakash Bapodara
Phone: +91 9819973583
Email: mail@biosolutions.co.in
biosolutions@rediffmail.com

Biomatiq Quality Solutions

Plot No.1386, Apuroopa Township, Pragathi Nagar
Medchal District, Hyderabad, India 500090
Contact: Omkaranath B.
Phone: +91 91000 50333/ +91 91210 10201
Email: contact@biomatiq.com / omkar@biomatiq.com
Website: www.biomatiq.com

IRAQ

Al-Banfsaj for Scientific Equipment

Building No. 82, Areyadeh Street
Arjan Amman 11196, Jordan
Contact: Mr. Mazin Mousa
Phone: +962 (06) 5565424
Mobile: +962 795514322
Fax: +962 (06) 5058273
Email: info@chromejo.com

Iraq office:

No 5/23 central Market St, Karada
Baghdad 10069-IRAQ
Contact: Eng Wajdi Khalaf
Phone: +9647500771483
Email: info@chromejo.com

ISRAEL

Eisenberg Bros. Ltd.

1 Sderot HaTa'asiyot,
Millennium Center, 4th FL
P.O. Box 184, Modi'in 7171101, Israel
Tel: +972-3-9777000
Fax: +972-3-9777001
Email: info@eb1.co.il
Website: www.eisenbros.co.il

ITALY

Bioclass s.r.l.

Via Rodolico, 152 – 51100
Pistoia, Italy
Contact: Gianluca Dami and Margherita Nannucci
Phone: +39 0573 365563
Fax: +39 0573 365610
Email: info@bioclass.it
Website: www.bioclass.it

IVORY COAST, GHANA, TOGO, MALI, SENEGAL, GUINEA,
ETHIOPIA, BURKINA FASO, AND BENIN

African Bio-expertise Services-CI Ltd

Cocody Angre Djibi, Boulevard 8e Tranche
28 B.P. 1080, Abidjan 28, Abidjan - Côte d'Ivoire
Contact: Stephane Damien
Phone: +225 22 42 89 02, +225 78 69 42 26
Email: stephane_damien@abs-ai.com
Website: www.abs-ai.com

JORDAN

Safe Way Medical Supply

Al-Bayader, Hosni Sober Street, Randa Complex 72
Amman, Jordan
Contact: Hamzeh Salameh
Phone: +962 6 5811 704
+962 797868687
Fax: +962 6 5811 704
Email: H.Salameh@sway-jo.com
info@sway-jo.com
Website: https://sway-jo.com

KAZAKHSTAN AND UZBEKISTAN

Algimed LLC

Auezova street, Building 8
Astana City, Kazakhstan
Contact: Anastassiya Sek
Phone: +7 700 613 13 07
Email: a.sek@algimed.kz
Website: www.algimed.kz

FUJIFILM WAKO LAL DISTRIBUTORS

KENYA, RWANDA, AND UGANDA

Aqualytic Laboratories Ltd

Ramco Court, Mombasa Road, Nairobi
Contact: John Kiiru
Phone: +254 722 54 73 44
Email: lab@aqualyticlab.com
Website: www.aqualyticlab.com

KYRGYZSTAN LLC “T&L”

720031 Bichkek, Karasaeva 84, Kyrgyzstan
Contact: Elrad Abdukadyrov
Phone: + 996 312 54 38 44 / + 996 777 47 90 14
Email: contact@tecdilog.kg
Website: www.tecdilog.kg

LEBANON

NUMELAB sarl. - Nuclear Medical Laboratory

457 New Naccache
Antelias, 70-410 Lebanon
Contact: Serge Daou
Phone: +961 1396677 ext. 120
Fax: +961 1396688
Email: serge.daou@numelab.com
Website: www.numelab.com

MACEDONIA, KOSOVO, ALBANIA

Biotek DOO

Zagrebska 28V 1000 Skopje
Macedonia
Contact: Milena Stojanova and Aleksandra Kasapovska
Phone: +389 02 3110106
Fax: +389 02 3220154
Email: mstojanova@biotek.com.mk
sales@biotek.com.mk
Website: www.biotek.com.mk

MEXICO

Bioprocesos Farmaceuticos Integral S.A. de C.V.

Cerro del Obispo 231 Lomas de Valle Dorado
Tlalnepantla de Baz, 54023, Mexico
Phone: +52 01 55 1509 1624 / +52 01 55 50128130 / +52 55 2869 0116
WhatsApp: +52 01 55 8398 6635
Email: hola@bpfintegral.com

City of Guadalajara
Rinconada del Camichin 1860 Int 4
Col. Real del Camichin, Zapopan Jalisco
Phone: +52 1 33 3311 8485
Contact: IBQ. Karla Gimena Velazquez Cansino (Commercial Director-Sales) / IBI. Luis Alberto Valderrama Lopez (General Director)
Email: suc_gdl@bpfintegral.com / direccion_comercial@bpfintegral.com / direccion_general@bpfintegral.com
Website: www.bpfintegral.com

MOROCCO

Labo City Agdal

Lot 3342 Secteur 3 AL WIFAK ERAQ - TEMARA - MAROC
Contact: Ezzahri Achraf
Phone: +212 (5) 37 64 68 17
Fax: +212 (5) 37 64 68 15
Email: labocity.ag@gmail.com / labocity3.lab@gmail.com

Website: https://labocityagdal.com

NEPAL

Rajshree Scientific Suppliers Pvt. Ltd.

Rajshree Scientific Suppliers Pvt. Ltd.
University Marga, Near Nabil Bank
Infront of Krishi Samagri Snasthan
Kuleshwor, Kathmandu, Nepal
Contact: Shankar Mainali
Phone: +977 1 5372428 / +977 9851195572
Email: rajshreesm@yahoo.com

NICARAGUA

Makol de Nicaragua S.A.

Plaza Eclipse
Next to Holiday Inn hotel
Managua, Nicaragua
Contact: María del Milagro Montero Rojas
Phone: +506 8872 4337
Email: mmontero@makolcr.com
Website: www.makolcr.com

NIGERIA AND CAMEROON

Etcon Analytical and Environmental Systems and Services Limited

445 Herbert Macaulay Way, Yaba
Lagos State, Nigeria
Contact: Osasogie Osazuwa
Phone: +234 8021422237
Email: osasogie.osazuwa@etconanalytical.com / Etconanalytical@gmail.com
Website: www.etconanalytical.com

PANAMA, HONDURAS, GUATEMALA, EL SALVADOR AND BELIZE

Allscience, LLC

8458 NW 70 St. Miami, FL 33166
Florida, USA
Contact: Roberto Shlesinger
Phone: +1 (305) 570-5730
Email: info@e-allscience.com
Website: www.allscience.com

PAKISTAN

Sind Trading Company

Sindh Trading Company
3/14 Arkay Square, Shahrah- e –Liaquat,
Karachi, 74000, Pakistan
Contact: Dr. Shujat Mohsin
Phone: +92 300 8240578 / +92 2132410550
Email: shujatmohsin@gmail.com / info@sindhtradingcompany.com / sind_trading_co@yahoo.com
Website: https://sindhtradingcompany.com/

PARAGUAY AND ECUADOR

Doma International LLC

Doma International LLC
5717 NW 125 Ave.
Coral Springs Florida, 33076.
Email: info@domaininternationalllc.com
Website: www.domaininternationalllc.com

FUJIFILM WAKO LAL DISTRIBUTORS

RUSSIA

Algimed LTD

Barklaya street, 6, Building 5
Moscow, Russian Federation
Contact: Anna Tutnova
Phone: +7 499 682 6109
Email: a.tutnova@algimed.ru
Website: www.algimed.ru

ROMANIA

Elta90 Medical Research

Industriilor Street no. 13, Ilfov County, Chiajna City
Contact: Dohotariu Dragos
Mobile Phone: +40 799767971
Phone: +40 212322694
Email: office@elta90mr.ro
Website: www.elta90mr.ro

SAUDI ARABIA

Fuad Abduljalil Alfadhli and Sons Trading Company

Prince Mashur Ibn Abdulaziz Street
Mursalat, Riyadh, Saudi Arabia
Contact: Mohamed Hussin
Phone: +966 920 002 306
Fax: +966 920 002 406
Email: info@fuadalfadhli.com
Website: www.fuadalfadhli.com

SERBIA, MONTENEGRO AND BOSNIA & HERZEGOVINA RTC d.o.o.

Vladete Kovacevica 14a, 11040 Belgrade, Serbia
Contact: Milica Kangrga
Phone: +381 11 3699 916
Email: office@rtc.rs
Website: www.rtc.rs

SE ASIA

FUJIFILM Wako Pure Chemical Corporation

1-2, Doshomachi 3-Chome,
Chuo-ku, Osaka 540-8605, Japan
Phone: +81 6 6203 3741
Fax: +81 6 6201 5964
Email: ffwk-cservice@fujifilm.com
Website: https://labchem-wako.fujifilm.com/asia

SPAIN

FONTLAB 2000, S.L.

P.I. Can Magre, C/ Carles Buigues, 76
08187 – Santa Eulalia de Ronçana
Barcelona, Spain
Contact: Pelayo Fontsaré
Phone: +34 93 844 60 42
Email: info@fontlab2000.com
Website: www.fontlab2000.com

TANZANIA

Total Med Lab Solutions Ltd.

P.O. Box 33374, Block 24
Plot No. 99 Kinondoni, Dar es Salaam
Tanzania
Contact: Shahida Ngenya
Phone: +255 22 266 7627 / +255 713 631 074
Email: info@tmls.co.tz

TUNISIA

TUNLAB

2 Rue Medenine. CP2000 Le Bardo- Tunisia
Contact: Dr. Bankaji Mahmoud
Phone: +216 71 220 299, +216 20 317 158
Email: mbancaji@gmail.com
Website: www.tunlab.com

TURKEY

Mikromed Medikal Paz. Ith. Ihr. San. Ve Tic. Ltd. Sti.

Mehmet Akif Mah, Elalmis Cad. Asya Sk,No.19 D.3
34774 Umraniye, Istanbul, Turkey
Contact: Esin Tor, Derya Kurfeyiz
Phone: +90 216 668 00 52/42
Fax: +90 216 526 29 43
Email: esin@mikromedltd.com
derya@mikromedltd.com
info@mikromedltd.com
Website: www.mikromedltd.com

UKRAINE

Friedman-Ukraine Ltd.

Pr. Peremogy, 68/1, r.62
Kyiv, 03113, Ukraine
Contact: Vitaliy Melnik
Phone: +38 067 504 20 79
Email: vitaliy@friedman.com.ua
Website: https://friedman.com.ua

UNITED ARAB EMIRATES

Microgen Medical Equipments Trading Est.

Office No. 10, New Industrial Area, Al Ain
United Arab Emirates
Contact: Mr. Labeeb Abdulla
Phone: +971 506956253
Email: sales@microgen-me.com
Website: www.microgen-me.com

URUGUAY, BOLIVIA AND PERU KOMPASS

Prudencio Vazquez y Vega 1112 C.P. 11300
Montevideo, Uruguay
Contact: Guillermo Gonzalez
Phone: +598 2 710 4373
+598 99 167879
Email: info@kompass.com.uy
Website: www.kompass.com.uy

UNITED STATES OF AMERICA

FUJIFILM Wako Chemicals U.S.A. Corp.

1600 Bellwood Road
North Chesterfield, VA 23237

Toll Free: 800-992-9256
Tel: 804-714-1919
Tel: 804-672-4655
Fax: 804-271-7791
Email: wkuspyrostarinfo@fujifilm.com
www.wakopyrostar.com

JAPAN

FUJIFILM Wako Pure Chemical Corp.

1-2 Doshomachi 3-Chome, Chuo-Ku, Osaka
540-8605, Japan

Telephone: + 81-6-6203-3741
Facsimile: + 81-6-6203-2029
<http://ffwk.fujifilm.co.jp/>

GERMANY

FUJIFILM Wako Chemicals Europe GmbH

Fuggerstraße 12
41468 Neuss, Germany

Telephone: + 49-2131-311-0
Facsimile: + 49-2131-311-100
www.wako-chemicals.de