ESC swab

Letheen Broth

(ISO 18593, ISO 17604)

DESCRIPTION

ESC (Easy Surface Checking) swab - Letheen Broth is a device used for the microbiological monitoring of surfaces by swab sampling method. Neutralizing agents included in the medium allow to assess the effectiveness of sanitization procedures in areas where residual disinfectants are presumably present. Letheen Broth is recommended by the AOAC for determining the phenol coefficient of cationic surfactants.

CONTENTS OF THE PACKAGES

The product consists of a transparent plastic tube containing the collection broth and a sterile swab attached to the cap.

Each package contains:

- 100 tubes with swab and liquid medium;
- 1 instructions sheet.

METHOD PRINCIPLE

The medium contains beef extract and peptic digest of animal tissue for supporting recovery of a variety of microorganisms. Neutralizing activity is due to lecithin that neutralizes quaternary ammonium compounds and polysorbate 80 which is effective against phenolic disinfectants and hexachlorophene. Sodium chloride is included to maintain the osmotic balance of the medium.

QUALITY CONTROL

Each batch of ESC swab - Letheen Broth is submitted to quality control using the following strains: *E. coli* ATCC® 25922, *S. aureus* ATCC® 6538, *S. Typhimurium* ATCC® 14028.

TEST PROCEDURE

1. Take a tube from the package and unscrew the cap with an attached swab.
2. Press the tip of the swab against the wall of the tube to remove excess liquid.
3. Streak the swab on a surface horizontally and vertically. A Sampling Template 10x10 (ref. 96762) can be used to delimit and sample an area of 100 cm².
4. Return the swab to the tube.
5. Screw the cap and record date and the sampling point.
6. Transport the device to the laboratory and examine according to methods in use.

STORAGE

10-25°C away from light, until the expiry date on the label. Eliminate if signs of deterioration or contamination are evident.

WARNING AND PRECAUTIONS

ESC Swab - Letheen Broth does not contain hazardous substances in concentrations exceeding the limits set by current legislation and therefore is not classified as dangerous. It is nevertheless recommended to consult the safety data sheet for its correct use. The product must be used only by properly trained operators.

BIBLIOGRAPHY

- ISO 18593:2004 - *Microbiology of food and animal feeding stuffs - Horizontal method for sampling techniques from surfaces using contact plates and swabs*. International Standards Organization, Geneva.
- ISO 17604:2003 - *Microbiology of food and animal feeding stuffs - Carcass sampling for microbiological analysis*. International Standards Organization, Geneva.
- Horwitz W. (2000) *Official Methods of Analysis*. Association of Official Analytical Chemists (AOAC) International. Gaithersburg, MD, USA.

PRESENTATION	Ref.	Contents
ESC swab - Letheen Broth	85607	100 x 10 ml tubes
ESC swab - Letheen Broth	85608	100 x 5 ml tubes
ESC swab - Letheen Broth	85621	100 x 1 ml tubes

TABLE OF SYMBOLS

LOT Batch code	Fragile, handle with care	Manufacturer	Use by
REF Catalogue number	Do not reuse	Contains sufficient for <n> tests	Caution, consult Instruction For Use

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BIBLIOGRAFIA

- ISO 18593:2004 - *Microbiology of food and animal feeding stuffs - Horizontal method for sampling techniques from surfaces using contact plates and swabs*. International Standards Organization, Geneva.
- ISO 17604:2003 - *Microbiology of food and animal feeding stuffs - Carcass sampling for microbiological analysis*. International Standards Organization, Geneva.
- Horwitz W. (2000) *Official Methods of Analysis*. Association of Official Analytical Chemists (AOAC) International. Gaithersburg. MD. USA.

PRESENTAZIONE	Ref	Contenuto
ESC swab - Letheen Broth	85607	Provette 100 x 10 ml
ESC swab - Letheen Broth	85608	Provette 100 x 5 ml
ESC swab - Letheen Broth	85621	Provette 100 x 1 ml

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