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### INSTRUCTION MANUAL

#### North American Catalog Number

US Cord(120V): 82007-202

US Cord(230V): 83007-210

#### European Catalogue Number(s):

Euro Plug: 700-0245

Swiss Plug: 700-0247

UK Plug: 700-0246

Version: 3

Issued: 06 / 20 / 2014

## Legal Address of Manufacturer

### **United States**

VWR International  
1310 Goshen Parkway  
West Chester, PA 19380  
800-932-5000  
<http://www.vwr.com>

### **Europe**

VWR International bvba/sprl  
Researchpark Hassrode 2020  
Geldenaaksebaan 464  
3001 Leuven  
+ 32 16 385011  
<http://www.be.vwr.com>

## Country of origin

**USA**

# Table of Contents

|                                                     |           |
|-----------------------------------------------------|-----------|
| <b>Safety information</b>                           | <b>4</b>  |
| <b>Package Contents</b>                             | <b>4</b>  |
| <b>Unpacking</b>                                    | <b>4</b>  |
| <b>Installation</b>                                 | <b>5</b>  |
| <b>Intended Use</b>                                 | <b>5</b>  |
| <b>Symbols and conventions</b>                      | <b>5</b>  |
| <b>Specifications</b>                               | <b>6</b>  |
| <b>Description of buttons and switches</b>          | <b>6</b>  |
| <b>Operation</b>                                    | <b>7</b>  |
| <b>Troubleshooting</b>                              | <b>7</b>  |
| <b>General maintenance</b>                          | <b>7</b>  |
| <b>User replaceable accessories and spare parts</b> | <b>8</b>  |
| <b>Technical Service</b>                            | <b>8</b>  |
| <b>Warranty</b>                                     | <b>8</b>  |
| <b>Disposal</b>                                     | <b>9</b>  |
| <b>Local VWR offices in Europe</b>                  | <b>10</b> |

## Safety Information

Please follow the guidelines below, and read this manual in its entirety to ensure safe operation of the unit.

- Hazardous samples should be mixed only in appropriate containment vessels.
- Do not mix flammable or explosive samples!

To avoid electrical shock:

- Do not use the mixer near water sources. Be sure that water doesn't get on the device, especially during the cleaning procedures!

Avoiding Damage to the Instrument

- Do not expose the mixer to aggressive vapors (e.g. organic solvents).

## Package Contents

| Description         | Quantity |
|---------------------|----------|
| VWR Mini Blot Mixer | 1        |
| Small Blot Box      | 1        |
| Large Blot Box      | 1        |
| Instruction Manual  | 1        |
| Warranty Card       | 1        |

## Unpacking

Before starting the installation, carefully examine the mixer.

- Open the box and carefully remove the padding. Check that the mixer has not been visibly damaged during transportation. Keep the packing material until you are sure that the mixer works properly.
- Carefully lift the device out of the box.
- Be sure that the following information is included on the serial label located on the bottom of the unit: Serial Number, Product number, Electrical rating.
- Check that the power cord plug has a pin configuration that matches a wall socket according to the local standard.

Should any kind of damage have occurred during transportation, contact the carrier immediately. The carrier is responsible for shipping damage.

## Installation

Place the mixer on a horizontal, level and stable surface. Connect the plug end of the cord to a grounded wall socket. Avoid lines on which powerful electric motors, refrigerators and similar devices are connected. Use the on/off switch to control power to the mixer.

**Use caution when plugging into a grounded wall socket!**

**Do not touch plug with wet hands!**

**Do not unplug by pulling on the cord!**

## Intended use

The VWR Mini Blot Mixer combines the actions of a rocker and a shaker to produce a three dimensional motion that is gentle enough for fragile gels and blots, yet provides thorough mixing. Designed for processing blots and staining applications, the pitch and speed of the mixer are optimally set to allow for use of a minimal amount of solution while preventing the membrane or gel from drying out. The flat rubber mat is ideal for holding blotting boxes or trays.

For research use only. Not intended for any animal or human therapeutic or diagnostic use.

## Symbols and conventions

The following chart is an illustrated glossary of the symbols that may be used in this manual or on the product.

|  |                                                                                                                                                                                                                                               |
|--|-----------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
|  | The electrical warning indicates the presence of a potential hazard which could result in electrical shock.                                                                                                                                   |
|  | <b>CAUTION</b> This symbol refers you to important operating and maintenance (servicing) instructions within the product Instruction Manual. Failure to heed this information may present a risk of damage or injury to persons or equipment. |
|  | This symbol identifies a Protective Earth (PE) terminal, which is provided for connection of the supply system's protective earth (green or green/yellow) conductor.                                                                          |

## Product Specifications

|                                 |                                                                                           |
|---------------------------------|-------------------------------------------------------------------------------------------|
| <b>Electrical requirements:</b> | 115V model 115V $\pm$ 10% - 50/60Hz, 0.03A<br>230V model 230V $\pm$ 10% - 50/60Hz, 0.015A |
| <b>Speed:</b>                   | 115V/230V model 18 rpm / 20 rpm                                                           |
| <b>Motion:</b>                  | 3-Dimensional rocking/rotating                                                            |
| <b>Platform Tilt angle:</b>     | 5°                                                                                        |
| <b>Platform Dimensions:</b>     | 20 x 16.5cm                                                                               |
| <b>Dimensions WxLxH:</b>        | 20 x 16.5 x 17cm                                                                          |
| <b>Platform Load Capacity:</b>  | 0.8kg                                                                                     |
| <b>Platform Capacity</b>        | 4 Small Blotting Boxes<br>2 Large Blotting Boxes                                          |
| <b>Weight:</b>                  | 0.85kg                                                                                    |
| <b>Operating conditions:</b>    | 4°C to 65°C, up to 85% RH, non-condensing                                                 |

## Description of Buttons and Switches

### Main Power Switch

## Operation

Blotting boxes or trays may be placed directly on the rubber mat. When placing boxes, it is important to arrange them evenly to distribute the weight across the platform.

Always load the platform in a uniform and balanced fashion.

Use the On / Off switch to turn the unit on. The direction of rotation in which the unit starts is random and does not affect mixing. If desired, the rotational direction can be reversed by holding the unit's base and very briefly touching the rotating platform on the bottom at one corner. This technique can also be used to select the quietest direction of rotation.

## Troubleshooting

Review the information in the table below to troubleshoot operating problems.

| Problem                                                | Cause            | Solution                                                    |
|--------------------------------------------------------|------------------|-------------------------------------------------------------|
| <b>Mixer does not start.</b>                           | Power connection | Check power connection                                      |
|                                                        | Platform blocked | Check for obstructions blocking the motion of the platform. |
|                                                        | Motor Problem    | Contact VWR Technical service                               |
| <b>Platform continually changes rotation direction</b> | Platform blocked | Check for obstructions blocking the motion of the platform. |

## Repair and maintenance of the VWR Mini Blot Mixer

- No routine maintenance or lubrication is required.
- Keeping the mixer clean will help to maintain it in proper working order

### Cleaning

The outside housing can be cleaned with a soft, damp cloth when needed.

Do not use any aggressive or abrasive cleaners (acetone, nitro, polish etc.) This may permanently damage the mixer finish.

**Before cleaning the mixer, always unplug the power cord from the wall socket!**

## User replaceable accessories and spare parts

The small and large blot boxes are available separately

| Description                | Part Number |
|----------------------------|-------------|
| Small Blot Box, Pack of 10 | 700-0248    |
| Large Blot Box, Pack of 10 | 700-0249    |

## Technical service

### Web Resources

Visit the VWR's website at [www.vwr.com](http://www.vwr.com) for:

- Complete technical service contact information
- Access to VWR's Online Catalogue, and information about accessories and related products
- Additional product information and special offers

**Contact us** For information or technical assistance contact your local VWR representative or visit. [www.vwr.com](http://www.vwr.com).

## Warranty

**VWR International** warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non-performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

## Equipment disposal

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste.

Instead it's your responsibility to correctly dispose of your equipment at lifecycle -end by handing it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

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## Ireland / Northern Ireland

VWR International Ltd /  
VWR International (Northern Ireland) Ltd  
Orion Business Campus  
Northwest Business Park  
Ballycoolin  
Dublin 15  
Tel.: 01 88 22 222  
Fax: 01 88 22 333  
E-mail: sales@ie.vwr.com

## Italy

VWR International PBI S.r.l.  
Via San Giusto 85  
20153 Milano (MI)  
Tel.: 02-3320311/02-487791  
Fax: 800 152999/02-40090010  
E-mail: info@it.vwr.com

## The Netherlands

VWR International B.V.  
Postbus 8198  
1005 AD Amsterdam  
Tel.: 020 4808 400  
Fax: 020 4808 480  
E-mail: info@nl.vwr.com

## Norway

VWR International AS  
Haavard Martinsens vei 30  
0978 Oslo  
Tel.: 02290  
Fax: 815 00 940  
E-mail: info@no.vwr.com

## Poland

VWR International Sp. z o.o.  
Limbowa 5  
80-175 Gdansk  
Tel.: 058 32 38 200 do 204  
Fax: 058 32 38 205  
E-mail: info@pl.vwr.com

## Portugal

VWR International -  
Material de Laboratório, Lda  
Edifício Neopark  
Av. Tomás Ribeiro, 43- 3 D  
2790-221 Carnaxide  
Tel.: 21 3600 770  
Fax: 21 3600 798/9  
E-mail: info@pt.vwr.com

## Spain

VWR International Eurolab S.L.  
C/ Tecnología 5-17  
A-7 Llinars Park  
08450 - Llinars del Vallès  
Barcelona  
Tel.: 902 222 897  
Fax: 902 430 657  
E-mail: info@es.vwr.com

## Sweden

VWR International AB  
Fagerstagatan 18a  
163 94 Stockholm  
Tel.: 08 621 34 00  
Fax: 08 621 34 66  
E-mail: kundservice@se.vwr.com

## Switzerland

VWR International GmbH  
Lerzenstrasse 16/18  
8953 Dietikon  
Tel.: 044 745 13 13  
Fax: 044 745 13 10  
E-mail: info@ch.vwr.com

## Turkey

Pro Lab Laboratuvar Teknolojileri Ltd.Şti.  
a VWR International Company  
Orta Mah. Cemal Gürsel Caddesi  
Ördekcioglu İşmerkezi No.32/1  
34896 Pendik - Istanbul  
Tel.: +90216 598 2900  
Fax: +90216 598 2907  
Email: info@pro-lab.com.tr

## UK

VWR International Ltd  
Customer Service Centre  
Hunter Boulevard - Magna Park  
Lutterworth  
Leicestershire  
LE17 4XN  
Tel.: 0800 22 33 44  
Fax: 01455 55 85 86  
E-mail: uksales@uk.vwr.com

## Australia

VWR International, Pty Ltd.  
Unit 1/31 Archimedes Place  
Murarrie, Queensland 4172  
Tel.: 1300 727 696  
Fax: 1300 135 123

## China

VWR International China Co., Ltd  
Rm.219, 2100 Dongming Road  
Pudong New District  
Shanghai 200123  
Tel.: +86-21-5898 6888  
Fax: +86-21-5855 8801  
E-mail: info\_china@vwr.com

## India

VWR Lab Products Private Limited  
135/12, Brigade Towers, 2nd Floor  
Front wing, Brigade Road,  
Bengaluru, India – 560 025  
Tel.: +91-80-41117125/26 (Bengaluru)  
Tel.: +91-2522-647911/922 (Mumbai)  
Fax: +91-80-41117120  
E-mail: vwr\_india@vwr.com

## New Zealand

Global Science - A VWR Company  
241 Bush Road  
Albany 0632, Auckland  
Tel.: 0800 734 100  
Fax: 0800 999 002  
E-mail: sales@globalscience.co.nz

## Singapore

VWR Singapore Pte Ltd  
18 Gul Drive  
Singapore 629468  
Tel.: +65 6505 0760  
Fax: +65 6264 3780  
E-mail: sales@sg.vwr.com

---

**GO TO VWR.COM FOR THE LATEST  
NEWS, SPECIAL OFFERS AND  
DETAILS OF YOUR LOCAL VWR  
DISTRIBUTOR**

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