VWR Reagenzglasschüttler

Test tube shaker

Agitateur de tubes

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Instruction manual (EN)

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Manual de instrucciones (ES)

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European Catalogue Number(s):

Euro: 444-1372

UK: 444-1378

Swiss: 444-1379

Version: 1

Issue: 07.01.2009

Legal Address of Manufacturer

United States

VWR International
1310 Goshen Parkway
West Chester, PA 19380
800-932-5000
<http://www.vwr.com>

Europe

VWR International bvba
Haasrode Researchpark Zone 3
Geldenaaksebaan 464
B-3001 Leuven
+ 32 16 385011
<http://be.vwr.com>

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Table of Contents

Warning	16
Safety Information	16
Package Contents	17
Unpacking	17
Installation	17
Brief description	18
Setting up	19
Symbols und conventions	20
Product Specifications (Technical Data)	20
Overview	21
Description of Buttons and Switches	22
Troubleshooting	23
Routine cleaning and maintenance	23
Dismantling, Transport and Storage	24
Declaration of Conformity	26
Equipment disposal	26
Technical service	63
Local VWR offices in Europe	64

Thank you for buying this appliance. You have acquired a product that was manufactured and tested in accordance with DIN EN ISO 61010.

We are confident it will enable you to carry out your work easily and to your complete satisfaction.

Warning

Unpack the appliance and parts with care. Compare the contents with the packing list. Check for damage and report any defects or deviations from the packing list to the supplier without delay.

Read the operating instructions carefully and in full and ensure that everyone who operates the appliance has read the instructions carefully before use.

Keep the operating instructions in a place which is accessible to everyone.

This appliance comes fitted with a Euro plug (DIN 49441 CEE 7/VII 10/16 A 250 V).

In Great Britain and Switzerland, this product is sold with the locally approved standard plug. In Great Britain with a BS1363 13 Amps 3-prong plug, in Switzerland with a SEV type 12 plug, approved by SEV1011.

If the appliance is to be connected up to a different grid, an approved adapter must be used or the plug supplied replaced by a qualified electrician with a 3-pin plug approved for the grid in question.

The appliance is earthed upon delivery. If the original plug is replaced, make sure the appliance's protective conductor is connected to the new plug.

Safety Information

Avoiding of Accidents and Damage to the Instrument

Please follow the guidelines below, and read this manual in its entirety to ensure safe operation of the unit.

VWR recommends against the use of:

Avoiding of Accidents

Observe all safety and accident prevention regulations in the laboratory!

Wear safety goggles and suitable working clothes when working with the shaker.

Extreme care is called for when handling toxic or other hazardous substances.

Avoiding Damage to the Instrument

Observe all safety and accident prevention regulations in the laboratory!

Avoiding Damage to the Instrument and Accidents

The motor works without sparks. Nevertheless, the manufacturer cannot guarantee that the appliance is explosion-safe.

Repairs must be carried out only by technicians authorized.

To prevent the shaker sliding off the table top as a result of the vibrations generated during operation, make sure the feet are securely attached to the table top.

Package Contents

Please make sure that the consignment contains the following parts:

Description	Quantity
Test tube shaker REAX top ECN 444-1372 (EU) ECN 444-1378 (UK) ECN 444-1379 (CH)	1
tray, small ECN 444-1382	1
Main lead (supplied loose) 14-007-003-81 (EU) 14-007-003-92 (CH) 14-007-005-45 (UK)	1
Instruction manual 01-005-004-06	1
Accessories only if ordered Tray, large ECN 444-1383	

Unpacking

Unpack the appliance and parts with care. Compare the contents with the packing list. Check for damage and report any defects or deviations from the packing list to the supplier without delay.

Installation

1. Setting up the appliance

Set up the test tube shaker on a smooth and solid table top. The weight of the appliance and the 4 parabolically shaped rubber feet make for stability.

The feet can adhere securely to the table top only if both the feet and the surface to which they are to be attached are clean. If necessary, clean the feet with a damp cloth and ethanol.

2. Electrical connection

Before connecting the appliance into the mains, make sure that the mains voltage matches that required on the type plate:


- (for other plug standards, see the General Information on page 16).
- the ON/OFF switch is switched off (in the “0” position).
- the speed setting knob is turned to the left as far as it will go.

Plug the shaker into the mains using the mains lead. The plug for the mains lead is at the rear of the appliance.

3. Assembly and accessories

Tray, small (accessory ECN 444-1382) and tray, large (accessory ECN 444-1383)

The tray is simply clipped onto and pulled off the appliance by hand. When clipping on the tray, make sure the slot in the coupling area of the tray is in line with the lug on the eccentric of the appliance.

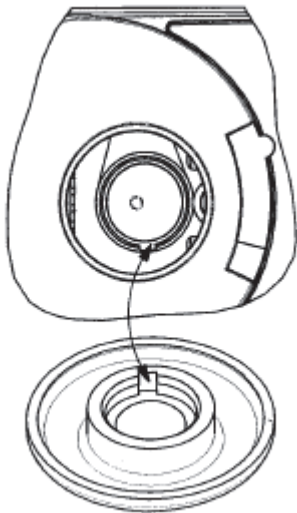Tray-Assembly

Brief description

Test tube shaker is equipped with speed control.

Shaking is actuated by the pressure of the test tube on the tray.

What is new is the practical switchover function from periodic to continuous operation.

Setting up

ON/OFF switch and setting speed

The appliance is switched on at the ON/OFF switch. The green control lamp in the ON/OFF switch indicates that the appliance is ready for operation.

The speed (shaking intensity) is selected using the speed setting knob. It can be increased by turning this knob to the right.

The test tube shaker has a dial.

Shaking with tilt operation

Before starting the shaker, select a low speed to prevent splashing when the appliance starts up.

The operating mode is selected with the slide switch. To select tilt operation, turn the slide switch to the right as far as it will go.

The symbol ▼ to the left of the slide switch indicates the operating mode.

Shaking is actuated by the pressure of the test tube on the tray and stops again automatically as soon as the test tube is removed.


Shaking in continuous operation


Before starting the shaker, select a low speed to prevent splashing when the appliance starts up.

The operating mode is selected with the slide switch. To select continuous operation, turn the slide switch to the left as far as it will go. The symbol ● to the right of the slide switch indicates the operating mode.

In this position, the tray is shaken continuously.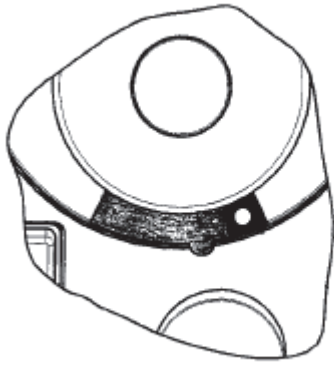

Operation: continuous operation

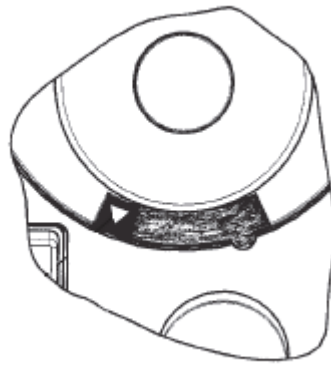tilt operation

Symbols and conventions


The following chart is an illustrated glossary of the symbols that are used in this manual.

	CAUTION This symbol indicates a potential risk and alerts you to proceed with caution
--	--

Product Specifications (Technical Data)

Order number		ECN 444-1372 (EU) ECN 444-1378 (UK) ECN 444-1379 (CH)
Type of motion		circular and vibrating
Shaking frequency	[1/min]	0-2400
Orbit [mm]		5
Speed display		dial
Control precision		uncontrolled
Switch		for periodic and continuous
Connected load	[W]	51
Unit weight	[kg]	2,8
Dimensions (B x H x D)	[mm]	134 x 105 x 172
Voltage/frequency		230V/50Hz (others also available)
Permissible duty factor		100%
Overheating protection prolonged		The motor switches off after overloading with self resetting function
Degree of protection		IP 22
Permissible ambient temperature		5-40 °C at 85% relative humidity, no condensation

Overview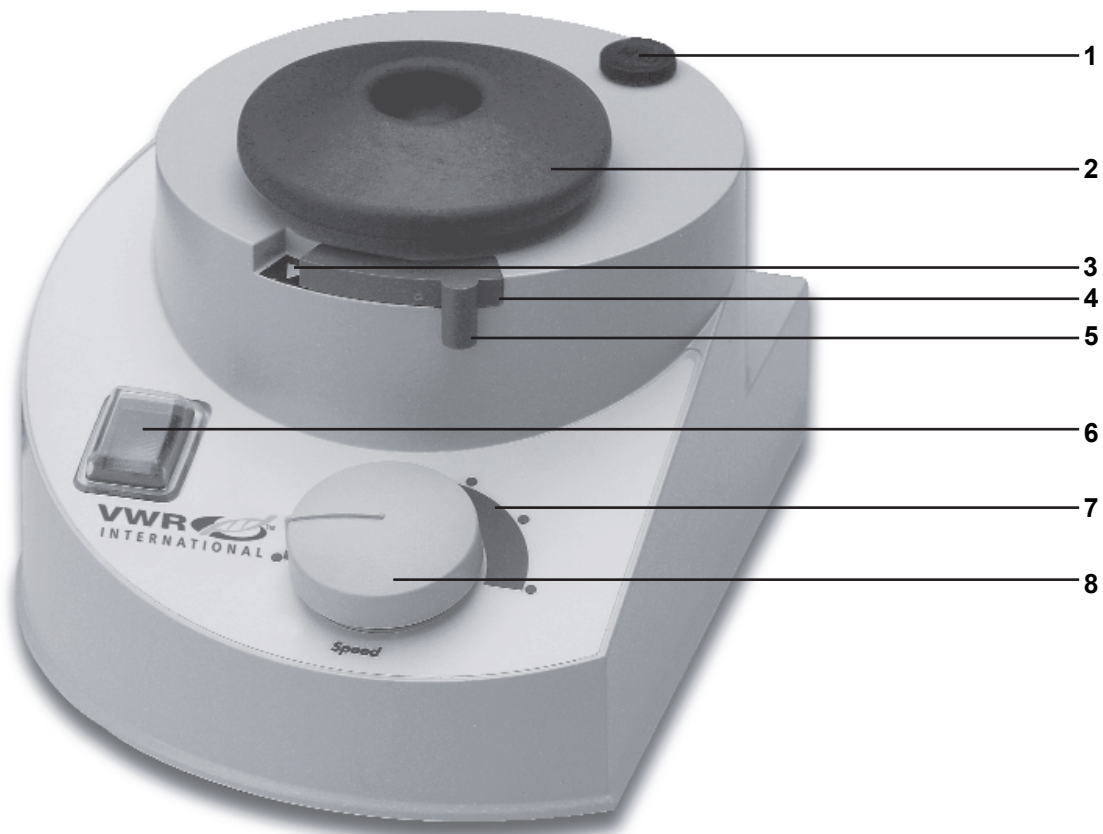

- 1: Befestigungsgewinde / attachment thread / filet de fixation / rosca de sujeción
- 2: Aufnahmeteller / tray / disque de logement / disco de alojamiento
- 3: Symbol ▼ / symbol ▼ / symbole ▼ / símbolo ▼
- 4: Symbol ● / symbol ● / symbole ● / símbolo ●
- 5: Schiebeschalter / slide switch / interrupteur à coulisse / interruptor deslizable
- 6: Netzschalter / ON/OFF switch / interrupteur du réseau / interruptor de red
- 7: Skala / Dial / échelle / escala
- 8: Drehzahleinstellknopf / speed setting knob / vitesse de rotation / botón de ajuste


Description of Buttons and Switches

Front View of ON/OFF switch

The appliance is switched on at the ON/OFF switch. The green control lamp in the ON/OFF switch indicates that the appliance is ready for operation.

Front View of operation switch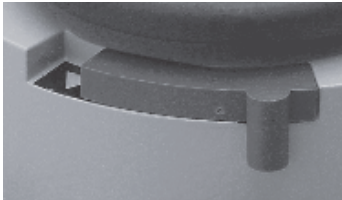

The operating mode is selected with the slide switch. To select tilt operation, turn the slide switch to the right as far as it will go. The symbol ▼ to the left of the slide switch indicates the operating mode.

To select continuous operation, turn the slide switch to the left as far as it will go. The symbol ● to the right of the slide switch indicates the operating mode.

Front View of setting speed switch

The speed (shaking intensity) is selected using the speed setting knob. It can be increased by turning this knob to the right. The test tube shaker has a dial.

Troubleshooting

Review the information in the table below to troubleshoot operating problems.

Problem	Cause	Solution
No shaking movement		Check the mains lead and power supply.
	The ON/OFF switch is not switched on (see Operation, page 19).	
	The slide switch is not in the right position (see Operation, page 19.).	
	The speed selected is too low.	increase speed
	In tilt operation, not enough pressure on the tray.	increase pressure by hand
	In tilt operation, the reaction vessel exerts pressure on the tray at an angle.	switch over to continuous operation, if necessary (see Operation, page 19.)
	The motor is defective.	(requires service)
	The switch is defective	(requires service)
	The control electronics are defective.	(require service)

Routine cleaning and maintenance

Use a damp cloth (mild soap solution) to wipe clean the housing and surfaces of the appliance.

Important: Never use chlorine bleach or chlorine-based detergents, scouring agents, ammonia, cleaning wool or cleaning agents with metallic components to clean the appliance, as these would damage the surfaces of the appliance.

The appliance is maintenance-free. Any repairs which may be necessary must be carried out by an authorized technician.

Please contact your local VWR International company.

Avoid of Accidents and Damage of the Instrument

To ensure the feet adhere securely to the table top, clean the feet at regular intervals using a damp cloth and ethanol.

Dismantling, Transport and Storage

Dismantling

1. Switch off the appliance and disconnect it from the mains (allow to cool).
2. Remove any apparatus around the shaker so that the appliance can be dismantled without difficulty.
Remove the glass set and dismantle the entire appliance and all the parts in the reverse order to assembly.

Transport and storage

1. Place the appliance and the parts in the original packing or in another suitable container in order to avoid damage during transit. Seal the packing with adhesive tape.
2. Keep the appliance in a dry place.

Caution: Avoid knocks and vibrations when transporting the appliance.

Warranty and Liability

Warranty

VWR International grants a three-year warranty on the products described here (except worn parts). This warranty covers faulty materials and workmanship.

Damage incurred in transit is not covered by the warranty.

In the event of a warranty claim, please notify VWR International. If the claim is for faulty materials or workmanship, VWR International will restore your appliance to good working order by repair or replacement. VWR International's warranty does not cover damage caused by improper use.

Any changes in this warranty declaration must be confirmed in writing by VWR International to be effective.

Exclusion of liability

VWR International cannot be held liable for any damage caused by the improper treatment and use of the appliance.

There shall be no liability for consequential damage.

If, after reading the operating instructions, you have any questions on the installation, operation or maintenance of the appliance, please contact:

VWR International nv/sa
Geldenaaksebaan 464
B-3001 Leuven / Belgium
Internet address:
www.vwr.com

If repairs are required, please contact your local VWR International company

Declaration of Conformity

We do hereby declare, that this product complies with the following standards or standardizing documents:

The German EMC Act:

EN 61326: 1997 + A1:1998 + A2:2001+ A3 2003

EN 61000-3-2: 2000

EN 61000-3-3: 1995 + 1997 + A1:2001

EN 61000-4-3:2002 + A1:2002

EN 61000-4-5:1995 + A1:2001

EN 61000-4-6:1996 + A1:2001

EN 61000-4-8: 1993

EN 61000-4-11:1994 + A1:2001

Low Voltage Act:

(73/23/EWG):

EN 61010-1 + EN 61010-2-051

Equipment disposal

Please make sure that any old or defective appliances are taken to a collection point and disposed of professionally. Collect the metal, glass and plastic parts separately.

Please also dispose of the packing materials in an environment-friendly manner (sorting).

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Technical service

Web Resources

Visit the VWR's website at www.vwr.com for:

- Complete technical service contact information
- Access to VWR's Online Catalogue, and information about accessories and related products
- Additional product information and special offers

Contact us For information or technical assistance contact your local VWR representative or visit.

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*14 Cent/Minute aus d. dt. Festnetz

Ireland

AGB Scientific Ltd - A VWR
International Company
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: 01 88 22 222
Fax: 01 88 22 333
e-mail info@agb.ie

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We reserve the right to make technical Changes without prior announcement

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Garantie / Responsabilité civile / Copyright
VWR International AB
VWR International AG
VWR International Ltd