

VWR® Gel Documentation system Basic Installation – Quick Guide

European Catalogue number: 1730-1470

■ System components

1. Camera
2. Lens
3. Hood
4. Bracket
5. UV filter
6. +4 dioptre filter
7. Screw and washers

■ Installation

To attach camera lens

1. To remove the lens cover press the black button next to the lens slot and rotate anti-clockwise. The lens cover also comes off by screwing the top anti-clockwise. The camera and the lens have corresponding white spots. Line these up and screw the lens on clockwise. Remove front lens cap

Camera set-up

1. To charge battery please refer to camera manual
2. Screw on the UV filter to the front of the lens
3. Install the camera software next. To do this refer to the instructions in the camera manual
4. It is recommended to have the camera set to take black and white images (this should have been already set), the camera should be in 'P' mode. To set B/W imaging turn on the camera making sure the settings indicator is on 'P'. Press the "Q" button and then move down the down arrow to Picture Style . Press Set then select monochrome. Then press Set to store
5. To use the camera in colour mode return the main selector to Auto
6. To use the camera in colour mode move the mode dial to the green square auto-mode



Hood assembly and camera set-up


1. Attach the camera on the bracket using the screw provided

2. To attach the camera to the bracket insert the screw through the bracket and washer between the screw and the bracket (2)
3. Align the camera with the hole of the hood (making sure there are no gaps between the black plinth and the camera lens).
4. The system is now ready for use, place the camera/hood assembly on the transilluminator over the gel you want to image.



Installing software

1. To install the software you need to be connected to the internet. From a computer access the internet and go to the following website www.canon.com/icpd
2. Select your country from the list
3. Select the support tab and enter model of camera e.g. EOS 4000D
4. Select the software tab and select to download EOS software
5. Enter your serial number which can be found on the bottom of the camera
6. Double click on the EOS camera software icon from your desktop
7. Select Remote Shooting from the menu

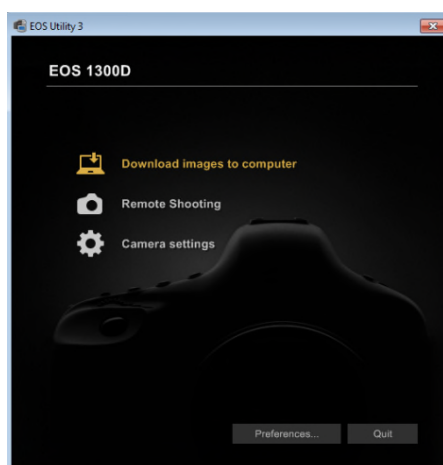

N.B. WiFi must be set to disabled to be able to connect to the PC with a USB cable. When you turn on the camera the software will start automatically

CONTACT VWR:
For further details , please consult the camera manual.
Visit www.vwr.com to find your local VWR contact

