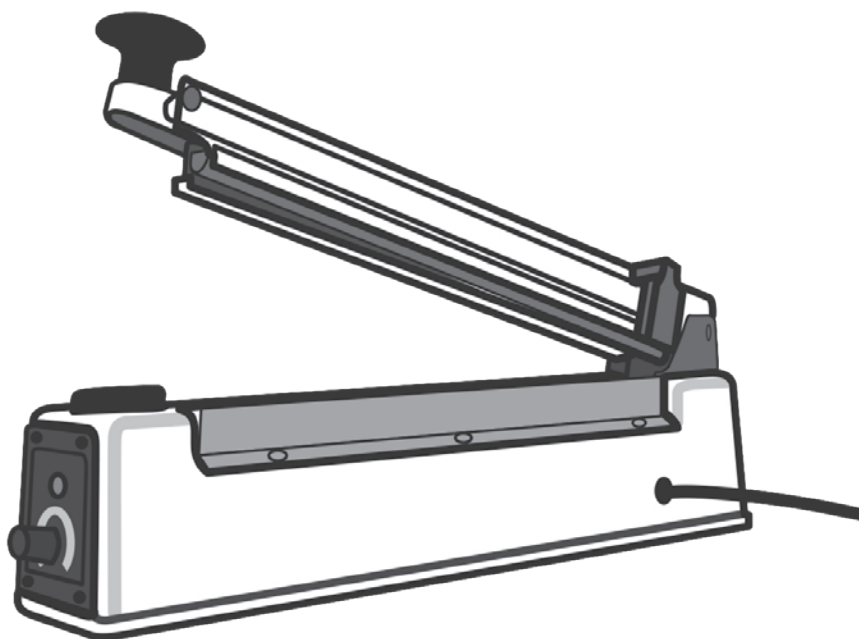
Electric Bag Sealer with Cutter

Instruction Manual

Instruction Manual.....	2
Manual de Instrucciones	9
Manuel d'instructions.....	17
Manuale d'istruzione.....	25
Bedienungsanleitung.....	33

Legal Address of Manufacturer		Country of Origin
United States	Europe	China
VWR International, LLC 100 Matsonford Rd Radnor, PA 19087 800-932-5000 http://www.vwr.com	VWR International bvba Researchpark Haasrode 2020 Geldenaaksebaan 464 B-3001 Leuven + 32 16 385011 http://be.vwr.com	

Catalog Number

Bag Sealer Type	Voltage	Catalog Number
U. S. Plug 	110/120 VAC	490007-638
European Plug 	220/230 VAC	129-1160
UK Plug 	220/230 VAC	129-1159
Replacement Parts	Blades (2) Heating element (2) PTFE cloths (2)	129-1177

Table of Contents

Warning/Safety Information	3
Package Contents	3
Unpacking & Installation	3
Components.....	4
Operating Instructions.....	4
Heating Adjustment Guide	5
Maintenance	5
Technical Service	6
Warranty	6
Compliance	7
Equipment Disposal	7
Appendix.....	8

Warning/Safety Information

This manual cannot address all safety hazards. It is the responsibility of the user to consult and observe all health and safety precautions and to assess the suitability of the sealer for the task.

1. Read the instruction manual before operating the sealer and cutter.
2. Observe general laboratory safety precautions and regulations.
3. Follow all precautions from the Material Safety Data Sheet (MSDS) for any reagent you use with this sealer.
4. Never operate the sealer in a hazardous or flammable environment.
5. To protect against risk of electrical shock, do not put sealer in water or other liquid.
6. Blade is extremely sharp. Keep hands, hair, clothing and other utensils away from blade during operation.
7. Always work in a manner which neither endangers the user, nor any other person.
8. Unplug the unit from the outlet before changing parts or cleaning.
9. Only use original manufacturer's accessories and spare parts.
10. Do not pull, twist or otherwise abuse the power cord. Never carry the sealer by the cord or pull from the cord to remove from an outlet. Always grasp the plug and pull to disconnect from an outlet.
11. Repairs are to be performed by trained and authorized service personnel only.

Package Contents

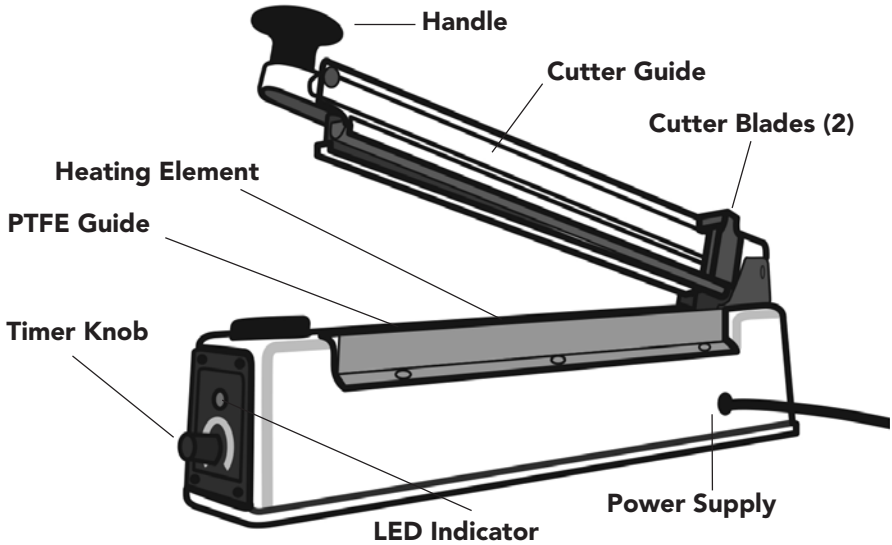
- 1 heat sealer
- 2 extra blades
- 2 heating elements
- 2 PTFE cloths

Unpacking and Installation

1. Remove all packing materials from the unit.
2. Plug the unit into the outlet.
3. See Operating Instructions for use.

Thank you for purchasing our 305mm (12") Electric Bag Sealer with Cutter. You will be able to quickly and efficiently seal polyethylene or polypropylene bags, sheets or flat tubing. A unique sliding blade removes excess material 19 mm from the seal.

Components

Operating Instructions

1. Set the timer based on the thickness of the material to be sealed.
(see Heating Adjustment Guide)
2. Sealing time is between 1 and 8 seconds. Use a higher number for thicker materials.
3. Place the area to be sealed across the bar. Press the handle down evenly.
4. Heating is only applied while the light is on.
5. When the light turns off either slide blade across to cut away excess material or lift handle to remove the item.
6. Adjust the timer if needed and repeat steps.

Heating Adjustment Guide

INDICATOR	POLYETHYLENE	POLYPROPYLENE OR HIGH-HEAT FILM
1	Under 0.06 mm	—
2	Under 0.1 mm	—
3	Under 0.14 mm	—
4	Under 0.2 mm	Under 0.03 mm
5	—	Under 0.04 mm
6	—	Under 0.06 mm
7	—	Under 0.08 mm

Maintenance

1. Unplug the unit.
2. On the base, loosen the 6 side screws that hold the PTFE cloth over the heating wire.
3. Remove the PTFE cloth; may need to replace.
4. The heating wire is held in place by two black plastic holders. Remove the screw on end with the timer knob.
5. Remove the wire on that end.
6. Pull the wire toward the timer knob end of the unit.
7. Place an extra screwdriver or pen in the opening behind the other back holder; on the end with hinge. This will keep that piece in place.
8. Remove the screw and the rest of the wire.
9. Replace the wire.
10. Screw the wire in place.
11. Remove the extra screwdriver or pen.
12. Replace or put on a new PTFE tape.
13. Tighten the 6 side screws.
14. Plug in the unit.
15. Check unit operation.

Technical Service

Web Resources

Visit the VWR website at www.vwr.com for:

- Complete technical service contact information
- Access to the VWR Online Catalogue, and information about accessories and related products
- Additional product information and special offers

Contact us for information or technical assistance contact your local VWR representative or visit. www.vwr.com.

Warranty

VWR warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non-performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

Compliance with local laws and regulations

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorizations necessary to run or use the Product in its local environment. VWR will not be held liable for any related omission or for not obtaining the required approval or authorization, unless any refusal is due to a defect of the product.

Equipment disposal

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste.

Instead it's your responsibility to correctly dispose of your equipment at lifecycle -end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

Thank you

Appendix

Local VWR offices in Europe and Asia

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India

VWR Lab Products Private Limited
No.139. BDA Industrial Suburb,
6th Main, Tumkur Road,
Peenya Post,
Bangalore, India – 560058
Tel.: +91-80-28078400
Email: vwr_india@vwr.com

Ireland / Northern Ireland

VWR International Ltd / VWR
International (Northern Ireland) Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: +353 01 88 22 222
Email sales.ie@vwr.com

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VWR International AB
VWR International GmbH