

SANI-SPONGE:

The SANI-SPONGE kit is designed to collect samples to detect the presence of microbiological contaminations such as *Listeria*, *Salmonella*, *E. coli*, and other food-borne pathogens on almost any surface.

SANI-SPONGE sampling kits are produced in accordance with USDA, ACIA and CFIA requirements.



STEP BY STEP INSTRUCTIONS

- 1 Remove the tear-off strip and carefully take the sponge out (sterile gloves are recommended). Do not touch the inside of the bag.
- 2 Swab the sampling surface (suggested 30x30 cm area (12"x12") with the sponge in a vertical (up and down) direction.
- 3 Turn the sponge over and swab the sample area again, this time in a horizontal (left to right) direction.
- 4 Return the sponge to the sterile sampling bag with care.
- 5 Identify the sample via the write-on area with a permanent marker.

