

VWR® Gel Documentation system Basic Image Capture - Quick Guide

European Catalogue number: 1730-1470

1 Turn on

Select




→ Turn on the camera (ensure the battery is charged)

→ The Canon control software launches automatically when you turn on the camera. If the software does not start automatically click on Start/All Programmes/Canonutilities/EOS Utility/EOS Utility

2 Capturing an image

Select

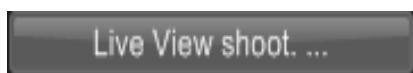

→ Select Live View shoot to open live window

→ Ensure the system has been installed correctly, as per the Installation Quick Guide

3 Settings

Select



Select



→ From the White Balance menu select Cloudy from the drop down menu

4 Turn on transilluminator


Select



→ If bands on gel are faint then increase the white balance (From the EOS 1200D window right click on the control to activate it). For very faint bands go to +2

5 Focus setting

Select

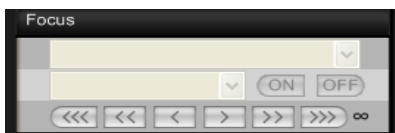

From the Remote Live View window

→ To adjust zoom turn the lens of the camera

→ To adjust focus use the arrows

6

Grab image

Press



From the EOS 1200D window

- ➔ The camera will set an exposure time relevant to the light levels coming from the sample
- ➔ Then turn off the camera

7

Save and export

- ➔ Images are saved as JPEG files
- ➔ Export as TIFF or BMP

CONTACT VWR:
For further details , please consult the camera Manual
Visit www.vwr.com to find your local VWR contact



BIC0616