Instruction for use

VWR® FACE MASK 3-PLY EARLOOPS BLUE

EU cat. No 113-0716

EN

Intended use

The VWR face masks are designed to reduce process and products contamination in controlled environments.
For use in industrial environment only

Read and Understand before Use.

- 1 This product is not a medical device in context of Regulation (EU) 2017/745/Directive 93/42/EEG nor personal protective equipment in context of Regulation (EU) 2016/425!
- 2 This product does not prevent Coronavirus transmission and is not intended for general public!
- 3 Before use, the wearer should always check the mask is in good condition - not dirt, no damage to head strap attachment etc.
- 4 Facial hair and certain facial characteristics will prohibit effective use of this product.
- 5 Please make sure the mask fit properly before entering working environment.
- 6 Those masks are for single use. Do not reuse nor clean!
- 7 Dispose the products according to applicable local legislations

General Specifications

- Made of non-woven fabric 20g PP/20g MB/25g PP
- Blue colour
- Ear loops
- Bacteria filtration rate: 99.7% in average (EN14683:2019+AC:2019)
- Breathability: 28.8Pa/cm² in average (EN14683:2019+AC:2019)
- Nature Rubber Latex Free

Packaging and Storage Specifications

Package/Quantity	50pcs/box, 20boxes/case
Best storage Conditions	0° - 50°C
Best storage Conditions	45% - 85% humidity
Shelf Life	without sunshine exposure
Country of Origin	3 years
	China

Fitting Instruction

- 1 Put loops around ears.
- 2 Press forefingers down on both sides of nose.
- 3 Adjust mask over nose and under chin.

WARNING!

For Industrial and clean room use only; do not use for protection against COVID-19!
This product is not a medical device for medical or clinical use within the meaning of Regulation (EU) 2017/745 / Directive 93/42/EEC!
This product is not a personal protective equipment (PPE) within the meaning of Regulation (EU) 2016/425!

Imported by VWR International bv
Researchpark Hazeleindree 2020, Gidselandsbosboom 464, B-3001 Leuven - vwr.be@vwr.com

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