VWR® Thermometer

INSTRUCTION MANUAL

North American Catalog Numbers: 61220-439 / 61161-310 /

23609-172 / 77776-720 / 37000-428 / 37000-426 / 62433-734

European Catalogue Numbers: 620-2020 / 620-2024 / 620-

2082 / 620-2095 / 620-1665

Version: 1

Issued: 27th April 2017

Legal Address of Manufacturer

United States

VWR International
Radnor Corporate Center
Building One, Suite 200
100 Matsonford Road, P.O. Box 6660
Radnor, PA 19087
800-932-5000
<http://www.vwr.com>

Europe

VWR International bvba
Researchpark Haasrode 2020
Geldenaaksebaan 464
B-3001 Leuven
+ 32 16 385011
<http://be.vwr.com>

Country of origin - China

Intended use: This item is intended for general laboratory use to monitor temperature.

61220-439 / 620-2020 Traceable® Digital Dial Thermometer

61161-310 / 620-2024 Traceable® Jumbo-Display Dial Thermometer

Specifications

Range: –58 to 302°F / –50 to 150°C

Accuracy:

Cat# 61161-310 / 620-2024- ±1°C between –20 and 100°C

Cat# 23609-172 / 620-2082- ±1°C between –20 and 100°C

Cat# 77776-720 / 620-2095- ±0.3°C between –20 and 100°C

Resolution:

0.1° above –20° and below 200°

1° below –20° and above 200°

Sampling Rate: 1 second

Ultra™ Thermometer Accuracy

Ultra™ thermometers are tested at selected test points to be within tighter than normal tolerances to assist in providing improved accuracy. Other points will not necessarily fall within the same accuracy as those found at the selected test points, but will be within an accuracy of ±1°C.

Operation

Slide the ON/OFF switch to ON to turn the unit on. Slide the °C/°F switch to °F to display the temperature in Fahrenheit or °C to display the temperature in Celsius. Insert the stem into the material to be measured. Slide the ON/OFF switch to OFF to turn the unit off. To conserve battery life, always turn the unit off when not in use.

All Operational Difficulties

If this unit does not function properly for any reason, replace the battery with a new high-quality battery (see “Battery Replacement” section). Low battery power can occasionally cause any number of “apparent” operational difficulties. Replacing the battery with a new fresh battery will solve most difficulties.

BATTERY REPLACEMENT

Erratic readings, a faint display, or no display are all indications that the battery must be replaced. To replace the battery, remove the battery cap screw. Remove the battery cap by turning it in the direction of the arrow (counter clockwise). Remove the exhausted battery and replace it with a new 1.5 volt size G-13 battery. Equivalent replacement batteries are: Ray-O-Vac RW42, Duracell D357, Eveready 357. Make certain that the positive (+) side is facing you. Replace the battery cap. Replace the battery cap screw and tighten securely.

Equipment disposal

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste. Instead it's your responsibility to correctly dispose of your equipment at lifecycle -end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment. For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

Compliance with local laws and regulations

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorizations necessary to run or use the product in its local environment. VWR will not be held liable for any related omission or for not obtaining the required approval or authorization, unless any refusal is due to a defect of the product.

Technical service

Web Resources

Visit the VWR website at vwr.com for:

- Complete technical service contact information
- Access to the VWR Online Catalogue, and information about accessories and related products
- Additional product information and special offers

Contact us For information or technical assistance contact your local VWR representative or visit vwr.com.

Warranty

VWR warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

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VWR International AB
VWR International AG
Teknolojileri Ltd.Şti.
VWR International Ltd