RFM700-M Series Refractometer

Thank you for purchasing this instrument. In order to ensure that this product provides many years of service please follow the guidance in this document. To register your instrument with the manufacturer and print a warranty certificate visit:

www.bellinghamandstanley.com

Unpacking

Carefully remove all of the packing material. It is recommended that the box and other packing materials are retained so that, should the need arise, the instrument can be safely returned to the supplier for service. Check that all parts listed below are present and that no transit damage has occurred. If any are damaged or missing, contact the supplier immediately.

Contents list

Quantity	Item	Code
1	RFM700-M Series Refractometer	See below
1	Mains leads	55-109
1	Power supply	55-105
1	Quick Start Guides	19-405

RFM700-M Refractometer code numbers:

Model	Complete refractometer	Refractometer module only
RFM712-M	19-00	19-712
RFM732-M	19-10	19-732
RFM742-M	19-20	19-742

Positioning the system

- Dry and indoors.
- Capable of holding the instrument's weight.
- Away from draughty or hot equipment like fans or heaters.
- Out of direct sunlight or strong ambient light.
- Away from potential sources of interference, such as RFI generating equipment.
- Within reach of a power point.
- Not using a power circuit that also has large motors or noise generating equipment connected to it.

Mains connection

The power supply adapter is supplied with a moulded mains cord and plug to suit one of several socket types. For UK leads, replace the fuse only with the type indicated on the plug.

- Voltage 110 to 230 V ~ ±10%
- Maximum current 2 A
- Frequency 50 to 60 Hz

RISK OF ELECTRIC SHOCK:

- For indoor use only.
- Must be kept dry.
- Disconnect the equipment from the mains supply before unplugging the mains lead from the instrument.
- Do not cover, designed to operate with free air convection.
- No cleaning required.
- Do not jet wash

QUICK START GUIDE

19-405'05 REV A, OCT-2023

Bellingham + Stanley, a Xylem brand, Xylem Analytics Germany GmbH, AM Achalaich 11, 82362 Weilheim, Germany

Email: sales.bs.uk@xylem.com

Switching on and off

Plug in the power supply and switch on the mains supply, after startup the instrument will show a loading screen indicating the software version. Once loaded, the instrument will be in standby mode. To switch on, press the green Power softkey.

To switch to standby mode press and hold the green Power softkey for several seconds and select Yes on the confirmation screen. It is then safe to turn off the mains supply and disconnect the power lead.

Navigating through menus

The softkey functions change depending upon what inputs are required at that time. If a softkey has no text then that button is disabled.

Menu items can be selected by either:


- Pressing the Up/Down softkeys until it is highlighted and then pressing Select.
- Pressing the item number key.

Alternatively, a USB keyboard, available from Bellingham + Stanley, can be used to select items and navigate menus.

If the menu item contains a list of options to choose from then either:

- Select the menu item and then choose the option from the displayed list.
- Press the Forward/Back arrow keys to step through the options.

Menu items that can be modified by using the arrow keys are identified by arrows being displayed next to the menu items value.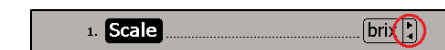

Return to the Measurement Display by pressing Quit to step back through the menus.

Setup Wizard

When the instrument is switched on for the first time, a Setup Wizard will step through a number of questions so that the system can be configured for a particular application. First select the desired language by pressing the associated number key. Press the Next softkey to access each setting screen:

Copy user guides and PC software to USB: Makes copies of the full User Guides and PC software compatible with the instrument to a removable USB storage device. Insert a USB removable storage device before selecting. Full User Guides provide detailed instructions on instrument operation.

Set the standard measurement conditions: Make changes to the Measurement Settings.

Set clock: Set the instrument to local time.


Set the printer options: Specify how reading results will be printed.

When the desired conditions are set in each screen press the Quit softkey to proceed.

Zero

The Zero calibration is displayed in °Brix with sucrose temperature compensation. Distilled water must be used. The instrument should be Zeroed regularly to ensure accurate readings.

- Press and hold the Zero softkey.
- Follow the onscreen instructions – ensure that the prism is thoroughly clean.
- Apply a sample of distilled water.
- The complete prism surface must be covered, 1ml is generally sufficient to achieve this.
- Press OK.
- The Zero point will then be set.

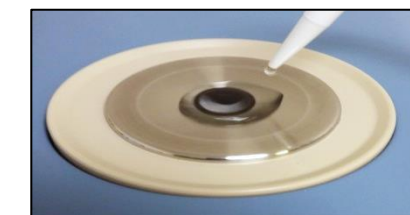Reading

- Ensure that the prism is thoroughly clean.
- Apply a sample to the prism, the complete surface must be covered, 1ml is generally sufficient to achieve this.
- A progress bar will appear on the Read softkey.
- A reading will be initiated once the progress bar is full.
- Push the read softkey to take further readings of the same application.

This symbol is an internationally agreed indicator that the product bearing it should not be disposed of as general waste or garbage which might end up in landfill sites, but should instead be sent for special processing and/or recycling in those countries where appropriate legislation and facilities are in place.

Full User Guides can be downloaded from the instrument's Help Menu or Setup Wizard

Caution or warning.

Electrical hazard.

RFM700-M Series Refractometer

Measurement Display
From the Measurement Display readings can be taken, viewed and printed.

- 1. Reading result:** After taking a reading, the final result will be displayed here.
- 2. Time:** The current time.
- 3. Temperature:** The temperature that the prism was at when the reading was recorded.
- 4. Quality:** The Quality figure is derived from the optical pattern that falls upon the instrument detection system. A high Quality value indicates that the optical pattern is well-defined, making the reading signal easy to resolve. A low Quality value, caused for instance by an opaque sample or poor sample application, means a less well defined optical pattern so the reading will be less reliable. A Quality value of 100 is set when zeroing the instrument, which can then be used as a reference with which to compare other measured samples.
- 5. Measurement settings:** The scale and temperature compensation used for the reading.
- 6. Softkeys:** Indicates the action that will happen when a softkey is pressed.

Read

Hold for 2 seconds to switch off the instrument.

Print

Press the softkey to print the currently displayed reading.
Hold for 2 seconds to change display brightness.

Zero

Hold the softkey for 2 seconds to initiate a Zero calibration. Follow onscreen instructions.

Settings

Hold the softkey for 2 seconds to change the Measurement Settings. The Supervisor Settings can also be accessed from here.

This symbol is an internationally agreed indicator that the product bearing it should not be disposed of as general waste or garbage which might end up in landfill sites, but should instead be sent for special processing and/or recycling in those countries where appropriate legislation and facilities are in place.

Full User Guides can be downloaded from the Instrument's Help Menu or Setup Wizard

Caution or warning.

Electrical hazard.

Menu Structure

	3. Supervisor Settings...	
	1. Scale	1. refractive index
	2. Temp Comp.	1. none 2. brix
1. Calibration		
1. Zero		
2. Span (Top Calibration)		
3. Undo		
3. Setup		
1. Setup Wizard		
2. Passwords		
1. operator PIN - 123		
2. supervisor PIN - 000		
3. Reading		
1. Recurrence		
1. single		
2. continuous		
3. auto single		
2. Resolution		
1. low		
2. medium		
4. Recording		
1. Mode		
1. none		
2. single		
3. save		
4. single + save		
2. Printer		
1. none 5. network		
2. serial		
3. usb		
4. pdf		
3. Record After Read		
1. no		
2. yes		
4. Traceability		
1. Batch Codes		
1. off		
2. number		
3. text		
4. dateinc		
2. Auto Increment		
1. off		
2. on		
3. Log Users		
1. off		
2. number		
3. text		
4. Remember Text		
1. none		
2. batch		
3. user		
4. batch and user		
5. Ports		
1. Serial Port		
1. Baud Rate		
2. Word Length		
1. DHCp		
2. Network		
1. IP Address		
3. Subnet Mask		
4. Default Gateway		
3. CSV (Lims)		
1. none		
2. serial		
3. network		
4. usb		
5. System		
1. Language		
2. Time / Date		
1. Date Format		
1. dd/mm/yy		
2. mm/dd/yy		
3. Summer Time		
1. no		
2. yes		
3. Display		
3. Set Clock...		
1. Reading		
2. Temperature		
3. Quality		
4. Configuration		
5. Softkeys		

4. Help
1. Information
2. Contact
3. User Guide
4. PC software
5. Update Software

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