

Catalog #507

DuoSource™ 75 Power Supply



Specifications:

Max Voltage:	75 Volts
Max Current:	300 Milliamps
Type Output:	Constant Voltage
Lead Inputs:	2 Sets, Recessed, Color Coded
Fuse:	1.0 Amp, 250 V Glass fuse
Input Power:	60 Hz, 110 Volts
Connection:	3-Wire Grounded Cord
Made in the USA	

EDVOTEK's DuoSource™ 75 Power Supply is designed to power any combination of horizontal or vertical electrophoresis units. It generates constant voltage of 75 volts and has a maximum current of 300 milliamps. Safety features include a fused UL-approved, three-pronged power plug and a non-conductive casing. The unit is equipped with an easy-to-replace 1 Amp slow burn fuse. The DuoSource™ 75 is backed by a limited warranty and free technical support.


TO OPERATE

Plug the unit into a grounded outlet and flip the switch to turn on the power. The lead inputs are color-coded. The black input is negative and the red input is positive.

MAINTENANCE

This power supply includes a fuse tray that is located directly under the power cord socket at the rear of the unit (Fig 1). A flat head screw driver can be used to remove the tray (Fig 2) and replace the blown fuse when necessary (Fig 3).

The unit includes one spare fuse in the fuse tray. If this spare has already been used, the blown fuse can be replaced with a standard 5 x 20 mm fuse.

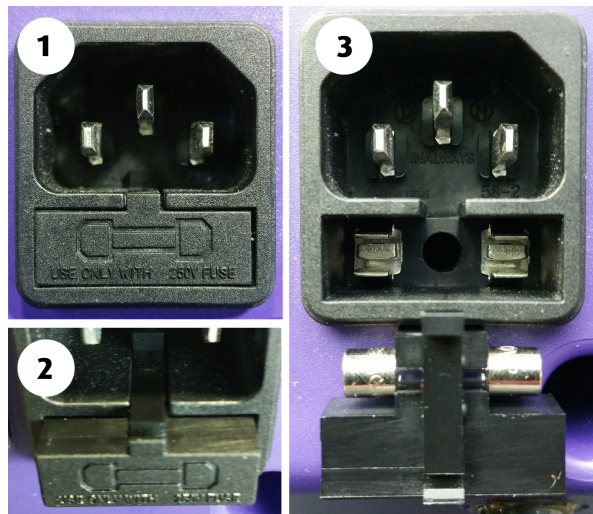

EDVOTEK®

1.800.EDVOTEK
www.edvotek.com