

Igientest Haccp

THE SIMPLEST QUICK TEST FOR HYGIENE CONTROL OF WORK SURFACE

Simple to use: it's enough to streak the wet swab on a surface and to insert the swab into the tube

Easy interpretation:
the comparison between the solution colour and the colorimetric scale confirms if the surface is clean or dirty

Minimal expiration date granted at the delivery time 36 months

Quick: results in only
10 minutes

Colorimetric scale printed on each single tube for easy interpretation

Application fields: food Industry, catering, dairy industry, health authority, HACCP consultants

Disposing procedure: at the end of the test the tube can be disposed as normal waste

Room temperature storage (from 10 °C to 30 °C)

HOW TO USE



take the swab out of the tube and streak on a surface measuring 100 cm² (10x10 cm)



insert the swab into the tube



break the glass vial contained in the tube by using the clamp cod. 91255



shake 5 times the tube grasping it by the tip



evaluate the result after 10 minutes evaluating the result by matching the colour with colorimetric scale



COLORE DOPO 10 MINUTI - COLOR AFTER 10 MINUTES

Luogo/Site



Igientest-HACCP

cod. 87883

THE
COLORIMETRIC SCALE
TO INTERPRET THE
RESULTS IS PRINTED
ON EACH TUBE

WHAT TO ORDER

CODE	DESCRIPTION	PACKAGING
87883	Igientest HACCP Surface hygiene monitoring test	50/box
Accessories		
91255	Clamp to break the glass vial of "Igientest HACCP"	1/box

Distributed By: