

Bone Clones®

800-914-0091



www.boneclones.com

OSTEOLOGICAL REPRODUCTIONS

Assembly Instructions:

SC-344-A

Articulated Large Dog Skeleton
(Bullmastiff)




Bone Clones, Inc. 9200 Eton Ave. Chatsworth, CA 91311

Phone: 818-709-7991 Fax: 818-709-7993

email: info@boneclones.com website: www.boneclones.com

Assembly Instructions
SC-344-A Articulated Large Dog Skeleton

All photos © Bone Clones, Inc.

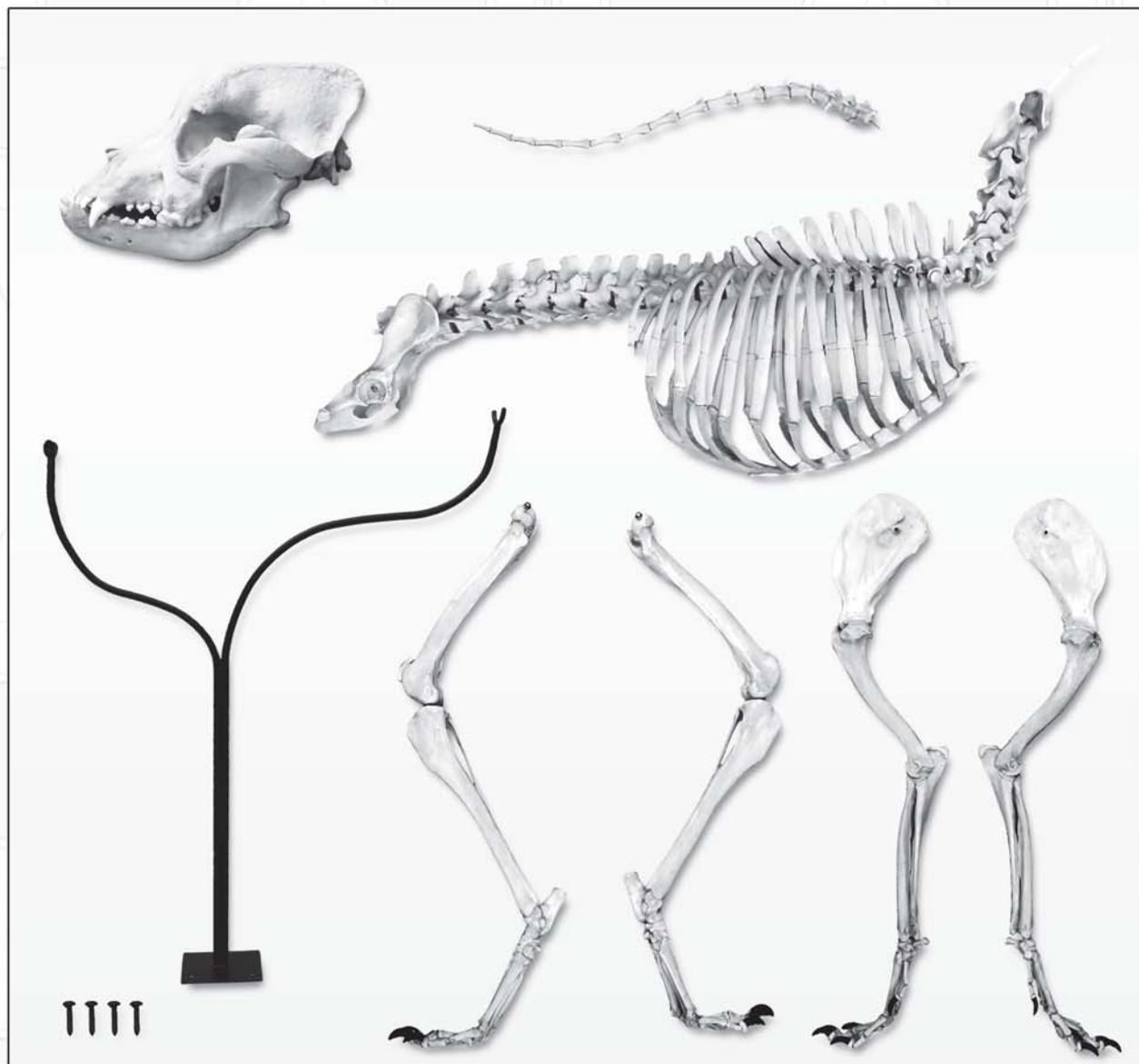

Parts Included:

- Skull
- Tail
- Torso
- Stand
- Screws (4)
- Front Legs (2)
- Back Legs (2)


***Note: Wooden display base S-SC-344 shown in these instructions is sold separately.**



Assembly Instructions

SC-344-A Articulated Large Dog Skeleton

All photos © Bone Clones, Inc.

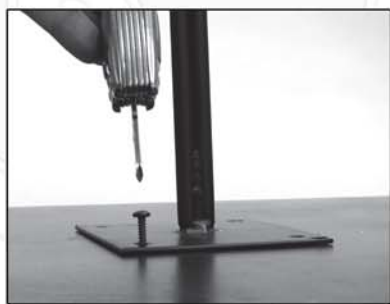

• Step 1


Attach the metal stand to the base of your choice.

Optional wooden base S-SC-344 is shown here.

Secure with the screws provided.

• Step 2

Place the torso on the stand as shown.

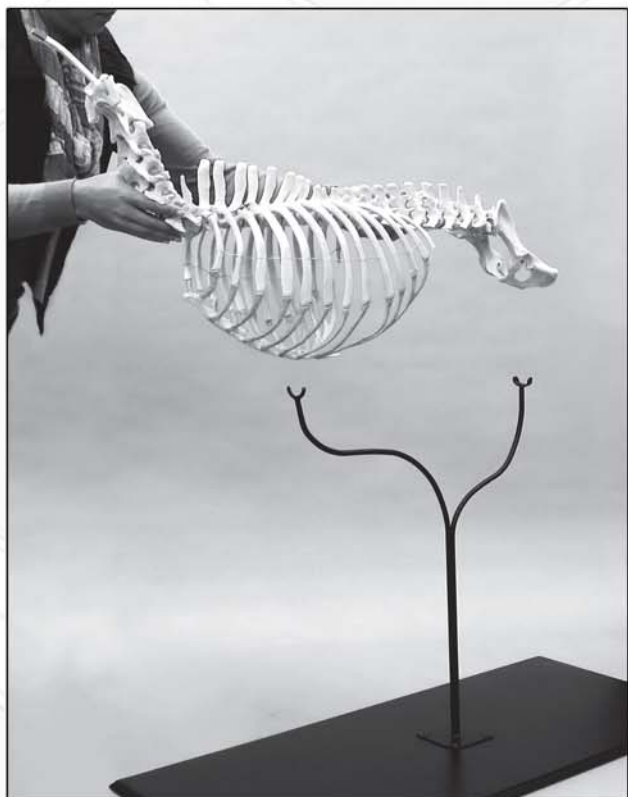

Assembly Instructions
SC-344-A Articulated Large Dog Skeleton

All photos © Bone Clones, Inc.



- Step 3

Attach one of the front legs to the threaded rod on the torso, and secure it with a cap nut.

- Step 4

Repeat for the other front leg.



Assembly Instructions
SC-344-A Articulated Large Dog Skeleton

All photos © Bone Clones, Inc.



- Step 5

Attach one of the back legs to the hip socket. Secure it with a cap nut.



- Step 6

Repeat for the other back leg.



Assembly Instructions
SC-344-A Articulated Large Dog Skeleton

All photos © Bone Clones, Inc.



- Step 7

Attach the tail
to the spine.

- Step 8

Attach the skull
to the skeleton.

