Instruction for use

VWR® FACE MASK PP 2-PLY EARLOOPS

EU cat. No 113-0713

EN

Intended use

The VWR face masks are designed to reduce process and products contamination in controlled environments.
For use in industrial environment only

Read and Understand before Use.

- 1 This product is not a medical device in context of Regulation (EU) 2017/745/Directive 93/42/EEC nor personal protective equipment in context of Regulation (EU) 2016/425!
- 2 This product does not prevent Coronavirus transmission and is not intended for general public!
- 3 Before use, the wearer should always check the mask is in good condition - not dirt, no damage to head strap attachment etc.
- 4 Facial hair and certain facial characteristics will prohibit effective use of this product.
- 5 Please make sure the mask fits properly before entering working environment.
- 6 Those masks are for single use. Do not reuse nor clean!
- 7 Dispose the products according to applicable local legislations

General Specifications

- Made of non-woven fabric 20g PP/25g PP
- White colour
- Ear loops
- Nature Rubber Latex Free

Packaging and Storage Specifications

Package/Quantity	50pcs/box, 20boxes/case
Best storage Conditions	0° - 50°C 45% - 85% humidity without sunshine exposure
Shelf Life	3 years
Country of Origin	China

Fitting instruction

- 1 Put loops around ears.
- 2 Press forefingers down on both sides of nose
- 3 Adjust mask over nose & under chin.

WARNUNG!

For industrial and clean room use only; do not use for protection against COVID-19!
This product is not a medical device for medical or clinical use within the meaning of Regulation (EU) 2017/745 / Directive 93/42/CEE!
This product is not a personal protective equipment (PPE) within the meaning of Regulation (EU) 2016/425!

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