Laboratory Equipment Manufacturer
www.mrclab.com

1 Litre Blender

Operating Manual

Blender 1 Litre - 230 volt

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Licuada de 1 litro - 230 voltios

Mode d'emploi

Mixer 1 Litre 230 volt

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Mixer 1 Liter 230 V

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Frullatore da 1 litro 230 volt

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1 Liter Blender 230 volt

3, Hagavish st. Israel 58817 Tel: 972 3 5595252, Fax: 972 3 5594529

mrc@mrclab.com MRC.10.17

IMPORTANT SAFEGUARDS

When using electrical appliances, basic safety precautions should always be followed including the following:

1. **READ ALL INSTRUCTIONS.**
2. To protect against electrical hazards, do not immerse the Blender Base in water or other liquid.
3. Close supervision is necessary when any appliance is used by or near children.
4. Unplug from outlet when not in use, before putting on or taking off parts, and before cleaning.
5. Avoid contacting moving parts.
6. Do not operate any appliance with a damaged cord or plug, or after the appliance malfunctions or is dropped or damaged in any manner. Return appliance to the nearest authorized service facility for examination, repair or electrical or mechanical adjustment.
7. The use of attachments, including canning jars, not recommended by the appliance manufacturer may cause hazards.
8. Do not use outdoors.
9. Do not let cord hang over edge of table or counter, or touch hot surfaces.
10. Keep hands and utensils out of container while blending to prevent the possibility of severe personal injury and/or damage to the blender. A rubber scraper may be used but must be used only when the blender is not running.
11. Blades are sharp. Handle carefully.
12. If your blender utilizes a removable Jar Base and Blade Assembly – to avoid injury, never place cutting assembly cutting blades on base without container properly attached.
13. Always operate blender with cover in place.
14. When blending hot liquids, remove center piece of two-piece cover on those units supplied with this construction. See Hot Blending Instructions.

SAVE THESE INSTRUCTIONS

SPECIFICATIONS

○ RATINGS	230 VAC – 50 Hz – SINGLE PHASE
○ MOTORS	PERMANENTLY LUBRICATED SERIES – WOUND UNIVERSAL TYPE
○ TIMER	SOME MODELS HAVE A TIMER, ADJUSTABLE FROM 1 TO 60 SECONDS FOR AUTOMATIC BLENDING, WITH A MANUAL SETTING FOR CONTINUOUS BLENDING.
○ BASES	ZINC DIE-CASTING WITH EPOXY ENAMEL FINISH OR POLISHED STAINLESS STEEL.
○ CONTAINERS	STAINLESS STEEL, BOROSILICATE GLASS, POLYCARBONATE OR ACRYLIC PLASTIC.
○ CORD	THREE CONDUCTOR GROUNDED CORD.
○ CE LISTED	
○ OPERATING CYCLE:	3 min ON / 1 min OFF

OPERATION

1. Before starting, make sure that the blender cord is unplugged and that the switch is set to the OFF (O) position.
2. Place container assembly on motor base so that it is firmly seated and not tilted to one side.
NOTE: If your blender has a container with a removable base and blade, the following points must be observed when assembling the container. Refer to Fig. 1:
 - Container gasket must be between the blade assembly and the container.
 - Container bottom must not be misthreaded onto the container.
 - Container bottom must be securely tightened to the container before placing food or liquid in the container.
 - Tighten the container bottom by holding the container upside down and rotating the container bottom **counterclockwise**, in the direction of the molded arrow on the underside of the container bottom.
 - Never place the container bottom or the blade assembly on the motor base unless they are completely attached to the container and securely tightened.
3. Be sure the switch is still in the OFF (O) position, and plug the power cord into a convenient outlet.
4. Put ingredients into container and cover it. Always add liquid ingredients first, unless otherwise specified in recipe. Do not fill above MAX line.
5. Set switch to the ON (I), LOW (1) or HIGH (2) POSITION. Always hold container with one hand when starting, and while blending. Never operate without jar lid in place.

When finished blending, set switch to OFF (O) position.

NOTE: For blenders with timers, turn the timer control knob clockwise to the desired time setting before starting the motor. If time setting is less than 10 seconds, turn knob past “10”, then back to desired setting.

When switch is set to the desired speed, motor will run. Motor will automatically stop at the end of the set time, and indicator lamp will light.

When switch is set to OFF (O) position, indicator lamp will go out.

Turning timer knob counterclockwise to the MANUAL position disables timer and indicator lamp, allowing blender to be turned on and off manually.

6. Allow motor to come to a complete stop before removing container. Never place the container on the motor base or remove it from the motor base while the motor is running.
7. Remove the container from the motor base and empty the contents.
8. Wash and dry container after each use. Containers with removable bottoms should be disassembled before washing.
 - The following parts are dishwasher safe: glass containers, stainless steel containers, removable container bottoms, blade assemblies, and gaskets.
 - **The following parts are not dishwasher safe and must be hand washed: plastic containers, container lids, and measuring cup lid inserts.**

After cleaning and drying, reassemble containers with removable bottoms.

Be sure gasket is between blade assembly and container as shown in Fig. 1.

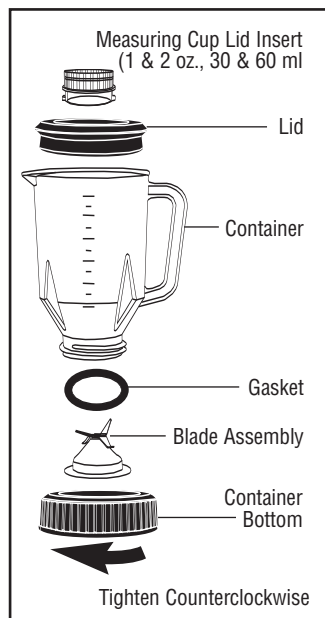
Fig. 1

Tighten container bottom securely before washing.

9. Not for waterjet cleaning.

HOT BLENDING INSTRUCTIONS: IMPORTANT!

To minimize the possibility of scalding when blending hot materials, the following precautions and procedures must be observed:

1. Limit the amount of material processed to 1/2 pint (235 ml).
2. Completely remove the measuring cup lid insert before turning on blender. This will allow venting of the rapidly heated and expanding air present inside container when blender is running.
3. Always start blending on lowest speed setting, then progress to desired speed.
4. Never stand directly over blender when it is being started or while it is running.

CIRCUIT BREAKER OPERATION

You will find a small, square, red plastic button located on the underside of the motor base of some models.

This is the reset button for a circuit breaker which prevents the motor from overheating in case of an overload.

If an overload occurs, the circuit breaker will shut off the motor and cause the button to protrude slightly from its normal position.

To reset the circuit breaker and resume operation, do the following:

- 1) SET THE POWER SWITCH TO ITS OFF (0) POSITION.
- 2) UNPLUG THE POWER CORD FROM ELECTRICAL OUTLET.
- 3) DETERMINE THE REASON FOR THE OVERLOAD, AND ELIMINATE THE CAUSE (FOR EXAMPLE: REDUCE THE AMOUNT OF MATERIAL BEING PROCESSED).
- 4) PUSH THE RED RESET BUTTON BACK INTO ITS NORMAL POSITION.
- 5) PLUG THE POWER CORD BACK INTO ELECTRICAL OUTLET.
- 6) RESUME NORMAL OPERATION.

MAINTENANCE

- **MOTOR**

Not for waterjet cleaning.

The motor requires no greasing or oiling. Sealed-in lubrication gives the unit lifetime smooth operation.

- **POWER SUPPLY CORD**

The power supply cord should not be replaced by the end user.

The power supply cord should only be replaced by the manufacturer's authorized importer (or, in the United States and Canada, by an authorized service center).

For repairs or further questions, please contact the manufacturer's authorized importer (or, in the United States and Canada, an authorized service center).

WARRANTY

This warranty is VOID if appliance is used on Direct Current (D.C.)

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3. DE REDEN VAN OVERBELASTING VASTSTELLEN EN DE OORZAAK ELIMINEREN (B.V.