Instruction manual

VWR® Digital High Speed Microplate Shaker

EU cat. no 444-7016

EN	- English		1
FR	- Français		11
ES	- Español		21
DE	- Deutsch		31
IT	- Italiano		41
PT	- Português		51
BG	- Български		61
CZ	- Čeština		64
DK	- Dansk		67
EE	- Eesti		70
FI	- Suomi		73
GR	- Ελληνικά		76
HR	- Hrvatska		79
HU	- Magyar		82
IS	- Íslenska		85
LT	- Lietuvių		88
LV	- Latviešu		91
NL	- Nederlands		94
NO	- Norsk		97
PL	- Polski		100
RO	- Română		103
SI	- Slovenščina		106
SK	- Slovenčina		109
SV	- Svenska		112

Legal address of Manufacturer

Europe

VWR International bv,
Researchpark Haasrode 2020
Geldenaaksebaan 464 B-3001
Leuven
+3216385011
be.vwr.com

UK Importer:

VWR International Ltd
Hunter Boulevard, Magna Park
Lutterworth, Leicestershire, LE17 4XN
uk.vwr.com

Country of Origin

United States

Table of Contents

Safety Information	4
Package Contents	4
Installation	4
Intended Use	5
Maintenance & Servicing	5
Environmental Conditions	5
Product Specifications	6
Control Panel	6
Operating Instructions	7
Troubleshooting	8
Technical Service	8
Warranty	8
Disposal	9
Distribution Contact Information	10

Safety Instructions

Please read the entire instruction manual before operating the unit.

WARNING! DO NOT use the **UNIT** in a hazardous atmosphere or with hazardous materials for which the unit was not designed. Also, the user should be aware that the protection provided by the equipment may be impaired if used with accessories not provided or recommended by the manufacturer, or used in a manner not specified by the manufacturer.

Always operate unit on a level surface for best performance and maximum safety.

DO NOT lift the High Speed Microplate Shaker by the tray.

CAUTION! To avoid electrical shock, completely cut off power to the unit by disconnecting the power cord from the *unit or unplug from the wall outlet*. Disconnect unit from the power supply prior to maintenance and servicing.

Spills should be removed promptly.

DO NOT immerse the unit for cleaning.

DO NOT operate the unit if it shows signs of electrical or mechanical damage.

Earth Ground - Protective Conductor Terminal

Alternating Current

Package Contents

Digital High Speed Microplate Shaker

Description	Quantity
234cm detachable power cord (EU, UK, & CH)	3

Installation

1. **Important:** Choose the space on your bench where the VWR High Speed Microplate Shaker will stay. Because of the amount of energy dissipated, the unit is held onto your lab bench with 12 rubber feet (suction cups). The unit is VERY difficult to remove from work surface and should be done with caution. Use paper under the feet to keep them from attaching to the bench while determining the final location. When final location is determined, remove the paper.
2. Be sure the power switch is in the off position.
3. The VWR High Speed Micro Plate Shaker comes with a connector plug that is inserted into the IEC connector on the back of the unit first, then it can be plugged into a properly grounded outlet. The 230v unit plugs into a 230 volt, 50/60 Hz source.
4. Your VWR High Speed Micro Plate Shaker is ready for use.

Intended use

This shaker is intended for general laboratory use.

Maintenance & Servicing

The VWR High Speed Microplate Shaker is built for long, trouble-free, dependable service. No lubrication or other technical user maintenance is required. It needs no user maintenance beyond keeping the surfaces clean. The unit should be given the care normally required for any electrical appliance. Avoid wetting or unnecessary exposure to fumes. Spills should be removed promptly. Do not use a cleaning agent or solvent on the front panel which is abrasive or harmful to plastics, nor one which is flammable. Always ensure the power is disconnected from the unit prior to any cleaning. If the unit ever requires service, contact your VWR representative.

Environmental Conditions

Operating Conditions: Indoor use only.

Temperature: 4 to 40°C

Humidity: 20% to 80% relative humidity, non-condensing

Altitude: 0 to 2000 M above sea level

Non-Operating Storage:

Temperature: -20 to 65°C

Humidity: 20% to 80% relative humidity, non-condensing

Installation Category II and Pollution Degree 2 in accordance with IEC 664.

Product Specifications

Overall dimensions (L x W x H): 39,4 x 30,5 x 32,5cm

Tray dimensions (L x W): 31,1 x 27,9cm

Electrical (50/60 Hz): 230 V - 1 amp, 120 watts

Fuses: 5mm x 20mm, 5 amp quick acting (2 required)

Speed range: 600 to 2500rpm

Timer: 1 second to 160 hrs

Orbit: 3,6mm


Controls: See page 6

Ship weight: 22,7kg

Capacity: 3,2kg

48 microplates in stacks up to 12,7cm high

Control Panel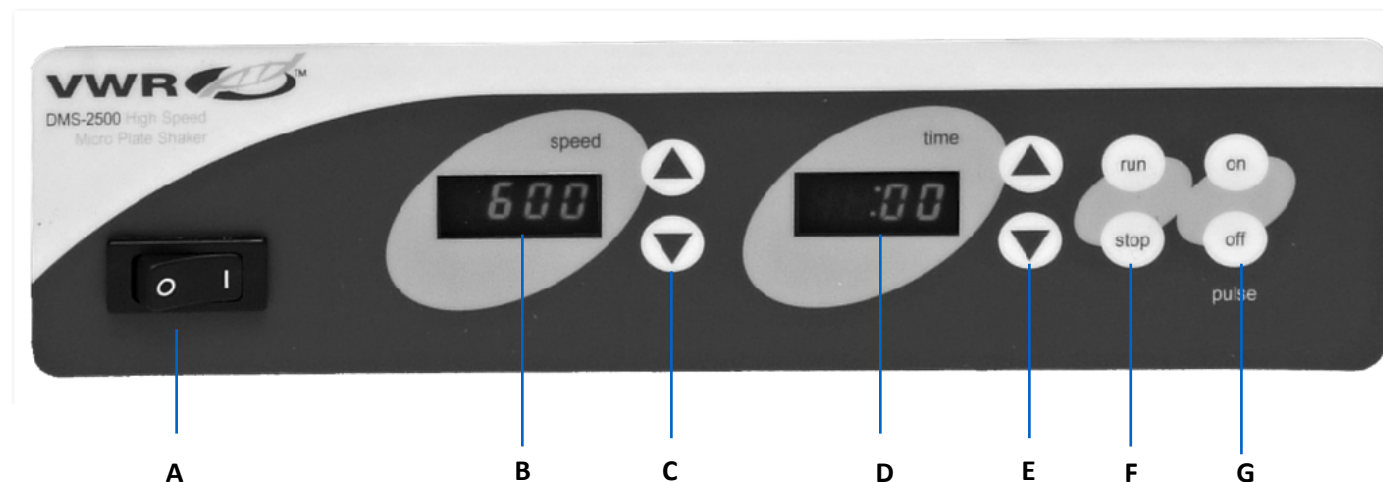

The front panel of the High Speed Microplate Shaker contains all the switches, controls and displays needed to operate the unit.

A. On/off rocker switch: Turns main power on/off.

B. Speed display: Displays the speed of the shaker.

C. Up/down arrows for set-point control.

D. Time display: Displays accumulated time (continuous mode) or how much time is remaining (timed mode). The display range is from 0 to 9999 minutes in one (1) second increments. The display will indicate minutes and seconds until the timer reaches 99 minutes and 59 seconds (99:59), then the display will automatically display minutes up to 9999.

E. Up/down arrows for set-point control.

F. Run/stop buttons: Activates shaking.

G. Pulse on/off buttons: Activates pulse mode

Operating Instructions

1. Flip up the four quick release latches which will release top tray from the 4 posts. DO NOT USE without both top and bottom tray foam in place.
2. Your Digital High Speed Micro Plate Shaker is designed to hold up to 3,2kg, which will accommodate deepwell blocks or stacked microplates up to 12,7cm high.
To avoid excessive noise and potential block damage, well blocks should always have space between them and should never touch each other.
3. Sample containers should be placed on bottom tray foam insert and centered on the tray to achieve optimal performance.
4. To place top tray on unit, place the short side to the front engaging the 4 posts. When top tray is lined up, press down on the center of the tray and engage by pushing the quick release latches down into the locked position. Repeat for all corners. When top plate is locked in place, tighten the knobs on the four adjustment blocks. The High Speed Micro Plate Shaker is ready to operate. When changing samples, disengage the four quick release latches to remove the top tray. Do not loosen the knobs on the adjustment blocks. Loosen and move the four adjustment blocks only when the sample heights have changed.
5. Turn the Power on by depressing the left side of the rocker switch on the bottom left side of the control panel. LED displays for Speed and Timer will light.
6. Set speed by pressing the up/down arrows to the right of the speed display until you reach the desired speed. Adjust Time using the up/down arrows to the right of the timer display. Press the run button. The unit will run for the programmed time. Timer LED will display time remaining.
7. To run in the un-timed mode, set the time to zero (0) using the up/down arrows to the right of the timer display. Press the run button and the unit will run until you press the stop button. Timer LED will display accumulated time.
8. To run in pulse mode unit must not be running. Press the "ON" button above the word "PULSE". To stop the pulse feature press the "OFF" button above the word "PULSE". Pulse mode is factory programmed for 2 seconds on 1 second off. The pulse "ON" and pulse "OFF" times can be adjusted between 1 and 59 seconds in 1 second intervals.

The Pulse time can be reprogrammed by following the directions below.

Pulse "ON" time

- a. Turn power off
- b. Turn power on while holding pulse "ON" button. The LED display for "SPEED" shows current pulse on time.
- c. Change the pulse on time using the up/ down arrows to the right of the "SPEED" display.
- d. When the power is turned off, the value on the display is stored as the new pulse on time.

Pulse "OFF" time

- a. Turn power off
- b. Turn power on while holding pulse "OFF" button. The LED display for "TIMER" shows current pulse off time.
- c. Change the pulse off time using the up/down arrows to the right of the "TIMER" display.
- d. When the power is turned off, the value on the display is stored as the new pulse off time.

Troubleshooting

Review the information in the table below to troubleshoot operating problems.

Problem	Cause	Solution
Unit fails to power on	Missing or blown fuse	Add or replace fuse as necessary. If problem persists, please contact your VWR representative for repair.
Unit is excessively noisy	Rattling or ticking sounds may indicate a loose screw in the tray	Ensure that tray is secured tightly. If problem persists, please contact your VWR representative for repair.
Unit vibrates abnormally	Suction cup feet	Firmly press down on the four (4) corners of the unit creating a strong suction to the work surface (DO NOT place on bench mat). If problem persists, please contact your VWR representative for repair.

Technical service

Web Resources

Visit the VWR website at www.vwr.com for:

- Complete technical service contact information
- Access to the VWR Online Catalogue, and information about accessories and related products
- Additional product information and special offers

Contact us For information or technical assistance contact your local VWR representative or visit. www.vwr.com.

Warranty


VWR warrants that this product will be free from defects in material and workmanship for a period of five (5) years from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non-performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

Compliance with local laws and regulations

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorisations necessary to run or use the Product in its local environment. VWR will not be held liable for any related omission or for not obtaining the required approval or authorisation, unless any refusal is due to a defect of the product.

Equipment disposal

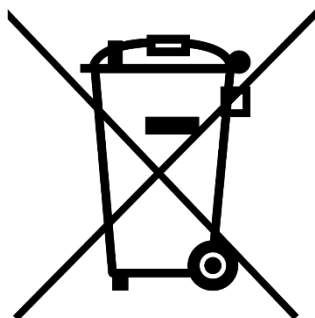This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste.

Instead it's your responsibility to correctly dispose of your equipment at lifecycle -end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

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India

VWR Lab Products Private Limited
No.139. BDA Industrial Suburb,
6th Main, Tumkur Road, Peenya Post,
Bangalore, India – 560058
Tel.: +91 80 28078400
Email: vwr_india@vwr.com

Ireland

VWR International Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: +353 01 88 22 222
Email: sales.ie@vwr.com

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