Instrucciones

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Instructions

Operation

Connect the sensor probe to the unit.
Open the battery cover; remove the battery strip to turn on the unit. The display will show full segment display for about 3 seconds and then display current temperature and MIN/MAX value.
Select the temperature unit by the [°C-°F] switch on the back of the unit.

To Measure Temperature

Insert sensor probe into test substance and take the reading when it becomes stable.

To Record an Event

The unit will automatically record the date and time of an out of alarm limit when the Low/High Alarm is turned on. It will record 10 events.

Change Display Mode

Press the [MODE] button once to change to display to the LO/HI alarm mode
Press the [MODE] button again to change to display to the DATE TIME mode
Press the [MODE] button again to change to display to the MIN/MAX mode

Note: When the unit is powered on, the default display is the MIN/MAX mode

Setting Date and Time

To set the date and time, press the [MODE] button to display the DATE TIME mode.
Press the [SET] button on the back of the unit to enter into date and time setting.
Press either the [▲] or [▼] button on the back of the unit to set the correct values.
The value to be set will flash.
Press the [EVENT] button to select either DD/MM or MM/DD display format.
Note, if the date and time change, memory will be reset.

Setting Low/High Alarm

To set the Low/High Alarm, press the [MODE] button to display the LO/HI Alarm mode.
Press the [SET] button on the back of the unit to enter into the low alarm setting, press either the [▲] or [▼] button on the back of the unit to set low alarm limited.
Press the [SET] button again to enter into the high alarm setting, press either the [▲] or [▼] button on the back of the unit to set high alarm limited.
Press the [SET] button three times to exit the alarm setting.

When the current temperature reading is either lower than low alarm value or higher than high alarm value, the unit will sound, the LO ALM or HI ALM icon will be flash and the LED in the upper right hand corner of the unit will flash.

The alarm will sound for 1 minute, then the alarm sound will be 2 "di" every 1 minute to save battery life. The sound will stop when the unit is back within the alarm limits but the LED will continue to flash until the memory is cleared.

Turn Low/High Alarm On or Off

In the LO/HI alarm display mode, press the [☒] button to turn on or off the alarm function. When the alarm is on, the [☒] icon will be display. The icon will be displayed in the alarm and Max/Min display modes.

Low/high Alarm Date Time Stamp

In the LO/HI alarm display mode, press the [EVENT] button to read the LO/HI alarm generated date time stamp.

Press either the [▲] or [▼] button on the back of the unit to scroll through the date time stamped events.
Press the [EVENT] button again to exit the reading mode.
Press the [CLEAR] button on the back of the unit to clear the memory.

MIN/MAX temperature

In the MIN/MAX display mode, the display can be read current temperature and MIN/MAX temperature memory.
Press the [EVENT] button to read the MAX temperature and the date and time of the reading.
Press the [EVENT] button to read the MIN temperature and the date and time of the reading.
Press the [CLEAR] button on the back of the unit to clear the memory.

NOTE:

Clean the probe immediately with cleaning wipes after each measurement to avoid cross contamination.
Do not expose the probe handle to temperatures that may exceed over 50°C/ 122 °F.

ERROR SYMBOLS:

Symbol	Description	Action required
	Low battery voltage	Replace the battery with the same type
	1) Sensor open circuit	1) Return the thermometer for repair
LLLL	2) The reading is out of low range (-50°C)	2) Keep the measurement above low range
	1) Sensor short circuit	1) Return the thermometer for repair
HHH.H	2) The reading is out of high range (200°C)	2) Keep the measurement below high range

Battery Handling/Packaging

- Remove batteries from equipment immediately after they fail to operate the equipment.
- If equipment uses two or more batteries, always replace batteries in complete sets.
- Do not attempt to recharge alkaline batteries, they are not rechargeable.
- Do not handle hot or warm batteries.
- Do not heat, incinerate, crush, puncture or mutilate batteries.
- Wear personal protective equipment if batteries show signs of leakage, bulging, swelling or deformity.

Warning: If potassium hydroxide electrolyte from an alkaline battery comes in contact with the skin, do not try to neutralize the electrolyte with vinegar or any other acidic solutions. Neutralization may trap electrolyte on the skin. Flush the affected skin area with copious amounts of water. If the battery electrolyte gets into your eyes, it can cause severe damage and or blindness.

Equipment disposal

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste.

Instead it's your responsibility to correctly dispose of your equipment at lifecycle -end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

Thank you

Technical service

Web Resources

- Visit the VWR website at **www.vwr.com** for:
- Complete technical service contact information
 - Access to the VWR Online Catalogue, and information about accessories and related products
 - Additional product information and special offers
- Contact us** For information or technical assistance contact your local VWR representative or visit **www.vwr.com**.

Warranty

VWR warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

Compliance with local laws and regulations

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorizations necessary to run or use the product in its local environment. VWR will not be held liable for any related omission or for not obtaining the required approval or authorization, unless any refusal is due to a defect of the product.

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