Instruction manual

- 471-1307 VWR Refrigerator +5C 118L sparkfree
- 471-1308 VWR Refrigerator +5C 118L
- 471-1309 VWR Refrigerator +5C 310L sparkfree
- 471-1310 VWR Refrigerator +5C 310L

Legal address of Manufacturer

Europe

VWR International bv,
Researchpark Haasrode 2020
Geldenaaksebaan 464 B-3001 Leuven
+3216385011
be.vwr.com

UK Importer:
VWR International Ltd
Hunter Boulevard, Magna Park
Lutterworth, Leicestershire, LE17 4XN
uk.vwr.com

Country of Origin

CHINA

Content

- Safety Precautions 1
 - Precautions for use.....4
- Product Installation5
- Component Names&Control Panel9
- Application method 12
- Alarm 13
- Cleaning and Maintenance.....14
- FAQ 15
- Refrigeration Schematic&Circuit Diagram 16
- Specification 17
- Technical service 19
- Warranty 19
- Compliance with local laws and regulations 19
- Equipment disposal20
- Local VWR offices in Europe and Asia Pacific21

Safety Precautions

Dear users:

Please make sure you have carefully read and observed the contents with following signs in the manual, for better understanding of this manual and better use of this product, so as to prevent personal injuries and refrigerator damage.

Actions or operations
which are prohibited

Actions or operations
which must be followed

Caution

-  Close down the gas leaking valve and open the door and windows for ventilation in case of leakage of inflammable gas like coal gas etc., do not insert or pull out the power plug of the refrigerator, otherwise it may trigger an explosion and fire.
-  Only the professional technicians or after-sale service men can install this refrigerator. Otherwise it may result in an electric shock or fire.
-  Make sure to place the refrigerator firmly on a solid and flat ground. If the ground is not stable or it is not properly placed, it will cause overturn of the refrigerator or personnel injury.
-  Please use the dedicated power supply indicated on the nameplate of the refrigerator, or otherwise, it may cause a fire or electric shock.
-  If it is required to lengthen the power cord, the section area of the lengthen cord must not be less than 2mm² and the length shall not be longer than 3m. Otherwise, it may cause a fire or electric shock.
-  The power cord of this refrigerator is provided with a three-wire (earthing) plug, complying with the standard 10A three-wire (earthing) socket. Never cut or remove the earthing pin of the power cord in any case. Make sure that the power plug and socket is inserted tightly and reliably, otherwise it may result in a fire.
-  Please use an earthed power supply socket for fear of electric shock.
-  If the power socket is not earthed, be sure to have the earth wire installed by a professional technician.
-  Do not use the refrigerator in the open air. When the refrigerator gets wet in the rainwater, it may cause leakage of electricity or electric shock.
-  Do not place the refrigerator in a damp site or a place where it may easily suffer from spraying water. Otherwise, it may cause leakage of electricity or electric shock due to degeneration of its insulation.
-  Do not pour water onto the refrigerator in a direct inclined way, otherwise it may cause an electric shock or a short circuit.
-  Do not place any vessel containing water or any heavy object on the refrigerator. If the object drops off, it may cause personal injury and the flowing water will degenerate the insulation and further cause leakage of electricity or electric shock.

-  Do not make the refrigerator earthed via a gas-pipe, power supply conduit, telephone line or lightning rod. The above earthing will cause an electric shock or other dangers.
-  Do not contact any electric parts or any switches such as power plug and so forth with a wet hand, or otherwise it may cause an electric shock.
-  When pulling out a plug from the power socket, hold the power plug firmly but do not pull the wire of the power plug. Pulling the wire with hand may cause an electric shock or lead to a fire due to short circuit.
-  If the refrigerator runs out of order, please pull out the power plug. Continuous running in an abnormal condition may cause an electric shock or fire.
-  User must not dismantle, repair or refit the refrigerator by itself. Otherwise, it may cause a fire or personal injury due to improper operation.
-  Before conducting any repair or maintenance of the refrigerator, be sure to disconnect the power of the refrigerator for fear of electric shock or personal injury.
-  Make sure not to suck in medicine or suspended particles inside and surrounding the refrigerator in time of repair and maintenance, otherwise, it may bring hurt to your health.
-  When storing toxic, hazardous or radioactive articles, please use the refrigerator in a safety zone.
-  Improper use may bring hurt to human health or environment.
-  If the refrigerator is left unused for a long time in an unattended area, make sure that children will not approach the refrigerator and the refrigerator door will not be closed completely.
-  Disposal of discarded refrigerator shall be conducted by the relevant personnel. The appliance contains reusable materials and should be disposed of properly - not simply with unsorted household refuse. Appliances which are no longer needed must be disposed of in a professional and appropriate way, in accordance with the current local regulations and laws.
-  When disposing of the appliance, ensure that the refrigerant circuit is not damaged to prevent uncontrolled escape of the refrigerant it contains (data on type plate) and oil.
-  Do not store corrosive articles such as acid, alkali etc. in the refrigerator. Otherwise, it may spoil the internal units or electric parts of the refrigerator.
-  Do not place the packing plastic bags at the locations where children can easily take as plastic bags may cause occurrence of suffocation accident.
-  Do not climb up to the refrigerator or place articles on the refrigerator, otherwise it may cause personal injury or damage of refrigerator due to overturn of the refrigerator.
-  Do not insert any metal objects like iron nails, iron-wire into any orifice and gap or any air vent used for internal air circulation of the refrigerator, otherwise, it may cause an electric shock or injury due to contact of the above objects with the moving parts.
-  If the supply cord is damaged, it must be replaced by the manufacturer, its service agent or similarly qualified persons in order to avoid a hazard.
-  Do not store explosive substances such as aerosol cans with a flammable propellant in this appliance.

-  Keep clear of obstruction all ventilation openings in the appliance enclosure or in the structure for building-in.
-  Do not use mechanical devices or other means to accelerate the defrosting process, other than those recommended by the manufacturer.
-  Do not damage the refrigerant circuit.
-  Do not use electrical appliances inside the food storage compartments of the appliance, unless they are of the type recommended by the manufacturer.
-  In order to reduce flammability hazards the installation of this appliance must only be carried out by a suitably qualified person.
-  For the product with non-conductive inner chamber(white), see the warning label of product.
-  Put explosive/flammable material in sealed container before storing into laboratory refrigerator.

 Caution

-  When restarting the refrigerator after a power failure or power-off, it is required to check the setting of the refrigerator. Change of setting may spoil the articles stored.
-  Once the power of the refrigerator is cut off, wait for over 5 minutes before it is reconnected so as not to damage the compressor or the system.
-  Put on your gloves at time of maintenance so as not to bump the sharp edges or corners, resulting in personal injury.
-  Close the door of the refrigerator with your hand holding the handle so that your fingers will not be pinched by the door.
-  The inclination angle shall be no greater than 45 ° when handling the refrigerator.
-  When handling the refrigerator, be careful not to trip over it, preventing damage to the refrigerator or personal injury.
-  Do not pull or carry the equipment by the door handle, preventing damage to the refrigerator or personal injury.
-  There shall be no obstacles around the refrigerator, and keep it well-ventilated and clear.
-  Do not spoil the refrigerating circuit.
-  Do not use any electric appliance in the store-room of the refrigerator except for the types that the manufacturer recommended for use.

Precautions for use

- Product structure and composition: The product adopts a refrigeration system attached to the thermal insulation box which is microcomputer controlled with air cooling design, and a rack type design that can be adjusted at discretion.
- Before putting the articles into the refrigerator, make sure firstly that the design temperature is reached inside the refrigerator, then put in the articles in batches. The articles put in every time shall not exceed 1/3 of the internal capacity so as to prevent temperature increase.
- The indicated temperature value of the refrigerator is the value indicated inside the refrigerator at the temperature sensing probe. Occasionally the indicated temperature has a certain difference with the actual temperature in the central part of the refrigerator, but it will gradually approach to the true temperature.
- Use the diluted mild detergent to clean the refrigerator. Do not use brushes, acid, gasoline, soap powder, polishing powder or hot water to clean the refrigerator as the above materials may damage the painting surface as well as the plastic and rubber parts and components. Take particular care not to use volatile solvent such as gasoline to wipe plastic and rubber parts and components.
- Cut off the power of the refrigerator when it is not in service for a long time.
- Every time when putting in and taking out the articles, please try to reduce the door opening time so as not to cause larger fluctuations of the temperature and humidity inside the refrigerator.
- After opening the door, the temperature inside the refrigerator will rise steeply in a short time, this is normal and the temperature will restore within one hour after the door is closed.
- The refrigerator is equipped with security door lock. Please keep the key properly. Once the key is lost, the door body of the refrigerator will not be opened.

Product Installation

■ Installation environment

- Environmental temperature: 16°C to 32°C , the optimal temperature is 18°C to 25°C , the air conditioning system shall be used when necessary.
- Ambient humidity: below 85% Rh.
- Avoid large quantities of dust.
- Avoid mechanical sway or vibration.
- The altitude of refrigerator's working position: lower than 2000m.
- Input voltage: within the range of 220V±10%.

Caution

- As the refrigerator is sensitive to the environmental temperature, it can not operate normally if it is installed in an environment other than the above. Please use it after improving the environment.
- It is prohibited to install the refrigerator in the open air environment. When the refrigerator gets wet in the rainwater, it may cause leakage of electricity or electric shock accident.

■ Installation site

For normal operation and optimum performance level of the refrigerator, its installation place shall conform to the following conditions:

- It can not be installed in a narrow, small and airtight space, the door frame of the room for placing it shall not be less than and lower than this product, and shall at least guarantee its normal passing in and out so as to avoid any maintenance difficulty and failure of in-time repair in occurrence of machine fault, causing spoilage to the stored articles.
- The ground at the installation place shall be solid and flat.
- It shall be well-ventilated without direct sunshine.
- Do not share a socket with other equipment, the plug and socket shall be well connected.
- Do not distort nor stress the power cord.
- If it is required to lengthen the power cord, the section area of the lengthened wire shall not be less than 2mm² and the length shall not be longer than 3m.
- Check the operating voltage before use, at the area where the voltage is not stable, use of a voltage regulator suitable for the motor load shall be considered for voltage stabilization. So that the normal input voltage is stabilized at 220V-240V~, the power of the voltage regulator is greater than 4000W.
- The refrigerator shall be guaranteed with reliable earthing.
- If the power socket has earth wire, check if the earthing is good before use.
- The interior is used for storing flammable and explosive reagents. Only applies Sparkfree/ATEX units;
- The appliance is not suitable for use in explosion-hazard areas.

Warning

- Do not make refrigerator earthing via gas piping, water supply pipe, telephone line or lightning rod that may easily cause an electric shock.
- After installation, the power plug must be accessible to facilitate unplugging of the power cord in emergency. No articles shall block out the air vent of the refrigerator.

Preparation before use

1. Take off the packing material and the packaging tape

Take off all packing materials and packaging tapes for transportation.

keep the plastic bag away from the children.

2. Check of accompanying accessories

Please check the box contents as per the packing list. If there is any mistake, please contact the aftersales department timely.

3. Storage condition

At least a 10cm gap shall be left around the storage box, to ensure the well air ventilation and to facilitate the heat radiation.


Sufficient space shall be kept at the engine-room grilling, to facilitate maintenance and ventilation.

4. Adjustment of leveling leg

Please use a wrench to turn the horizontal leveling leg clockwise so that the horizontal leveling leg are extended outwards and supported on the ground. Make sure that the refrigerator is not movable in time of use.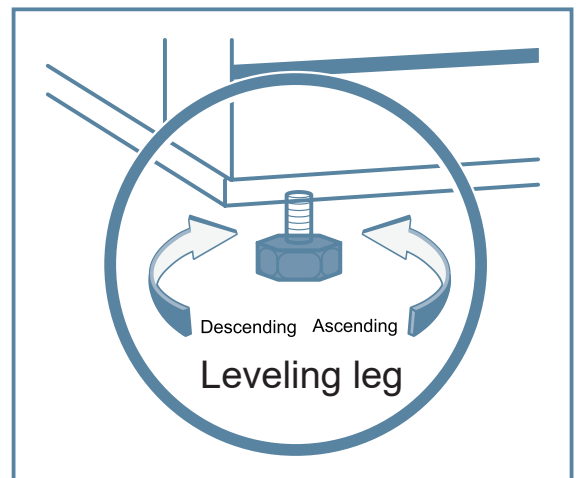

5. Standing

Do not power on the refrigerator firstly after leveling and cleaning. Let it stand still for over 24 hours before power it on so as to ensure normal running of the refrigerator.

6. Water collection tray

There is a water tray to collect defrosting water of evaporator.

For model 471-1307 and 471-1309: the water tray is placed below the evaporator, it need to be dumped manually to prevent the water from spilling out.


For model 471-1308 and 471-1310: the water tray is placed on the bottom of the unit, and the water will be evaporated automatically.

Changing over door hinge

Change over the door hinge (e.g from right hinged to left hinged)

Caution: at least 2 trained staffs are needed to proceed this operation

• For Model 471-1307 & 471-1308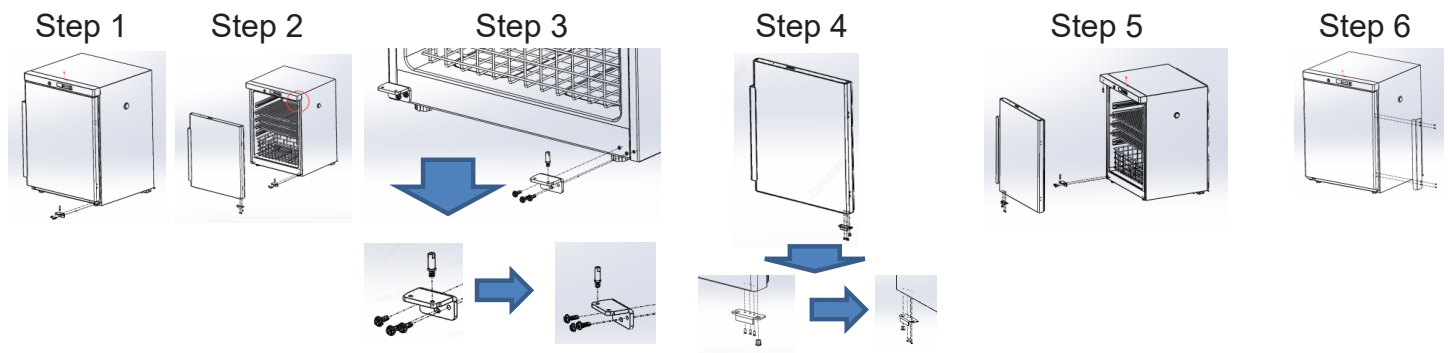

Step 1. Unscrew the right lower hinge (Caution: the door must be hold during this process) and then remove the door

Step 2. Move the pin from the right upper hinge to the left one


Step 3. Install the pin onto the left hole of left lower hinge

Step 4. Move the door stopper at the right bottom of the door to the left; meanwhile move the sleeve onto the left hole

Step 5. Hold the door and insert the left upper hinge into the shaft hole at the upper left side, then insert the left lower hinge into the shaft hole at the lower left side, and then screw the left lower hinge

Step 6. Move the handle from the left side to the right side

• For Model 471-1309 & 471-1310

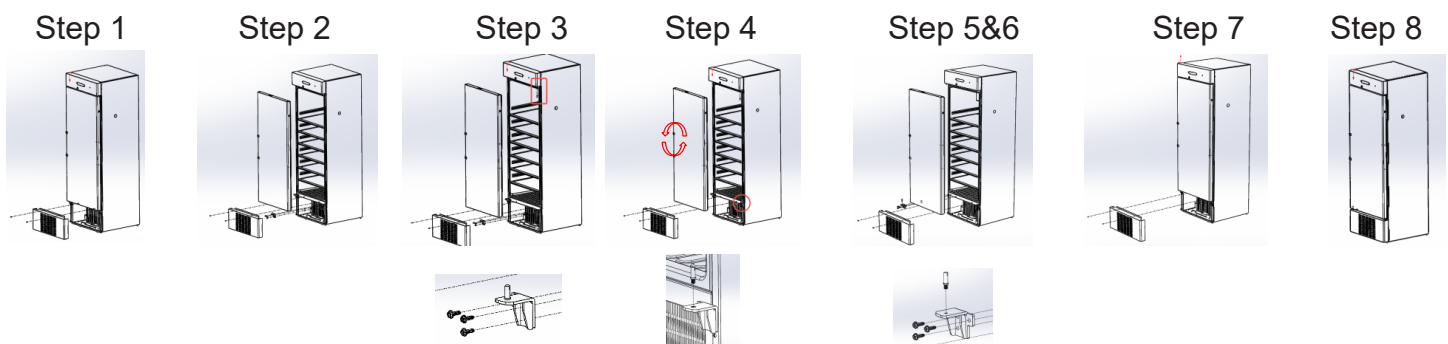Step 1. Unscrew the front grid at the bottom

Step 2. Unscrew the right lower hinge and then remove the door (Caution: the door must be hold during this process)

Step 3. Move the pin from the right upper hinge to the left one

Step 4. Install the right lower hinge back on its original place, then remove the pin on it then rotate the door up side down.

Step 5. Unscrew the left lower hinge, and install the pin on it

Step 6. Hold the door and insert the left upper hinge into the shaft hole at the upper left side, then insert the left lower hinge into the shaft hole at the lower left side, and then screw the left lower hinge

Step 7. Install the front grid at the bottom

Step 8. Move the handle from the left side to the right side

■ Initial Power-on

Please observe the following rules when conducting initial start-up and continuous running:

Connect the power cord to the dedicated socket of proper gauge when the refrigerator is empty, make sure the power supply is 220V-240V~/50Hz.

Make sure you switch on the battery switch.

This refrigerator is preset at below 5°C before factory delivery.

The sound alarm will ring after power-on and this is normal. At this moment, press the "Alarm " key on the control panel to cancel the alarm beep. The alarm indicator keeps flickering till the temperature sensor reaches the temperature range required.

The temperature can be stabilized at the set temperature after the refrigerator has been running for a period of time.

Start putting articles gradually into the refrigerator after completing a thorough inspection of the refrigerator.

- ! A person shall be specially assigned for daily check and record of the product operating conditions, in case the temperature inside the refrigerator is on the high side or on the low side, just transfer the articles to the other storage boxes till the fault is removed and then replace back the articles.
- ! The indicated temperature of this product may have a certain difference with the actual temperature inside the boxes due to refrigeration inertia, this is normal.
- ! Do not put any articles in the space between the bottom of the refrigerator's internal bladder and the bottom rack for fear of air duct clogging and affecting the temperature uniformity inside the boxes.
- ⊘ All the refrigerators are of storage equipment, and it is strictly forbidden to put excessive articles that are relatively too hot in one go, this will cause long time of running of the compressor without stop, slow drop of temperature, thus affecting service life of the compressor, the articles must be put in batches so that the temperature drops in steps till the required temperature is reached.
- ⊘ Do not spoil the refrigerating circuit.
- ⊘ Inside the refrigerator, do not use the electric appliance without production permission.

■ Operation of over-current protector

1. Mounting position: This device is mounted on the backside of the product, and there is a special warning sticker to mark it.

2. Working principle: In occurrence of any abnormalities found with the interior circuit of the product that exceed the protection limit of the over-current protector, this device will automatically spring up, cut off the connection with the external

circuit and play the role of protecting the entire equipment circuit; at time of reuse, it is required to manually push down the device and the product is powered on for running.

■ Operation after power failure

The refrigerator has memory function to the setting values, after re-electrified following a power failure, the refrigerator will continue running as per the parameters set up before the last power failure.

Caution

- Once the power of the refrigerator is cut off, wait for over 5 minutes before it is reconnected so as not to damage the compressor or the system.
- Pull out the power plug when the refrigerator is not in service for a long time so as to prevent electric shock, leakage of electricity or fire due to aging of the power cord.
- If the refrigerator is left unused for a long time in an unattended area, make sure that children will not approach the refrigerator and the refrigerator door can not be closed completely.

Component Names&Control Panel

■ Component Names

• 471-1307

• 471-1308

• 471-1309


• 471-1310


■ Control panel

•471-1307 / 471-1308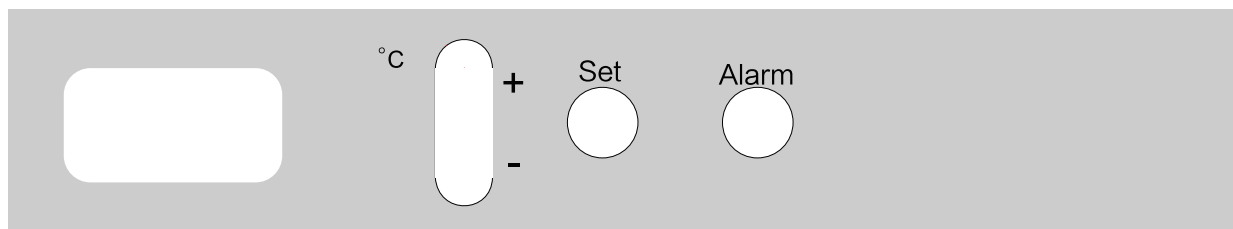

•471-1309 / 471-1310

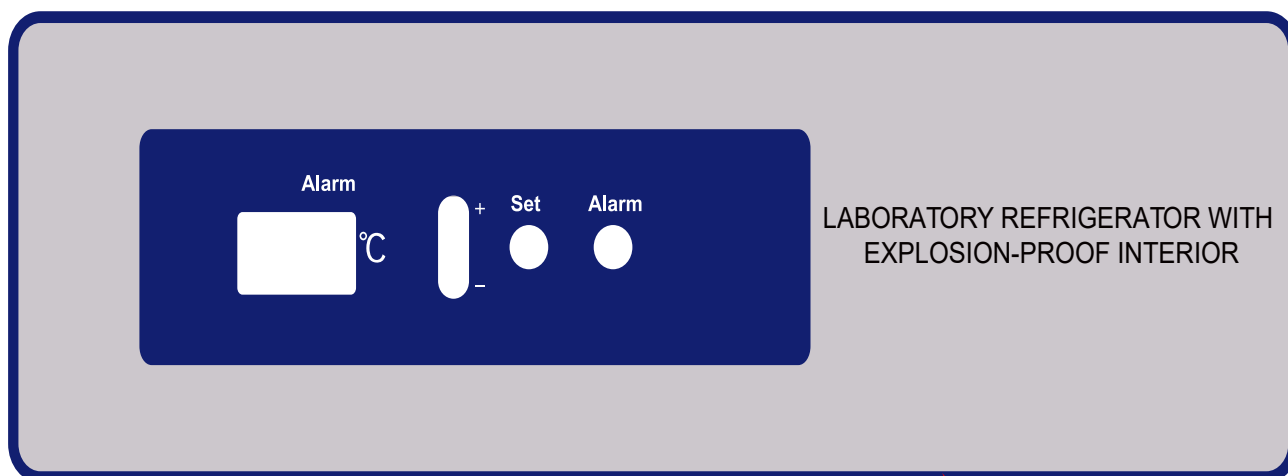
Application method

■ Temperature indicator

The front control panel of the refrigerator is equipped with a temperature indicator that facilitates adjustment of the box temperature inside the refrigerator.

For first use, the refrigerator shall run at least for over 30 minutes before it reaches the stable state, do not make any adjustment before it is stabilized.

Schematic diagram of temperature indicator :

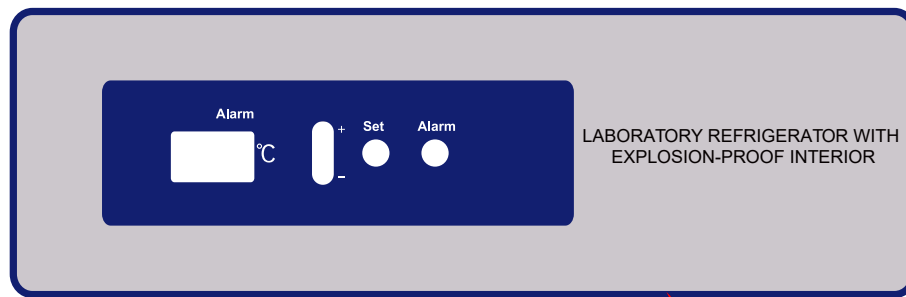- Temperature adjustment

Factory pre-set to 5°C .

Set inner temperature: Press “Set” button for 5 seconds, LED display “TS”. Press “Set” again, the current set temperature value flashes. Make adjustment to specified value through Up Key “+” and Down Key “-”. Press “Set” to save and return.

The set temperature will be reached after running for a period of time.

- Error display

When there is error of sensor, the LED displays “EE” and alarms. The refrigerator would still cool in error safety mode in a short period of time. Please move the stored things to other proper refrigerator ASAP to avoid loss. Then call for service.

- High/low temperature alarm

When the inner temperature exceeds the range of high and low alarm temperature limits, alarm light flashed and buzzer sounds. Press “Alarm” button to silent the alarm but the alarm light keeps flashing until inner chamber temperature within alarm limits.

- Check recent highest and lowest inner temperature

Press both “Alarm” and “+” buttons to check the recent highest inner temperature. Release the buttons to return to the inner cabinet temperature display.

Press both “Alarm” and “-” buttons to check the recent lowest inner temperature. Release the buttons to return to the inner cabinet temperature display.

- Set values of high temperature alarm “ALH” and low temperature alarm “ALL”

Press “Set” button for 5 seconds. LED display “TS”. Press “+” button to show “ALH”. Press “Set” again, the current high alarm temperature value flashes. Make adjustment to specified value through Up Key “+” and Down Key “-”. Press “Alarm” to return.

Press “+” button or “-” to show “ALL”. Set low temperature alarm value in the same way.

Alarm

■ Alarm function

The refrigerator has the alarm function in the table below:

Alarm	Code	Meaning	Alarm method
High temperature alarm	E00	The display temperature exceeds the set value for 15 minutes	Beep + alarm light flashing
Low temperature alarm	E01	The display temperature exceeds the set value for 15 minutes	Beep + alarm light flashing
Power off alarm	E02	Main power off or fault	Alarm lamp flashing + power supply lamp flashing + beep+Screen alternates temperature and code + Remote
Communication failure alarm	E03	Display board and power board can't send and receive data for 5 times in a row	Beep + alarm light flashing + code
Control sensor failure alarm	E04	Control sensor fault (open circuit or short circuit)	Alarm lamp flashing + beep + screen alternate display temperature and code
Open door alarm	E06	The opening time is more than 2 minutes	Alarm light flashing + beep

Cleaning and Maintenance

Component Cleaning

Warning

- In order to prevent an electric shock or personal injury, be sure to cut off the power of the refrigerator before conducting any repair to the refrigerator.
- Make sure not to suck medicine or suspended particles around the storage box when maintaining the refrigerator, otherwise, it will hurt your health.

Inner Cleaning

For model 471-1307 and 471-1309: use a damp cloth only, a dry cloth is not allowed to be used.

For model 471-1308 and 471-1310: both a damp cloth or a dry cloth could be used.

EXPLOSION-PROOF INTERIOR

Interior Temperature Range : +3 ~ +16 °C

Internal  II3/- G Ex ic nA IIC T6 Gc

Serial No.:In the Barcode

Anti-static operating instruction

WARNING

1. clean plastic parts with a damp cloth only.
2. Do not use a dry cloth to clean plastic parts.
3. Danger of sparks due to electrostatic discharge.
4. Warning - potential electrostatic charging hazard - see instructions.

Fan Cleaning

Danger of sparks produced by friction due to dust on the fan blades. Clean the ventilation slots on the re-circulated air fan every month using a vacuum cleaner.

Replacing the illuminating lamp (For model 471-1308 and 471-1310:)

The LED lamp is mounted in the upper front part of the refrigerator, exercise cautions when replacing the lamp:

- Unplug power plug;
- Remove the safety cover of the LED lamp;
- Take out the LED lamp;
- After changing with a new LED lamp, put it back to its original position;
- Turn on the power.

Stopping Refrigerator


- Cut off the refrigerator power if the refrigerator is not in service for a long time.
- Open the refrigerator's door and take out the inner plate glass shelf.
- Make thorough cleaning of the refrigerator.
- Clean the plate glass shelves.
- Wait till the Refrigerator and the plate glass shelves are dried in the air, and replace the plate glass shelves in the refrigerator.
- Close the door of the refrigerator and cover it with a plastic bag for sealing up.

FAQ

Fault	Troubleshooting
The refrigerator does not work	Inspect the power supply to make sure it meets the requirement
	Inspect the connection between the power plug and the socket
	Check the power cord for any obvious damage
	Whether the voltage is too low
Refrigeration effects are not apparent, temperature exceeds the standard	The unit is loaded with too much warm sample and product
	There is a lack of space between stored samples
	There might be a direct sunlight or other heat radiation energy on the refrigerator
	The door is opened too frequently
	Ambient temperature is too high
	Air duct is blocked
Unit seems to emit too much noise.	Check the installation of the unit to make sure the unit does sit on a solid floor without vibration
	Part of the refrigerator leans against a wall
	As a reference, the published sound data is obtained at specific laboratory condition. It is normal that the actual sound level in the field differs due to subtle installation and operation conditions
Alarm light flashes and alarm buzzer sounds	Warm products have been placed into the unit recently and the refrigerator has not stabilized yet. The alarm condition will be automatically corrected once the temperature reaches the set value
	The door is not closed properly
	The unit lost power recently and it is in the process of cooling down
	Ambient temperature might exceed the designed range

Refrigeration Schematic&Circuit Diagram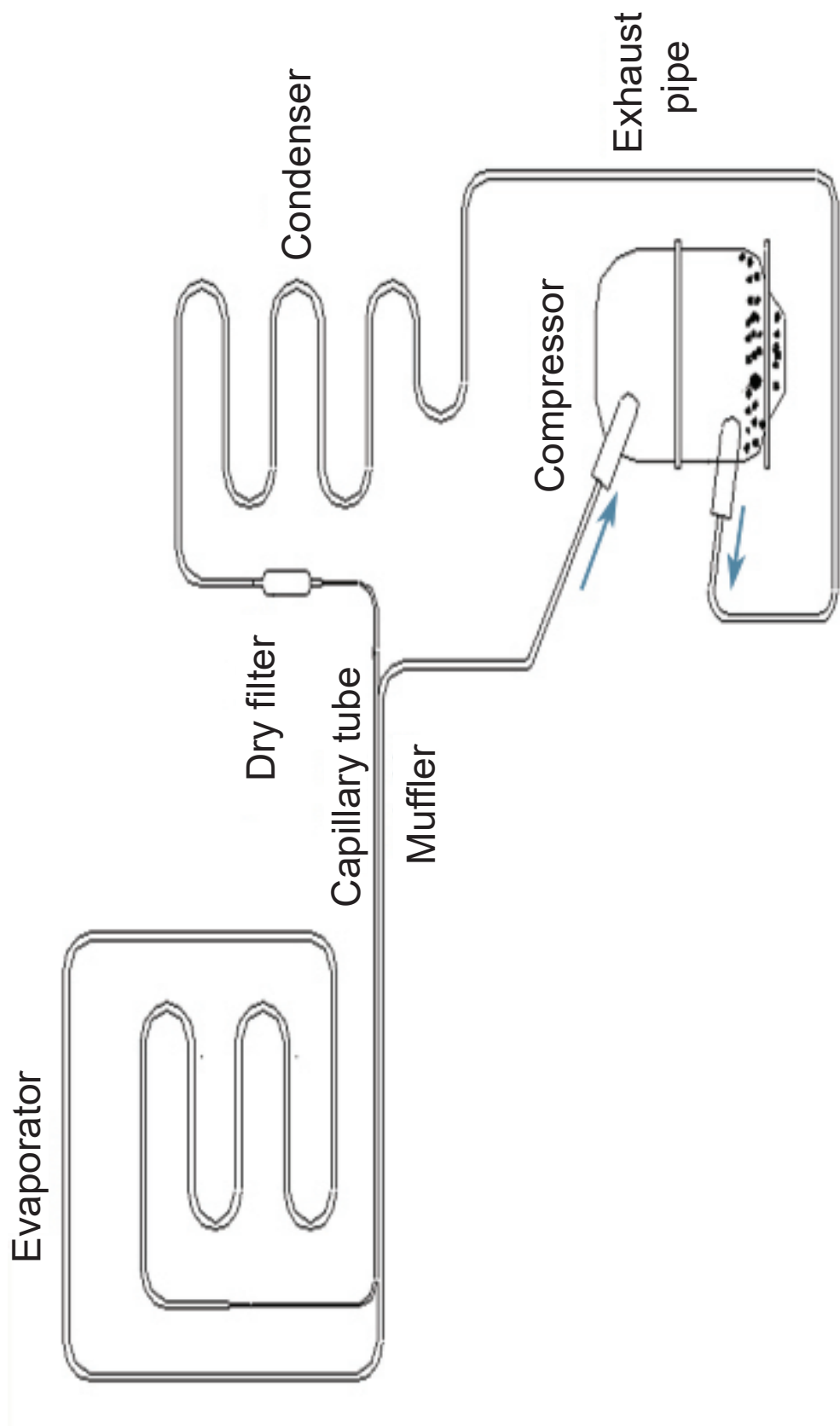

Specification

■ Specification

Model	471-1307	471-1308	471-1309	471-1310
Cabinet Type	Under bench		Upright	
Climate Class	N			
Cooling Type	Air cooling			
Defrost Mode	Auto			
Refrigerant	R600a			
Internal Temperature Uniformity(℃)	±3.0℃			
Sound Level(dB(A))	40		41	
Temperature Range(℃)	2~8℃			
Controller	Microprocessor			
Display	Digital			
Power Supply(V/Hz)	220~240/50			
Capacity(L)	118		310	
ext.dimensions (HxWxD cm)	81x60×64		184x60×60	
int.dimensions (HxWxD cm)	63×52×42		129×56×50	
Net weight (kg)	41		72	77
Gross Weight (kg)	49	46	83	87

Packing List

Model	471-1307	471-1308	471-1309	471-1310
Operation Manual	1	1	1	1
Factory inspection report	1	1	1	1
Key	2	2	2	2
Basket	1	1	1	1
Plastic bag	1	1	1	1
Shelf	2	2	4	4
Water tray	1	/	1	/
Handle	/	/	1	1
Handle Cover	/	/	1	1
Power cord	3	3	3	3

Technical service

• Web Resources

Visit the VWR website at www.vwr.com for:

- Complete technical service contact information
- Access to the VWR Online Catalogue, and information about accessories and related products
- Additional product information and special offers

Contact us For information or technical assistance contact your local VWR representative or visit. www.vwr.com.

Warranty


VWR warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non-performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

Compliance with local laws and regulations

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorisations necessary to run or use the Product in its local environment. VWR will not be held liable for any related omission or for not obtaining the required approval or authorisation, unless any refusal is due to a defect of the product.

Equipment disposal

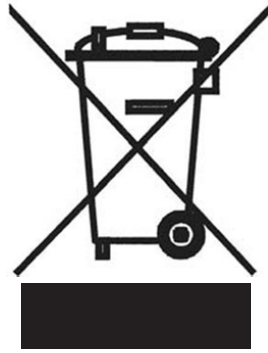This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste.

Instead it's your responsibility to correctly dispose of your equipment at lifecycle -end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment. For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

Thank you

Local VWR offices in Europe and Asia Pacific

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India

VWR Lab Products Private Limited
No.139. BDA Industrial Suburb,
6th Main, Tumkur Road,
Peenya Post,
Bangalore, India – 560058
Tel.: +91-80-28078400
Email: vwr_india@vwr.com

Ireland

VWR International Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: +353 01 88 22 222
Email sales.ie@vwr.com

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VWR International AB Esbogatán 16
VWR International GmbH Lerzenstrasse 16/18 8953 Dietikon