Multilanguage short instruction manual & Troubleshooting

VWR[®] pH METER PH-130

EU Catalogue Number: 662-2351

Version: 1.0
Issued: 01, December, 2020

LANGUAGES


English.....	3
Français.....	5
Deutsch.....	7
Italiano.....	9
Español.....	11
Português.....	13
Hrvatski.....	15
Český.....	17
Dansk.....	19
Nederlandse.....	21
Suomen.....	23
Magyar.....	25
Latviešu.....	27
Lietuvių.....	29
Norsk.....	31
Polski.....	33
Română.....	35
Slovenský.....	37
Slovenski.....	39
Svenska.....	41

FOR MORE DETAILS, PLEASE CONSULT THE FULL MANUAL

English


The PH-130 meter allows you to measure pH, mV and temperature. Turn the device on with the  key. Other operating activities are performed using the touch screen.

Sockets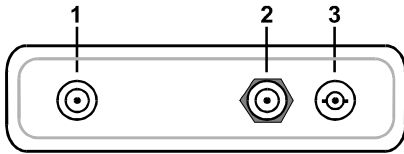


1	Power adapter socket
2	Temperature sensor socket
3	BNC-50 socket of the pH electrode

Measurement screen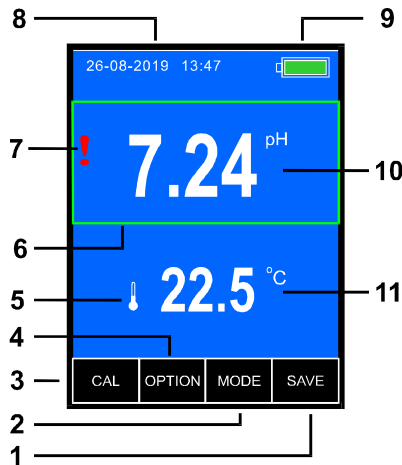


1	Collect results key
2	Parameter selection key
3	Key to enter the calibration mode (hold down for 2s)
4	Key for selecting the measurement function
5	Symbol of the connected temperature sensor
6	Fixed measurement frame
7	pH electrode warning symbol
8	Date and time
9	Battery status
10	Measurement result with unit
11	Temperature measurement result with unit

Options screen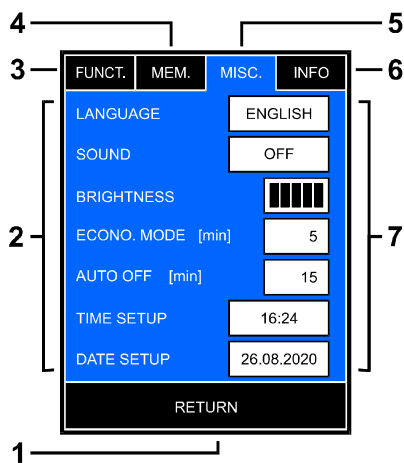

1	Key to return to the previous screen
2	Option parameters
3	Function selection tab key
4	Memory tab key
5	Miscellaneous tab key
6	Instrument and manufacturer information
7	Value input and selection windows

Calibration screen

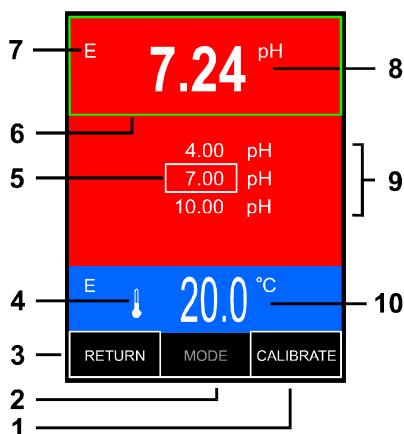1	Calibration result confirmation key
2	Parameter selection key
3	Key to return to the previous screen
4	Symbol of the connected temperature sensor
5	Confirmation of calibration in point
6	Stabilised measurement frame
7	pH electrode warning symbol
8	pH measurement result
9	Calibration points
10	Temperature measurement result with unit

TROUBLESHOOTING

Problem	Cause	Solution
Meter not switching on / No power	Discharged batteries	Charge batteries
	Incorrect battery polarity	Check the polarity
	Battery failure	Replace batteries with AA (R6) Rechargeable type
	Power adapter failure	Check with other 5V USB power adapter (for mobile)
Unstable reading	Poor connection	Check electrode connection to the unit
	pH electrode junction / membrane contaminated	Clean the electrode according to its manual
	pH electrode defective	Replace electrode
Meter is not calibrating	Incorrect value buffer used	Use buffers 4.00, 7.00 10.00 pH
	Buffer solution contaminated	Use new buffer solutions
	Calibration mode not entered properly	Touch the result to show frame around it. Press and hold the CAL button till displaying red background colour.
	Intermittent or no connection of pH electrode	Check electrode connection to the unit
	pH electrode contaminated	Clean the electrode according to its manual
	pH electrode defective	Replace electrode
Result different than expected	Electrode not calibrated	Calibrate electrode
	Buffer solution contaminated	Use new buffer solutions
	Calibration not correct	Make 2 or 3 points calibration
Temp. value incorrect  or  symbol displayed	Temperature sensor not connected or poor connection	Connect the temp. sensor / check the connection
	Breakage in temp sensor cable or connector / defective sensor	Check connection / replace sensor
Information: pH calibration readout error!	pH electrode calibration data loss	Recalibrate the pH electrode
	If the situation appears after each switching on	Contact distributor or service agent
Information: Manufacturer's settings readout error!	Internal hardware or software problem / Losing factory settings	Contact distributor or service agent

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