Instruction manual

VWR® Analog & Digital Multi-Tube Vortexer

EU cat. no

Analog Multi-Tube Vortexer:

444-7063

Digital Multi-Tube Vortexer:

444-7061

EN	- English		1
FR	- Français		13
ES	- Español		25
DE	- Deutsch		37
IT	- Italiano		49
PT	- Português		61
BG	- Български		73
CZ	- Čeština		76
DK	- Dansk		81
EE	- Eesti		85
FI	- Suomi		89
GR	- Ελληνικά		93
HR	- Hrvatska		97
HU	- Magyar		101
IS	- Íslenska		105
LT	- Lietuvių		109
LV	- Latviešu		113
NL	- Nederlands		117
NO	- Norsk		121
PL	- Polski		125
RO	- Română		129
SI	- Slovenščina		133
SK	- Slovenčina		137
SV	- Svenska		140

Legal address of Manufacturer

Europe

VWR International bv,
Researchpark Haasrode 2020
Geldenaaksebaan 464 B-3001
Leuven
+3216385011
be.vwr.com

UK Importer:

VWR International Ltd
Hunter Boulevard, Magna Park
Lutterworth, Leicestershire, LE17 4XN
uk.vwr.com

Country of Origin

United States

Table of Contents

Introduction	4
Safety Information	4
Package Contents	4
Installation	4
Intended Use	5
Maintenance & Servicing	5
Environmental Conditions	5
Product Specifications	
<i>Analog Multi-Tube Vortexer</i>	5
<i>Digital Multi-Tube Vortexer</i>	7
Operating Instructions	
<i>Analog Multi-Tube Vortexer</i>	6
<i>Digital Multi-Tube Vortexer</i>	7
Troubleshooting	9
Technical Service	9
Warranty	9
Compliance with Local Laws and Regulations	9
Disposal	10
Accessories	11
Distributor Contact Information	12

Introduction

This manual contains installation, operation and maintenance instructions for the VWR Advanced Series. Please read the manual completely before using.

Safety Information

Please read the entire instruction manual before operating the Multi-Tube Vortexer.

WARNING! DO NOT use the Multi-Tube Vortexer in a hazardous atmosphere or with hazardous materials for which the unit was not designed. Also, the user should be aware that the protection provided by the equipment may be impaired if used with accessories not provided or recommended by the manufacturer or used in a manner not specified by the manufacturer.

Always operate unit on a level surface for best performance and maximum safety.

DO NOT lift the Multi-Tube Vortexer by the support plate or bottom tray. The support plate is removable.

CAUTION! To avoid electrical shock, completely cut off power to the unit by disconnecting the power cord from the unit or unplug from the wall outlet. Disconnect unit from the power supply prior to maintenance and servicing.

Spills should be removed promptly. **DO NOT** immerse the unit for cleaning.

DO NOT operate the unit if it shows signs of electrical or mechanical damage.

Earth Ground - Protective Conductor Terminal

Alternating Current

Package Contents

Description	Quantity
Multi-Tube Vortexer	1
12mm test tube foam rack	1
Tray pad set (Top & Lower pad)	1 set
234cm detachable power cord	1

Installation

Upon receiving the VWR Multi-Tube Vortexer, check to ensure that no damage has occurred in shipment. It is important that any damage that occurred in transport is detected at the time of unpacking. If you do find such damage the carrier must be notified immediately.

After unpacking, place the Multi-Tube Vortexer on a level bench or table, away from explosive vapors. Ensure that the surface on which the unit is placed will withstand typical heat produced by the unit and place the unit a minimum of 15cm from vertical surfaces. Always place the unit on a sturdy work surface.

The Multi-Tube Vortexer is supplied with a power cord that is inserted into the IEC connector on the back of the unit first, then it can be plugged into a properly grounded outlet. The 230V unit plugs into a 230-volt, 50/60 Hz source.

Intended Use

The Multi-Tube Vortexers are intended for use in research laboratories and for general laboratory use.

Maintenance & Servicing

The Multi-Tube Vortexer is built for long, trouble-free, dependable service. No lubrication or other technical user maintenance is required. It needs no user maintenance beyond keeping the surfaces clean. The unit should be given the care normally required for any electrical appliance. Avoid wetting or unnecessary exposure to fumes. Spills should be removed promptly. **DO NOT** use a cleaning agent or solvent on the front panel which is abrasive or harmful to plastics, nor one which is flammable. Always ensure the power is disconnected from the unit prior to any cleaning. If the unit ever requires service, contact your VWR representative.

Environmental Conditions

Operating Conditions: Indoor use only.

Temperature: 4 to 40°C

Humidity: 20% to 85% relative humidity, non-condensing

Altitude: 0 to 2000 M above sea level

Non-Operating Storage:

Temperature: -20 to 65°C

Humidity: 20% to 85% relative humidity, non-condensing

Installation Category II and Pollution Degree 2 in accordance with IEC 664.

Product Specifications

Analog Multi-Tube Vortexer

Overall dimensions (L x W x H): 24,1 x 38,4 x 40,6cm

Electrical (50/60 Hz): 230 V - 1 amp, 75 watts

Fuses: 230 V - 5mm x 20mm, 1 amp quick acting

Speed range: 1200 to 2400rpm

Accuracy: +/- 5%

Orbit: 3,6mm


Weight capacity: 4,5kg

Timer: mechanical timer runs 0 to 60 seconds

Controls: rocker switch: run, standby, timer, speed knob, variable 1 to 10 dial markings, time knob, variable 1 to 10 dial markings

Ship weight: 20,4kg

European Catalogue Numbers: 444-7063

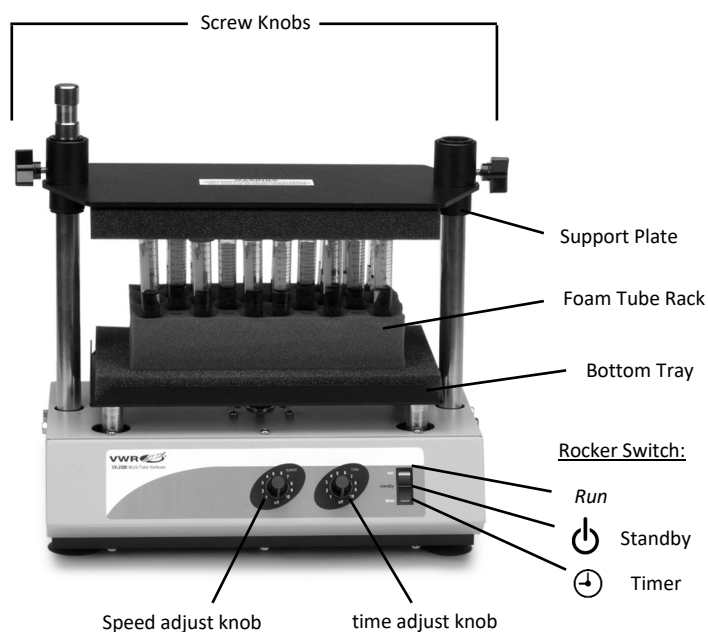Operating Instructions

Analog Multi-Tube Vortexer

The Multi-Tube Vortexer is used to prepare samples by mixing tubes just prior to testing.

IMPORTANT: This unit is designed for intermittent use only. **DO NOT** operate this unit for more than two hours at any one time.

1. Getting ready:

- a. The speed and time knobs have a built-in on/off switch at the extreme counterclockwise position. Turn the speed and time knobs to the off position.
- b. Press the rocker switch to the standby position, in the center.
- c. Plug the power cord into a properly grounded outlet.

2. Loading bottom tray:

- a. Loosen both screw knobs on the support plate. Lift the plate up. The screw knobs may be tightened to hold it in place, or if desired it may be swung out of the way. To do this, lift until the plate comes off the short right post and the left knob is up into the circular groove around the top of the left post. Screw the left screw knob in the circular groove but not quite tight. In this condition the support plate may be swung around the left post and out of the way.
- b. Fill the tubes to be vortexed and place them in the correct size test tube foam rack. Best results will be obtained if all tubes are filled to the same level, and not exceeding half of the tube capacity. Distribute tubes uniformly throughout the test tube foam rack. Always be sure that a tube is in each of the four corner holes on the test tube foam rack to properly hold the support plate. Empty tubes may be used for this purpose.
- c. Lower the support plate onto both posts and down on top of the tubes. Let the support plates weight rest on the tubes but **DO NOT** press down. Tighten both screw knobs securely.

3. Setting speed:

- a. Turn the speed knob clockwise to the 1 position.
- b. Press the rocker switch up to the continuous run position. The Multi-Tube Vortexer will start.
- c. Increase the speed by turning the speed knob clockwise until the desired speed is reached. Always use the slowest suitable speed.
- d. To stop vortexing, press the rocker switch to the standby position, in the center.

4. Setting time:

- a. For automatic timed operation, adjust the speed as in step 3, then set the time knob to the desired time, clockwise to increase and counterclockwise to decrease time.
- b. Press the rocker switch briefly to the timer position, downward, and release. The Multi-Tube Vortexer will run for the selected time and shut off automatically.
- c. To repeat for the same time, simply press the timer switch again. To interrupt an automatic timing cycle before it is complete, turn the time knob to off, extreme counterclockwise position.

When operating the Multi-Tube Vortexer, use the rocker switch to start and stop operations, leaving the speed knob on. When the unit is not being used turn the speed knob to off, extreme counterclockwise position.

Product Specifications

Digital Multi-Tube Vortexer

Overall dimensions (L x W x H): 24,1 x 38,4 x 40,6cm

Electrical (50/60 Hz): 230 V - 1 amp, 75 watts

Fuses: 230 V - 5mm x 20mm, 1 amp quick acting

Speed range: 500 to 2500rpm

Accuracy: +/- 5%

Orbit: 3,6mm


Weight capacity: 4,5kg

Timer: digital, 1 second to 9999 minutes
(increased in 1 second increments)


Controls: see page 7

Ship weight: 20,4kg

European Catalogue Number: 444-7061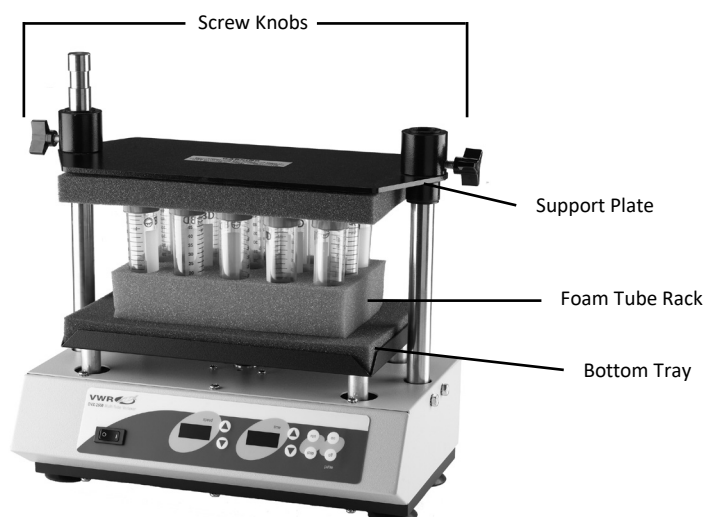

Control Panel

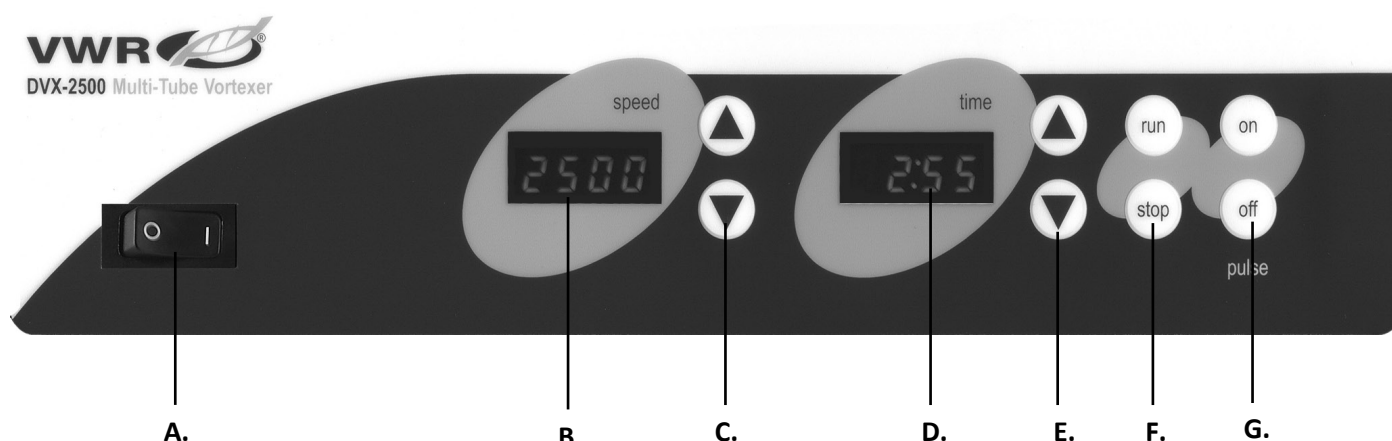Digital Multi-Tube Vortexer

The front panel of the Multi-Tube Vortexer contains all the switches, controls and displays needed to operate the unit.

A. On/off rocker switch: Turns main power on/off.

B. Speed display: Displays the speed of the vortexer.

C. Up/down arrows for set-point control.

D. Time display: Displays accumulated time (continuous mode) or how much time is remaining (timed mode). The display range is from 0 to 9999 minutes in one (1) second increments. The display will indicate minutes and seconds until the timer reaches 99 minutes and 59 seconds (99:59), then the display will automatically display minutes up to 9999.

E. Up/down arrows for set-point control.

F. Run/stop buttons: Activates vortexing.

G. Pulse on/off buttons: Activates pulse mode

Operating Instructions

Digital Multi-Tube Vortexer

The Multi-Tube Vortexer is used to prepare samples by mixing tubes just prior to testing.

IMPORTANT: This unit is designed for intermittent use only. **DO NOT** operate this unit for more than two hours at any one time.

1. Getting ready:

- By sure the rocker switch is in the off position, depressed to the right.
- Plug the power cord into a properly grounded outlet.
- Press the rocker switch to the on position, depressed to the left. The speed and time displays will illuminate.

2. Loading bottom tray:

- Loosen both screw knobs on the support plate. Lift the plate up. The screw knobs may be tightened to hold it in place, or if desired it may be swung out of the way. To do this, lift until the plate comes off the short right post and the left knob is up into the circular groove around the top of the left post. Screw the left screw knob in the circular groove but not quite tight. In this condition the support plate may be swung around the left post and out of the way.
- Fill the tubes to be vortexed and place them in the correct size test tube foam rack. Best results will be obtained if all tubes are filled to the same level, and not exceeding half of the tube capacity. Distribute tubes uniformly throughout the test tube foam rack. Always be sure that a tube is in each of the four corner holes on the test tube foam rack to properly hold the support plate. Empty tubes may be used for this purpose.
- Lower the support plate onto both posts and down on top of the tubes. Let the support plates weight rest on the tubes but **DO NOT** press down. Tighten both screw knobs securely.

3. Setting speed:

Press the up/down arrows to the right of the speed display until you reach the desired speed.

4. Setting timed mode:

To run in timed mode, press the up/down arrows to the right of the time display until you reach the desired time. Press the run button. The unit will run for the programmed time. Time LED will display time remaining.

5. Setting continuous mode:

To run in continuous mode, set the time to zero (0) using the up/down arrows to the right of the time display. Press the run button and the unit will run until you press the stop button. Time LED will display accumulated time.

6. Setting pulse mode:

- Make sure the unit is not running.
- Press the on button above the word pulse. The unit must be stopped before pressing the pulse-on button.
- To stop the pulse feature press the off button above the word pulse. Pulse mode is factory programmed for 2 seconds on, 1 second off. The pulse-on and pulse-off times can be adjusted between 1 and 59 seconds in 1 second intervals.

7. Reprogram the pulse-on and/or pulse-off time as follows:

Pulse-on time

- Press rocker switch to the off position, to the right.
- Press rocker switch to the on position, to the left, while holding pulse-on button. The LED display for speed displays current pulse-on time. The pulse-on button can be released, if necessary, once the display is illuminated.
- Change the pulse-on time using the up/down arrows to the right of the speed display.
- When the rocker is pressed off, the value on the speed display is stored as the new pulse-on time.

Pulse-off time

- Press rocker switch to the off position, to the right.
- Press rocker switch to the on position, to the left, while holding pulse-off button. The LED display for time displays current pulse-off time. The pulse-off button can be released, if necessary, once the display is illuminated.
- Change the pulse-off time using the up/down arrows to the right of the time display.
- When the rocker is pressed off, the value on the time display is stored as the new pulse-off time.

Troubleshooting

Review the information in the table below to troubleshoot operating problems.

Problem	Cause	Solution
Unit fails to power on	Missing or blown fuse	Add or replace fuse as necessary. If problem persists, please contact your VWR representative for repair.
Unit not reaching set speed	Top plate too tight	Re-seat top plate not to exceed 11,3 kg of force. If problem persists, please contact your VWR representative for repair.
Unit is excessively noisy	Rattling or ticking sounds may indicate a loose screw on the tray	Ensure that tray is secured tightly. If problem persists, please contact your VWR representative for repair.

Technical service

Web Resources

Visit the VWR website at www.vwr.com for:

- Complete technical service contact information
- Access to the VWR Online Catalogue, and information about accessories and related products
- Additional product information and special offers

Contact us For information or technical assistance contact your local VWR representative or visit. www.vwr.com.

Warranty


VWR warrants that this product will be free from defects in material and workmanship for a period of five (5) years from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non-performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

Compliance with local laws and regulations

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorizations necessary to run or use the Product in its local environment. VWR will not be held liable for any related omission or for not obtaining the required approval or authorization unless any refusal is due to a defect of the product.

Equipment disposal

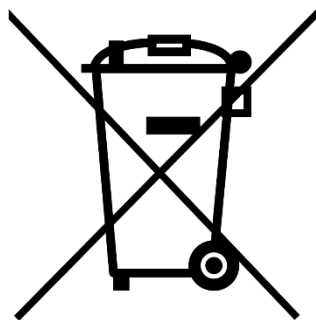This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste.

Instead it's your responsibility to correctly dispose of your equipment at lifecycle -end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

Thank you

Accessories

Description	Tube Capacity	Color	Cat. No.
10mm test tube foam rack	50	Gray	444-7065
12mm test tube foam rack	50	Blue	444-7066
13mm test tube foam rack	50	Yellow	444-7067
16mm test tube foam rack (for 15mL centrifuge tubes)	50	Pink	444-7068
25mm centrifuge tube foam rack	28	White	444-7091
29mm centrifuge tube foam rack (for 50mL centrifuge tubes)	15	Red	444-7092
Replacement tray pad set (top & lower pad)	N/A	Gray	444-7064
Post extension kit	N/A	N/A	444-7062

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India

VWR Lab Products Private Limited
No.139. BDA Industrial Suburb,
6th Main, Tumkur Road, Peenya Post,
Bangalore, India – 560058
Tel.: +91 80 28078400
Email: vwr_india@vwr.com

Ireland

VWR International Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: +353 01 88 22 222
Email: sales.ie@vwr.com

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