VWR

Microscope  **VisiScope®** series 100

INSTRUCTION MANUAL

Model	European Catalogue Number
ML103	630-1550
ML114	630-1551
ML124	630-1552
BL114	630-1553
BL124	630-1554

Version: 1
Issued: 01, 09, 2012

Legal Address of Manufacturer

United States

VWR International
Radnor Corporate Center
Building One, Suite 200
100 Matsonford Road, P.O. Box 6660
Radnor, PA 19087
(610) 386-170

Europe

VWR International bvba
Researchpark Haasrode 2020
Geldenaaksebaan 464
B-3001 Leuven
+ 32 16 385011
<http://be.vwr.com>

Country of origin: **ITALY**

Table of Contents

Warning

Safety information

Package contents

Unpacking

Intended use

Symbols and conventions

Specifications

Instructions for use

Overview

Description of buttons and switches

Operation

Troubleshooting

Repair and maintenance

Accessories and spares

Technical service

Warranty

Disposal

Warning

This microscope is a scientific precision instrument designed to last for many years with a minimum of maintenance. It is built to high optical and mechanical standards and to withstand daily use.

We remind you that this manual contains important information on safety and maintenance, and that it must therefore be made accessible to the instrument users.

We decline any responsibility deriving from incorrect instrument use that does not comply with this manual.

Safety Information

Avoiding Electrical Shock

Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off position.

Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users have full responsibility to use this equipment safely.

Please follow the guidelines below, and read this manual in its entirety to ensure safe operation of the unit.

Package Contents

DESCRIPTION	QUANTITY
Microscope stand with 4 position nosepiece, stage, condenser	1
Optical head (ML103, ML114, ML124: monocular; BL114, BL124: binocular)	1
Objective 4x	1
Objective 10x	1
Objective 40x	1
Objective 100x (or Objective 60x for mod. ML114, BL114. NOT included for mod. ML103)	1
Eyepiece WF10x/18mm	1 (MONOCULAR) 2 (BINOCULAR)
Green filter 30mm diameter	1
Immersion oil	1
Dust cover	1
Power supply output 5Vdc	1

Unpacking

The microscope is housed in a moulded Styrofoam container. Remove the tape from the edge of the container and lift the top half of the container. Take care to ensure that the optical items (objectives and eyepieces) do not fall out and get damaged. Using both hands (one around the neck and one around the base), lift the microscope from the container and put it on a stable desk.

Place the observation head onto the top of the neck and tighten the lock-screw. Insert the eyepieces into the eye tubes.

Connect the provided power supply to the power supply input jack on the rear of the microscope.

Intended use

For research and teaching and teaching use only. Not intended for any animal or human therapeutic or diagnostic use.

Symbols and conventions


The following chart is an illustrated glossary of the symbols that are used in this manual.


	CAUTION This symbol indicates a potential risk and alerts you to proceed with caution
---	---

PRODUCT SPECIFICATIONS

Model	Head	Eyepiece	Objectives	Stage	Condenser	Focusing system	Illuminator
ML103	Monocular	Widefield 10x/18mm	Achromatic 4X/10X/40X	Plain stage, 125x120mm with sample clips	0.65 N.A. with diaphragm	Coaxial coarse and fine focusing mechanism with limit stop	1W LED with external power supply
ML114	Monocular	Widefield 10x/18mm	Achromatic 4X/10X/40X/60X	Mechanical stage, 125x116mm with specimen holder	1.2 N.A. with diaphragm	Coaxial coarse and fine focusing mechanism with limit stop	1W LED with external power supply
ML124	Monocular	Widefield 10x/18mm	Achromatic 4X/10X/40X/100X	Mechanical stage, 125x116mm with specimen holder	1.2 N.A. with diaphragm	Coaxial coarse and fine focusing mechanism with limit stop	1W LED with external power supply
BL114	Binocular	Widefield 10x/18mm	Achromatic 4X/10X/40X/60X	Mechanical stage, 125x116mm with specimen holder	1.2 N.A. with diaphragm	Coaxial coarse and fine focusing mechanism with limit stop	1W LED with external power supply
BL124	Binocular	Widefield 10x/18mm	Achromatic 4X/10X/40X/100X	Mechanical stage, 125x116mm with specimen holder	1.2 N.A. with diaphragm	Coaxial coarse and fine focusing mechanism with limit stop	1W LED with external power supply

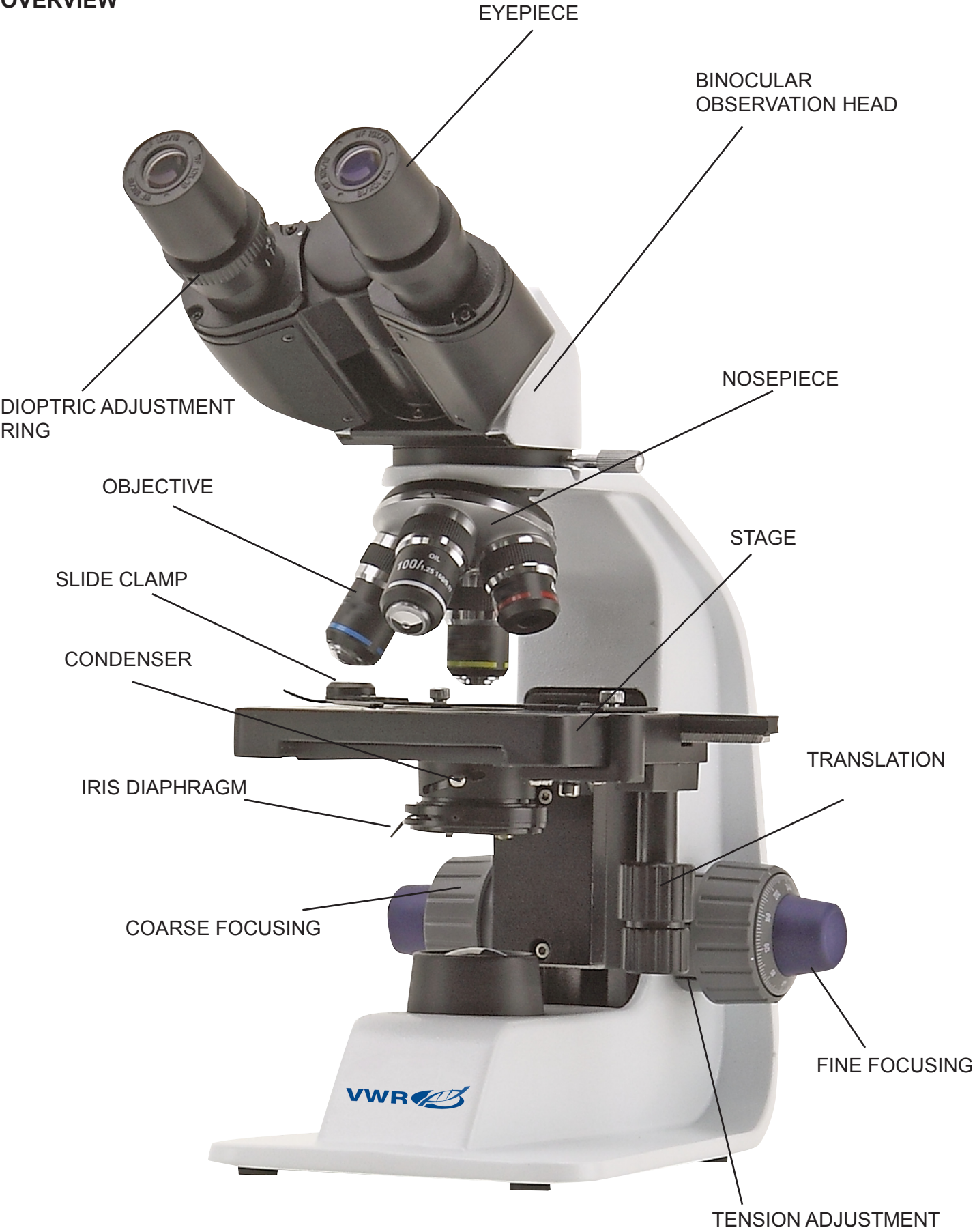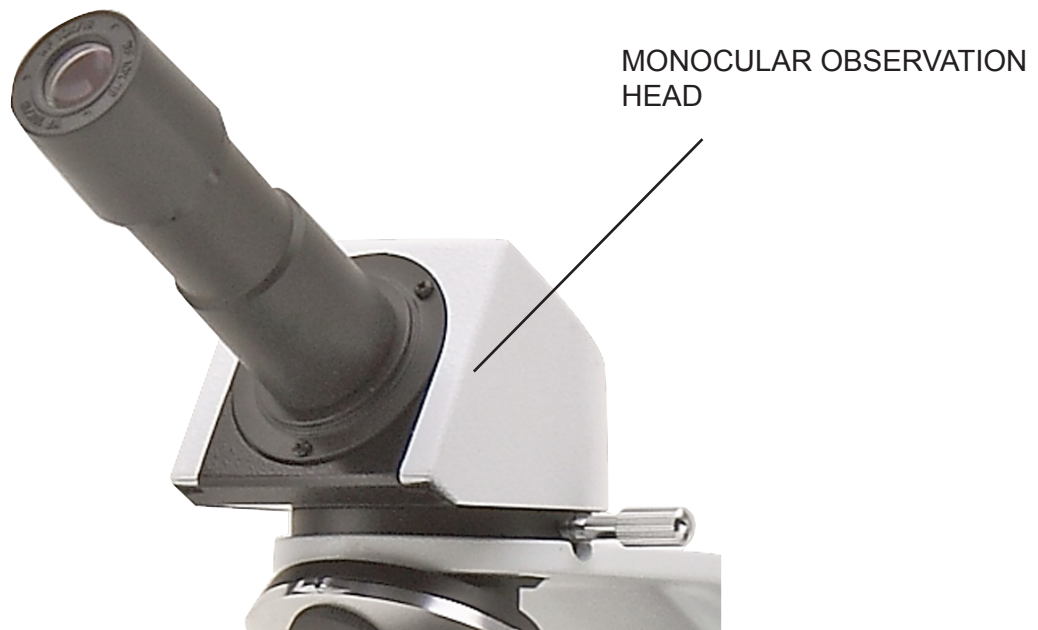OVERVIEW

DESCRIPTION OF BUTTONS AND SWITCHES

Rear View

Operation

Adjust the observation head

Loosen the lock-screw, turn the observation head to a comfortable position for observation, and then lock the lock-screw.

Place the specimen on the stage

Lock the specimen slide on the mechanical stage using the slide holder. Ensure that the specimen is centred over the stage opening by adjusting the coaxial knobs of the stage.

Lamp settings

The microscope is fitted with a power white LED illuminator. Before turning on the light, read the section about electrical safety precautions. Insert the plug of the 7,5Vdc power supply into the jack on the rear and turn on the ON/OFF switch. Turn the brightness adjustment knob to a brightness suitable for observation.

Adjust interpupillary distance (binocular models)

Hold the right and left side of the observation head with both hands and adjust the interpupillary distance by turning the two parts until one circle of light can be seen, when looking down the microscope with both eyes.

Focus and diopter adjustment (binocular models)

Rotate the coarse focusing knob to bring the slide into focus using a low magnification objective. Adjust the fine focusing knob to get the image sharp and clear while observing with your stronger eye, then turn the dioptic adjustment ring on the opposite eyepiece until a sharp and clear image can be observed also with the other eye. Turn the tension adjust knob to achieve a suitable focus tension without the specimen dropping out of focus. When the image is in focus, turn the nosepiece to choose the objective you need.

Condenser adjustment

Hold the notched ring at the base of the condenser and rotate it up or down until you reach a clear and uniform sample illumination.

Set the numerical aperture

Adjust the aperture of the iris diaphragm to set the numerical aperture of the illuminator, thus controlling image contrast and resolution.

Troubleshooting

Review the information in the table below to troubleshoot operating problems.

PROBLEM	CAUSE	SOLUTION
LIGHT DOESN'T TURN ON	Power supply not connected	Check that the power supply jack is well inserted on the rear of the microscope
	Connected	Rotate the brightness adjustment control and check if there is an increase in the light output
IMAGE CANNOT BE SEEN OR IS DARK	The iris diaphragm aperture is not completely opened	Completely open the iris diaphragm aperture
	Brightness level is low	Rotate the brightness adjustment potentiometer
	Objective is not aligned with the optical axis	Rotate the nosepiece until an objective is well inserted in the optical path (it "clicks")
IMAGE IS UNCLEAR, BLURRED OR HAS INSUFFICIENT CONTRAST	Objectives or filters are dirty	Wipe them clean
	The iris diaphragm aperture is not opened correctly	Open the iris diaphragm aperture completely
	Condenser at wrong height	Hold the condenser and rotate it until you see a uniform illumination

Repair and maintenance

Microscopy environment

This microscope is recommended for use in a clean, dry and shock free environment with a temperature of 0-40°C and a maximum relative humidity of 85 % (non condensing). Use a dehumidifier if needed.

After using the microscope

The microscope should always be kept vertical when moving it so that no moving parts, such as the eyepieces, fall out.

Never mishandle or impose unnecessary force on the microscope.

Never attempt to service the microscope yourself.

After use, turn down the illumination intensity control and turn the light off. Cover the microscope with included the dust cover, and keep it in a dry and clean place.

Electrical safety precautions

Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off-position.

Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users do have full responsibility to use this equipment safely.

Cleaning the optics

If the optical parts require cleaning first use compressed air.

If that is not sufficient use a soft lint-free cloth with water and a mild detergent.

And as a final option use the piece of cloth moistened with a 3:7 mixture of ethanol and ether.

Note: ethanol and ether are highly flammable liquids. Do not use them near a heat source, near sparks or near electric equipment. Use these chemicals in a well ventilated room.

Remember to never wipe the surface of any optical items with your hands. Fingerprints can damage the optics. Do not disassemble objectives or eyepieces in an attempt to clean them.

For the best results, use the VWR cleaning kit (see catalogue number below).

If you need to send the microscope to manufacturer for maintenance, please use the packaging if possible.

User replaceable accessories and spare parts

DESCRIPTION	QUANTITY	CAT. NO.
H5x eyepiece	2	VWRI630-1624
WF10x/18mm eyepiece	2	VWRI630-1625
WF16x/12mm eyepiece	2	VWRI630-1626
WF10x/18mm micrometric eyepiece	1	VWRI630-1627
Achromatic objective 4x/0.10	1	VWRI630-1628
Achromatic objective 10x/0.25	1	VWRI630-1629
Achromatic objective 20x/0.40	1	VWRI630-1630
Achromatic objective 40x/0.65	1	VWRI630-1631
Achromatic objective 60x/0.85	1	VWRI630-1632
Achromatic objective 100x/1.25 (oil)	1	VWRI630-1633
Polarising set (filters only)	1	VWRI630-1635
Dust cover, type 3	1	VWRI630-1636
Solar battery pack SBP 10	1	VWRI630-1637
Attachable mechanical stage	1	VWRI630-1634
Cleaning kit	1	VWRI630-1803

Technical service

Web Resources

Visit the VWR's website at **www.vwr.com** for:

- Complete technical service contact information
- Access to VWR's Online Catalogue, and information about accessories and related products
- Additional product information and special offers

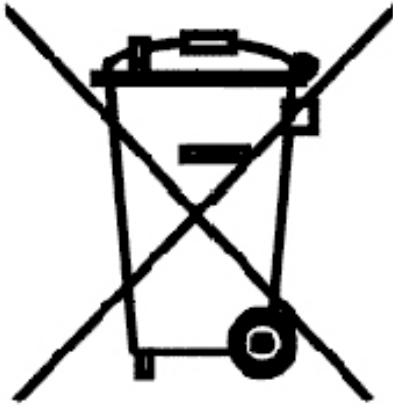
Contact us for information or technical assistance contact your local VWR representative or visit www.vwr.com.

Warranty

VWR International warrants that this product will be free from defects in material and workmanship for a period of five (5) years from date of purchase. If a defect is present, VWR will, at its option, repair, replace, or refund the purchase price of this product at no charge to you, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear.

For your protection, items being returned must be insured against possible damage or loss. This warranty shall be limited to the replacement of defective products. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

Equipment disposal

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste.

Instead it is your responsibility to correctly dispose of your equipment the end of its life cycle by handling it over to an authorized facility for separate collection and recycling. It is also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

Thank you

Local VWR offices in Europe and Asia Pacific

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VWR International China Co., Ltd
VWR Lab Products Pvt. Ltd
VWR Singapore Pte Ltd