

HIGH-DENSITY STORAGE RACK INSTALLATION INSTRUCTIONS

PRODUCT SERIES: SU780XLE

NO TOOLS REQUIRED

PROCEDURE:

1. Unpackage rack system components.



2. Components should include the following:

Tower (4 pieces)




Numbered Trays (140 pieces)



Blank Box (1 piece)





3. Remove and dispose of any shelves and shelving clips from the freezer. Insert lower left three-column tower (two spacers on back). Push against left inner side and back wall of cabinet.

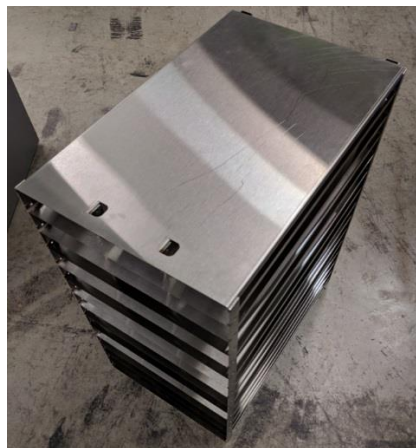

4. Insert lower right two-column tower (two spacers on back and right side). Push against right side of lower left three-column tower and back wall. Tower should be oriented so that the two small slots are on top.


Spacers

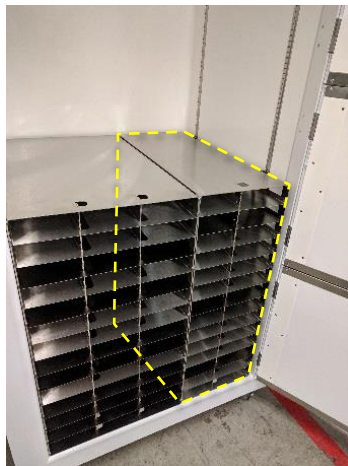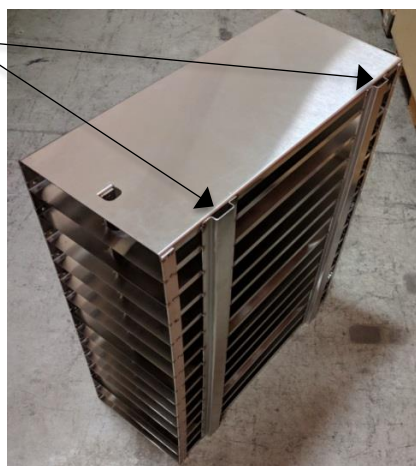

5. Insert top left two-column tower (two spacers on back, cutout in top front left corner). Push against left inner side and back wall of cabinet.



6. Insert top right three-column tower (two spacers on back and right side). Push tower against right side of top left two-column tower and back wall.

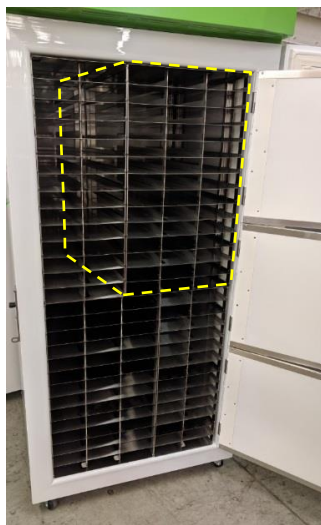

7. Place blank box in the first slot of tray #1. This is to prevent a sample box from blocking the freezer ports.



8. Insert tray #1 into the top most left slot of the top left three-column tower. Continue inserting numbered trays sequentially in a vertical manner until column is filled.



9. Insert number tray "29" into the first slot of the second column. Continue inserting numbered trays sequentially in a vertical manner until the column is filled. Repeat with remaining columns until all slots are filled with "140" numbered trays.

