

SANI-STICK

The SANI-STICK is designed to collect samples to detect the presence of microbiological contaminations on any surface. The SANI-STICK allows swabbing of hard to reach areas. These sponges on sticks, are widely used in food industry applications, such as quality assurance procedures (QA/QC).



STEP BY STEP INSTRUCTIONS

- 1 Remove the tear-off strip and carefully take the SANI-STICK by the handle.
- 2 The wet sponge is already on the handle; swipe it on the determined surface.
- 3 Press on the bottom of the handle to gently release the sponge into the sterile sampling bag.
- 4 Twirl the bag three to four times and close it with the tabs.
- 5 Identify the sample via the write-on area with a permanent marker.

