

Short user's manual

Gel Documentation Basic Image Capture

EU cat. No: 730-1470

Manufactured in UK

1 Turn on

Select



➔ Turn on the camera (ensure the battery is charged)

➔ The Canon control software launches automatically when you turn on the camera. If the software does not start automatically click on Start and in the search box type EOS Utility.

2 Capturing an image

Select



➔ Select Live View shoot to open live window

➔ Ensure the system has been installed correctly, as per the Installation Quick Guide

3 Settings

Select




Select



➔ The AWB button and from the White Balance menu select Cloudy from the menu

4 Focus setting

Select

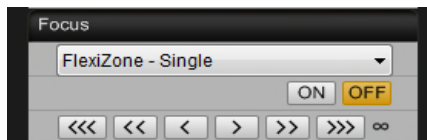

➔ **From the Remote Live View window**

To adjust zoom turn the lens of the camera

➔ To adjust focus use the arrows

5 Grab image

Press



From the EOS 2000D window

➔ The camera will set an exposure time relevant to the light levels coming from the sample

➔ Then turn off the camera

6

Save and export

- ➔ Images are saved as JPEG files
- ➔ Export a TIFF or BMP